



General Purpose Cabinet Installation and Maintenance Instructions

Model No. 6075

Please unpack these cabinets carefully. If any damage is evident, immediately notify the delivering freight carrier and the supplier of your SciMatCo cabinet.

You have just purchased the finest cabinet designed solely for the safe storage of hazardous chemicals. Please read the manual thoroughly before using this product. Only store chemicals in the compartments which are designed to hold these chemicals. Please keep this manual on hand for future reference.

Caution: Always use eye, face, hand, and body protection when using this cabinet.

I. Unpacking the Cabinets

The best method of unpacking the cabinets is as follows.

- A) Remove staples, tape and banding material from the top and bottom of the carton.
- B) Open the top of the box and remove the four (4) packing corners. Gently lift the carton off the product. Be careful not to scratch the cabinet with the box.
- C) Move the cabinet to its appropriate location. Make sure the surface on which you would like to set the cabinet can support the weight of the cabinet and its contents.

Caution: Do not locate your cabinet near open flame, spark or heat or any source of ignition.

II. Assembly and Set-Up

Your General Purpose cabinet comes fully assembled so no assembly is needed. Once you have moved your cabinet into its proper location, you may need to adjust the leveling feet to insure the proper stability.

Warning: Always close and secure cabinet doors after removing your chemicals.

III. Maintenance

Your SciMatCo General Purpose cabinet is constructed of 1" thick, high density plywood, painted with several coats of a lacquer paint and then painted with three (3) coats of clear epoxy. This cabinet was designed to give you years of trouble free service with minimal maintenance.

Caution: Always use eye, face, hand and body protection when cleaning this cabinet.

The inside of the cabinet should occasionally be wiped out using a solution of any household cleaner. Having a good cleaning program and following it will assure you that you will get many years of service out of your cabinet.

Any scratches in the painted surface of the cabinet should be repainted at once. You can order an epoxy touch up paint kit from us. Please refer to the replacement parts list at the end of this manual for the appropriate catalog number.

IV. Materials to be stored

This General Purpose cabinet is designed for the storage of flammable liquids, poisons and some oxidizers.

Warning: Absolutely no corrosive materials should ever be stored in this cabinet.

The metal hinges and lock used in this cabinet, while required for use as a flammables cabinet, will not hold up to strong oxidizers such as calcium hypochlorite, iodine or corrosive materials such as acids.

Never store acids or any material which will corrode the metal components in this cabinet.

Warning: If this cabinet is to be used for the storage of **Flammable Materials** it must have the proper label. Please contact SciMatCo for the proper label.

V. Replacement Parts List

Description	Part No.
Paint Touch-up Kit Gray	2060-1
Extra Keys General Purpose	7021-1
New Locks General Purpose	7021-4
Miscellaneous Parts Dispensing Tray	8041-6

If you have any questions or need any replacement parts, please contact your supplier:

SciMatCo Product Warranty

Scientific Materials Company warrants its products to be free from defects in materials and workmanship for a period of one (1) year from the date of purchase. This warranty excludes damage caused by abuse, misuse, accidents or the failure to provide proper maintenance.

SciMatCo

Division of Flinn Scientific, Inc.
P.O. Box 305 Batavia, IL 60510
(630) 879-7620 FAX (630) 879-6962