**SPDC Data Collection  
Instruction Manual**

**SPDC Data Collection  
Manual de instrucciones**

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## 1. INTRODUCTION

This software automatically detects data from the instrument and enables the user to record the received data into a Text, Word, Excel or Access file.

## 2. INSTALLATION

Execute the “setup.exe” file to install the software.


## 3. START-UP

After installation has been completed, start the software by clicking on either of the following:

- The “SPDC DataCollection” icon on desktop.
- The “SPDC DataCollection” icon in programs menu.
- The “SPDC DataCollection” file in installation folder.

## 4. MAIN WINDOW


When the software has been launched, the main window will open.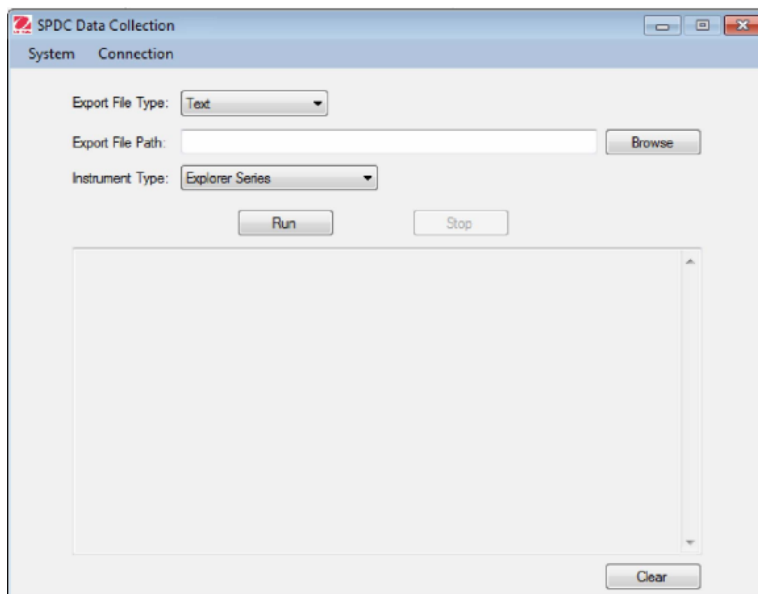

## 5. SERIAL PORT CONFIGURATION

Click the “Config” link in “Connection” menu, the <Connection Configuration> window will appear.

Click the “RS232”. The default configuration is 9600-8-N-1-N.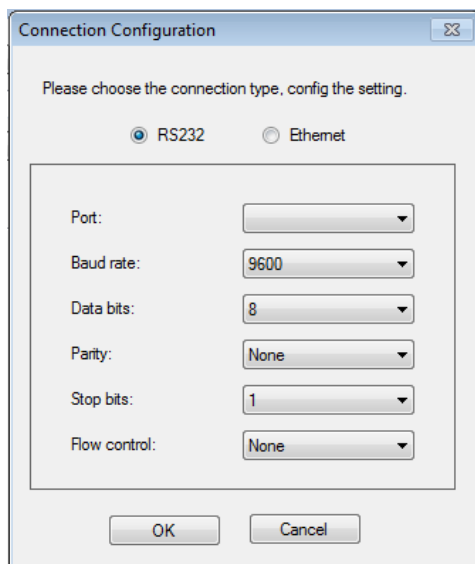

The following can be set:


- The serial port number to which the instrument is connected.
- The baud rate.
- The number of data bits.
- The parity.
- The number of stop bits.
- The flow control.

**Warning:** The baud rate, the number of bits, the parity and the number of stop bits selected in the software must correspond exactly to those of the instrument settings. Otherwise, the data will not be received or will be misinterpreted.

## 6. ETHERNET CONFIGURATION

Click the “Config” link in “Connection” menu, the <Connection Configuration> window will appear.

Click the “Ethernet”. Then the IP address and Port can be set.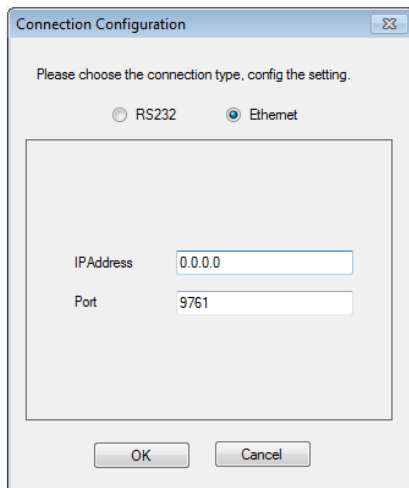


**Warning:** The IP address and Port filling in the software must correspond exactly to those of the instrument settings. Otherwise, the data will not be received or will be misinterpreted.

## 7. EXPORT FILE TYPE

Click the dropdown list, and choose the type of file that you want to export the received data to from the instrument.

- Text File
- Word File
- Excel File
- Access File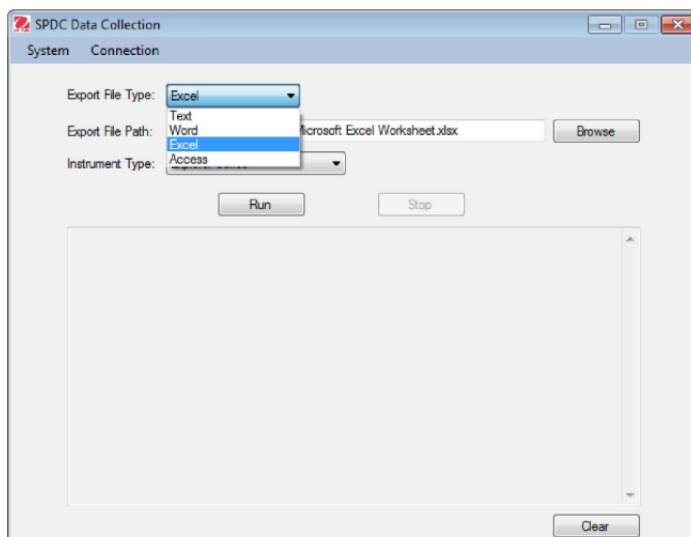


## 8. EXPORT FILE PATH

According to the Export File Type, you may choose a file to which you want to export the received data.

- Input file path manually or choose file by clicking “Browse” button.
- The export file can be an existing file, or you may create a new file.

## 9. EXPORT INSTRUMENT TYPE

Click the dropdown list and choose the type of instrument that you want to link.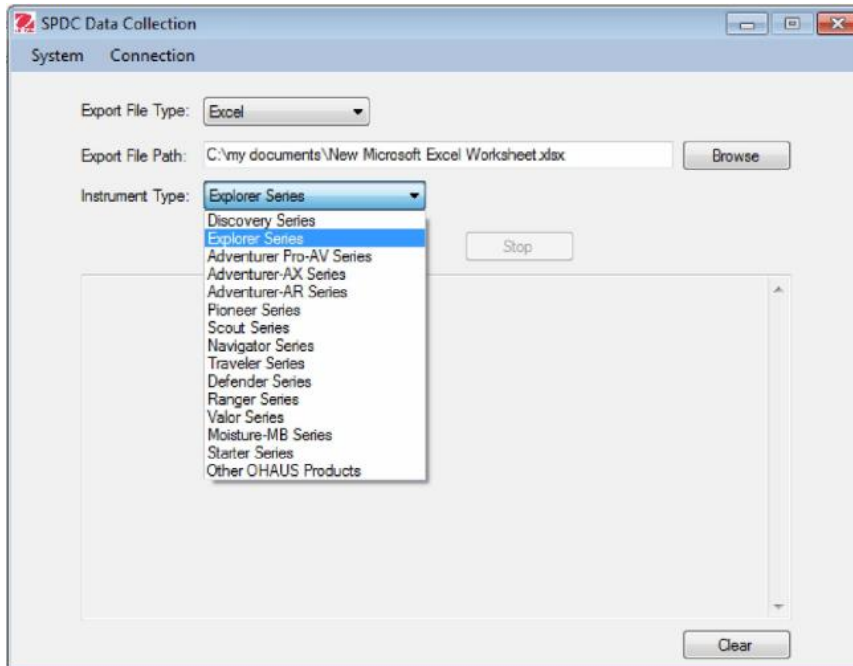

### Data Collection

#### 1. Prerequisite:

- The serial port settings must correspond exactly to instrument settings.
- The chosen export file must exist and be writable.

#### 2. Start:

Click the “Run” button.

The software will now detect the data received from the balance.

When there are data detected from instrument, the software will export the received data to both file and screen automatically.

**Note:** For Word and Excel files, the file will be opened and the received data will display synchronously.

#### 3. Field input:

For Excel files, the input field could be customized by point the cursor to certain grid.

#### 4. Stop:

Click the “Stop” button.


The software will stop detecting data received from the instrument.

**Note:** For Word and Excel files, if the file is closed the detection will also stop.


#### 5. Clear data:

Clear data in the screen by clicking the “Clear” button.

Examples: When instrument transfer to Word file.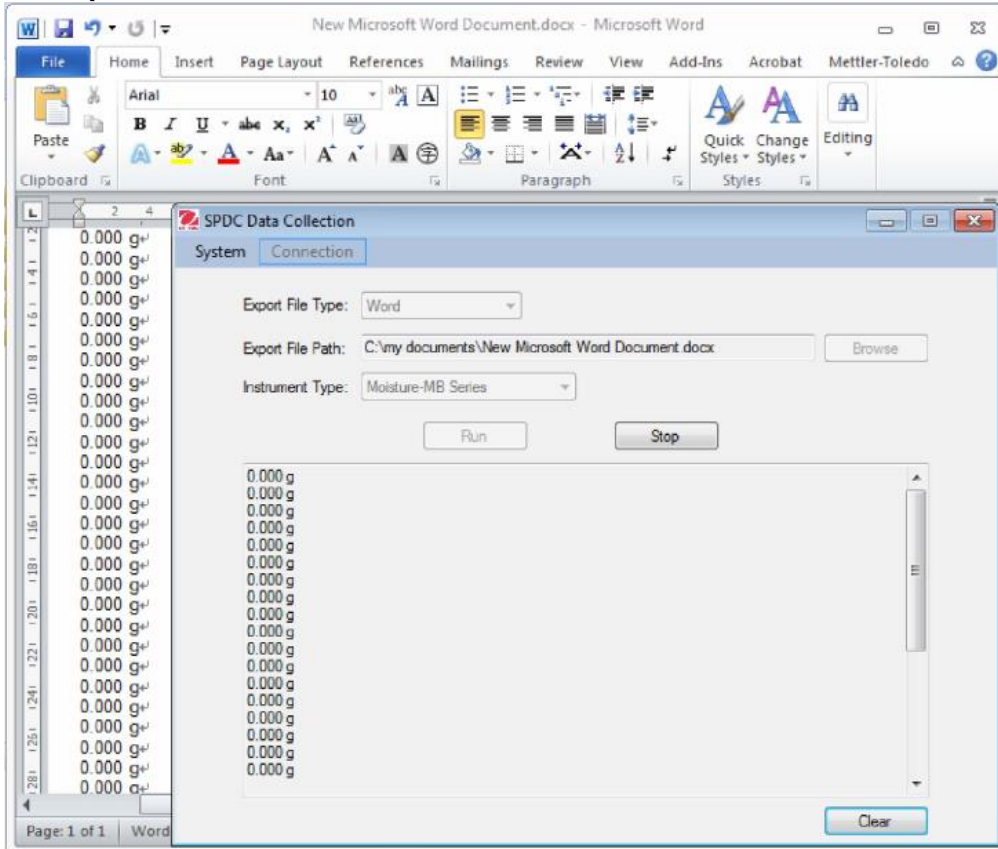

Examples: When instrument transfer to Excel file.

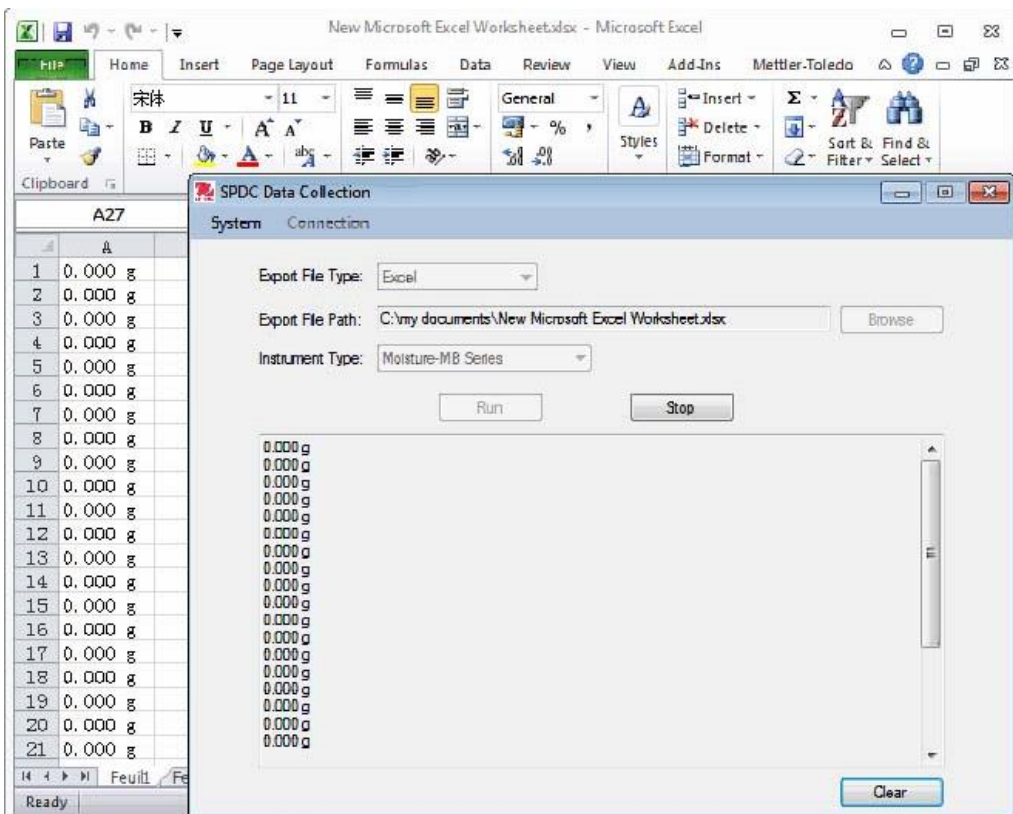**10. EXIT**

Exit the software by clicking:

- \_ The "Exit" link in "System" menu.
- \_ The "X" button in software main window.

**11. UNINSTALL**

Uninstall the software by clicking:

- \_ The "Uninstall" icon in programs menu.
- \_ The "Uninstall.exe" file in installation folder.

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P/N 30240594 | © 2018 OHAUS Corporation, all rights reserved / todos los derechos reservados / tous droits réservés / Alle Rechte vorbehalten / tutti i diritti riservati
Ohaus Corporation