prac-tec®

USER GUIDE

G-202 Digital Scale

GB

Thanks for purchasing the G-202 Gram Scale! This User Guide is designed to get you up and running quickly with your scale. Scales are precision instruments and require proper care and skill to get the best results.

What's In The Box

User Guide

100g
Calibration
Weight

G-202 Digital Scale
& Zippered Case

Key Descriptions

R1/R2 – Changes the Capacity and Resolution
(See Capacity Setting).


MODE – Changes the weighing unit.

ZERO/TARE – Zeroes the scale. Also subtracts the weight
of an empty container (See Tare Function).

ON/OFF – Turns the scale ON or OFF.

Weighing Your First Item

1. Remove protective cover/tray.
2. Press the ON/OFF key to turn the scale on.
3. Place the item on the centre of the weighing platform.
4. Take your reading from the display. You can press the MODE key to convert between weighing units.
5. The scale/s auto-off function will turn the scale off after 3 minutes of inactivity or you can press the ON/OFF key to turn the scale off.

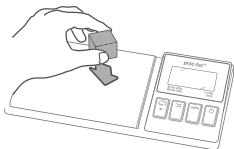Things To Avoid

Unlevel Surfaces

**Radio Frequency
Interference**

High Humidity

Air Currents


Vibration

Extreme Temps

Tare Function

The Tare function allows you to subtract the weight of a container so that the display only shows the weight of the contents.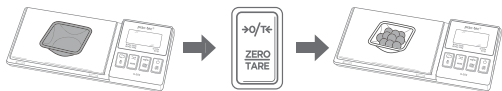

Place an empty tray or container on the platform.

Press the **ZERO/TARE** key to subtract the container's weight.

Add items to the container to see their weight on the display.

NOTE: The weight of the container will subtract from the available capacity and will leave a negative value on the display once the weight is removed. Press **ZERO/TARE** again to return the display to zero.

Capacity Setting

Your scale has two capacities and the ability to change between R1 (200g/0.01g) and R2 (250g/0.1g). The default capacity is set to R1 (200g/0.01g). To change the capacity setting:

1. Turn the scale on and press the **R1/R2** key. The display shows 200.00.
2. Press **R1/R2** to toggle between **R1** (200g/0.01g) and **R2** (250g/0.1g).
3. To confirm your capacity selection, press **ON/OFF**.
4. After 2-3 seconds the display shows 'PASS' then returns to normal weighing mode.

Auto-Off

By default, the scale will auto-off after 180 seconds of inactivity. To change or disable the auto-off timer:

1. Turn the scale off and hold down the ZERO/TARE key: At the same time quickly press the ON/OFF key 6 or 7 times until the display shows the current auto-off setting (eg180), then release.
2. Press the MODE key to change the auto-off setting between 180 sec, 120 sec, 60 sec, or 0 sec (auto-off disabled).
3. Press ON/OFF to confirm.
4. The display shows „PASS“ and setting is complete.

Accuracy

Your scale comes calibrated from the factory and most users won't need to calibrate their scale for a long time. You should check your scale's accuracy periodically using the included calibration weight and perform calibration if necessary. **NOTE:** Only use trusted calibration weights.

Calibration

Calibration is permanent and should only be performed using accurate calibration weights that match the requirements for your scale. Do not use coins or other household items to calibrate.

1. Prepare the required weight for your scale (see chart on back).
2. Find a flat and level surface to perform calibration and let the scale adjust to room temperature.
3. Press **ON/OFF** to turn the scale on and wait for the scale to stabilise.
4. Press and hold the **MODE** key until the display shows 'CAL'.
5. Make sure nothing is on the platform then press **MODE** to confirm the zero point.
6. The display will then flash the required calibration weight.
7. Place the required calibration weight on the centre of the platform.
8. The display will show 'PASS' then return to normal weighing mode. Calibration is complete. You may now turn the scale off or continue weighing.

Specifications

G-202

Max Capacity	R1 200g R2 250g
---------------------	----------------------------------

Readability	R1 0.01g R2 0.1g
--------------------	-----------------------------------

Weighing Units	g, oz, ozt, dwt, ct, gn
-----------------------	-------------------------

Dimensions	Overall:	163 mm x 74 mm x 23 mm
	Platform:	98 mm x 66 mm
	Display:	35 mm x 17 mm

Calibration Weight	100g (included)
---------------------------	-----------------

Power	2 x AAA Batteries (Sold Separately) 5V/1000mA Micro-USB Adapter (Sold Separately)
--------------	---

Warranty	1 Year Limited
-----------------	----------------

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