

★ Storage

Store at Room Temperature.

Contents

- Product Manual
- Ponceau S Solution

ALL PRODUCTS SOLD BY GenDEPOT ARE INTENDED FOR RESEARCH USE ONLY UNLESS OTHERWISE INDICATED. THIS PRODUCT IS NOT INTENDED FOR DIAGNOSTIC OR DRUG PURPOSE

★ Shipping Condition

Ship at ambient.

★ Procedure

1. Wash the blot once with Ponceau S Solution.
2. Add fresh Ponceau S Solution and incubate for 5-15 min at RT with agitation.
3. Destain the blot in multiple changes of PBS.

Alternatively the blot can be destained in 5% acetic acid if destaining is difficult in PBS.

★ Related Products

| Product Name | Cat No |
|---|--------|
| Albumin, Ultra pure Bovine Serum Albumin | A0100 |
| Xpert Protease Inhibitor Cocktail Solution (100X) | P3100 |
| Xpert 2 Prestained protein marker | P8502 |
| Tween 20, Molecular Biology Grade | P9100 |
| West-Ez Blocking Buffer, 3% BSA | W3710 |
| West-Q Chemiluminescent Substrate Kit | W3650 |
| West-Q Chemiluminescent Substrate kit, Plus | W3651 |
| West-Ez Stripping Buffer | S2100 |