
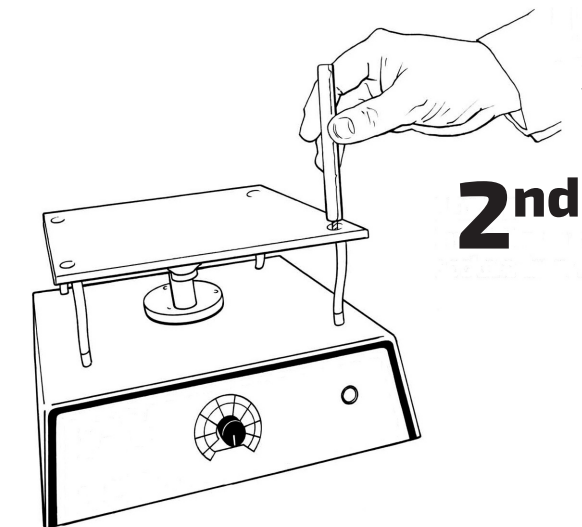



The Belly Dancer[®] 2nd Story Platform Installation

Warning: Prior to starting, ensure unit is unplugged from electrical source

First, remove the four screws attaching The Belly Dancer[®] platform to the flexible stabilizer legs at each corner. Use the provided hex wrench enclosed in the kit, save the detached screws for later use. Do NOT remove threaded standoffs from bottom of acrylic platform.

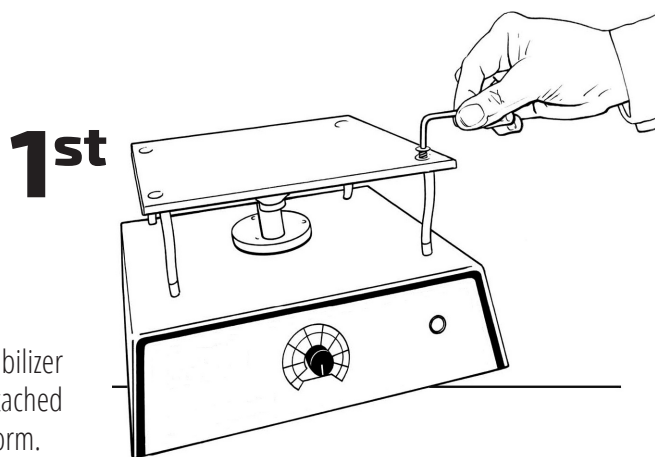

Second, install the aluminum spacing rods by threading them into the existing standoffs (through the acrylic platform) in which the screws were just removed. Hand-tight is fine for these items.

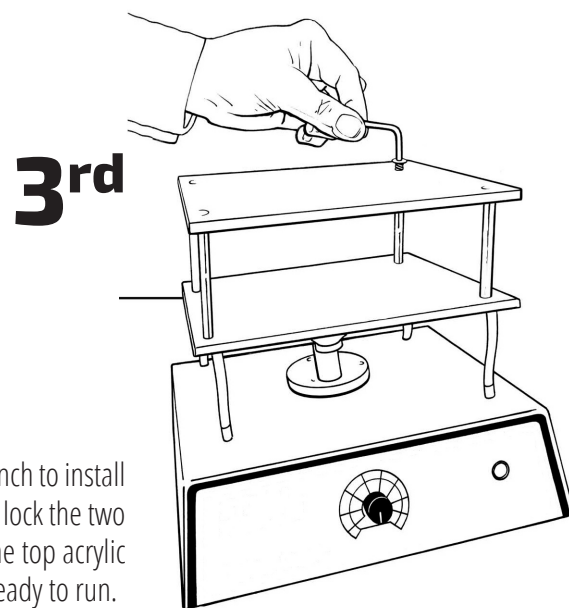

Finally, align the new 2nd story platform with the four spacer legs. Use the provided hex wrench to install the screws, previously removed in step 1. Tighten all four screws, in all four corners, this will lock the two platforms together. Replace the bottom silicone mat, and place the new supplied mat on the top acrylic platform. Reconnect your Belly Dancer[®] unit to an electrical power source and you will be ready to run.



SCIENTIFIC

7445 Chavenelle Road • Dubuque, IA 52002
(800) 253-4942 • (563) 690-0484

ibisci.com

Building
Quality
Scientific.