


ASSEMBLY INSTRUCTIONS SIGNATURE SCIENCE LAB TABLE (SLT4 / SLT5 SERIES)


TABLE COMPONENTS

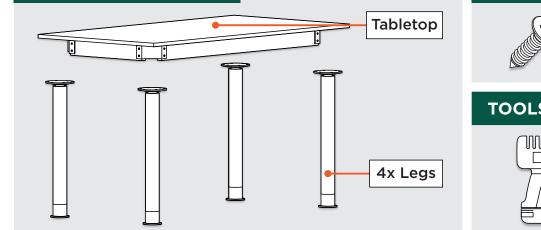






24x Screws

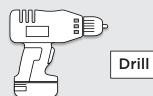

LEG INSTALLATION


Secure the legs to tabletop by installing 6 screws through the pre-drilled holes on the tabletop. Ensure that the numbers on the leg is facing outwards.

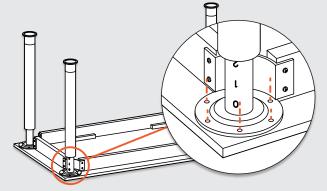

If a screw is difficult to secure tightly, uninstall the screw and carefully deepen the predrilled hole slightly, using the correct size drill bit.

HEIGHT ADJUSTMENT

Adjust the height by raising the table and it will automatically lock into place. To lower the table, raise it to the top and after hearing the click, lower it to the bottom.

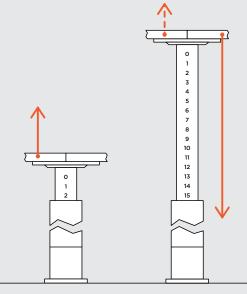



When moving the table - carry it. Do not drag or push it across the floor.

Congratulations, your table is now assembled and ready to be used!


Dear Customer, Thank you for your purchase. Although extreme care goes into manufacturing & packaging of this product, mistakes are possible. If you encounter any missing parts or difficulty in assembly, please contact our customer service Toll Free at **1-800-261-4112** or **info@nationalpublicseating.com** for courteous & immediate resolution to your problem.

BOOK COMPARTMENTS INSTALLATION INSTRUCTIONS

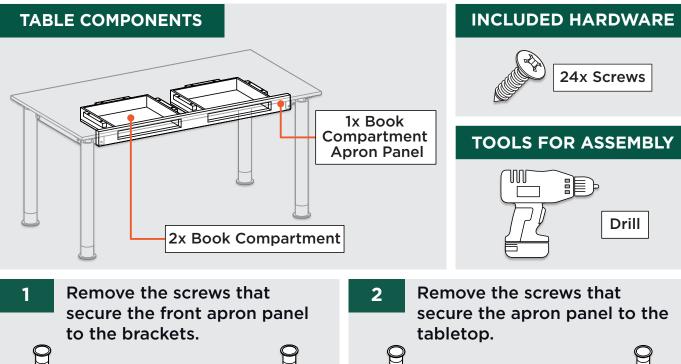


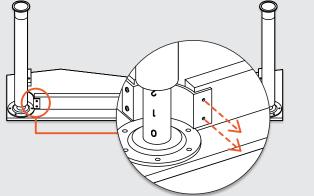

3 Place the book compartment apron panel in the location where the original apron was located and attach the brackets with two screws on each side.

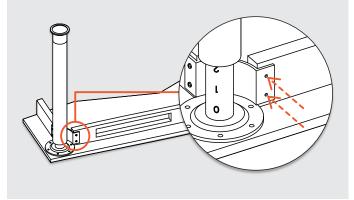

4 Secure the book compartment to tabletop by installing 6 screws through the holes of the book compartment reinforcement blocks.

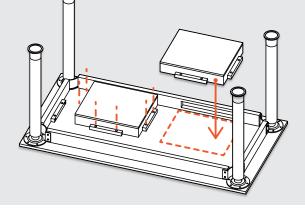

Congratulations, your book compartments are now assembled and ready to be used!