### WATER TEST KIT - TOTAL BACTERIA+ E. COLI

EN

AquaVial™ Plus is designed for testing well and ground water and consists of two tests: AquaVial™ Total Bacteria and AquaVial™ E. coli. AquaVial™ Plus is a quick and convenient screening test and does not replace mandatory water testing through an accredited laboratory.

AquaVial™ Total Bacteria detects bacteria and fungi concentrations as low as 500 CFU/ml (500 bacteria per ml) in 15 minutes at 35-40 °C (95-104 °F), or in 30 minutes at 20-25 °C (68-77 °F).

AquaVial™ E. coli detects E.coli and coliform concentrations as low as 1 CFU/ml (one bacteria per ml) of E. coli and coliform bacteria within 24 hours when incubated at 35-40 °C (95-104 °F), or 48 hours at 20-25 °C (68-77 °F).

#### KIT CONTENT:

- 1 T-Cup unit
- 1 AquaVial™ Total Bacteria vial with pink solution and red cap
- 1 AquaVial™ E. coli vial with yellow dry media and yellow cap
- 1 saline solution pouch
- 1 set of Testing Instructions

#### IMPORTANT SAFETY NOTES:

- The solutions in the vials are non-toxic and; therefore pose no significant hazard provided that the directions for use are followed. **DO NOT INGEST.**
- Not recommended for children (small parts)
- Store at temperatures above freezing. For best results, test before the expiration date indicated on the box

#### TIP:

Have a white background (i.e. a piece of paper) and a camera ready before starting the tests. This will help observing the colour change.

#### DISPOSAL:

T-Cup components, are recyclable. Please recycle whenever possible. Vials can be disposed of in the trash. Warning: Do not attempt to open the vials after test completion or dispose of their contents into your sink.

#### IMPORTANT:

Please read the directions carefully and watch the instructional video, before starting the test. Follow the steps below in the exact order to obtain an accurate reading of your water quality. Any deviation may cause cross contamination or ruin the sample, possibly resulting in false positive or false negative results.

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
##### Step 1: Collect and Concentrate Water Sample

**OPEN T-CUP:** Separate the top and bottom halves of the T-Cup by twisting the top half counter clockwise.

**ADD WATER SAMPLE:** Fill the bottom of the T-Cup halfway with the sample water.

**CONCENTRATE:** Twist top of the T-Cup clockwise into bottom section until it stops, forcing water through the integrated T-Cup filter. Water will pour through and spill over the sides of the plunger. This is normal. The T-Cup is intended to sample this exact amount of water.

**DISCARD:** Drain excess water from the top and sides of the T-Cup.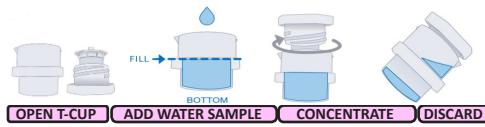

##### Step 2: Add Saline Solution


This step is required to ensure the water sample collected and concentrated in Step 1 is washed of chemicals, like chlorine, used to treat the water, which may affect your test results. The addition of saline solution will not dilute or interfere with the water sample to be tested. The T-Cup system has a built-in filter that retains the bacteria at the bottom, while eliminating chemicals through the filter.

**OPEN T-CUP:** Open the T-Cup by twisting the top half counter-clockwise.

**ADD SALINE SOLUTION:** Carefully tear the top section of the saline solution pouch and pour saline solution into the bottom half of the T-Cup.

**CONCENTRATE:** Twist the top of the T-Cup into the bottom clockwise to close the T-Cup.

**DISCARD:** Pour out any excess solution from the top of the T-Cup.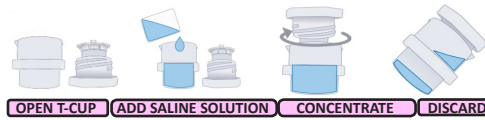


##### Step 3: Test Water Sample

**OPEN T-CUP:** Open the T-Cup by twisting the top half counter-clockwise.


**OPEN VIAL:** Hold AquaVial™ Total Bacteria vial (pink solution & red cap) into your hand and twist the red cap counter clockwise. **BE CAREFUL NOT TO TILT THE VIAL OR SPIL AQUAVIAL™ SOLUTION.**

**FILL AQUAVIAL™:** Pour the unfiltered water from the bottom of the T-Cup into the AquaVial™ Total Bacteria vial. In most cases the mixed liquid will fill to the 4 ML line (+/- 1ML). Sample water should not exceed the 5 ML line; or stop filling water in the vial after this line. Close AquaVial™ Total Bacteria with its original red cap. Tip: Take a picture of the vial against a white background.

**INCUBATE:** Let the AquaVial™ Total Bacteria vial incubate in an upright position for 30 minutes and check for changes in colour. If you are not sure if the vial changed color, let the vial incubate another 30 minutes and check color again.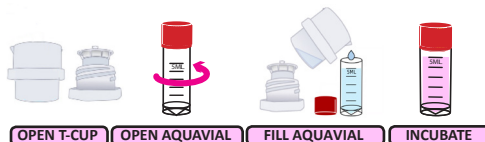

Compare the color of the reagent in the test vial against the color diagram below: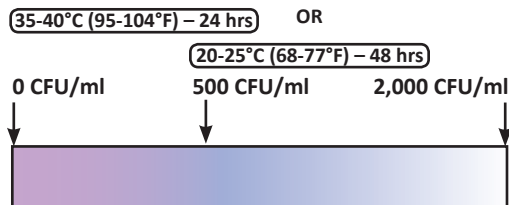

#### RESULTS INTERPRETATION:

If the solution changes its colour to clear, extreme caution should be exercised, and a sample should be taken to a laboratory for further testing. This test is intended for rapid screening only.

#### E. COLI AND COLIFORM TEST

STEP-BY-STEP VIDEO INSTRUCTIONS:  
[www.aquavial.ca/AquavialPLUS](http://www.aquavial.ca/AquavialPLUS)

##### Step 1: Test Water Sample

**OPEN AQUAVIAL™:** Twist the AquaVial™ yellow cap counterclockwise to open the vial.


**FILL AQUAVIAL™:** Fill AquaVial™ E. coli vial with sample water. Sample water should not exceed the 5 ml fill line; or stop filling water in the vial after this line. Twist clockwise to close the vial. **DO NOT LEAVE THE VIAL OPEN.**

**SHAKE:** Shake the vial well to dissolve the dry media deposited on the side wall of the vial. **Tip:** Take a picture of the vial against a white background.


**INCUBATE:** Let the AquaVial™ E. coli vial incubate in an upright position for 24 hours at 35 - 40°C (95 - 104°F) or 48 hours at 20-25°C (68 - 77°F) and check for changes in colour. Compare the color of the reagent in the test vial against the colour diagram below:

#### WARNING:

Do not attempt to open the vials or dispose of their contents into your sink after test completion.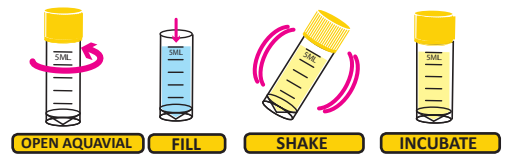

35-40°C (95-104°F) – 24 hrs      OR      20-25°C (68-77°F) – 48 hrs

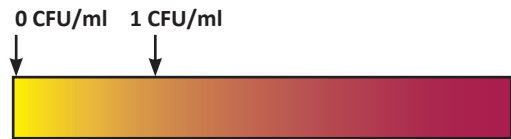#### RESULTS INTERPRETATION:

If the solution changes its colour to orange or purple, extreme caution should be exercised, and you should not come into contact, ingest or shower with this water, until further testing has been conducted and additional analysis performed. This test is intended for rapid screening only.

**IMPORTANT: Any colour change after 48 hours should be disregarded.**

The test is highly sensitive and even dead (unharmful) bacteria can change the colour after more than 48 hours.

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