

SMART Spectro® 2. QUICK START GUIDE

1. Press ON . The LaMotte logo screen will appear for about 2 seconds. After initialization the MAIN MENU will appear.	MAIN MENU 12:00:00
	PROGRAMMED TESTS %T/ABS PC LINK SYSTEM SETUP

2. Press ▲ or ▼ to scroll to PROGRAMMED TESTS .	MAIN MENU 12:00:00
	PROGRAMMED TESTS %T/ABS PC LINK SYSTEM SETUP

3. Press *ENTER to select PROGRAMMED TESTS .	PROGRAMMED TESTS 546 NM
	SEQUENCE 1 SEQUENCE 2 SEQUENCE 3 ALL TESTS EDIT

4. Press ▲ or ▼ to scroll to ALL TESTS .	PROGRAMMED TESTS 546 NM
	SEQUENCE 1 SEQUENCE 2 SEQUENCE 3 ALL TESTS EDIT

5. Press *ENTER to select ALL TESTS .	ALL TESTS 546 NM
	2 ALUMINUM 3 AMMONIA-N L F 4 AMMONIA-N L S 5 AMMONIA-N H 7 BIGUANIDE 8 BORON 9 BROMINE-LR 12 CADMIUM

6. Press ▲ or ▼ to scroll to the desired test.	ALL TESTS	546 NM
	1 ALUMINUM	
	2 ALKALINITY-UDV	
	3 AMMONIA-N L F	
	4 AMMONIA-N L S	
	5 AMMONIA-N H	
	6 ARSENIC	
	7 BIGUANIDE	
	8 BORON	

7. Press *ENTER to select the test.	2 ALUMINUM	535 NM
	SCAN BLANK	
	SCAN SAMPLE	
	TEST RESULT LIST	

8. Insert the blank. Close lid. Press *ENTER to scan the blank. BLANKING will momentarily appear on the top line of the display. The blank has been stored.	2 ALUMINUM	535 NM
	SCAN BLANK	
	SCAN SAMPLE	
	TEST RESULT LIST	

9. Insert the reacted sample. Close lid. Press *ENTER to scan the sample. The result will be displayed.	2 ALUMINUM	535 NM
	<p>T = 13.5%T</p> <p>A = 0.8707A</p> <p>C = 0.26 ppm</p> <p>STORE PRESS [ENTER]</p>	

After obtaining test results, scroll with ▲ or ▼, and make another selection with *ENTER. Press *ENTER to store the result to the Test Result List. If Store Method is automatic test results will be stored automatically. To print the test results, scroll to Test Result List. Press *ENTER to print.

After obtaining test results, scroll with ▲ or ▼, and make another selection with *ENTER. Press EXIT to escape to previous menu.



802 Washington Ave Chestertown MD 21620
800-344-3100 410-778-3100 f 410-778-6394
www.lamotte.com