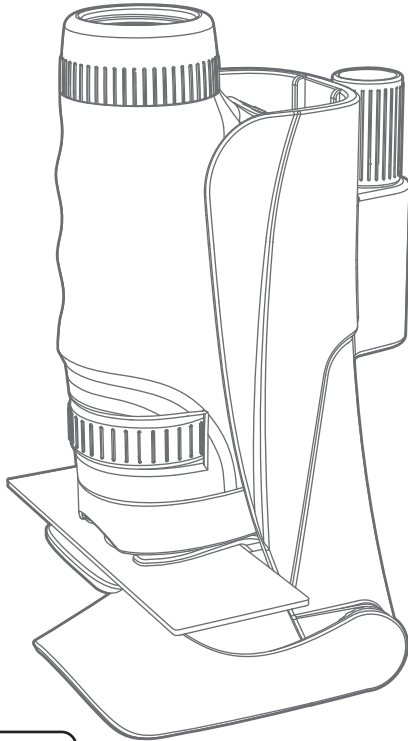
PORTABLE SCIENCE MICROSCOPE

STEAM
EDUCATIONAL PRODUCT

8+
Year

 **WARNING:** CHOKING HAZARD-SMALL PARTS,
NOT SUITABLE FOR CHILDREN UNDER 3 YEARS.

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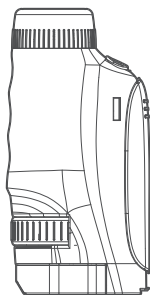
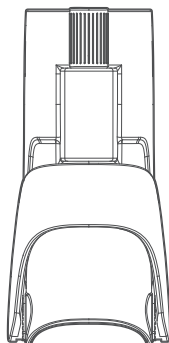
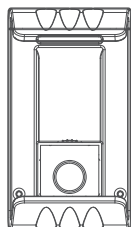
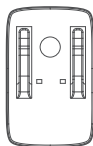
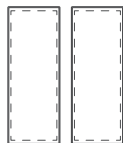
EN:

WARNING:

All packaging materials, such as tapes, plastic sheets, binding wires and labels, are not part of the product and are not suitable for children to play with. For the safety of your child, please handle packaging materials properly. The user manual contains important information, please keep it for later use.

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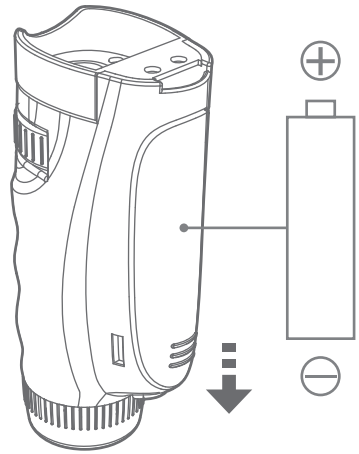
EN:

BATTERY INSTALLATION:

1. Confirm the machine is turned off.
2. Uncover the battery box as shown in the diagram.
3. Install 1 new 1.5V AAA (LR03/7) battery with correct polarity (+, -), as shown in the diagram on the right.
4. Cover the battery box.

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EN:

NOTE FOR BATTERY USAGE:

1. Alkaline, carbon and zinc battery can be used. Do not mix alkaline standard(carbon-zinc), or rechargeable(nickel-cadmium)battery.
2. Non-rechargeable battery are not to be recharged;
3. For electric toys using rechargeable battery, the battery should be charged under adult supervision.
4. The rechargeable battery should be removed from the toy before charging.
5. Different types of battery or new and used battery are not to be mixed;
6. battery are to be inserted with the correct polarity(+and-);
7. Exhausted battery are to be removed from the toy;
8. The supply terminals are not to be short-circuited.
9. Remove the battery when not in use for a long time.
10. Don't put the battery in the fire.

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EN: 1. Hand-held Observation

Step 1: Turn on the LED lamp, and look through the eyepiece viewing port while aiming the object lens at the object of observation;

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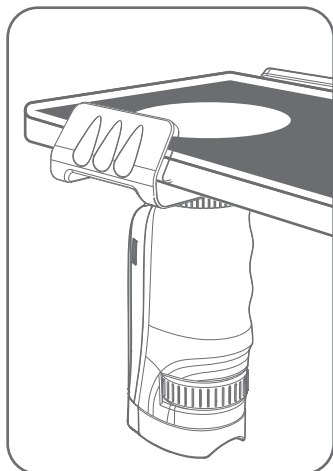
EN: Step 2: Put the microscope close to the observed object, turn on the mobile camera, and manually adjust the zoom wheel until the screen picture becomes clear. Take pictures.

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EN: 3. Desktop Observation

Step 1: Open the folding support, and attach the microscope to the support;

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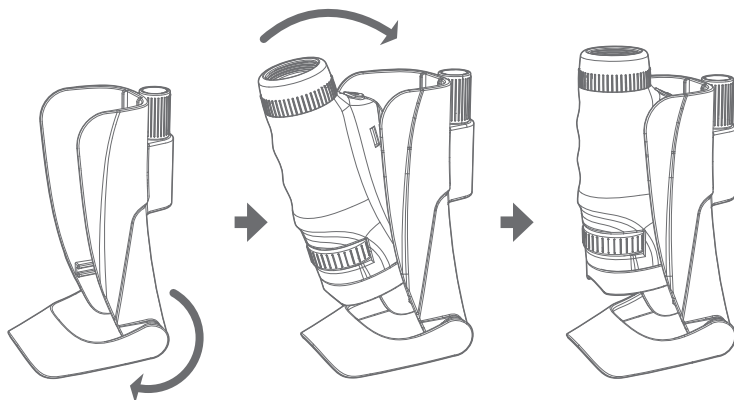
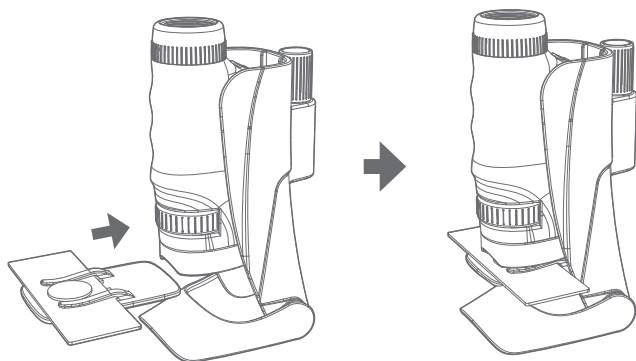
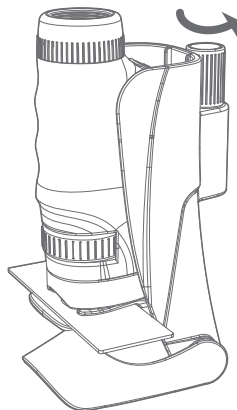
EN: Step 3: Adjust the knob on the back of the support so that the microscope reaches the bottom, and turn on the LED lamp;

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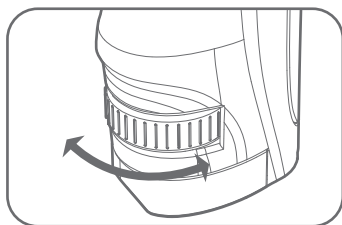
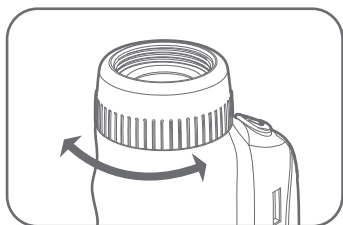
EN: Step 4: Manually adjust the zoom wheel until the screen picture becomes clear.

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EN:

MAINTENANCE:

1. Wipe the surface of the main unit gently with soft dry cloth to keep it clean.
2. Keep away from direct sunlight or heat source.
3. Please take out the battery if the device is to be left unused for a long time.
4. Do not bump against a hard object or try to disassemble the main unit.
5. Do not immerse the main unit in water.

COMMON TROUBLE SHOOTING:

If the main unit stops responding for some reason, please:

1. Shut down the main unit;
2. Take out the battery and disconnect power supply;
3. Install the battery again after a few minutes;
4. Restart the device for use.
5. Replace with new battery if still not working.

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©2020 TOPBRIGHT ANIMATION all rights reserved. MADE IN CHINA
Manufacturer: Jiangsu TopBright Wisdom Information Technology Co.Ltd TEL: +(86)0571-86879391
TOPBRIGHT GmbH, Brunnengasse 65, 90402 Nürnberg, Germany