

Receiver Plates

1 Contents

Receiver Plates 10 µm	
REF	740989.4
Receiver Plates 10 µm	4
Receiver Plates 35 µm	
REF	740512.4
Receiver Plates 35 µm	4
Receiver Plates 35 µm hydrophilized	
REF	740513.4
Receiver Plates 35 µm hydrophilized	4
Receiver Plates 50 µm	
REF	740688.4
Receiver Plates 50 µm	4
Receiver Plates 50 µm hydrophilized	
REF	740689.4
Receiver Plates 50 µm hydrophilized	4

2 Product description

Receiver Plates are 96-well plates available with inserted filter frits of 10 µm, 35 µm, or 50 µm pore size. They can be used for general filtration purposes as well as for retaining chromatographic resins (e.g., Protino® Ni-NTA Agarose, Protino® Glutathione Agarose 4B, Sephadex® G25, G50, or Sephacryl® S200). The Receiver Plates can be used under centrifugation using standard benchtop centrifuges as well as under vacuum. The plates are also offered with hydrophilized filter frits which enable the plates to be used additionally with gravity flow.

3 Specifications

Plate material:	Polypropylene
Filter frit material:	Polyethylene untreated (for use under centrifugation and vacuum)
	Polyethylene hydrophilized (for use under gravity, centrifugation, and vacuum)
Filter frit pore size:	approx. 10 µm, 35 µm, or 50 µm
Filter frit diameter:	8.3 mm
Effective filtration area:	54.1 mm ²
Volume capacity per well:	1.5 mL

4 Safety instructions

Receiver Plates do not contain hazardous contents.

5 Product use restriction / warranty

Receiver Plates are developed, designed, and sold for research purpose only. All goods have to be checked immediately upon receipt. Damage claims are only acceptable in writing within 8 working days of receipt of all the goods. In case of legitimate claims the buyer can only require replacement of the goods. If replacement is not possible, the buyer has the right to choose between alternative products with same value or refund. The buyer cannot claim further compensation. All returns must first be authorized by us in writing. We offer a one year warranty from date of delivery that our products will conform to applicable specifications set forth in the product specifications if not sold to persons set forth in § 13 of German Civil Code. In such a case, the provisions of the German Civil Code shall be valid. This warranty does not cover defects in or damage to the products which are due to improper installation or maintenance, misuse, neglect or any use other than ordinary commercial application.

Any discharge from liability will be void if a defect results from a heavily negligent or intentional breach of contract or if the buyer will be bodily injured as a result from a negligent or intentional breach.

The seller shall not be liable for the goods being fit for a particular purpose unless otherwise agreed upon, to which the buyer intends to put them.

This warranty is strictly exclusive. Any further damage compensation is impossible.

MACHEREY-NAGEL MAKES no other warranty of any kind whatsoever, and SPECIFICALLY DISCLAIMS AND EXCLUDES ALL OTHER WARRANTIES OF ANY KIND OR NATURE WHATSOEVER; DIRECTLY OR INDIRECTLY; EXPRESS OR IMPLIED; INCLUDING; WITHOUT LIMITATION; AS TO THE SUITABILITY; REPRODUCIBILITY; DURABILITY; FITNESS FOR A PARTICULAR PURPOSE OR USE; MERCHANTABILITY; CONDITION; OR ANY OTHER MATTER WITH RESPECT TO MACHEREY-NAGEL PRODUCTS OR THE SALE OF MACHEREY-NAGEL PRODUCTS.

If not superseded by these terms the current terms and conditions for the sale of goods (export version) are additionally valid.

Trademarks:

Protino is a registered trademark of MACHEREY-NAGEL GmbH & Co. KG

Sephadex and Sephacryl are registered trademarks of GE Healthcare companies

MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6–8 · 52355 Düren · Germany
Tel.: +49 24 21 969-0 · Fax: +49 24 21 969-199 · tech-bio@mn-net.com · www.mn-net.com