



# AccuClean<sup>®</sup> Advanced Rapid Protein Residue Test

## AccuClean<sup>®</sup> Advanced Rapid Protein Residue Test

AccuClean Advanced delivers industry-leading speed and sensitivity for determining the cleanliness of food contact surfaces and equipment. AccuClean Advanced is a simple, 10-second visual test that reveals detected protein residue through an easy-to-interpret color change. With a protein detection limit of 10  $\mu\text{g}$ , the test provides sensitive, reliable results.

### Features

- ✓ **Fast:** 10 seconds to result
- ✓ **Sensitive:** As low as 10  $\mu\text{g}$  protein per sample
- ✓ **Simple:** No equipment required



# AccuClean® Advanced

Note: Please read kit instructions completely before performing test.


## How it works

AccuClean Advanced determines cleanliness by detecting any protein residues left behind from previously processed food and liquids. If no protein is detected, the solution in the bottom of the cup will remain copper-colored to indicate the surface is clean. A color change to gray/green or blue indicates protein has been detected from the surface sample



### Sample

Remove the plunger portion of the device and place the end with the pad on the surface to be tested. To sample, crisscross a 4 x 4 inch (10 x 10 cm) portion of the surface. For best results, sample in a back-and-forth motion within the sampling area as shown.

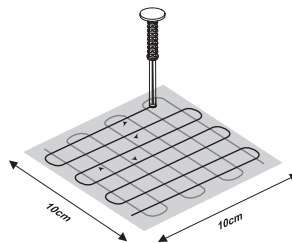

### Activate

Return the plunger to the cartridge and activate by pressing it fully into the cartridge. Gently swirl the cartridge for ten seconds.



### Interpret

Read by examining the color of the remaining solution using the provided color comparator.



Call **800/234-5333** to order or visit us online at  
[foodsafety.neogen.com](http://foodsafety.neogen.com)

#### #9960 AccuClean Advanced Rapid Protein Residue Test

Materials Provided:  
100 AccuClean Advanced Samplers  
1 Kit Insert  
1 Color Comparator



800-234-5333 (USA/Canada) or 517-372-9200  
[foodsafety@neogen.com](mailto:foodsafety@neogen.com) • [foodsafety.neogen.com](http://foodsafety.neogen.com)

