

SE572

Installation, Maintenance & Operation Instructions

Manufactured in Australia to comply with AS4775-2007.

DESCRIPTION

Deck mounted EMERGENCY AERATED EYE WASH. Supplied with 15mm stay-open full flow ball valve, activated by stainless steel push hand flag handle with (2) spray outlets with flow restrictor, and flow off dust covers. Unit swivels 90° for easy access to sink when not required, and mounts on counter tops up to 1-1/4" thick.

SPECIFICATIONS

SPRAY HEADS:Aerated Stainless Steel.VALVE:15mm 316 Stainless SteelACTIVATOR:Stainless Steel Push Handle.

INSTALLATION

At desired location drill a 25mm (1") diameter hole through countertop (1-1/4" thick max.) Be sure the hole is drilled as straight as possible. Attach eyewash to countertop using supplied hardware, and connect a water supply to unit that is a 12mm minimum uninterrupted source of potable water (use pipe sealant on connection). Failure to use potable water can result in emergency units producing impure or contaminated water causing possible further injury.


TESTING PROGRAM


The AS 4775 - 2007 states that all safety emergency equipment shall be activated on a weekly basis to flush the line and verify proper operation. All units furnishes a testing tag (SE Tag). On this tag the date of the inspection and the inspectors initials should be noted.

<u>NOTE</u>: This unit should be connected to an uninterrupted source of potable water, with a minimum Flowing pressure of 30 Kpa and a maximum static pressure of 125 PSI.

OPERATION


Be sure unit is over the sink before activating. The eyewash is activated by pushing the push handle backwards 90°. Once the unit is activated, water will discharge from the spray outlets until the push handle is manually pulled forward to the off position.

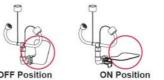

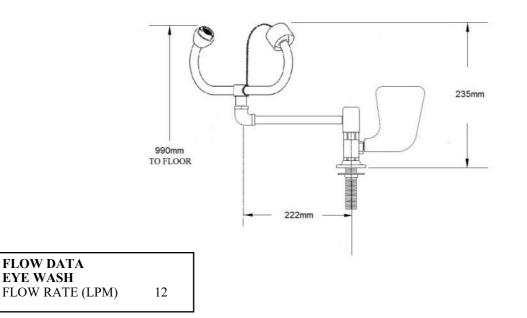

WARNING : Use only genuine Speakman parts when replacing or repairing components. When ordering quote correct part numbers listed.

BIG Safety PO Box 2117 Normanhurst NSW 2076 12A/10 Pioneer Ave, Thornleigh NSW 2120 P: 1300 73 22 74 E: sales@bigsafety.com.au www.bigsafety.com.au

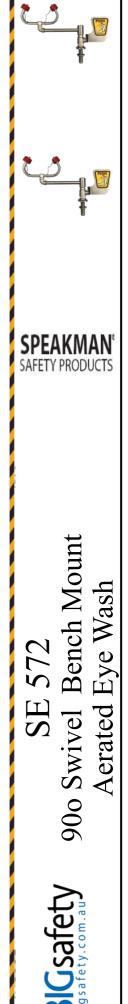

Installation, Maintenance & Operation Instructions

MAINTENANCE

1

3

Maintenance should be carried out weekly and yearly as detailed in the schedule. Should you need to repair this unit, please reference the spare parts list for the correct part numbers. Before any maintenance is done be sure the water supply is shut off.

MAINTENANCE SCHEDULE

| WEEKLY | Activate eyewash unit using push handle | |
|--------|---|--|
| | Allow water to flow for 30 seconds to flush line. | |
| | Ensure water flows evenly from both outlets. | |
| | Ensure water is clean. | |
| YEARLY | Ensure unit complies with AS 4775-2007 | |

Part No: SA 01003




Complete Bull Horn Assembly

Part No. RPG05 1177 Aerator Screen & Cap Assembly



Sold as a packet of 6.

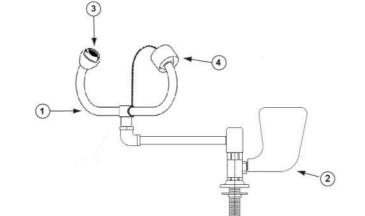

SPARE PARTS LIST

| Item | Description |
|------|-------------|
| | |

- 1. 2. 3. Complete Aerated Bullhorn Assy.
- 15mm Ball Valve & Push Handle Aerator Screen & Cap Assembly
- 4. Red Dust Cap & Chain

| Part No |
|------------|
| SA 01003 |
| SE 910V |
| RPG05 1177 |
| SA 01008 |
| |

WARNING : Use only genuine Speakman parts when replacing or repairing components. When ordering quote correct part numbers listed.

BIG Safety PO Box 2117 Normanhurst NSW 2076 12A/10 Pioneer Ave, Thornleigh NSW 2120 P: 1300 73 22 74 E: sales@bigsafety.com.au www.bigsafety.com.au

