
Digital Microscope AD1605 User Manual

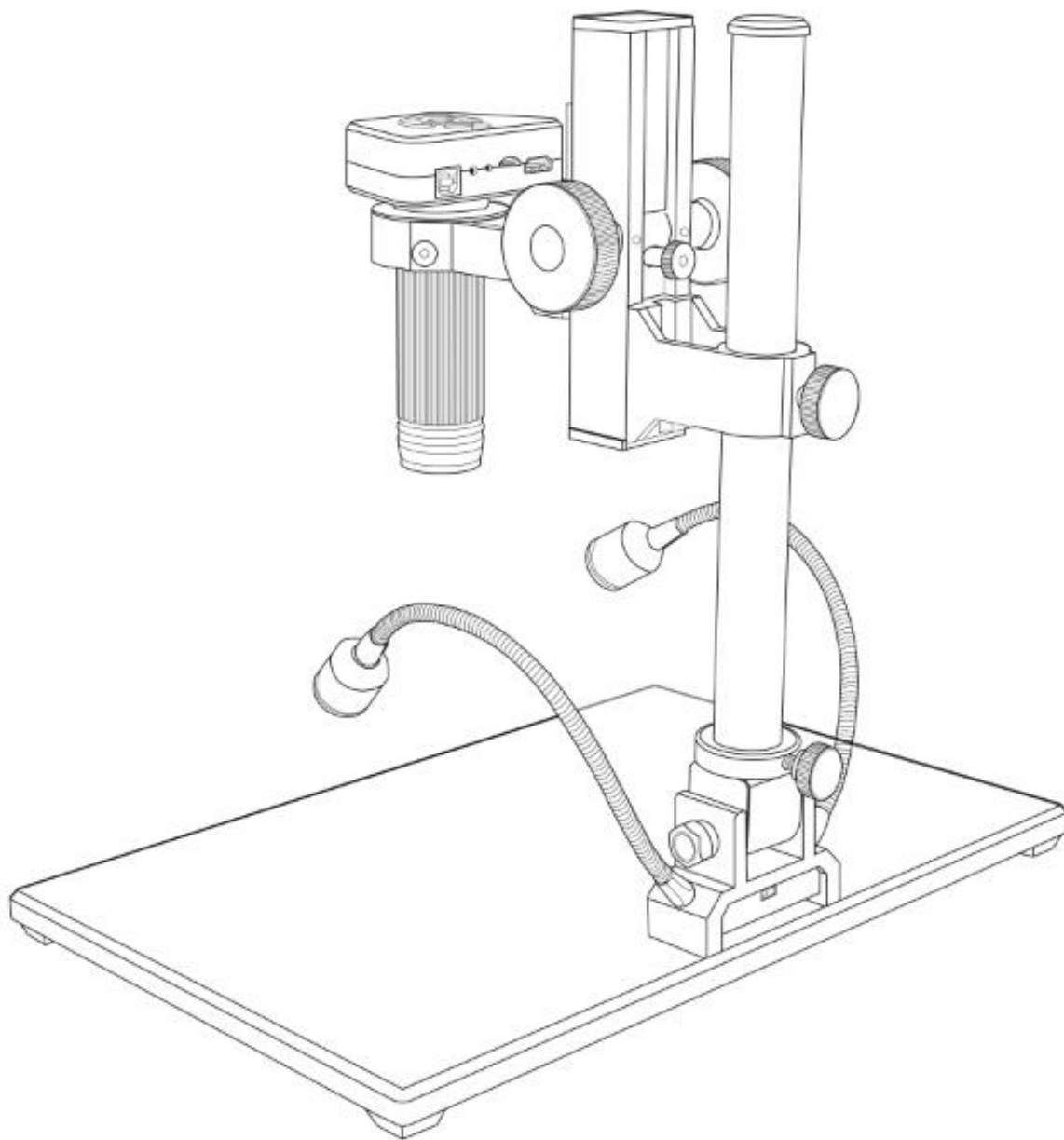Digital Microscope AD1605 User Manual-----1-9

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LIST OF CONTENTS

Preface	1
Basic Parameters	2
What's in the box	2
Assembly	3
Button Instructions	4
Remote controller	5
Connection and output	6
Quick Start	7
◆ Independent use	7
◆ HDMI output	8
FAQ	8

Preface

AD1605 is a multifunctional digital industrial microscope. It integrates 4K high-definition output, high object distance, multi-function and convenient operation, which fully meets the various needs of industrial maintenance, teaching demonstration and other fields. Its exquisite appearance design took ergonomics into consideration, and the high-quality materials provide users with a more comfortable experience. We hope this product can truly free your eyes and bring improvements to your work and life. Thank you for your support of this product.

◆In order to use this product better, we hope you can keep this manual forever. Any damage caused by non-compliance with this manual will be invalidated.

◆In order to ensure a better experience for customers, we may make small changes to the product in the process of perfecting. Where the description differs from the product, please refer to the actual product. We hope that consumers will understand.


◆Safety and Maintenance Instructions:

1. Keep the equipment in a clean and dry environment. Keep away from oil, water, gas and other materials that may cause corrosion. Storage and operation can refer to general electronic equipment, such as laptop computers.
2. The product contains small parts and the overall volume is relatively large, please keep children away from this product without a guardian to avoid accidental ingestion of parts or accidental shaking.
3. When the image quality has declined, the protective lens under the main body lens should be cleaned. Use a soft clean cotton cloth (such as glasses cloth) dipped in a proper amount of alcohol and wipe it carefully.
4. If you need to wipe the microscope with a cotton cloth dampened with water, please be sure to cut off the power first. Please be sure to check whether the machine has returned to dryness before turning on the power again.
5. People with disabilities, mental illnesses or people with allergic symptoms are not suitable to use this product.
6. Don't overdo it. The adjustments between the various parts of the product are limited. To ensure the best working condition, please keep the overall balance of the equipment.

Basic Parameters

Image sensor	14 Mega Pixels HD Sensor
Video output	1920x1080 60 FPS; 1920x1080 30 FPS 2560x1440 30 FPS; 3840x2160 30 FPS
Video format	MP4
Magnification Ratio	Up to 150 times (output to 29-inch high-definition display)
Photo resolution	MAX 14M (4320*3248)
Photo format	JPG
Focus distance	5CM-29CM
Frame Rate	Max 60f/s
Storage	Micro-SD card, up to 32G (not included)
Video output interface	HDMI
Power source	5V DC
Stand size	37.5cm*25cm*47cm
Package Included	Microscope ×1 Bracket ×1 Adapter ×1 Power cable ×1 HDMI cable ×1 User manual ×1 UV filter×1 IR remote×1
Packing data	40×30×16cm/ 5.5kg

What's in the box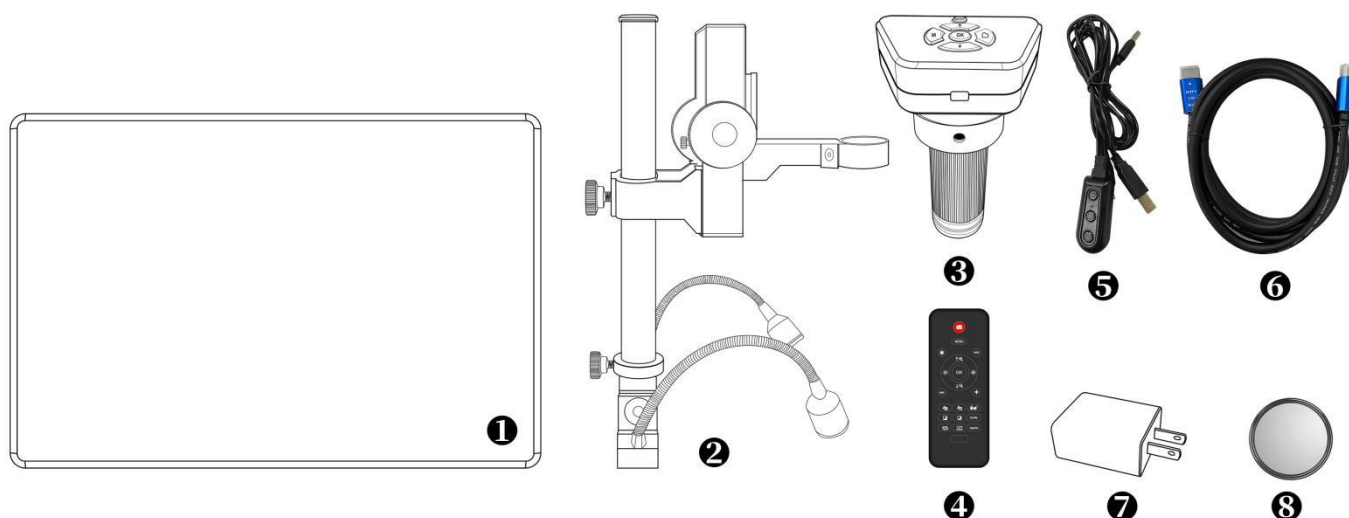

- ① Metal base
- ② Bracket
(including: universal LED light*2, beam, buffer ring)
- ③ Digital microscope host
- ④ Remote control

- ⑤ Power cable
- ⑥ HDMI cable
- ⑦ Power adapter
- ⑧ UV filter

Assembly

1. Install the bracket to the metal base (see Figure 2 Installation 1).
2. Raise the bottom white buffer ring (see Figure 2①) to the below of the beam (see Figure 2②), leaving some gaps, and tighten the screws to fix it (see Figure 2⑩). This step plays a buffering role for the beam to fall when it is not fixed properly, avoiding direct fall to damage the host.
3. Unscrew the top cover of the bracket (see Figure 2③), insert the black beam part (see Figure 2②) into the bracket (see Figure 2④), align the direction of the host hole (see Figure 2⑥), and tighten the rear fixing screws. After completing this step, turn the top cover of the bracket back and tighten it.
4. Insert the host into the host hole, adjust the position and direction of the host, and tighten the host fixing screw (see Figure 2⑦). Be careful not to jam the screw on the threaded focusing tube. After completion, try to rotate the focusing tube left and right to ensure that there is no jam.

After installation, if the operation is correct, the angle between the focusing tube, the bracket and the metal base can be adjusted freely. If there are any problems during the installation process, please contact us in time to avoid damage to the product and affecting normal use due to improper installation.

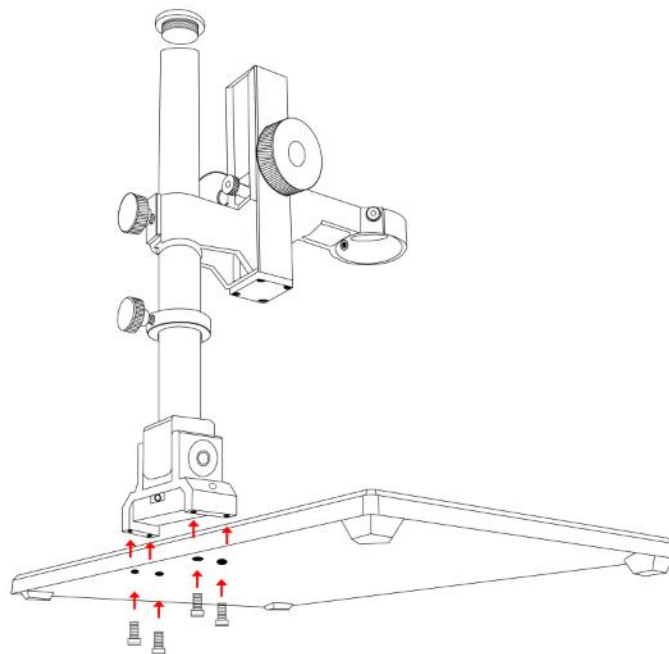
Figure 2 Installation 1



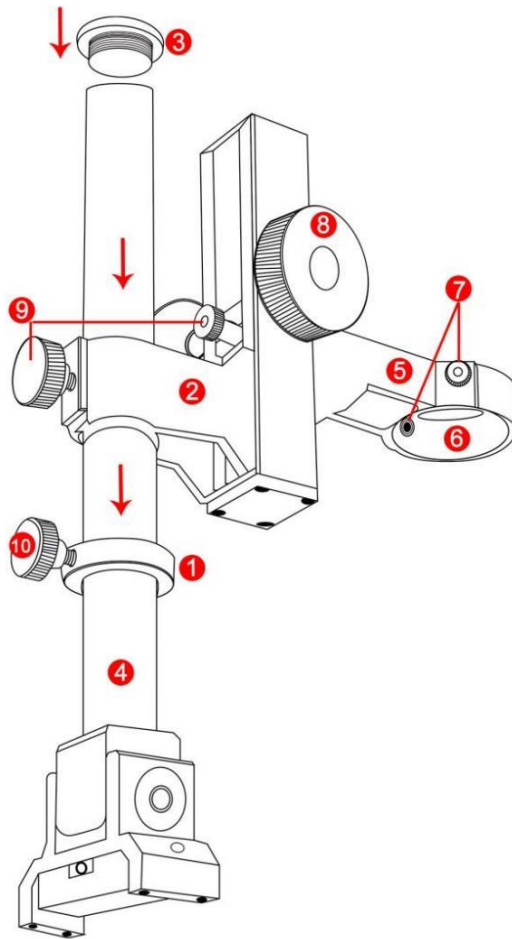
Figure 2 Installation 2

- ① Buffer ring ② Large beam ③ Bracket top cover ④ Bracket ⑤ Small beam
- ⑥ Host hole ⑦ Host fixing screw ⑧ Object distance adjusting pulley
- ⑨ Beam fixing screw ⑩ Buffer ring fixing screw

Button Instructions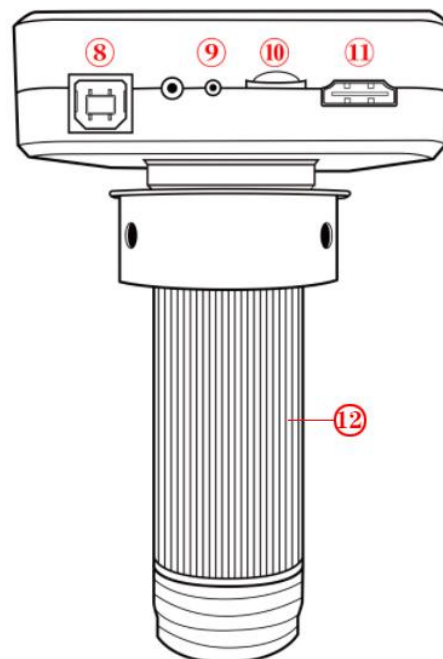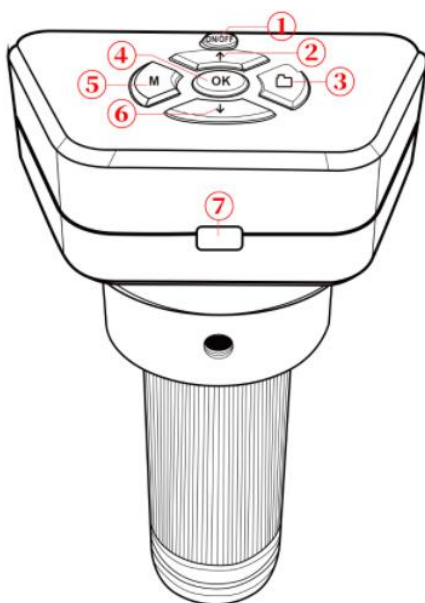

- ① Power on/off
- ② Up
- ③ Picture capture
- ④ OK

- ⑤ Menu/Mode
- ⑥ Down
- ⑦ Infrared sensor
- ⑧ Power interface


- ⑨ Power Indicator
- ⑩ Micro-SDcard slot
- ⑪ HDMI port
- ⑫ Focusing tube

Remote controller


- *Frozen: Keep the image on the screen. It works at standby, record and capture modes.
- *Brightness up&down: exposure compensation from +2.0 to -2.0.
- *Sharpness up&down: there are three levels of sharpness: Soft/Normal/Strong. Those two buttons are both used to switch the levels, but in different orders.
- *Contrast up&down: same principle as Sharpness.
- *Inverse: refer to pictures in below:

Normal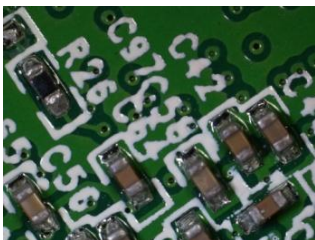



Inverse

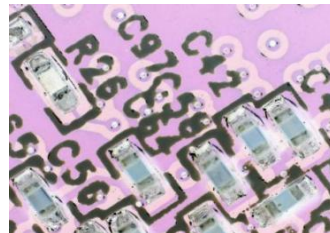- *Black&white: refer to pictures in below:

Normal



Black&white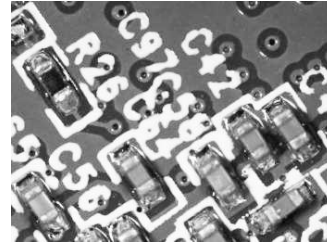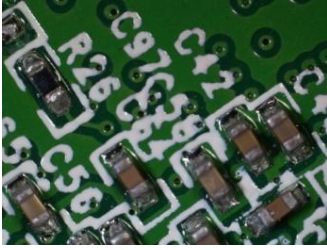

Connection and output

AD1605 has two output methods: HDMI and memory card.

The picture below shows the different ports on the host. The details of the HDMI output method will be detailed in the following chapters.

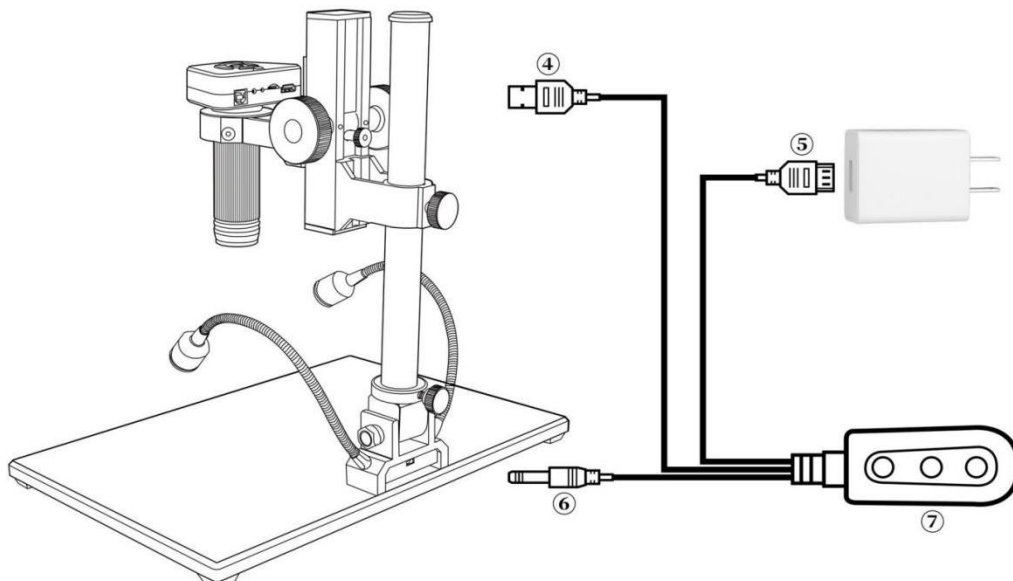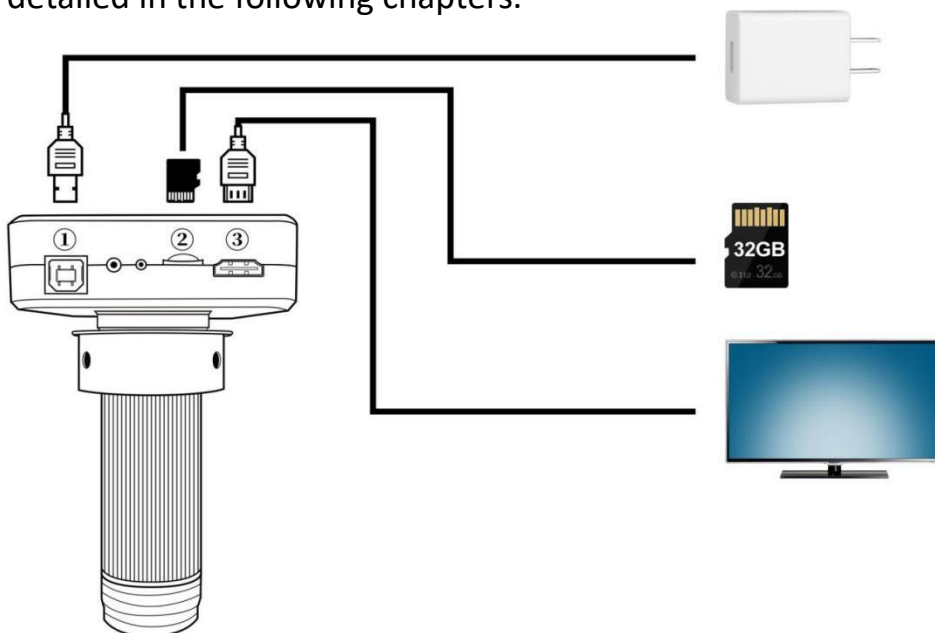①: Power interface.

- ②: Memory card. (Used to store photos or videos)
- ③: HDMI output. (Output to HDMI display screen through HDMI cable)
- ④: Microscope power cable.
- ⑤: USB power plug.
- ⑥: Power supply cable for LED lights on the base.
- ⑦: Switch. There are three buttons: the left is the "on/off" button to disconnect the device or connect the power; the other two buttons are the base lights brightness adjustment buttons, the middle is "Brightness +", and the right is "Brightness -".

Quick Start

◆Independent use

*How to turn on

1. Connect the three cables to the positions according to the diagram in the "Connection and Output" chapter ④⑤⑥, connect the HDMI cable to output the host image to the HDMI high-definition display, and keep the power supply in the power-on state.
2. Press and hold the Power button on the host (After connecting the power supply, you need to manually press the power button to turn on the microscope.)
3. Adjust the brightness and angle of the LED lights on the base, and adjust the object distance and the focus manually to obtain the desired image.

* Mode switch

1. Short press the "M" key to switch between different modes. (Or press the "MODE" button on the remote control)
2. Interpretation of icons in different modes:
: Video Standby mode;  : Photo mode;  ● : Video mode;
 :Playback mode (when the file is a video, press "OK" to play the video file.During video playback, press the down button to fast forward, press the up button to enter slow playback)

*Video settings menu

1. In the Video Standby mode, press and hold the "M" key on the main unit until the menu is displayed on the screen. (Or press the "MENU" button on the remote control)
2. Press the "up/down" key to select the parameter to be set, and press the "OK" key to enter the content to be set.
3. After completing the settings, press the "M" key to return to the Video Standby mode.

*Photo setting menu

1. In the Video Standby mode, press and hold the "M" key on the main unit until the

menu is displayed on the screen, and then press the "M" key again.

(Or use the remote control to operate, press the "MENU" button twice).

2. Press the "up/down" key to select the parameter to be set, and press the "OK" key to enter the resolution setting.

3. After completing the settings, press "OK" to return to the video setting menu, then press the "M" key to return to the Video Standby mode.

* Take pictures, 2 ways

1. In the Video Standby mode, press the "Capture" button to take a photo.

2. Using the remote control, press the "Capture" button.

*Video recording steps

1. In the Video Standby mode, press "OK" to start recording.

2. Press the "OK" button to stop recording.

◆ HDMI output

• First of all, make sure that the monitor connected to the microscope is a 4k HD monitor, otherwise the image effect will not reach 4k.

• Then, power up the microscope and connect it to the high-definition display. The microscope does not contain a built-in battery, so please make sure that the microscope is always energized during use.

• Third, press the "On/Off" button to turn on the microscope, place the observed object on the stage and adjust the focus and object distance. Note that the closer the lens is to the object, the greater the magnification.

• After inserting the SD card, you can take a video.

• If the display screen only supports 1080P, you can press and hold "OK" to automatically switch to 1080P mode to make the microscope output signal consistent with the display screen.

FAQ

• Why there is no image on the display?

Please check whether the cables are correctly connected and whether the power switch is turned on. If the connected monitor only supports 1080P, you need to press and hold the "OK" button to automatically switch to 1080P mode to make the microscope output signal consistent with the display on the screen.

• Why the image is not clear?

Please adjust the focus or focus pulley, the angle of light irradiation, and check whether the lens is dirty.

• Why the remote control does not work?

The remote control is not equipped with batteries at the factory, please check

whether a valid battery is installed.

- Why there is no image on the display screen when the playback mode is selected?
It is possible that the video is not saved completely, the shooting is not successful, or the video file of the video is too large, and the system needs running time.(It takes about 1 minute to display the image after switching from standby mode to playback mode)

- Why the photo is not clear?
Please check whether the photo setting is the highest pixel to ensure high-quality photo output.

- Prompt memory card error?
First make sure that the memory card is a genuine memory card. Then choose to format the memory card in the system and try to use it again.
Due to the higher performance requirements of the memory card, when recording video, if it still cannot be used normally after formatting, please replace the high-quality TF card with class 10 or above.

Warranty Card	
User's name:	Address:
Phone number:	Post code: email:
Model name:	Purchase date:
Fault feedback:	
Date:	Fault:
Notice:	
Our warranty doesn't cover:	
1	If users can not provide the purchase proof or the Warranty.
2	If users use it in an improper environment, such as incompatible power supply, high temperature environment etc.
3	If the fault caused by accident, oversight, misoperation or natural disasters etc.
4	If the fault caused by people who are not belong to authorized organizations of our company during repacking, repair, dismantle, Or if users repack, repair or dismantle the device not following our advice.
5	Out of warranty period.

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