

# 3-D ORBITAL SHAKER ATTACHMENT ACCESSORY FOR THE ROTO-SHAKE GENIE® Catalog No. SI-1150



1. Remove magnetic platform from the Roto-Shake Genie (see Roto-Shake Genie manual).
2. Place 3-D Orbital Shaker Attachment on the Roto-Shake Genie base, aligning the hex shaft with the hex on the pulley of the 3-D Orbital Shaker Attachment. Make sure the two feet are in contact with the Roto-Shake Genie base.
3. Use the Roto-Shake Genie speed control knob for controlling the speed of your 3-D Orbital Shaker Attachment. Speed range is 10-70RPM.
4. Place Non-Skid Mat on the 3-D Orbital Shaker platform. Do Not use attachment without Non-Skid Mat.
5. Place all loads centered on 3-D Orbital Shaker platform. Please note, platform capacity is 1 Kg (2.2 lb).



70 Orville Drive, Bohemia, NY 11716 U.S.A.  
Phone: (631) 567-4700 • Fax: (631) 567-5896 Toll Free: 888-850-6208  
Customer Service: [custsvc@scientificindustries.com](mailto:custsvc@scientificindustries.com)  
[www.scientificindustries.com](http://www.scientificindustries.com)