

EnSURE™ Touch



Quick Test



Locations



Plans



Users



Results &




Search

EnSURE™ Touch Features


- **Quick Test** – Results in as little as 2 taps
- **Location Test** – Easily find any location with details
- **Users** – Add and manage users all with administrative control on the device
- **Plans** – Quickly create, manage, and view plans directly on the device
- **Retest** – Swipe through failed test locations and instantly run re-tests to show corrective action
- **Reports** – Testing results and trend analysis are just a tap away
- **Sync** – Sync program information and results to the cloud, results send real time as tests are taken
- **Remote Support** – Built-in screen sharing capabilities for support right on your device
- **RLU Conversions** – RLU converter to make it easy to convert to Hygiena without changing threshold limits




Common Functions

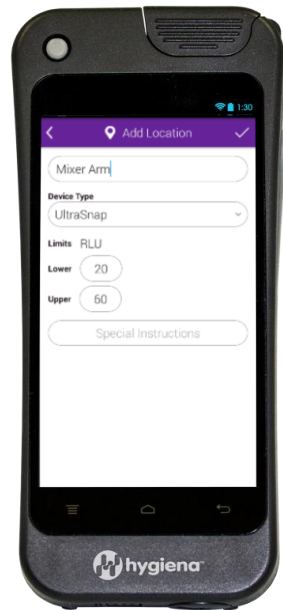


Manage Test Point Locations

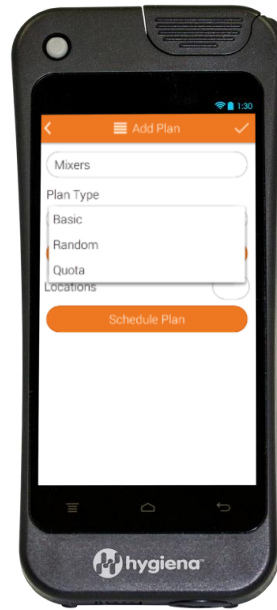


Manage Test Plans

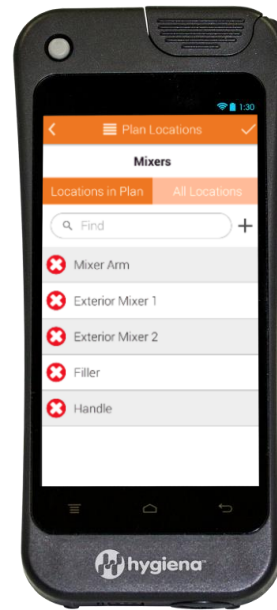


Configure Test Plan Location

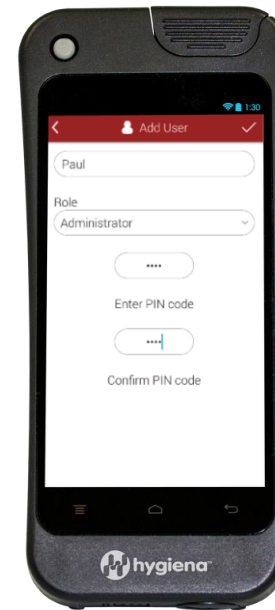

Manage Users

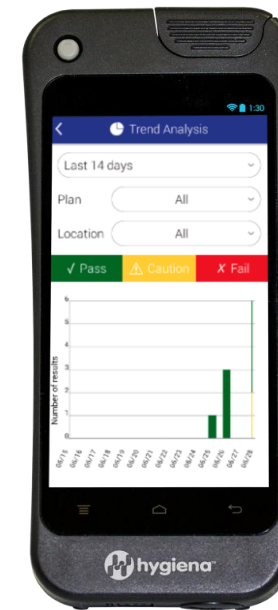


View Reports

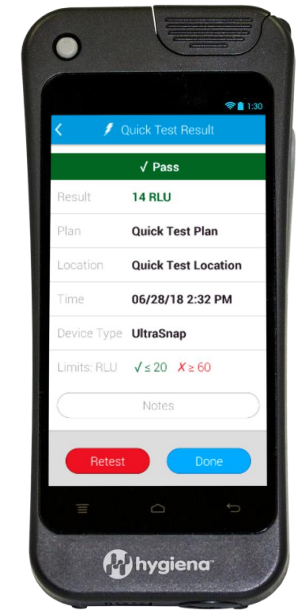


Perform Tests


Key Usability Advantages

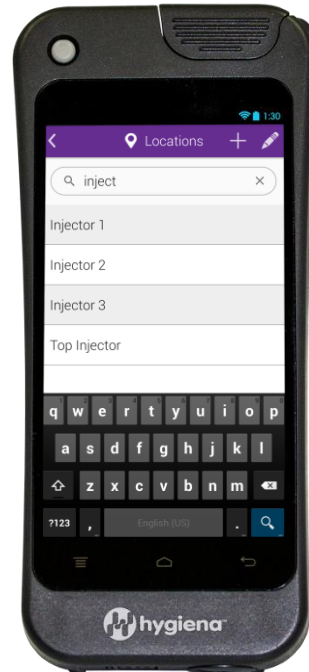

Finding Data like Locations, Plans, Results

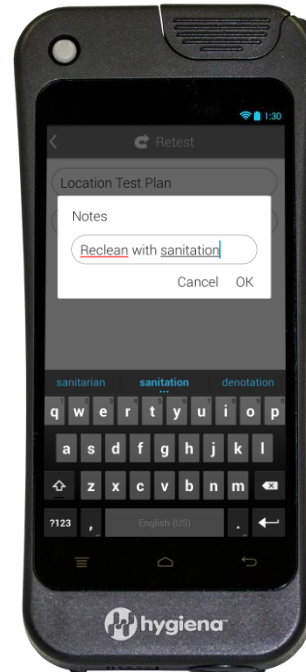

Full Query keyboard for data collection

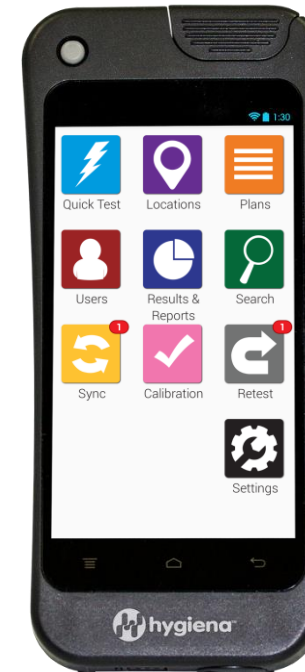


Notifications and Badges




Unicode support for any Language



Remote Support

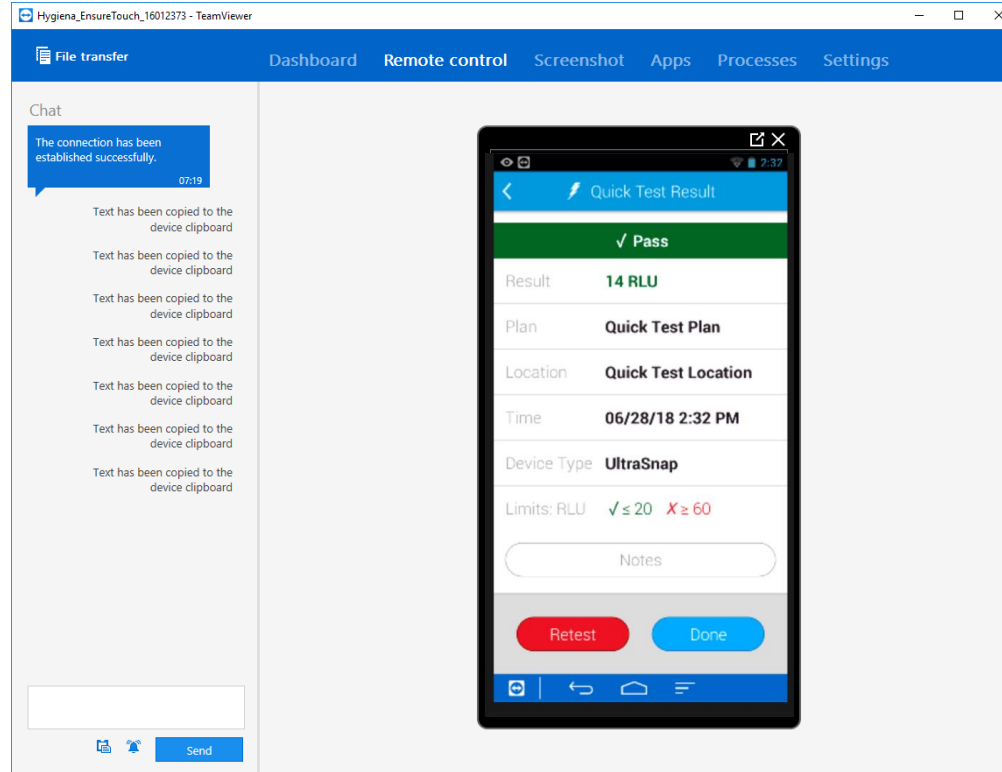


EnSURE™ Touch

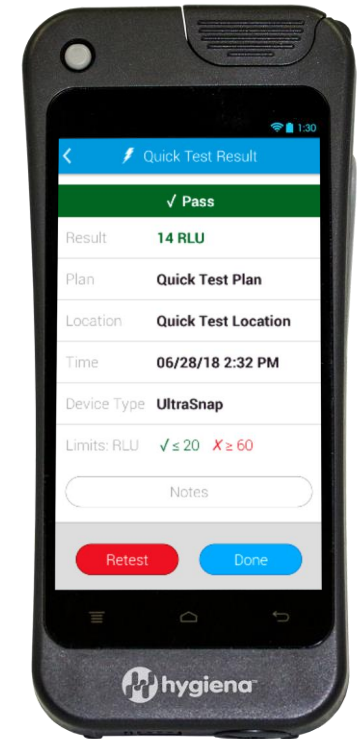



SureTrend Features

- Microsoft Azure Cloud
- Multitenant WebApp with IoT features for instrument
- Easily configure and manage instruments
- Consolidate results from multiple instruments
- Multiple language support

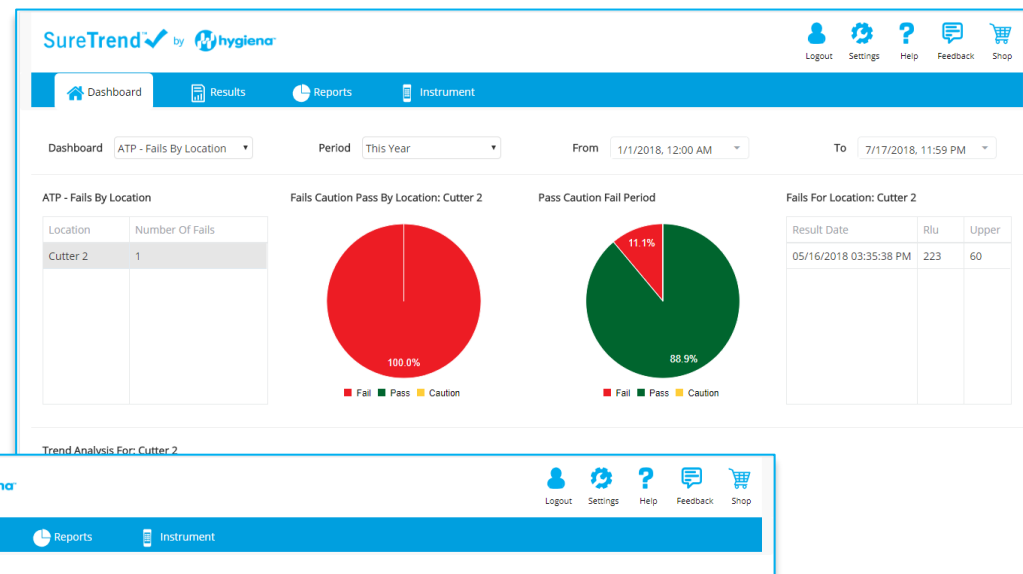

The Results table displays the following data:

Result State	Site Name	Is Retest	Result Date	Location Name	Unit No	Unit Name	Plan Name	Username
Pass	Default Site		05/30/2018 2:06:20 PM	Quick Test Location	16012373	Johns Unit	Quick Test Plan	John
Pass	Default Site		05/29/2018 10:09:10 /	Quick Test Location	16012373	Johns Unit	Quick Test Plan	John
Pass	Default Site		05/29/2018 10:07:53 /	Quick Test Location	16012373	Johns Unit	Quick Test Plan	John
Pass	Default Site		05/29/2018 10:07:30 /	Quick Test Location	16012373	Johns Unit	Quick Test Plan	John
Pass	Default Site		05/22/2018 8:14:15 AM	Quick Test Location	16012373	Johns Unit	Quick Test Plan	John
Pass	Default Site		05/16/2018 11:16:27 /	Exit door knob	16012373	Johns Unit	Location Test Plan	
Fail	Default Site		05/16/2018 8:35:38 AM	Cutter 2	16012373	Johns Unit	Location Test Plan	
Pass	Default Site		05/16/2018 8:35:02 AM	Cutter 1	16012373	Johns Unit	Location Test Plan	
Pass	Default Site		05/15/2018 4:40:50 PM	Quick Test Location	16012373	Johns Unit	Quick Test Plan	




SureTrend Cross Platform

- Cloud, private cloud, local server, or desktop
- Microsoft ASP.Net MVC Core (Any Platform)
- Cloud version has some special features and value
- Data in the cloud is very valuable


“Through digital technology, we are increasing our ability to offer real-time, actionable insights and smarter solutions . . . and deliver a superior customer experience and better outcomes.”

–Doug Baker

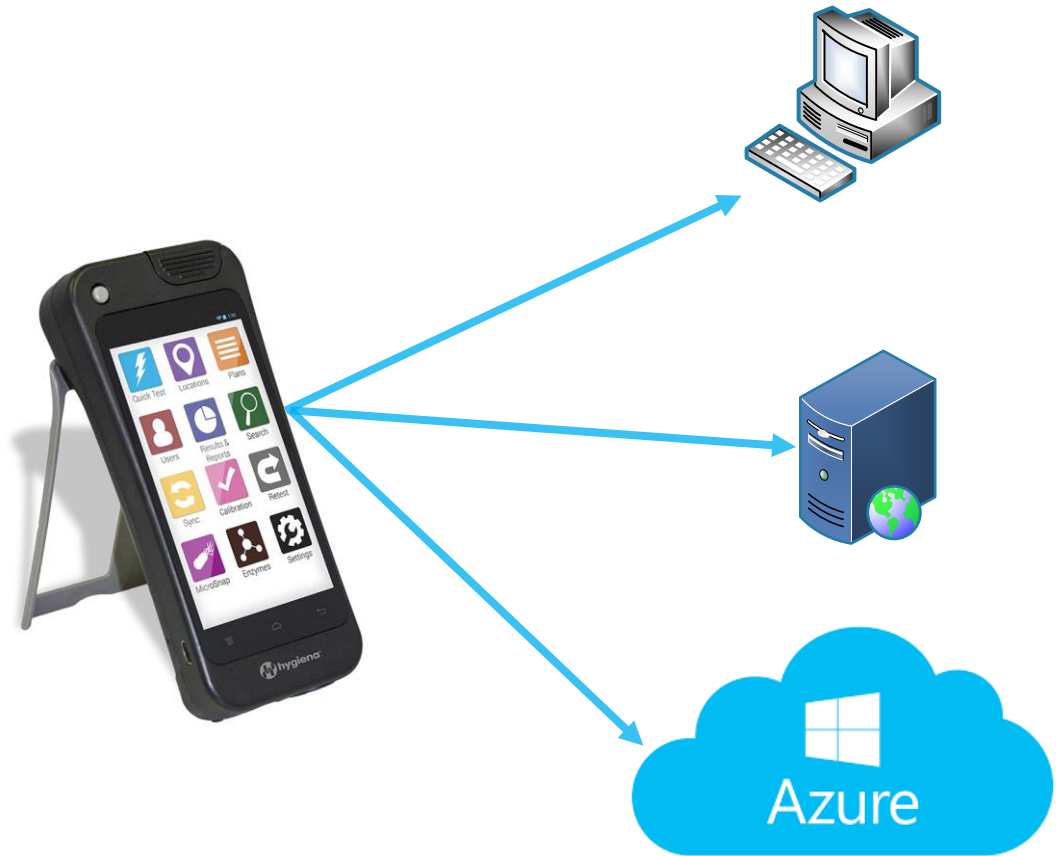

Bioluminescent Chemistry

- Liquid stable chemistry up to 15 Month 4°C
 - Efficient and inexpensive manufacturing for reproducible results
 - Removes variation from rehydrating lyophilized reagents at point of use
- Premoistened swab with lysing agent

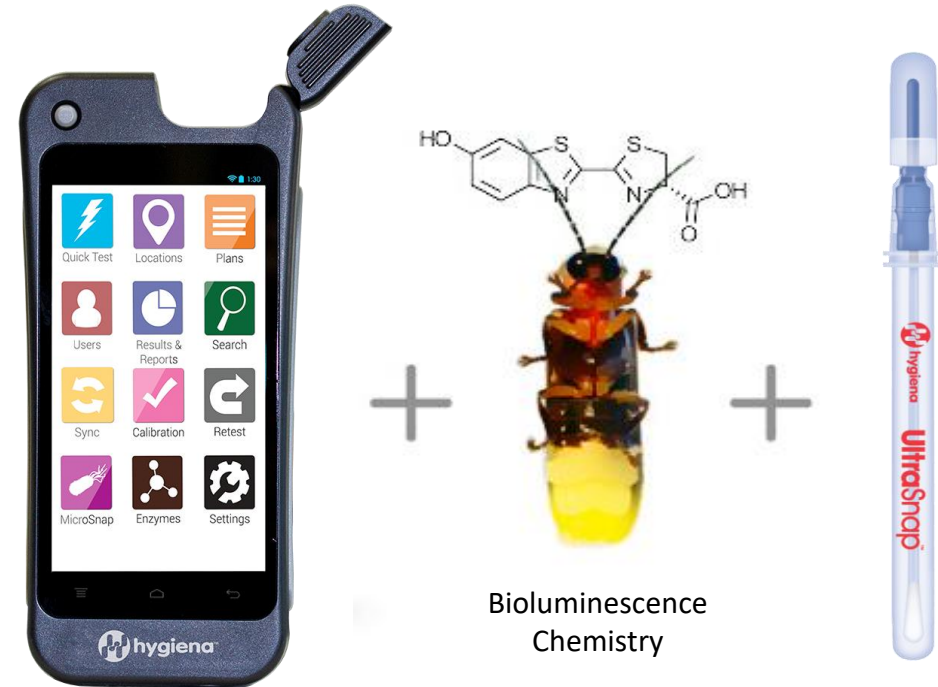

Test Device Format

- Easy-to-use Snap-Valve technology delivers the reagent directly to the collected sample
 - No need to break multiple foil barriers with swab potentially losing some of your sample
- Recyclable & non-toxic plastics



RLU Conversion

- **RLU Conversions** – RLU converter to make it easy to convert to Hygiena without changing threshold limits
- EnSURE TOUCH with UltraSnap ATP swabs can accurately emulate a competitor's system across a large dynamic range.
- RLU scale will be displayed in the competitor format
 - Current threshold limits in SOPs do not need to be changed

