

Upgrading to a New Version of QRlogix Software

For V1.6 and newer. *For earlier versions, please contact support@QRlogix.com


1. Export any current templates to the Templates folder or another safe location.
 - a. C:\Users\Public\Documents\QRlogix\Templates
2. Take screenshots of your template form in case they need to be rebuilt.

The screenshot shows the 'My template' form in the QRlogix Software. The form is titled 'My template' and has a sidebar on the left with navigation options: Dashboard, Template 1, Cassette DS1, My template (selected), EB DS1, and File Drop 1. The main form area contains fields for: Prefix (SL23), Accession / Quantity (400), Specimen / Quantity (A), Block / Quantity (1), Name (PAN, PETER), and Stain (H&E). There is a 'PRINT' button at the bottom right of the form. On the right side of the window, there is a preview of the label output, which reads: 'SL23-400-A1', 'Lev 1 H&E', 'PAN, PETER', and '03/14/2024'. Below the preview is a 'Print History' section with a 'Show All' button and a list of print records, each showing 'Printed by: Administrator' and a timestamp.


3. Take a screenshot of Label Builder

The screenshot shows the 'Label Builder' window. The window has a title bar 'Label Builder' and a sidebar on the left with 'Properties' and 'Field' sections. The 'Field' section has 'Position X' and 'Position Y' both set to 120. The main area is a grid where the label text is being built. The text on the grid is: 'B22-1001-A1', 'Lev 1 H&E', 'PAN, PETER', and '03/14/2024'. At the bottom of the grid, it says 'Showing sample data for preview' and 'Get support and watch tutorials'. There are 'Cancel' and 'Save' buttons at the bottom. On the right side, there is an 'Elements Added' section with a list of elements: 'Aa Concatenate B22-1001-A1', 'Aa Stain H&E', 'Aa Concatenate Lev 1', 'Current date 03/14/2024', and 'Aa Name PAN, PETER'.

4. Backup contents of your Sqlite folder located at: C:\ProgramData\QRlogixLMS\Sqlite (or it could be located at C:\ProgramData\QRlogixSoftware\Sqlite)
 - a. Create new folder = Old
 - b. Place all contents of Sqlite folder into “Old” folder

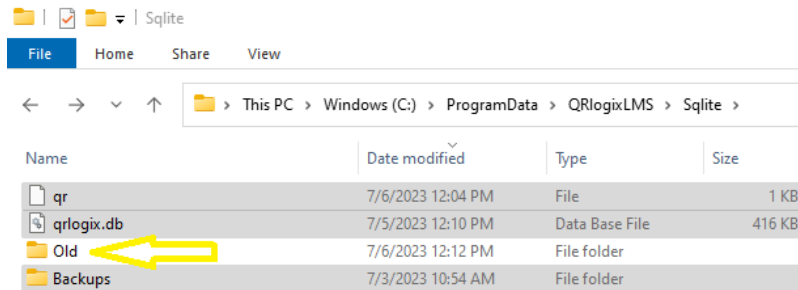


5. Uninstall software in Control Panel

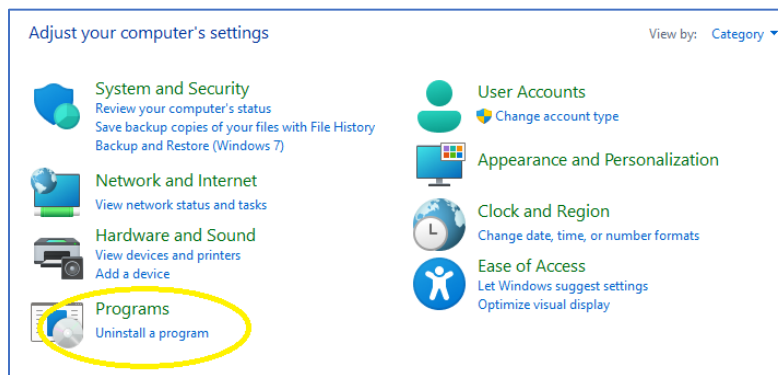

6. If PrintShare (File Drop) is installed, uninstall the AutomationService (File Drop Service) application in Control Panel.

Confirm the following folder was removed during uninstallation:
C:\AutomationService or C:\FileDropService

7. Download the new version of the software from the link provided.
8. Download the new version of File Drop if you are printing from CSV/TXT or connecting to LIS.
9. Follow the software installation instructions found in the User Manual.
10. If connecting to File Drop (formerly PrintShare), follow the installation instructions found in the User Manual.
11. Use the Import Template button to load previous templates.

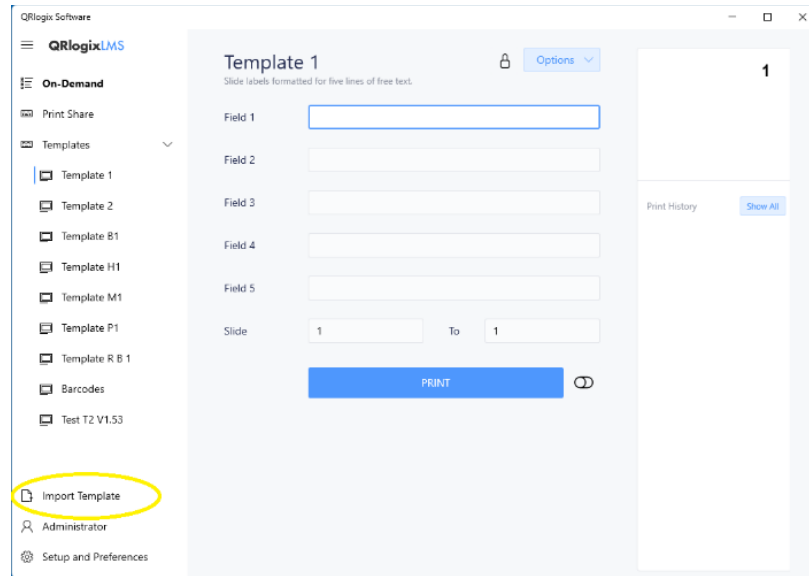

12. If you have any problems loading old templates, email the exported template files, along with the form screenshot and label designer code to support@QRlogix.com and we'll rebuild the templates in the current version and send you the files.

Support:

Phone: 800-650-0632

Email: Support@QRlogix.com

<https://qrlogix.com/pages/qrlogix-software-support>