***Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung***

MST DIGITAL Magnetic Stirrer

F203A0450

General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise

Before using the unit, please read the following instruction manual carefully.

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.

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This unit must be used for laboratory applications indoor only. The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a not specified way by the manufacturer or with non specified accessories, product's safety may be compromised.

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This unit has been designed and manufactured in compliance with the following standards:

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**IEC/EN 61010-1
IEC/EN 61010-2-051**

Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it.

It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres, in bain-marie and to stir combustible liquids that have a low combustion temperature. The minimum fire point of flammable solution is 200 °C. Only small amounts (< 50 ml) of flammable liquid can be used with the device.

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In case of black-out, once the power is back the device will restart automatically.

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The MST magnetic stirrer is a small, simple and effective stirrer.

Thanks to the new brushless motor, the stirrer stays cold even after several days of continuous operation.

Thanks to the white cover surface it is specially designed for microtitrations, and it is extremely useful where reliable and small instrument is needed.

The stirring speed can be set using the frontal knob within a range of 80 - 1500 rpm

The stirrer display shows the speed set point.

The instrument is equipped with advanced functions as auto-reverse of the stirring direction, intermittent mode programs and integrated timer to meet users' workflow and safety lock feature to avoid possible alteration of the working conditions.

NOTE: *it's important to choose the most suitable magnetic stirring bar related to the quantity and the liquid to be stirred as well as to the type of the beaker. The magnetic stirring bar that satisfies most of the application is A00001056 (6x35mm).*

2. Installation

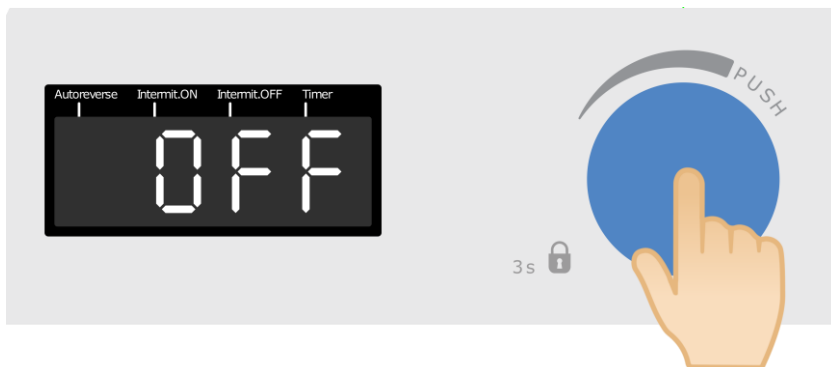
- Unpacking
 - Check the integrity of the unit after unpacking

- The box includes
 - MST Digital magnetic stirrer
 - Instruction manual
 - Plug for power supply
 - Switching power supply 100-240 V / 12 V


- Installation
 - Place the unit on a non-flammable surface
 - Make sure that the technical specification of the unit and the power supply line are the same
 - Place the flasks containing the sample and a suitable magnetic stirring bar on the stirring plate
 - Connect the instrument to an easily accessible socket (compliant with the current safety norms), using only the provided power supply and dedicated plug

3. Operation

- Push the knob to start stirring.

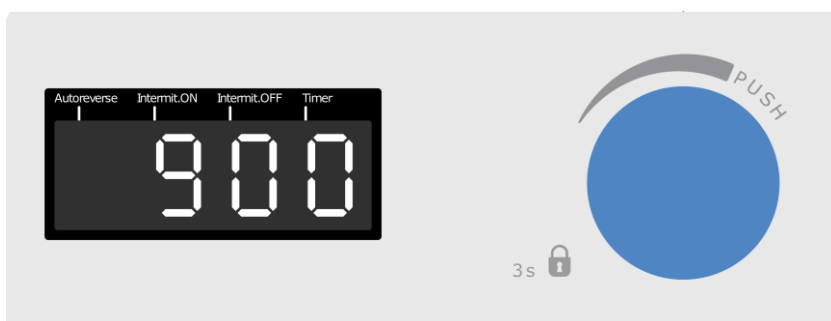
- Rotate the knob to set the stirring speed.


Note: push the knob to stop stirring.

Autoreverse, intermittent mode and timer functions

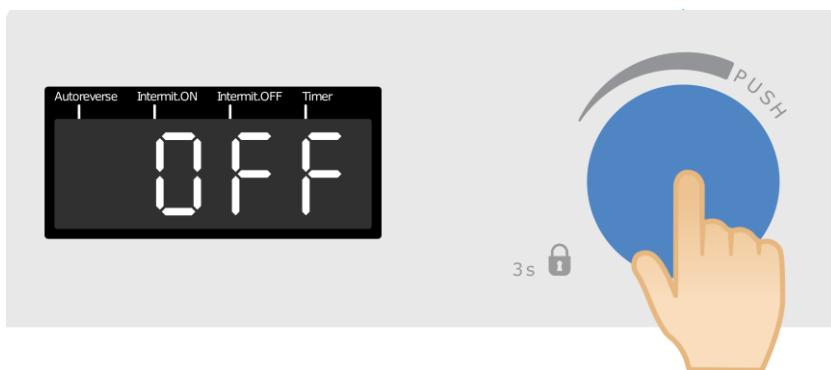- While the display shows OFF, push the knob for 3 seconds.

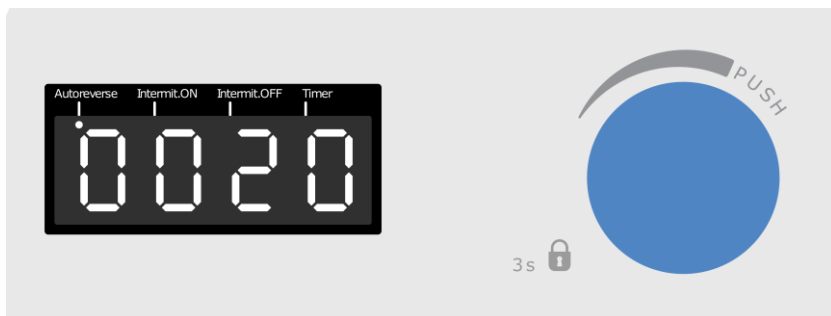
- Rotate the knob to set the autoreverse time (minutes/seconds).

The rotation direction will change after each set interval time.


Note: select *OFF* to disable the function.

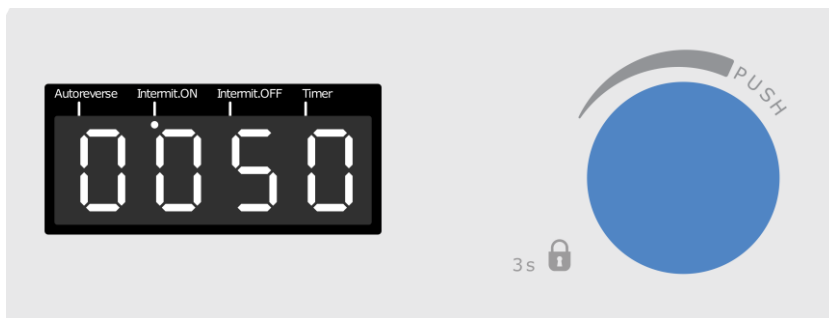- Push the knob. Then rotate it to set the run time of the intermittent mode (minutes/seconds).

Note: select *7755* to disable the function.

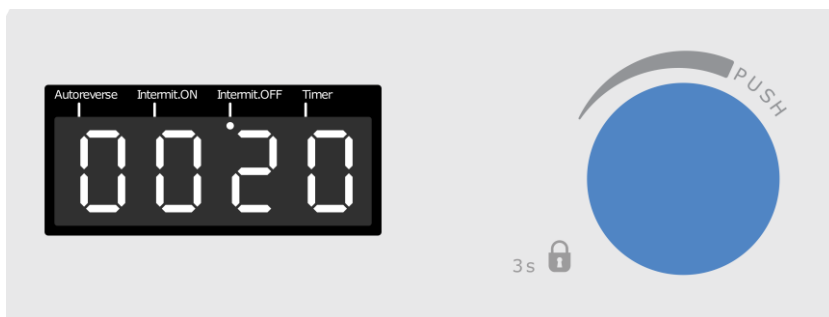
- Push the knob. Then rotate it to set the pause time of the intermittent mode (minutes/seconds).

The stirring function will be interrupted cyclically. The run time and the pause time should both be set.

Note: select *7755* to disable the function.

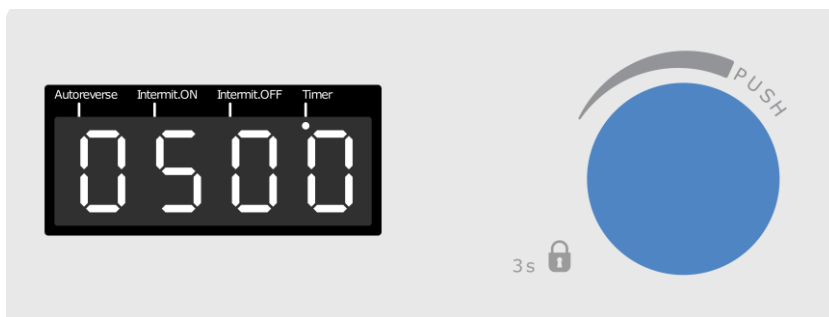- Push the knob. Then rotate it to set the timer (hours/seconds).

After the timer has elapsed, the stirring will be stopped and the display will show *End*.

Note: select *7777* to disable the function.

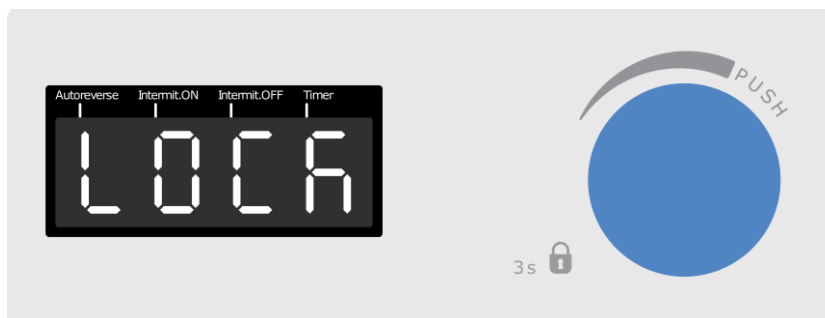
- Push the knob or wait 10 seconds to confirm.

Dedicated blinking leds show the active functions.

Note: each function can be set individually or in combination with the others.

Lock function

- While the unit is working, push the knob for 3 seconds to enable the lock function.

If the lock function is enabled, it is not possible to change the work conditions. The display shows *Lock* if the knob is pressed or rotated.

Note: push the knob for 3 seconds to disable the function.

4. Maintenance

No routine or extraordinary maintenance is necessary except periodically cleaning.

Cleaning

Disconnect the unit from the power supply and use a cloth dampened with a non-flammable non-aggressive detergent.

Repair

Repairs must be carried out by authorized Velp personnel only.

The instrument must be moved using the original shockproof packaging when freight forwarders, couriers or other are taking care of the unit transport. Follow the indications on the original packaging (e.g. place the instrument on a pallet). It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

5. Technical data

General features	Model	F203A0450
	Voltage	100 – 240 V / 50 – 60 Hz
	Power input	12 VDC 2.5 W
	Current consumption	0.2 A
	Dimensions (WxHxD)	130x50x150 mm
	Weight	0.4 kg
	Construction material	PA
	Timer	1 min ÷ 99 h 59 min
	Autoreverse / Intermittent mode	5 s ÷ 99 min 55s
	Working in continuous	Admitted
	Noisiness	< 35 dBa
	Environmental temperature admitted	+5...+40 °C
	Storage temperature admitted	-10...+60 °C
	Max humidity	80%
	Overvoltage category	II
	Pollution degree CEI EN61010-1	2
Max altitude	2000 m	
Stir	Stirring capacity	5 L H ₂ O
	Speed range	0 ÷ 1500 rpm
	Motors type	BLDC
	Speed selection	10 rpm

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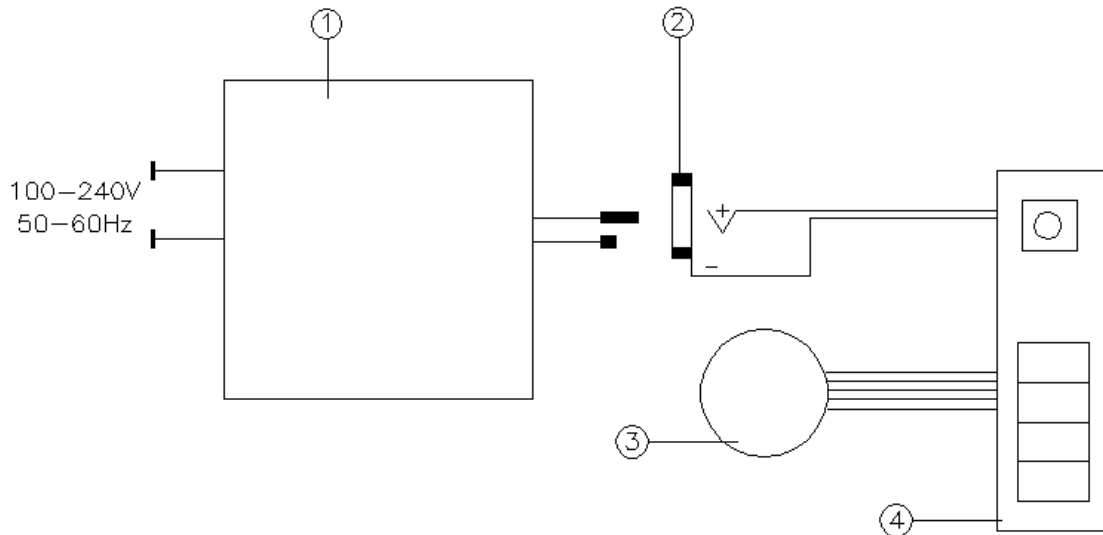
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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2011/65/EU (RoHS)

2012/19/UE (RAEE)

and satisfies the essential requirements of the following directives:

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Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Homogenizers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Radiation Detector
Open Circulating Baths
Pumps

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VELP Scientifica Srl