# QUICK START GUIDE

**MOTICAM**S SERIES

## EN Quick Start Guide

Thank you for purchasing this Moticam. This camera is a delicate instrument and should be handled with care. When not in use, please store the camera in a safe place with the dust cap attached. Do not let the camera get in contact with water or expose it to extreme temperatures (below 0 or above 50 degrees Celsius). In order to receive the highest level of service, please retain your receipt as proof of purchase and **register your camera online at [www.motic.com/As\\_Support](http://www.motic.com/As_Support) today.**

Should you have any questions please do not hesitate to **contact our customer service.**

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#### EN Using the camera on a Trinocular Microscope

1. Install the Motic Images Plus application program from the USB stick included.
2. Remove the Dust Cap (B) from the camera (A) and attach the microscope's C-Mount adapter (\*Please note that the C-Mount adapter is specific to your microscope. For Motic microscopes, please consult the relevant Motic Accessory list and choose an appropriate adapter for your Moticam. Depending on the type of adapter and microscope, the C-Ring (C) may be necessary between the camera and the C-mount adapter).
3. Connect the USB cable (D) onto the camera and then plug it into the computer
4. Install the driver if necessary by following the instructions on screen. The driver for MoticamS and MoticamProS is located in the Motic USB Driver Subfolder on the USB stick.

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