



# Liquinox™

## Critical-Cleaning Liquid Detergent

- Concentrated to save you money
- Replaces corrosive acids and hazardous solvents
- Phosphate free, biodegradable and readily disposable
- Free rinsing to give you reliable results and no interfering residues
- Use to pass your cleaning validation tests for lab accreditation and plant inspection approval

## Used to clean

Healthcare instruments, laboratory ware, vacuum equipment, tissue culture ware, personal protective equipment, sampling apparatus, catheters, tubing, disk drives, clean rooms, medical devices, optical parts, electronic components, pharmaceutical apparatus, cosmetics manufacturing equipment, metal castings, forgings and stampings, industrial parts, pipes, tanks and reactors. Authorized by USDA for use in federally inspected meat and poultry plants. Passes inhibitory residue test for water analysis. Used for phosphate sensitive analysis ware. FDAcertified.

## Used to remove

Soil, grit, grime, slime, grease, oils, blood, tissue, particulates, deposits, chemical and solvents.

## Surfaces cleaned

Corrosion inhibited formulation recommended for glass, metal, stainless steel, porcelain, ceramic, plastic, cement and fiberglass. Can be used on soft metals such as copper, aluminum, zinc and magnesium if rinsed promptly. Used for art restoration. Corrosion testing may be advisable.



## Cleaning method

Soak, brush, sponge, cloth, ultrasonic, flow through clean-in-place. Will foam—not for spray or machine use.

## Directions

Make a fresh 1% solution (2 1/2 Tbsp. per gal., 1 1/4 oz. per gal. or 10 ml per liter) in cold, warm or hot water. If available, use warm water. Use cold water for blood stains. For difficult soils, raise water temperature and use more detergent. Clean by soak, circulate, wipe or ultrasonic method. Not for spray machines, will foam. RINSE THOROUGHLY preferably with running water. For critical cleaning, do final or all rinsing in distilled, deionized or purified water. For food contact surfaces, rinse with potable water. Used on a wide range of glass, ceramic, plastic and metal surfaces. Corrosion testing may be advisable.



Convenient sizes	Alconox cat#
Case 12 x 1 Quarts	AL-1232
Case 4 x 1 Gallons	AL-1201
15 Gal. Drum	AL-1215
30 Gal. Drum	AL-1230
55 Gal. Drum	AL-1255

1 gal. makes 100 gal. cleaning solution

### Chemical Description

Liquinox consists primarily of a homogeneous blend of sodium linear alkylaryl sulfonate, sodium xylene sulfonate, alkanolamide, and ethoxylated alcohol. Liquinox is anionic in nature.

### Cleaning Validation Methods

Test a parameter of rinse water before and after rinsing the cleaned surface, or test the clean surface. No significant change in the parameter indicates no detectable residue. Parameters measured include: pH, conductivity, UV, TOC, HPLC, sodium concentration, phosphorus concentration, anionic surfactant concentration using inexpensive detergent water testing kits, surface tension, and surface analysis. For details see the Pharmaceutical Cleaning Validation References at [www.alconox.com](http://www.alconox.com)

### Health Safety Information

**OSHA Hazardous Ingredients:** None  
**RCRA Hazard Class:** Non-hazardous  
**RoHS:** No RoHS hazardous ingredients  
**EPA Priority Pollutants:** None  
**DOT Hazard Class:** Non-hazardous  
**Flammability:** Non-flammable  
**Latex Content:** None in detergent, packaging materials or adhesives.  
**Oral Toxicity:** LD50 >5000 mg per kg Rat, Oral)  
**Eye Irritation:** Mild to Moderate eye irritant if not rinsed  
**Inhalation Toxicity:** Non-irritating  
**VOC Content:** 0%  
**Carcinogenicity:** NTP = No IARC = No OSHA = No  
 All ingredients in Liquinox are listed in TSCA inventory

### Physical data

Physical data	Typical value
pH as is	8.5
Specific gravity (g/ml)	1.07
Density (lbs./gal.)	8.9
Vapor pressure (mm Hg)	10.5
Flash Point (degrees F)	None
Phosphate Content (as Phosphorus)	0%
Organic Carbon (% calculated w/w)	21%
Fragrance Content	0%
Surface Tension 1% Sol'n (Dyne/cm)	32

Color: Pale Yellow

Form: Liquid

**Solubility in Water:** Completely soluble in all proportions

**Hard Water Effectiveness:** Highly Effective

**Biodegradability:** Biodegradable

**Foam Tendency:** High Foaming

**Shelf Life:** Two years from the date of manufacture

### Precautions

No special precautions other than good industrial hygiene and safety practices employed with any industrial chemical (see Directions). A Material Safety Data Sheet is available at [www.alconox.com](http://www.alconox.com) or by calling fax-on-demand at 914-948-4040 and following the prompts. The Liquinox catalog number starts with 12. Contact Alconox, Inc. for purchase specifications. Information presented is typical. Not to be taken as specifications. Typical data is not a specification While the information in this report should not be considered to be a product warranty, we urge you to investigate, test and verify the suitability of Alconox detergents for your specific application. We, of course, can not give permission to use, or recommend the use of, our detergents where they infringe patents. No representation or warranty is made as to the safety of products or materials mentioned under the Federal Food Additives Amendment of 1958.

#### AL-1201-1

1 gallon plastic bottle

