

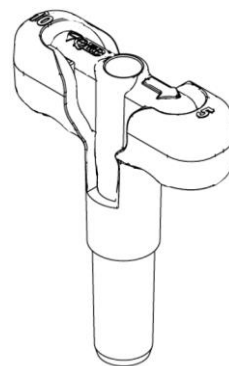
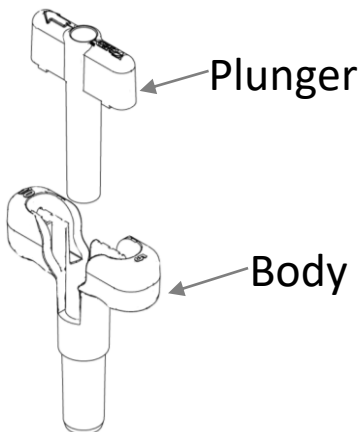
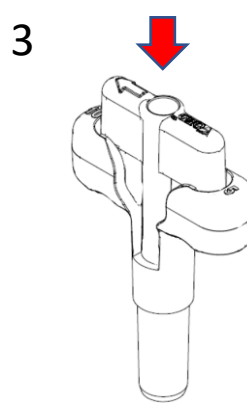
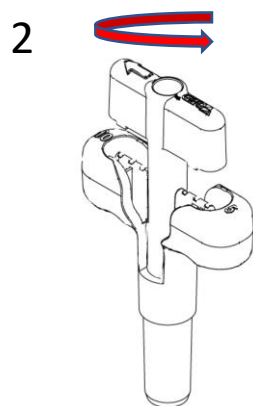





# Combo Core™ Sampler Instructions

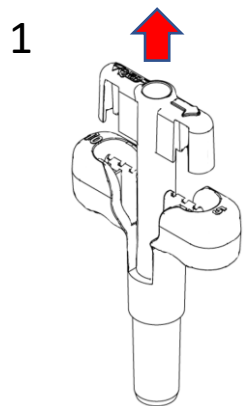


**Collecting 5 gram samples:**  
The sampler is shipped in the 5 gram sample configuration. The top of the plunger is flush with the top of the body handle, and the arrow on the top of the plunger points to the "5" marking on the top of the handle.

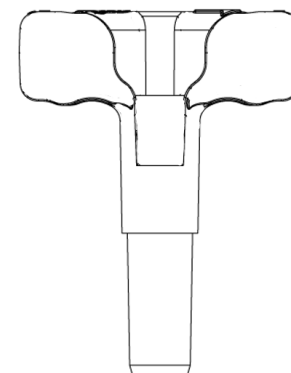

**Collecting 10 gram samples:**  
To switch to the 10 gram configuration, lift the plunger (1) and rotate it 180° (2), then lower the plunger to engage the positioning tabs (3). The plunger sits above the top of the handle in this position, and the arrow on the top of the plunger points to the "10" marking on the top of the handle.

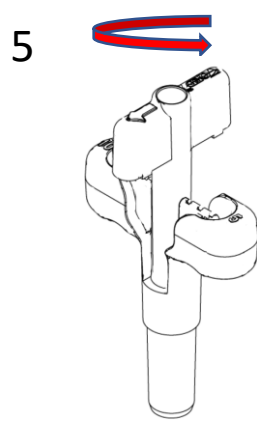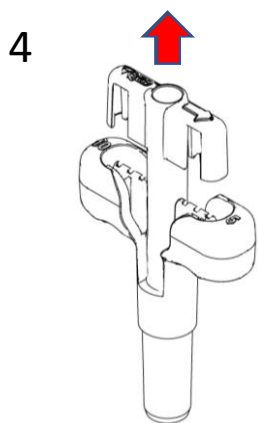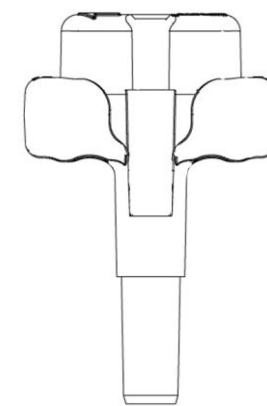

**Sample ejection:**  
Insert the barrel into your sample vial. Lift the plunger (4) and rotate 90° in either direction to line up the plunger with the slot in the body (5), then push the plunger down until it bottoms out in the slot (6), ejecting the sample.

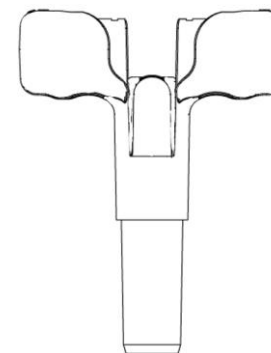

Scan here to view the Combo Core user video

