

# Product Data Sheet





Product Name *3-D Life* RGD Peptide

Catalog Number 09-P-001

**Description** RGD sequences are motifs on extracellular matrix proteins promoting adhesion to cellular integrin receptors. The *3-D Life* RGD Peptide contains a RGD motif and a thiol group on the N-terminus. The thiol group is used to immobilize the peptide on thiol-reactive polymers in *3-D Life* Hydrogels. The immobilized peptide promotes adhesion of cells carrying the appropriate integrin receptor. For instructions, please consult General Protocols GP-1 "Preparation of *3-D Life* Fast Gelling Hydrogels" or GP-2 "Preparation of *3-D Life* Slow Gelling Hydrogels" and the *3-D Life* Hydrogels User Guide on [www.cellendes.com](http://www.cellendes.com).

**Quantity** 1 µmol

| Material   | Quantity | Concentration of reactive groups | Storage   |
|--|----------|----------------------------------|---|
|  RGD Peptide, lyophilized | 1 µmol   | 20 mmol/L*                       | Lyophilisate and after reconstitution: -20°C or lower |
|  Water                  | 600 µl   | n/a                              | Room temperature or lower                             |

All materials are filter-sterilized.

\*Volume/concentration after reconstitution of lyophilisate.

**Reconstitution** RGD Peptide:

1. Briefly centrifuge vial containing the RGD Peptide lyophilisate to make sure that the entire product is at the bottom of the reaction tube.
2. Add 48 µl *3-D Life* Water per tube of RGD Peptide to dissolve the lyophilisate at a concentration of 20 mmol/L thiol groups.
3. Close tube and briefly vortex.
4. Incubate for 5 min.
5. Briefly vortex and centrifuge again.
6. The RGD Peptide is now ready for use.

**Pept. Sequence** Acetyl-Cys-Doa\*-Doa-Gly-Arg-Gly-Asp-Ser-Pro-NH<sub>2</sub>

\*: Doa:8-amino-3,6-dioxaoctanoic acid

**Molar Mass** 1022 g/mol

**Purity** >90% (HPLC)

*Intended for research use only. Not for use in human therapeutic or diagnostic applications.*



Distributed by:  
Ilex Life Sciences LLC  
Tel: (828) 531-9949  
Email: [info@ilexlife.com](mailto:info@ilexlife.com)  
Web: <https://ilexlife.com/>