Cellpuri™

Disposable cell enrichment chip

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⚠ Warnings & Precautions

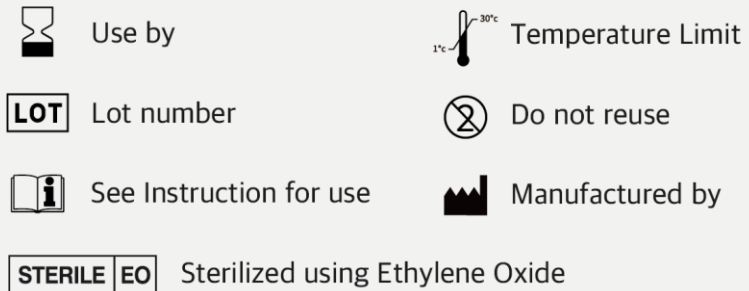
» Storing Cellpuri

- Do not store at extreme temperature and humidity.
- Keep the product at a room temperature between 1 ~ 30 c and relative humidity.
- Keep the product in the original box to protect from light and physical damage.
- Cellpuri can be easily damaged by external shock. Be careful not to drop the product.

» Using Cellpuri

- Cellpuri is a single-use, sterilized product. Do not reuse.
- Do not use if the package is open or damaged.
- Read the instructions carefully before use.
- Use immediately after opening the package.
- Make sure that the product is facing upward and placed on a flat rigid surface.
- Do not add any solution other than cell suspension.
- Be careful of product's sharp edges.

Safety symbols

» Using Syringe pump

- Operate the device following the operating condition.
- Ensure the voltage matches with the device's power supply voltage.
- Check if the power cable is properly grounded to avoid potential electric shock.
- Use only the components provided and authorized by Curiosis Inc.
- Do not disassemble the device and contact an authorized service center if required.

Specifications

Cellpuri	
Dimension	76 x 25 x 23 mm
Weight	15.3 g
Sample size	7~15 μ m size cells
Sample concentration	$\sim 5 \times 10^5$ cells/ml
Enrichment	20-fold \sim
Yield	90~100%
Loading volume	2~10 ml

Syringe pump	
Dimension	17 x 26 x 28 cm
Weight	4 kg
Volume range	2~10 ml
Flow rate	1 ml/min
Power	DC 24V, 2A
Operating condition	Temperature 20~35 $^{\circ}$ C Relative humidity < 95 %

Ordering information

Catalog no.	Product Name	Description
CRCPR-CE10	Cellpuri	Cell Enrichment chip (10 pc/box)
CRCPR-SP01	Syringe pump	Motorized vertical syringe pump

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