

BUG TONGS

SCIENCE UNDERSTANDING NATURE

PRODUCT GUIDE

IP 040259

THIS PRODUCT GUIDE MAY BE PHOTOCOPIED FOR CLASS USE ONLY

CONTENTS

1 x set of metal Bug Tongs 1 x clear pot 1 x clear lid The Invicta Bug Tongs are used with a scissor action to capture insects and small organisms.

If a creature is found on a leaf (or similar thin object) then the lid of the open Bug Tongs should be gently placed over it, taking care

ASSEMBLY

The Bug Tongs are supplied assembled with the clear pot inserted into the large loop of the tongs and the clear lid attached onto the small loop of the tongs.

CLEANING

The Bug Tongs can be wiped with a damp cloth or washed in soapy water after use. Ensure they are completely dry prior to storage.

not to damage fragile legs or wings, and the Bug Tongs should be closed. Once closed the creature can be carefully slid off its leaf, gently dropping it into the pot.

For larger creatures the Bug Tongs can be used the other way up with the pot placed over the creature and then closed so that the specimen is enclosed into the sealed pot and lid, which can then be turned the correct way up for viewing.

The specimens can be observed in the Bug Tongs or they can be carefully transferred into Bug Pots or Magnifying Pots for further detailed examination.

Always ensure that children release their specimens carefully where they were found.

Invicta Education Leicester, England

Tel: +44(0)116 281 7164 email: sales@invictaeducation.com web site: www.invictaeducation.com WARNING: NOT SUITABLE FOR CHILDREN UNDER 36
MONTHS BECAUSE SMALL PARTS MAY CAUSE
A CHOKING HAZARD. TO BE USED ONLY UNDER ADULT SUPERVISION.
Please retain the information on this pack for future reference. We reserve
the right to alter designs and specifications (including colours and materials) if and when

such changes are unavoidable. This product conforms to the safety requirements of EN71, ASTM, 16 CFR, CCC and The Canadian Hazardous Products (Toys) Regulations. MADE IN CHINA





