
**BT100S/BT300S/BT600S / BT100S-1**

**Variable-Speed Peristaltic Pump**

**Operating Manual**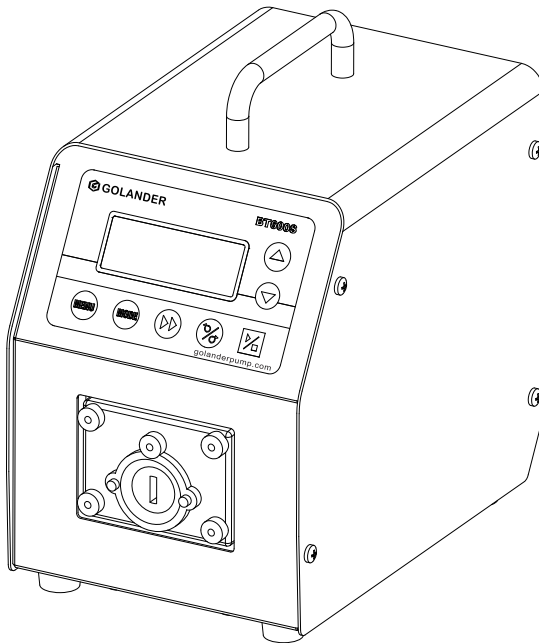

# Contents

Safety Precautions .....	1
1 Description .....	6
2 Functions and Features.....	6
3 Components and Connectors.....	7
4 Display Panel and Operating Keypads .....	8
4.1 Keypad.....	8
4.2 LCD Screen Display .....	9
5 Parameter Settings .....	11
6 System Advanced Parameter Settings .....	13
7 External Control Interface.....	16
8 Operating Instructions .....	16
8.1 Before Operation.....	16
8.2 Power Connection.....	17
8.3 Mode Change.....	17
8.4 Internal Control Mode.....	17
8.5 External Control Mode.....	18
8.6 Time Dispense Mode.....	20
8.7 Logic Level 1 Control Mode (footswitch).....	21
8.8 Logic Level 2 Control Mode (footswitch and direction switch).....	22
8.9 Communication Mode .....	24
8.10 Speed Setting.....	25
9 Maintenance .....	25
9.1 Warranty.....	25
9.2 Regular Maintenance .....	25
9.3 Malfunction Solutions .....	25
10 Dimensions .....	27
11 Naming Rule.....	28
12 Specifications .....	29

## Safety Precautions

### Danger

- Use the correct voltage indicated on the rating plate label of the pump to avoid any damage.
- Do not make any unauthorized dismantling, changes or modifications to the pump which could result in malfunctions or even potential accidents.
- Turn off the pump drive before installing or removing tubing, attempting any maintenance, cleaning or repair of the drive, connecting or disconnecting external control devices or a communication interface. Fingers or loose clothing could get caught in the drive mechanism.

### Warning

- Ensure no chemical reactions occur between the handled fluid with the material of the pump head and tubing before use.
- Tubing should be checked regularly to avoid breakage. Tubing breakage may result in fluid being sprayed from the pump. Use appropriate measures to protect the operator and equipment. The operator is solely liable for damages resulting from tubing breakage, particularly the leakage of toxic or valuable liquids.
- The pump is provided with a grounded plug which must be well grounded at all times.
- This device is not designed for nor intended for usage in patient connected applications, including but not limited to medical and dental use.
- Observe all other applicable regulations concerning working safety, operational safety, environmental protection and relevant local regulations.

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# 1 Description

BT100S/300S/600S variable-speed peristaltic pump provides not only basic functions such as start/stop, reversible direction and adjustable speed, but also the Time Dispense Mode and Anti-Drip function. With the MODBUS RS485 interface, the pump is easy to communicate with external devices, such as PC, HMI or PLC. This pump series includes:

- **BT100S/BT100S-1**, flow rate 0.0001-720 mL/min, speed 0.1-150 rpm
- **BT300S**, flow rate 0.006-1600 mL/min, speed 0.1-350 rpm
- **BT600S**, flow rate 0.006-2900 mL/min, speed 0.1-600 rpm

# 2 Functions and Features

Peristaltic pumps can handle fluid that is particularly abrasive, corrosive or viscous. There are no seals in contact with the medium pumped and no valves to clog. The inner surfaces are smooth and easy to clean. The fluid contacts only the tubing or tube material. Suction lift and priming can be up to 8m water column at sea level. It can handle the most shear sensitive of fluids like latex or firefighting foam with low shearing. It is capable of running dry and pumping fluids with high quantities of entrained air, such as black liquor soap. The high volumetric efficiency allows operation in metering or dosing applications where high accuracy is required. Tubing and tube materials are available for food and pharmaceutical use.

- LCDs displays speed and working mode.
- Membrane keypad.
- Reversible direction, start/stop control and adjustable speed.
- Time dispense mode: automatically dispense by setting the duration for each dose, lag time between doses and the number of cycles.
- 0.5% high precision rotating speed control with 0.1 rpm speed resolution when speed is between 0-100 rpm, 1 rpm speed resolution when speed is between 100-600 rpm.

- An external logic level signal can control start/stop, direction and dispense functions; an external analogue signal can adjust the rotating speed. The signal is optically isolated.
- The RS485 MODBUS interface allows easy communication with external devices.
- Internal double-layer isolation structure; circuit board with conformal coating makes the pump dust- and moisture-proof.
- Anti-electromagnetic interference feature; wide input voltage range for complex power environment.
- Stainless steel enclosure, easy to clean, resistant to the corrosion of acid, alkali, sodium and organic solvents.

### 3 Components and Connectors

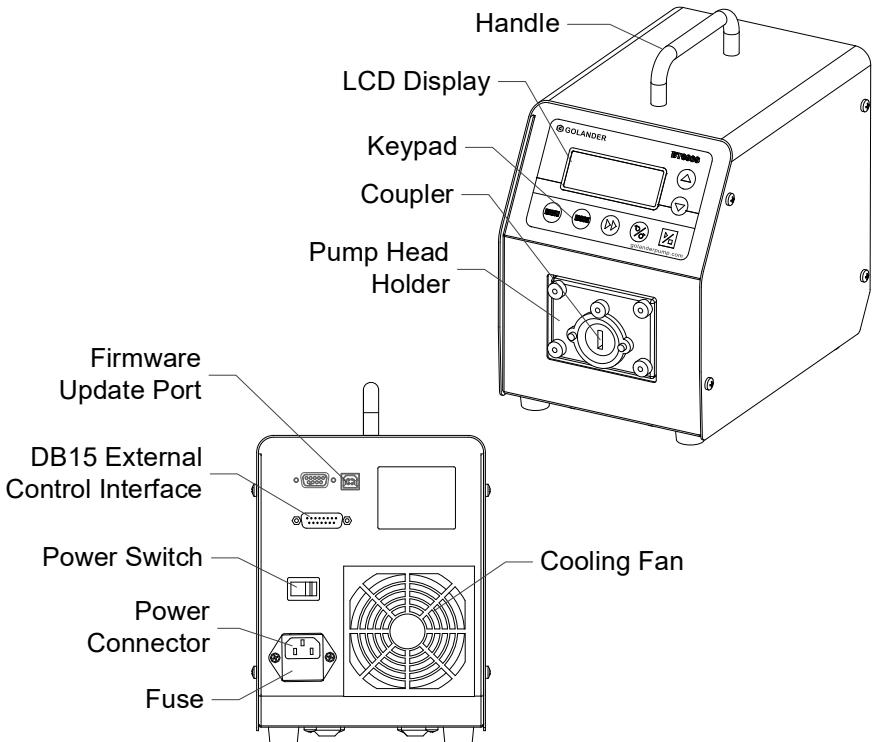
Figure 1. Components and Connectors

## 4 Display Panel and Operating Keypads

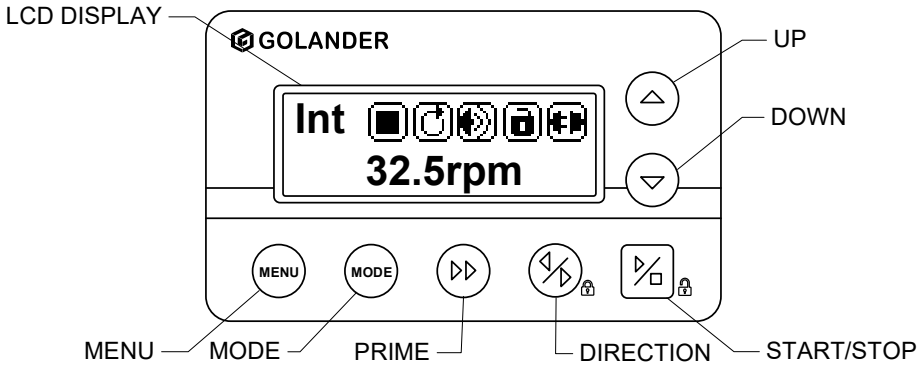
Figure 2. Display Panel

### 4.1 Keypad

**UP Key.** When pressed shortly, the last digit of the value will increase by 1. Hold the key to increase the value quickly.

**DOWN Key.** When pressed shortly, the last digit of the value will decrease by 1. Hold the key to decrease the value quickly.

**MENU Key.** When on the main screen, press the **MENU** key to enter the setting menu. When on the setting menu, press the **MENU** key to switch between the different menus. Press and hold it to return to the main screen. When the drive is running, this key is disabled.

**DIRECTION Key.** Press to key to change the drive rotating direction, clockwise or counterclockwise.

**MODE key.** When the drive is not running, use the **MODE** key to change the working mode: Internal Control mode, External Control mode, Time Dispense mode, Logic Level 1 control mode or Logic Level 2 Control mode.

**PRIME key.** Press the key to run the pump at the maximum allowed speed in the direction shown on the display. Press again to return to the previous state.

**START/STOP key.** Press to start or stop the drive.

## 4.2 LCD Screen Display

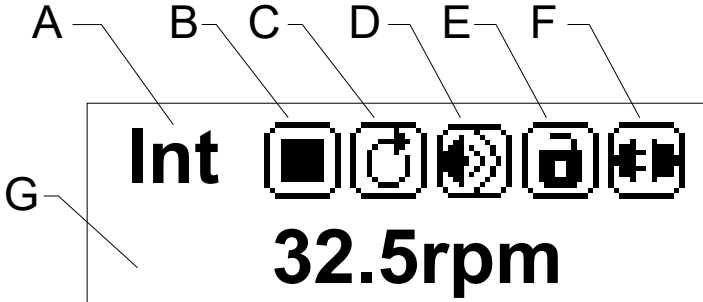
Figure 3. Display screen

### 4.2.1 A. Control Mode

It displays the current control mode. Press the **MODE** key to switch the control mode: Internal Control Mode, External Control Mode, Time Dispense Mode, Logic Level 1 Control Mode or Logic Level 2 Control Mode.

- Internal Control Mode: Use the keypad to operate the pump. Use an optional external pulse signal to control the start and stop.
- External Control Mode: Use an external analogue signal to control rotation speed. Use external logic level signal to control direction, start and stop. The keypad is disabled.
- Time Dispense Mode: Dispense fluid automatically by setting the duration for each dose, the time between doses and the number of cycles.
- Logic Level 1 Control Mode (footswitch): Use an external logic-level signal to control the start and stop. Use the keypad to control direction and speed.
- Logic Level 2 Control Mode (footswitch and direction switch): Use external logic level signal to control start, stop and direction. Use the keypad to adjust speed.

## 4.2.2 B. Running State

It shows current drive running state.

Stop

Running

Pause

Figure 4. Running State

## 4.2.3 C. Direction State

It shows current drive direction setting.

Clockwise

Counterclockwise

Figure 5. Direction

## 4.2.4 Key Tone Sate / Wi-Fi Signal Strength

When Wi-Fi is enabled in **System Advanced Parameter Settings** , it shows the state of the key tone, on or off.

Tone on

Tone off

When Wi-Fi is disabled, it shows the state of the key tone, on or off.

Note: Wi-Fi function is not available for US/European markets.

Excellent

Good

Poor

No Connection

Figure 6. Key Tone State / Wi-Fi Signal Strength

### 4.2.5 E. Keypad Lock State

When the keypad is locked, only the **START/STOP** key will work. On the main screen, press and hold the **DIRECTION** key to lock the keypad; press and hold the **START/STOP** key to unlock the keypad.

Keypad unlocked

Keypad locked

*Figure 7. Keypad Lock*

### 4.2.6 F. Communication State

It shows the current RS485 communication state.

Communication  
disconnected

#1

Connected, pump  
number is set to 1

*Figure 8. Communication state*

### 4.2.7 G. Rotating Speed State

It shows the current rotating speed, i.e revolutions per minute. When the drive is running at full speed, the display will be >>>>>>.

## 5 Parameter Settings

On the main screen, press the **MENU** key to enter the setting menu, press the **UP** or **DOWN** key to choose the parameter to adjust, press the **START/STOP** key to show the value of the parameter, and use the **UP** or **DOWN** key to adjust the value. Press the **START/STOP** key to confirm and return to the setting menu. To return to the main screen, press and hold the **MENU** key or press the **MENU** key a few times until it returns to the main screen.

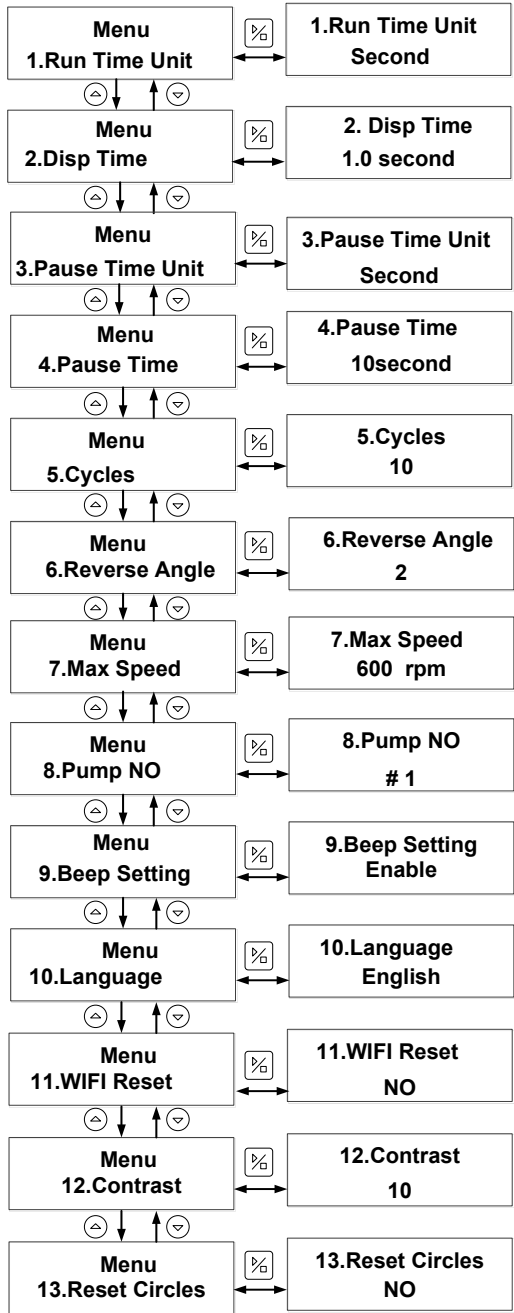
Figure 9. Parameter Setting Flow Chart

1. Run Time Unit: Time unit for the dispense duration of Time Dispense mode. The unit can be in days, hours, minutes or seconds.
2. Disp Time: Dispense time for Time Dispense mode. It is the dispense duration for each dose. The range is 0.1-999 seconds/minutes/hours/days.
3. Pause Time Unit: Time unit for the interval time of Time Dispense mode. The unit can be in days, hours, minutes or seconds.
4. Pause Time: Pause time for dispense mode. It is the lag time between doses when the number of cycles set is more than 1. The range is 0.1-999 seconds/minutes/hours/days.
5. Cycle: Number of cycles for dispense mode. The range is 0-999 cycles. When it is set to 0, the dispensing process will keep running until **START/STOP** is pressed. When it is set to the other value, the pump will stop when the set number of dispensing cycles is completed (*Figure 17*).
6. Reverse Angle: To minimize the drip after a dispense, the drive can reverse direction to draw the fluid back at the end of the tubing. The reverse angle range is 0-720 degrees. When it is set to 0, the anti-drip function is disabled.
7. Max Speed: Maximum speed for the External Control Mode. It is the maximum speed that the external analogue signal can control.
8. Pump No: It is the pump communication address for the communication mode. A restart of the drive is required for this to take effect.
9. Beep Setting: Set the key tone on or off.
10. Language: System language setting, English or Chinese.
11. Wifi resetting.
12. Contrast: LCD backlight contrast setting.
13. Reset Circles: Recount tubing life after a tubing is installed or replaced.

## 6 System Advanced Parameter Settings

In the main interface, press **MENU** and **UP** keys to enter the system parameter setting interface. In this interface, press **UP** and **DOWN** keys for advanced parameter selection, press **START/STOP** key to enter the submenu, press **UP** or **DOWN** key to adjust parameters. To return to the previous menu, press

**START/STOP** key. To return to the main interface, press **MENU** key.


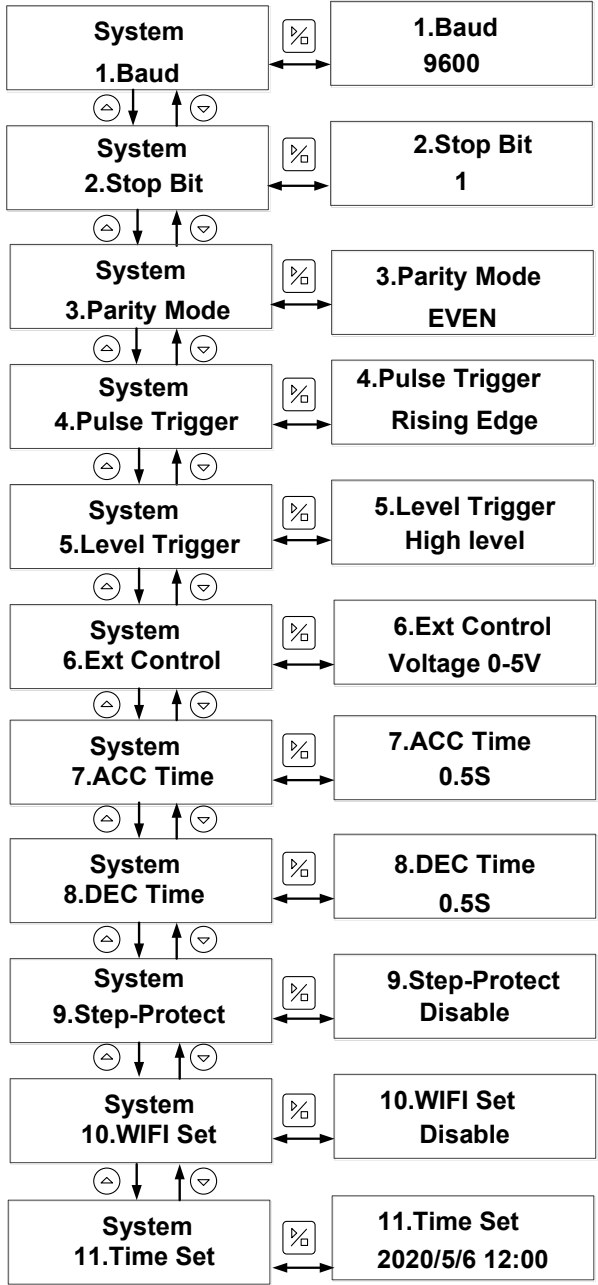
Figure 10 System advanced parameter setting flow chart

- 1. Baud: In the communication mode, set the baud to 4800, 9600, 19200, 38400 (the default is 9600).
- 2. Stop Bit: In the communication mode, set the size of the stop bit, which can be set to 1 and 2 (the default is 1).
- 3. Parity Mode: In the communication mode, set the parity type, which can be set as odd parity, even parity and none parity (the default is even parity).
- 4. Pulse Trigger: In Internal Control or Time Dispense mode, set the pulse trigger type of external control start / stop signal of the driver, which can be set as rising edge trigger and falling edge trigger (the default is rising edge trigger).
- 5. Level trigger mode: in External Control or Level mode, set the level trigger type when the driver is controlled externally, which can be set as high-level trigger and low-level trigger (the default is high-level trigger).
- 6. External Control mode: In the External Control mode, set the parameter type of analog signal to control the driver speed, which can be set to voltage mode 0-5V, voltage mode 0-10V, and current mode 4-20mA (the driver speed will change linearly with the change of external control analog signal).
- 7. Acceleration time: Set the acceleration time when the drive is started and running. The unit is seconds (the default value is 0.5 seconds).
- 8. Deceleration time: Set the deceleration time when the drive stops running. The unit is seconds (the default value is 0.5 seconds).
- 9. Step-Protect: The setting of step protection is off by default (WIFI version has this function).
- 10. WIFI setting: The WIFI switch of the driver is off by default (WIFI version has this function).
- 11. Time setting: Set the current date and time. Press **PRIME** key to change the set parameters, press **UP** or **DOWN** key to adjust the time, and press **START/STOP** key to save and exit.

## 7 External Control Interface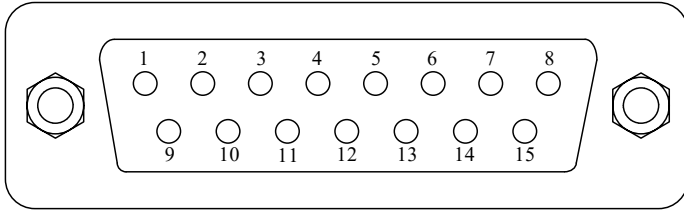

DB15	Mark	Note
1	ADC_W	Positive of external analogue input
2	B	Communication interface, B pole of RS485
3	A	Communication interface, A pole of RS485
4	VCC_W	External DC power input
5	-	-
6	CW_W	External input signal to control the direction
7	PWM	Pulse output
8	COM	Ground of external power
9	AGND	Negative of analogue signal input
10	+12V	Positive of internal +12V power source
11	GND	Ground of Internal power source
12	CW	Internal direction signal output
13	RS_W	External start/stop signal input terminal
14	PWM_W	External pulse signal input
15	RS	Internal start/stop signal output

## 8 Operating Instructions

### 8.1 Before Operation

- 1) Please check the packing slip to ensure that all parts are included and intact in the package. If there is a problem, please contact the

manufacturer or distributor.

- 2) Read the instruction.
- 3) A minimum of 200 mm space from the back of the pump should be maintained when it is in operation.

## 8.2 Power Connection

The voltage of the power supply should match what is indicated on the rating plate label of the pump. Please plug the power cord into the IEC Power Connector on the rear of the pump and plug the opposite end of the power cord into an electrical outlet. Flip the power switch located on the rear of the pump.

## 8.3 Mode Change

Turn on the power switch. The display will show a welcome message then the main screen. Press the **MODE** key to change the working mode (*Figure 11*).


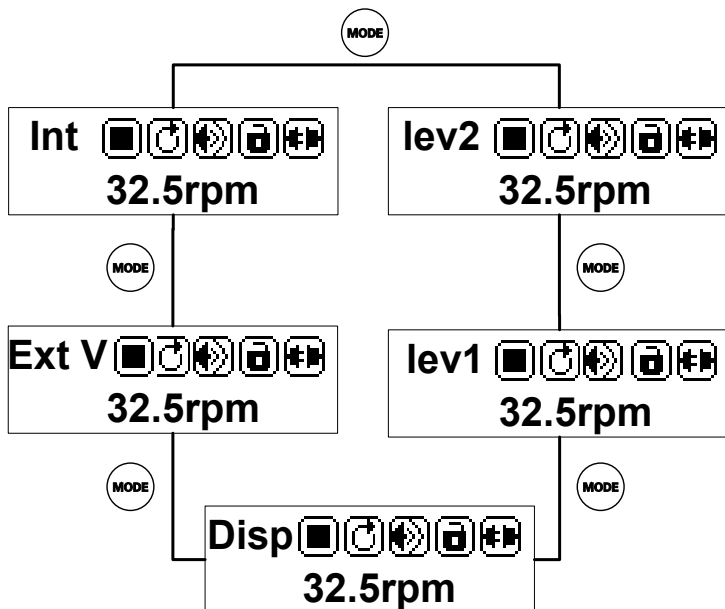
Figure 11. Change Working Mode

## 8.4 Internal Control Mode


Use the keypad to operate the pump. Use an optional external pulse signal to

control the start/stop.

- 1) Turn on the power switch. The pump will display the main screen.
- 2) Press the **MODE** key to change the mode to Internal Control mode (Int shown on the screen).
- 3) Press the **UP** or **DOWN** key to adjust the speed.
- 4) Press the **DIRECTION** key to change the rotating direction.
- 5) Press the **START/STOP** key to start or stop the pump.
- 6) Press the **PRIME** key, the pump will run at the maximum allowed speed.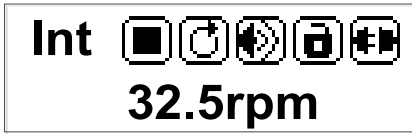

Normal Speed

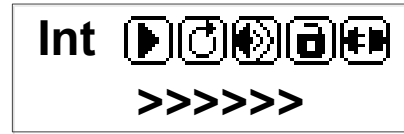Priming

Figure 12. Normal vs Priming

## 8.5 External Control Mode

In this mode, use an external logic level signal to control direction, start and stop; use an external analog signal to control rotation speed. The keypad is disabled.

For External Voltage Signal  
0-5V/0-10V

For External Current Signal  
4-20mA

Figure 13. External Control Mode

To control the pump by an external signal:

- 1) Switch the power off. Wire the DB15 connector as shown in [Figure 14](#) or [Figure 15](#). Connect it to the DB15 port on the rear of the pump.

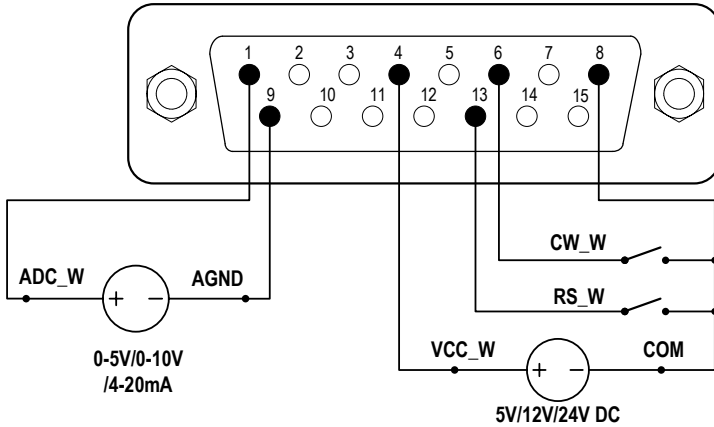
Figure 14. DB15 Wiring with an External Power Source

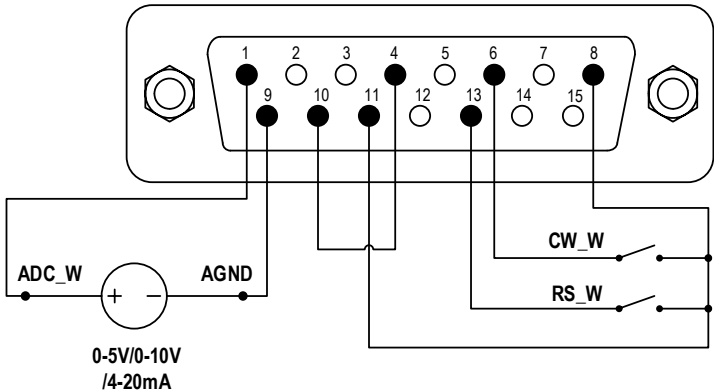
Figure 15. DB15 Wiring with the Internal 12VDC Power Source

- 2) Turn on the power switch. The pump will display the main screen.
- 3) Press the **MODE** key to change the mode to External Control mode (ExtV or ExtI shown on the screen).
- 4) Close the external RS\_W switch, and turn on the external analog signal power source. The speed will change according to the intensity of the input signal. Open the RS\_W to stop the drive.
- 5) Opening the CW\_W switch, the pump will run in the clockwise direction. Closing the CW\_W switch, the pump will run in the counterclockwise direction.

## 8.6 Time Dispense Mode

The pump will dispense fluid automatically by setting the duration for each dose, pause time between doses and the number of cycles. When dispensing, the display will show the dispensing time or lag time (1.2s in *Figure 16*), and the total cycles (4 in *Figure 16*) that has dispensed.

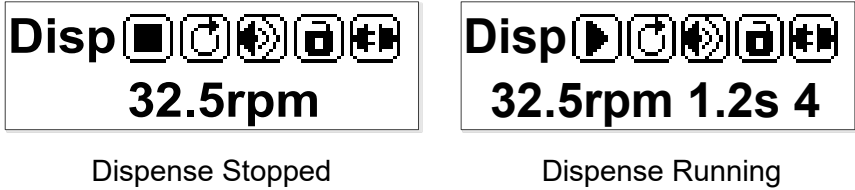
Figure 16. Time Dispense Mode

To set the Time Dispense Mode:

- 1) Turn on the power switch. The pump will display the main screen.
- 2) Press the **MODE** key to change the mode to Time Dispense mode (Disp shown on the screen).
- 3) Press the **MENU** key to enter the setting menu.
- 4) Set duration for each dose, pause time between doses and number of cycles.
- 5) Return to the main screen.

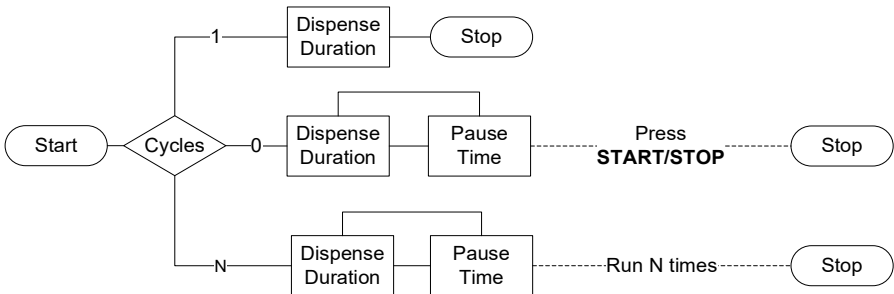
Figure 17. Dispense Cycle Setting

To run the Time Dispense mode:

- 1) Press the **DIRECTION** key to change the running direction, clockwise or counterclockwise.
- 2) Press the **START/STOP** key to start dispensing.
- 3) When the drive is running, press the **START/STOP** key to stop the drive at any time.
- 4) A footswitch can be used to start/stop the pump.

## 8.7 Logic Level 1 Control Mode (footswitch)

Use an external logic-level signal to control the start and stop. Use the keypad to control direction and speed.

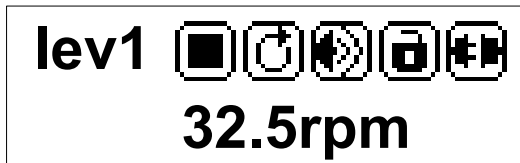
Figure 18. Logic Level 1 Control Mode

- 1) Switch the power off. Wire the DB15 connector as shown in [Figure 19](#) or [Figure 20](#), and connect it to the DB15 port on the rear of the pump.

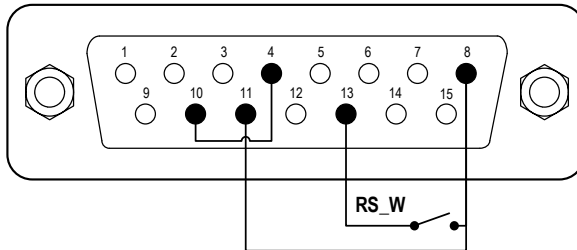
Figure 19. Logic Level 1 Control with Internal 12V Power Source

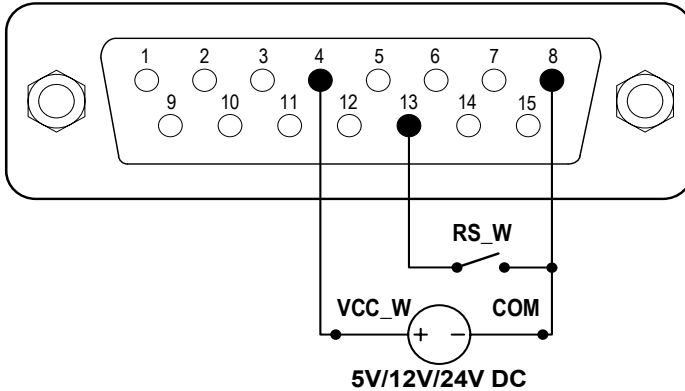
Figure 20. Logic Level 1 Control with an External Power Source

- 2) Turn on the power switch. The pump will display the main screen.
- 3) Press the **MODE** key to change the mode to Logic Level 1 control mode (shows lev1 on the screen).
- 4) Press the **UP** or **DOWN** key to adjust the speed.
- 5) Press the **DIRECTION** key to change the rotating direction.
- 6) When the switch is closed, the drive will start running; when the switch is open, the drive will stop.

Note: Use this mode to work with a TIME CONTROLLER.

## 8.8 Logic Level 2 Control Mode (footswitch and direction switch)

Use an external logic-level signal to control start, stop and direction. Use the keypad to control speed.

Figure 21. Logic Level 2 Control Mode

- 1) Switch the power off. Wire the DB15 connector as shown in [Figure 22](#)

or Figure 23, and connect it to the DB15 port on the rear of the pump.

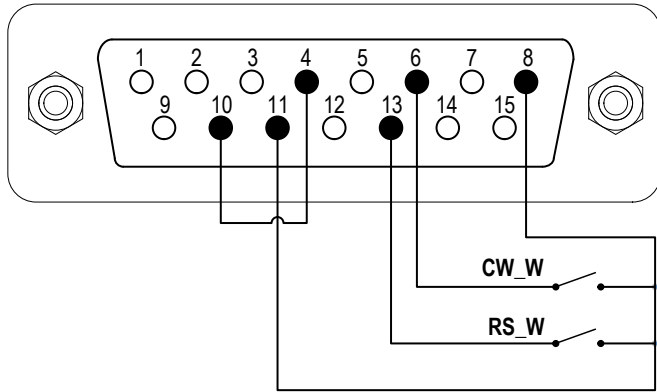
Figure 22. Logic Level 2 Control with the Internal 12V Power Source

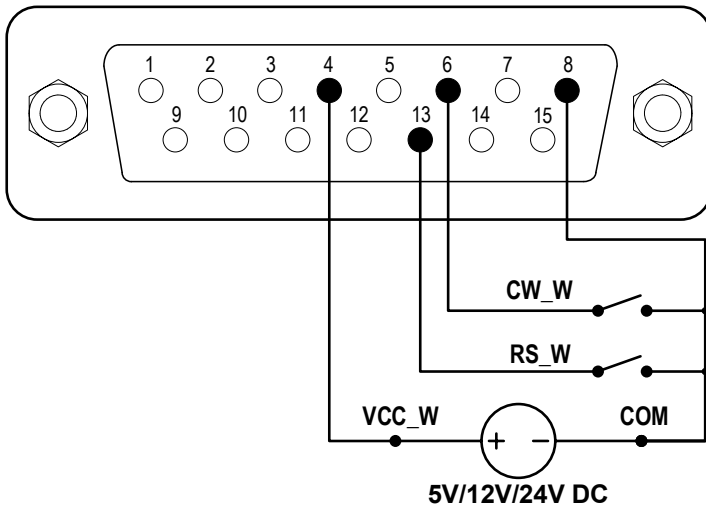
Figure 23. Logic Level 2 Control with an External Power Source

- 2) Turn on the power switch. The pump will display the main screen.
- 3) Press the **MODE** key to change the mode to Logic Level 2 control mode (shows lev2 on the screen).
- 4) Press the **UP** or **DOWN** key to adjust the speed.
- 5) When the switch RS\_W is closed, the drive will be running at the set

speed; when the switch RS\_W is opened, the drive will stop.

- 6) When the switch CW\_W is open, the drive will be running clockwise. When the switch CW\_W is closed, the drive will be running counterclockwise.

## 8.9 Communication Mode

The RS485 interface supports the standard MODBUS protocol. The pump can communicate with external devices via the communication port. Please refer to the [Communication Instruction manual](#) for the parameters and supported commands.

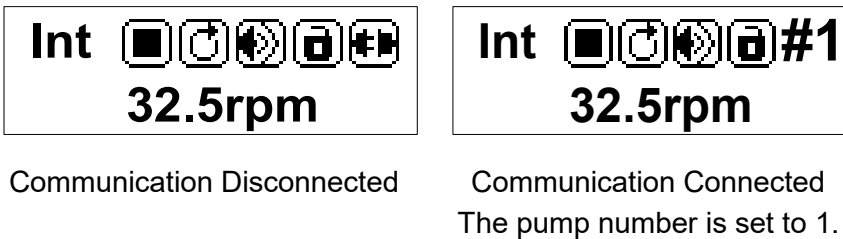
Figure 24. Communication Mode

- 1) When the power is off, wire the DB15 connector as shown in [Figure 25](#) and connect it to the DB15 port on the rear of the pump. An external DC power source is recommended to avoid electrical interference.

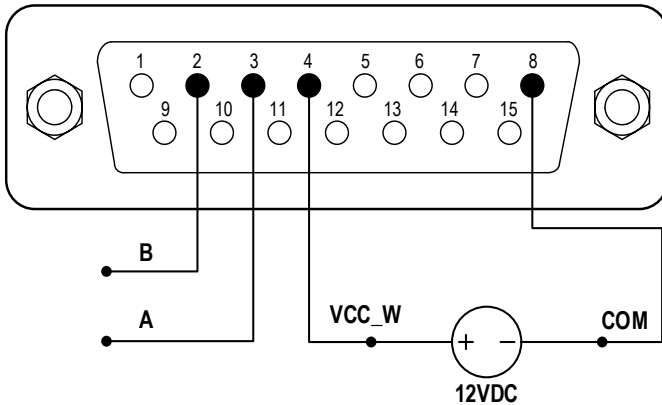
Figure 25. RS485 MODBUS Wiring

- 2) Turn on the power switch. The pump will display the main screen.

- 3) Press the **MODE** key to change the mode to Internal Control mode or Time Dispense mode.
- 4) Control pump with the communication interface.
- 5) Press the **START/STOP** key to stop the drive at any time.

## 8.10 Speed Setting

The speed resolution is 0.1 rpm when the speed is between 0.1 and 100 rpm; the speed resolution is 1 rpm when the speed is between 100 and 600 rpm. On the main screen, the speed can be set by pressing the **UP** or **DOWN** key. Pressing the **UP** or **DOWN** key shortly, the last digit of the value will change by 1. Hold the **UP** or **DOWN** key to change the value fast.

Hold the **PRIME** key and press the **UP** key to set the speed to the maximum directly. Hold the **PRIME** key and press the **DOWN** key to set the speed to the minimum (0.1 rpm) directly.

## 9 Maintenance

### 9.1 Warranty

The product comes with one-year labor and parts warranty. The limited warranty does not cover any damage that is caused by improper usage and handling.

### 9.2 Regular Maintenance

- 1) Always check the tubing and connections to avoid leakage.
- 2) Do not cover the fan on the rear of the pump.
- 3) Do not use water to wash the pump. Keep the pump head dry.
- 4) Do not use chemical solvents to clean the pump and pump head.

### 9.3 Malfunction Solutions

No.	Mal-function	Description	Solution
1	Hardware	No display	1. Check the power cord 2. Check the fuse. If it was blown, replace it with a 1A slow-blow fuse

BT100S/BT300S/BT600S/ BT100S-1 Peristaltic Pump

			3. Check the internal power cord connection inside the pump.
2	Hardware	Motor does not work	<ol style="list-style-type: none"> <li>1. Check the indicator of the driver board.</li> <li>2. Check the wire connection between the motor and the driver board.</li> <li>3. Check the wire connection between the driver and the mainboard.</li> <li>4. Check the power voltage for the pump.</li> </ol>
3	Hardware	Motor is trembling	<ol style="list-style-type: none"> <li>1. Check the wire connection between the motor and the driver board.</li> <li>2. The motor is overloaded. Check the mechanical connection.</li> </ol>
4	Hardware	Motor only runs in one direction	Check the connection between the drive board and the main control board.
5	Hardware	Keypad does not work	<ol style="list-style-type: none"> <li>1. Check the wire connection between the keypad and the mainboard.</li> <li>2. Check if the key is broken.</li> </ol>
6	Hardware	External control does not work	<ol style="list-style-type: none"> <li>1. Check the wiring of the connector.</li> <li>2. Check if the external control power voltage is provided.</li> <li>3. Check the connections of the external control board.</li> </ol>
7	Hardware	RS485 com does not work	<ol style="list-style-type: none"> <li>1. Check the wiring of the connector.</li> <li>2. Check if the external control power voltage is provided.</li> <li>3. Check the connections of the communication board.</li> </ol>
8	Hardware	Noisy when running	Check the screws and level on the pump head to make sure they are secure.
9	Software	External control does not work	Check if the pump is on External Control Mode.

10	Software	RS485 does not work properly	<ol style="list-style-type: none"> <li>1. Check if the display shows that the communication is ready.</li> <li>2. Reset the address of the pump.</li> <li>3. Check whether on the bus there are two pumps using the same address</li> </ol>
----	----------	------------------------------	---

If the problem cannot be solved, please contact the manufacturer or distributor.

## 10 Dimensions

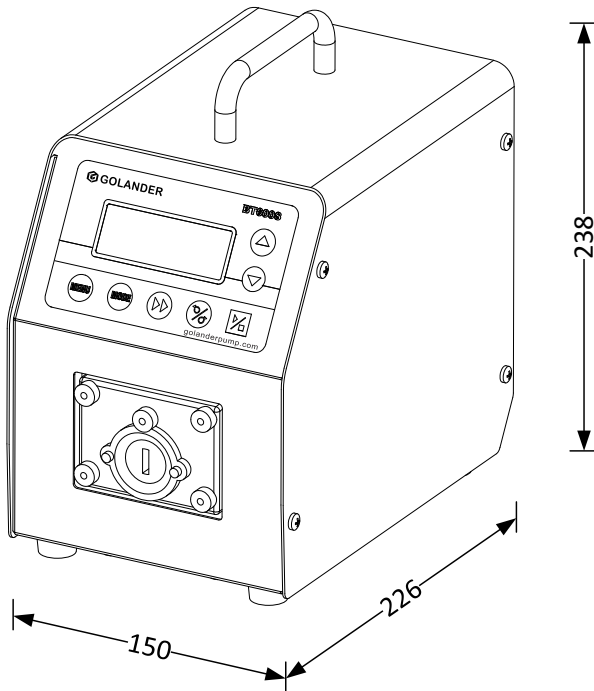
Figure 26. BT100S/BT300S/BT600S Dimensions (mm)


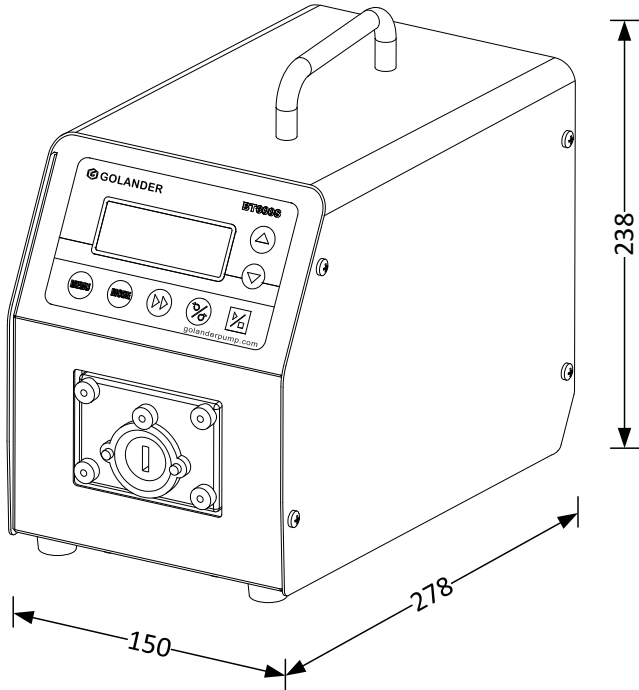
Figure 27. BT100S-1 Dimensions (mm)

## 11 Naming Rule
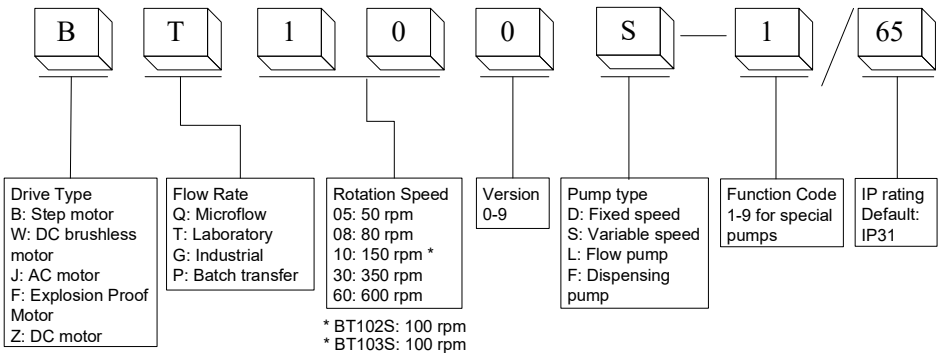
## 12 Specifications

Speed resolution	0.1-99.9 rpm, 0.1 rpm resolution; 100-600 rpm, 1 rpm resolution
Speed accuracy	0.5%
Power supply	AC 100-240V, 50/60Hz
Power consumption	BT100S: < 40W BT300S:< 50W BT600S: < 60W BT100S-1: <40W
External logic level control signal	5V, 12V (standard), 24V (optional)
External analog control signal	0-5V (standard); 0-10V, 4-20mA (optional)
Communication interface	RS485 MODBUS
Operating condition	Temperature 0-40°C Relative humidity <80%
IP grade	IP31
Display	132x32 LCD
Compliance (For CE Marking)	2014/35/EU (Low Voltage) 2014/30/EU (Electromagnetic Compatibility)
Dimensions (LxWxH)	BT100S/BT300S/BT600S: 238 x 150 x 226 mm (9.37 x 5.91 x 8.90 inch) BT100S-1: 278 x 150 x 238 mm (10.94 x 5.91 x 9.37 inch)
Weight	BT100S: 4.7 kg (10.4 lbs) BT300S: 4.9 kg (10.8 lbs) BT600S: 5.2 kg (11.5 lbs) BT100S-1: 4.9 kg (10.8 lbs)

**BT100S Suitable Pump Heads and Tubing, Flow Parameters**

<b>Pump head</b>	<b>No. of Channels</b>	<b>Tubing size (mm)</b>	<b>Flow rate per channel (mL/min)</b>
DG6-1 (6rollers)	1	Wall:0.8~1, ID:≤3.17	0.0002-49
DG10-1 (10rollers)	1	Wall:0.8~1, ID:≤ 3.17	0.0001-41
DG6-2 (6rollers)	2	Wall:0.8~1, ID:≤ 3.17	0.0002-49
DG10-2 (10rollers)	2	Wall:0.8~1, ID:≤ 3.17	0.0001-41
DG6-4 (6rollers)	4	Wall:0.8~1, ID:≤ 3.17	0.0002-49
DG10-4 (10rollers)	4	Wall:0.8~1, ID:≤ 3.17	0.0001-41
DT10-18	1	13# 14#, Wall: 0.8~1, ID:≤3.17	0.0002-82
DT10-28	2	13# 14#, Wall: 0.8~1, ID:≤3.17	0.0002-82
DT10-48	4	13# 14#, Wall: 0.8~1, ID:≤3.17	0.0002-82
YZ15	1	13# 14# 16# 19# 25# 17#	0.006-420
YZ25	1	15# 24#	0.16-420
2 x YZ15	2	13# 14# 16# 19# 25# 17#	0.006-420
2 x YZ25	2	15# 24#	0.16-420
YT15	1	13# 14# 16# 19# 25# 17# 18#	0.006-570
YT25	1	15# 24# 35# 36#	0.17-720
2 x YT15	2	13# 14# 16# 19# 25# 17# 18#	0.006-570
2 x YT25	2	15# 24# 35# 36#	0.17~720
DT15-14	1	19# 16# 25# 17#	0.07-560
DT15-24	2	19# 16# 25# 17#	0.07-560
DT15-44	4	19# 16# 25#	0.07-330

**BT300S Suitable Pump Heads and Tubing, Flow Parameters**

<b>Pump Head</b>	<b>No. of Channels</b>	<b>Tubing size</b>	<b>Flow rate per channel (mL/min)</b>
YZ15	1	13# 14# 16# 19# 25# 17#	0.006~990
YZ25	1	15# 24#	0.17~990
2 x YZ15	2	13# 14# 16# 19# 25# 17#	0.006~990
2 x YZ25	2	15# 24#	0.17~990
DT15-14	1	19# 16# 25# 17#	0.07~1300
DT15-24	2	19# 16# 25# 17#	0.05~930
YT15	1	13# 14# 16# 19# 25# 17# 18#	0.006~1300
YT25	1	15# 24# 35# 36#	0.17~1690
2 x YT15	2	13# 14# 16# 19# 25# 17# 18#	0.006~1300
2 x YT25	2	15# 24# 35# 36#	0.17~1690

**BT600S Suitable Pump Heads and Tubing, Flow Parameters**

<b>Pump head</b>	<b>No. of Channels</b>	<b>Tubing size</b>	<b>Flow rate per channel (mL/min)</b>
YZ15	1	13# 14# 16# 19# 25# 17#	0.006~1700
YZ25	1	15# 24#	0.17~1700
2 x YZ15	2	13# 14# 16# 19# 25# 17#	0.006~1700
2 x YZ25	2	15# 24#	0.17~1700
DT15-14	1	19# 16# 25# 17#	0.07~2240
DT15-24	2	19# 16# 25# 17#	0.07~2240
YT15	1	13# 14# 16# 19# 25# 17# 18#	0.006~2300
YT25	1	15# 24# 35# 36#	0.17~2900
2 x YT15	2	13# 14# 16# 19# 25# 17# 18#	0.006~2300

**BT100S-1 Suitable Pump Heads and Tubing, Flow Parameters**

<b>Pump head</b>	<b>No. of Channels</b>	<b>Tubing size (mm)</b>	<b>Flow rate per channel (mL/min)</b>
DG6-8 (6 rollers)	8	wall thickness 0.8 ~ 1, ID≤3.17	0.00016 ~ 49
DG10-8 (10 rollers)	8	wall thickness 0.8 ~ 1, ID≤3.17	0.00011 ~ 41
DG6-12 (6 rollers)	12	wall thickness 0.8 ~ 1, ID≤3.17	0.00016 ~ 49
DG10-12 (10 rollers)	12	wall thickness 0.8 ~ 1, ID≤3.17	0.00011 ~ 41
DG6-16 (6 rollers)	16	wall thickness 0.8 ~ 1, ID≤3.17	0.00016 ~ 49
DG10-16 (10 rollers)	16	wall thickness 0.8 ~ 1, ID≤3.17	0.00011 ~ 41
DG6-24 (6 rollers)	24	wall thickness 0.8 ~ 1, ID≤3.17	0.00016 ~ 49
DG10-24 (10 rollers)	24	wall thickness 0.8 ~ 1, ID≤3.17	0.00011 ~ 41
DT10-88	8	wall thickness 0.8 ~ 1, ID≤3.17, 13# 14#	0.0002 ~ 82
YZ15	1	13# 14# 16# 19# 25# 17#	0.006 ~ 420
YZ25	1	15# 24#	0.16 ~ 420
2×YZ15	2	13# 14# 16# 19# 25# 17#	0.006 ~ 420
2×YZ25	2	15# 24#	0.16 ~ 420
4×YZ15	4	13# 14# 16# 19# 25# 17#	0.006 ~ 420
4×YZ25	4	15# 24#	0.16 ~ 420

BT100S/BT300S/BT600S/ BT100S-1 Peristaltic Pump

KZ25	1	15# 24# 35# 36#	0.25 ~ 1500
2×KZ25	2	15# 24# 35# 36#	0.25 ~ 1500
DT15-44	4	19# 16# 25# 17#	0.05 ~ 400
YT15	1	13# 14# 16# 19# 25# 17# 18#	0.006 ~ 570
YT25	1	15# 24# 35# 36#	0.17 ~ 720
2×YT15	2	13# 14# 16# 19# 25# 17# 18#	0.006 ~ 570
2×YT25	2	15# 24# 35# 36#	0.17 ~ 720
4×YT15	4	13# 14# 16# 19# 25# 17# 18#	0.006 ~ 570
4×YT25	4	15# 24# 35# 36#	0.17 ~ 720

**Golander LLC**

4405 International Blvd  
Ste B117, Norcross, GA 30093  
USA

Tel: +1 678-587-8806

[info@golanderpump.com](mailto:info@golanderpump.com)

[www.golanderpump.com](http://www.golanderpump.com)

**Golander GmbH**

Dechant-Heimbach-Str. 29  
53177 Bonn  
Germany

Tel: +49 228 50446952

[info@golander.de](mailto:info@golander.de)

[www.golander.de](http://www.golander.de)