
WT600S-65

Variable-Speed Peristaltic Pump

Operating Manual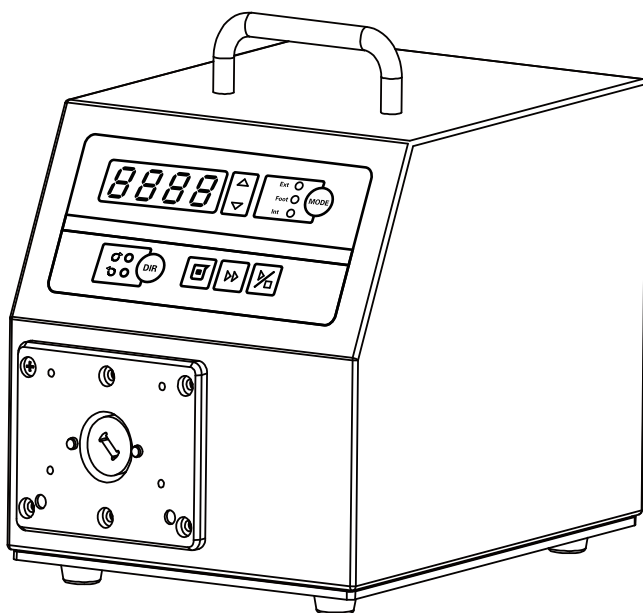

Contents

| | |
|---|----|
| Safety Precautions | 1 |
| 1 Description | 6 |
| 2 Functions and Features..... | 6 |
| 3 Components and Connectors..... | 8 |
| 4 Display Panel and Operating Keypads | 9 |
| 4.1 Keypad..... | 9 |
| 4.2 Digital LED and Indicators..... | 10 |
| 5 External Control Interface..... | 12 |
| 6 Operation Instructions | 13 |
| 6.1 Prepare before operating | 13 |
| 6.2 Power Connection..... | 13 |
| 6.3 Mode Change..... | 13 |
| 6.4 Internal Control Mode..... | 13 |
| 6.5 External Control Mode..... | 14 |
| 6.6 Footswitch Control Mode..... | 16 |
| 6.7 Logic Level Control Mode..... | 18 |
| 6.8 Communication Mode | 19 |
| 6.9 Footswitch Control | 19 |
| 7 Maintenance | 21 |
| 7.1 Warranty..... | 21 |
| 7.2 Regular Maintenance | 21 |
| 7.3 Malfunction Solutions | 21 |
| 8 Dimensions | 23 |
| 9 Naming Rule | 23 |
| 10 Specifications..... | 24 |

Safety Precautions

Danger

- To avoid damage, use the correct voltage indicated on the rating plate label of the pump.
- Make no unauthorized dismantling, changes or modifications to the pump. This may result in malfunctions or potential accidents.
- Turn off the pump drive before installing or removing tubing, attempting any maintenance, cleaning or repair of the drive, or connecting or disconnecting external control devices or a communication interface. Fingers or loose clothing may get caught in the drive mechanism.

Warning

- Ensure no chemical reactions occur between the handled fluid and pump head/tubing material before use.
- Check tubing regularly to avoid breakage. Breakage may result in fluid being sprayed from the pump. Use appropriate measures to protect the operator and equipment. The operator is solely liable for damages resulting from tubing breakage, particularly the leakage of toxic or valuable liquids.
- The pump is equipped with a plug which must be well grounded.
- This device is not designed nor intended for usage in patient-connected applications, including but not limited to medical and dental use.
- Observe all other applicable regulations concerning working and operational safety, environmental protection, and relevant local guidelines.

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1 Description

The WT600S-65 variable-speed peristaltic pump offers a flow range from 3 to 6000 mL/min and a speed range of 30-600 rpm. The IP65-rated enclosure ensures protection against dust and resistance to water. In addition to fundamental functionalities like reversible direction, start/stop, and adjustable speed, it includes the *Time Dispense* feature. The brushless drive provides substantial torque, enabling compatibility with multiple pump heads. Equipped with a MODBUS RS485 interface, the pump facilitates easy connection to external devices such as computers, human-machine interfaces, or PLCs.

2 Functions and Features

Peristaltic pumps excel in managing abrasive, corrosive, or viscous fluids. By eliminating valves that could clog and ensuring no seals directly contact the pumped medium, these pumps provide sleek inner surfaces that are easy to clean, with the fluid only interacting with the tubing or tube material. With a suction lift and priming capacity of up to 8 meters of water column at sea level, the pump demonstrates versatility. It handles shear-sensitive fluids, like latex or firefighting foam, and can operate in a dry state while efficiently pumping fluids with high levels of entrained air, such as black liquor soap. The pump's elevated volumetric efficiency renders it suitable for precise metering or dosing applications. Moreover, tubing options are available to meet the specific requirements for applications in the food and pharmaceutical industries.

- Four-digit LED displays indicate speed and working mode.
- LED indicator reflects the current working state.
- Membrane keypad for user input.
- Reversible direction, start/stop control, and adjustable speed.
- Time Dispense Mode allows setting dose duration for one-press dispensing.
- Brushless drive with large torque for multiple pump heads; no maintenance needed.

- 3% precision rotating speed control with 1 rpm resolution.
- External signals control start/stop, direction, and dispense; optically isolated.
- RS485 MODBUS interface for easy external control.
- Internal double-deck isolation, conformal coating for dust and moisture protection.
- Anti-electromagnetic interference, wide input voltage range.
- Stainless steel enclosure for easy cleaning and resistance to corrosion.

3 Components and Connectors

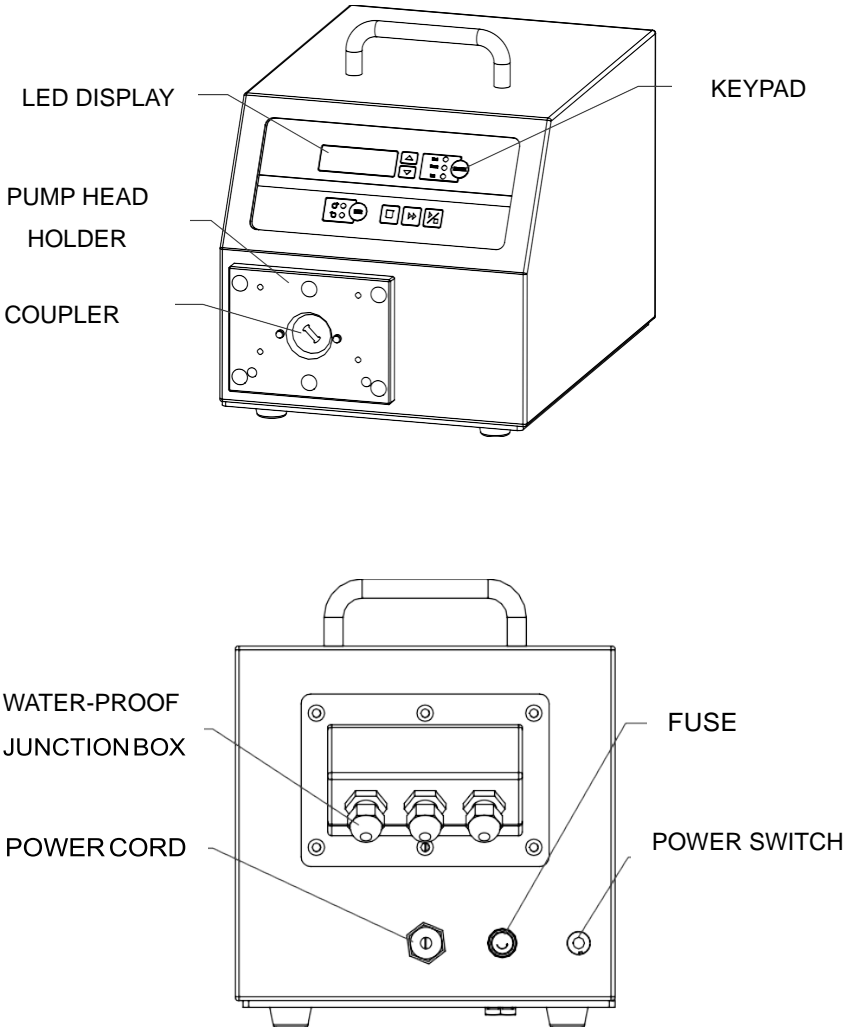
Figure 1. Components and Connectors

4 Display Panel and Operating Keypads



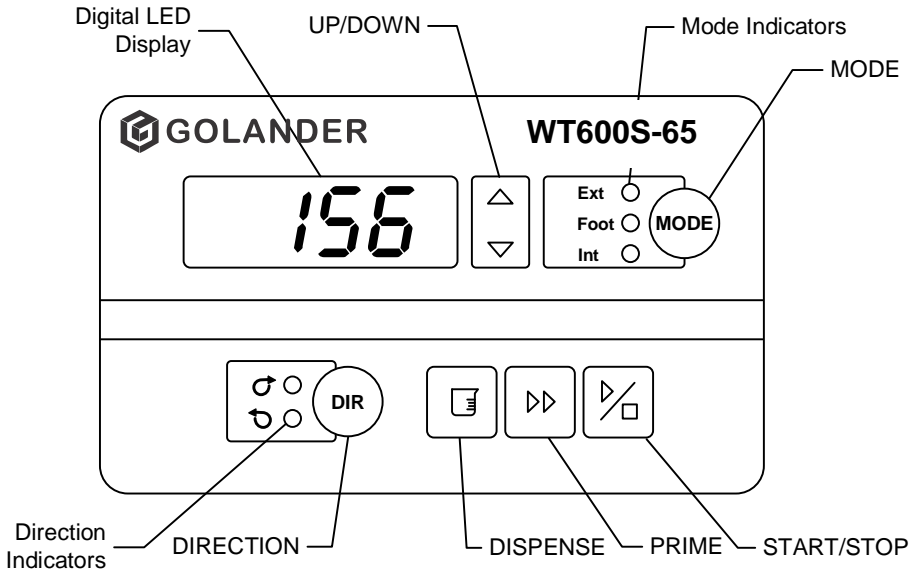
Figure 2. Display panel

4.1 Keypad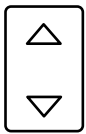

UP Key. When held momentarily, the last digit of the value will increase by 1. Hold the key to quickly increase the value.

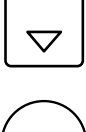DOWN Key. When held momentarily, the last digit of the value will decrease by 1. Hold the key to decrease the value quickly.

MODE key. When the drive is not running, use the MODE key to change the working mode: Internal Control mode, External Control mode, Footswitch Control Mode or Logic Level Control Mode.

DIRECTION Key. Press the direction key to change the rotating direction between clockwise and counterclockwise

DISPENSE key. When the drive is not running, use the DISPENSE key to dispense. Press and hold the key to enter the Time Dispense Setup Mode.


PRIME key. Press the key to run the pump at the maximum speed in the direction on the display. Press again to return to the previous state.


START/STOP key. Press the key to start or stop the pump.

4.2 Digital LED and Indicators


4.2.1 Indicators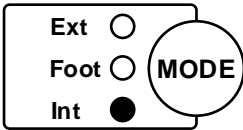


Int: Internal Control Mode. Use the keypad to operate the pump. A footswitch can be used to control start /stop.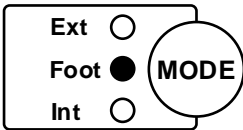


Foot: Footswitch Control Mode. Use a footswitch or the DISPENSE key to start or stop dispensing. Use the keypad for other operations.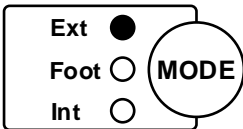


Ext: External Control Mode. Use an external analog signal to control rotation speed. Use an external logic level signal to control direction, start and stop. The keypad is disabled.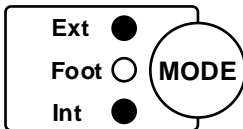

Int+Ext: Logic Level Control Mode. Use a logic-level signal to control the start and stop. Use the keypad to control direction and speed.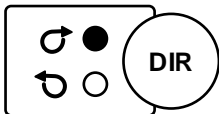

Clockwise Direction: The motor runs clockwise.

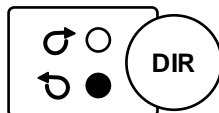Counterclockwise Direction: The motor runs counterclockwise.

4.2.2 Digital LED Display

The digital LED displays the rotating speed and working mode.

Figure 3. Internal Control Mode/Footswitch control Mode, 215 rpm

Figure 4. External Control Mode, 355 rpm

Figure 5. Running Time Dispense at 115 rpm

Figure 6. When flashing, the pump is in Time Dispense Setup Mode

Dispense Duration 8.5 seconds


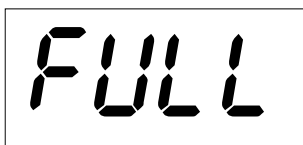
Figure 7. Pump priming, the drive is running at full speed

5 External Control Interface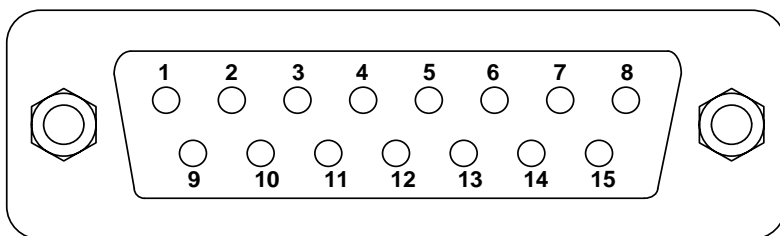

| DB15 | Mark | Note |
|------|-------|--|
| 1 | ADC_W | Positive of external analog input |
| 2 | B | Communication interface, B pole of RS485 |
| 3 | A | Communication interface, A pole of RS485 |
| 4 | VCC_W | External DC power input |
| 5 | | |
| 6 | CW_W | External input signal to control the direction |
| 7 | | |
| 8 | COM | Ground of external power |
| 9 | AGND | Negative of analog signal input |
| 10 | +12V | Positive of internal +12V power source |
| 11 | GND | Ground of Internal power source |
| 12 | | |
| 13 | RS_W | External start/stop signal input terminal |
| 14 | | |
| 15 | | |

6 Operation Instructions

6.1 Prepare before operating

- 1) Please check the packing slip to ensure all parts are included and intact. If there is a problem, please contact the manufacturer or distributor.
- 2) Read the instructions.
- 3) At least 200mm of space from the back of the pump should be maintained when in operation.

6.2 Power Connection

The voltage of the power supply should match what is indicated on the rating plate label of the pump. Plug the power cord into the IEC Power Connector on the rear of the pump and plug the opposite end of the power cord into an electrical outlet. Flip the power switch on the rear of the pump.

6.3 Mode Change

Turn on the power switch. The display will be on. Press the MODE key to change the working mode.


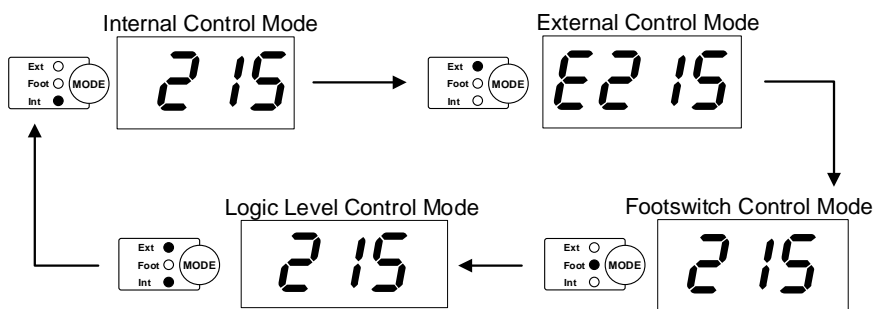
Figure 8. Change Working Mode

6.4 Internal Control Mode

In this mode, use the keypad to operate the pump.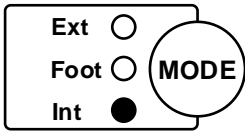

- 1) Turn on the power. The display will be on.
- 2) Press the MODE key to change the mode to Internal Control Mode (the Int indicator is on).
- 3) Press the UP or DOWN key to set the speed.
- 4) Press the DIRECTION key to change the rotating direction.
- 5) Press the START/STOP key to start or stop the drive.
- 6) Press the PRIME key, the drive will run at full speed.

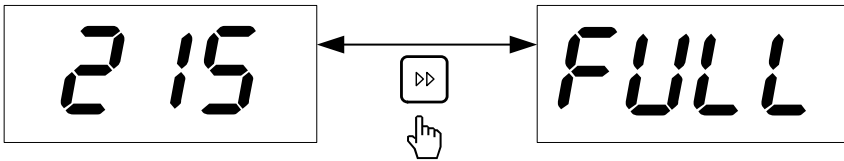
Figure 9. Prime the Pump

When in the Internal Control Mode, an optional footswitch can be used to start/stop the pump. The drive will be running when the

6.5 External Control Mode

Use the external logic level signal to control direction, start and stop. Use an external analog signal to control rotation speed. The keypad is disabled. The analog signal could be 0-5V, 0-10V or 4-20mA. The signal is 0-5V by default. For 0-10V or 4-20mA, the dip switch setting on the analog signal control board must be changed.

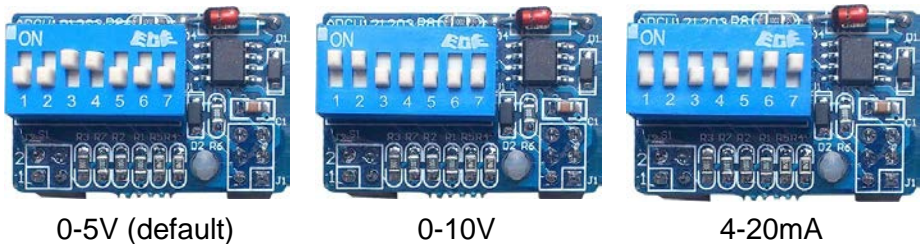
Figure 10. Analog Signal Control Board Setting

To control the pump by an external signal

- 1) Switch the power off. Wire the DB15 connector as shown in Figure

11 or Figure 12. Connect it to the DB15 port on the rear of the pump.

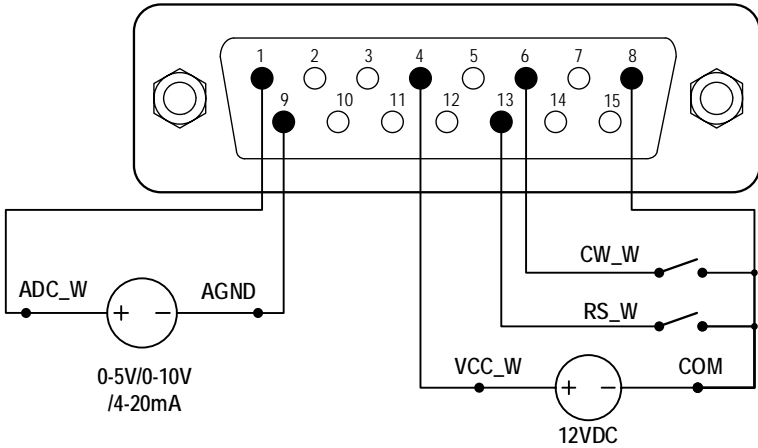
Figure 11. DB15 Wiring with External 12VDC Power Source


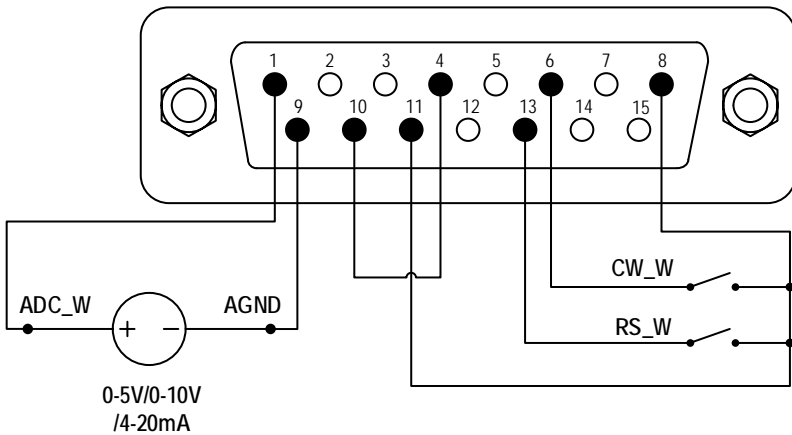
Figure 12. DB15 Wiring with Internal 12VDC Power Source

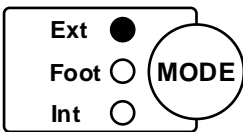- 2) Turn on the power switch. The LED display will be on.
- 3) Press the MODE key to change the mode to External Control Mode.
- 4) Close the external RS_W switch, and turn on the external analog signal power source. The speed will change according to the

intensity of the input signal. Open the RS_W to stop the drive.

- 5) Opening the CW_W switch, the drive will run clockwise. Closing the CW_W switch, the drive will run counterclockwise.

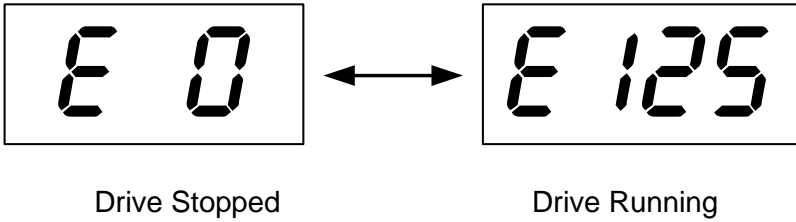
Figure 13. External Control Mode

Note: The external DC power source can be 5V and 12V. If it is 24V, a 1.5K resistor is needed to protect the internal circuit.


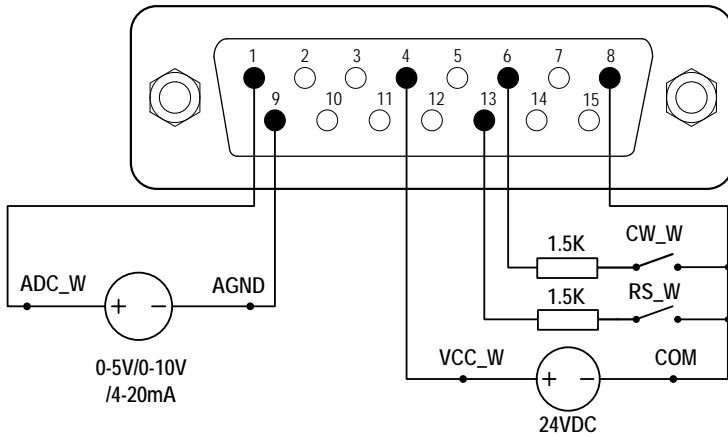
Figure 14. DB15 Wiring with External 24VDC Power Source

6.6 Footswitch Control Mode

The pump will dispense fluid by setting the duration time for each dose. The drive will stop automatically when dispensing is completed.

To set the duration time for each dose

- 1) Turn on the power switch. The display will be on.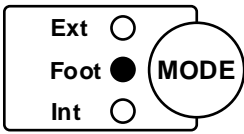

2) Press the MODE key to change the mode to Internal Control Mode or Footswitch Control Mode.

3) Press and hold the DISPENSE key for 3 seconds until the four-digit LED display starts to flash. The pump is in the Time Dispense Setup Mode.

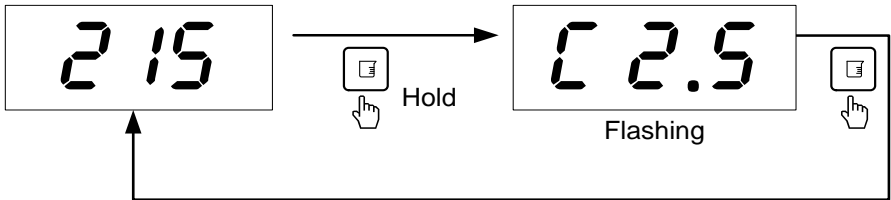
Figure 15. Time Dispense Setup Mode

- 4) In the Time Dispense Setup Mode, the duration time can be set by pressing the UP or DOWN key. The time range is 0.1 to 999 seconds.
- 5) Press the DISPENSE key again to exit the Time Dispense Setup Mode.

To Dispense Fluid

- 1) Press the DIRECTION key to adjust the running direction, clockwise or counterclockwise.
- 2) In the Internal Control Mode or Footswitch Control Mode, press the DISPENSE key to dispense one dose. Press the DISPENSE key again to repeat dispensing.
- 3) When in the Footswitch Control Mode, an optional footswitch can be used to dispense. Close then open the footswitch to dispense one dose. Close then open the footswitch again to repeat dispensing (see section [6.9](#)).
- 4) When the drive is running, press the START/STOP key to stop the process anytime.

6.7 Logic Level Control Mode

In this mode, use a logic-level signal to control the start and stop. The keys on the front panel can control direction and speed.

- 1) Switch the power off. Wire the DB15 connector as shown in [Figure 16](#) or [Figure 17](#), and connect it to the DB15 port on the rear of the pump.

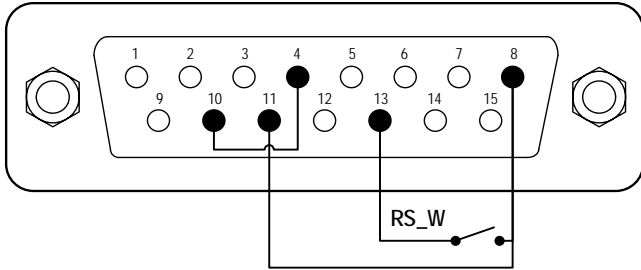
Figure 16. Logic Level Control with Internal 12V Power Source


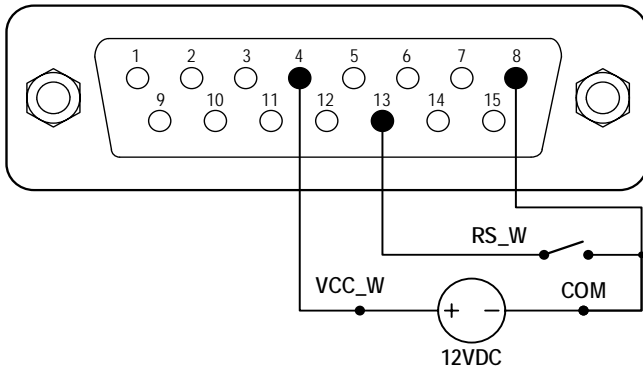
Figure 17. Logic Level Control with External 12V Power Source

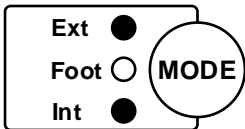- 2) Turn on the power switch. The display will be on.
- 3) Press the MODE key to change the mode to Logic Level Control Mode.
- 4) Press the UP or DOWN key to set the speed.
- 5) Press the DIRECTION key to change the rotating direction.
- 6) When the RS_W switch is closed, the drive will be running; when

the RS_W switch is open, the drive will stop.

Note: Use this mode to work with a TIME CONTROLLER.

6.8 Communication Mode

The RS485 interface supports the standard MODBUS protocol and allows the pump to communicate with external devices via the communication port. Please refer to the [Communication Instruction Manual](#) for the parameters and supported commands.

- 1) When the power is off, wire the DB15 connector as shown in [Figure 18](#), and connect it to the DB15 port on the rear of the pump. An external DC power source is recommended to avoid electrical interference.

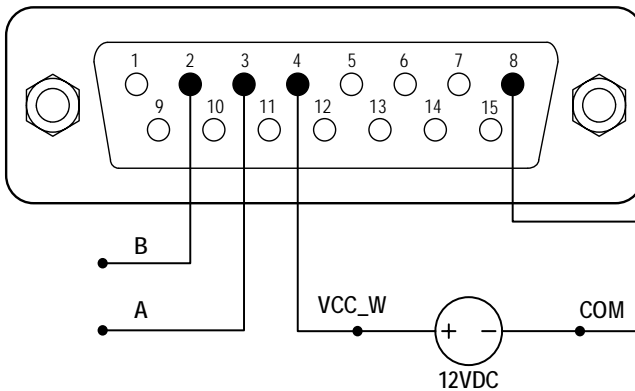
Figure 18. RS485 MODBUS Wiring

- 2) Turn on the power switch. The LED display will be on.
- 3) Press the MODE key to change the mode to Internal Control Mode or Footswitch Control Mode.
- 4) Control pump with the communication interface.
- 5) Press the START/STOP key to stop the drive anytime.

6.9 Footswitch Control

- 1) Switch the power off. Wire the DB15 connector as shown in [Figure 19](#) or [Figure 20](#), and connect it to the DB15 port on the rear of the pump.

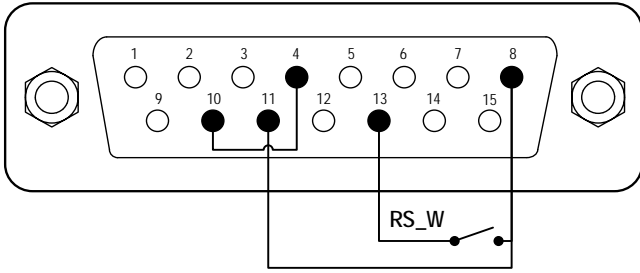
Figure 19. Footswitch with internal 12V Power Source

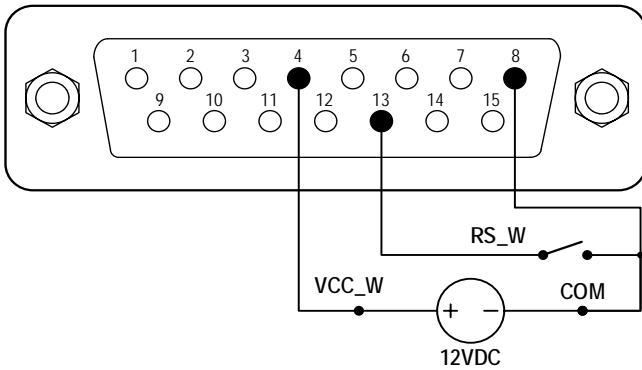
Figure 20. Footswitch with External 12V Power Source

- 2) Turn on the power switch. The LED display will be on.
 - 3) Press the MODE key to change the working mode.
- In the Internal Control Mode: When the switch RS_W is closed and then opened, the drive will start. When the switch RS_W is closed and then opened again, the drive will stop.
 - In the Footswitch Control Mode: When the switch RS_W is closed and then opened, the pump will dispense one dose and then stop.
 - In Logic Level Control Mode: Close the switch RS_W to start the drive. Open the switch to stop the drive.

7 Maintenance

7.1 Warranty

The product comes with one-year labor and parts warranty. The limited warranty covers no damage that is caused by improper usage and handling.

7.2 Regular Maintenance

- 1) To avoid leakage, always check the tubing and connections.
- 2) Do not cover the fan on the rear of the pump.
- 3) Do not use water to wash the pump. Keep the pump head dry.
- 4) Do not use chemical solvents to clean the pump and pump head.

7.3 Malfunction Solutions

| No. | Malfunction | Description | Solution |
|-----|-------------|---------------------|--|
| 1 | Hardware | No display | <ol style="list-style-type: none"> 1. Check the power cord 2. Check the fuse. If it was blown, replace it with a 1A slow-blow fuse 3. Check the internal power cord connection inside the pump. |
| 2 | Hardware | Motor does not work | <ol style="list-style-type: none"> 1. Check the indicator of the driver board. 2. Check the wire connection between the motor and the driver board. 3. Check the wire connection between the driver and the main board. 4. Check the power voltage for the pump. |
| 3 | Hardware | Motor is trembling | <ol style="list-style-type: none"> 1. Check the wire connection between the motor and the driver board. |

WT600S-65 Variable-Speed Peristaltic Pump

| | | | |
|----|----------|----------------------------------|--|
| | | | 2. The motor is overloaded. Check the mechanical connection. |
| 4 | Hardware | Motor only runs in one direction | Check the connection between the drive board and the main control board. |
| 5 | Hardware | Keypad does not work | 1. Check the wire connection between the keypad and the main board. 2. Check if the key is broken. |
| 6 | Hardware | External control does not work | 1. Check the wiring of the connector. 2. Check if the external control power voltage is provided. 3. Check the connections of the external control board. |
| 7 | Hardware | RS485 com does not work | 1. Check the wiring of the connector. 2. Check if the external control power voltage is provided. 3. Check the connections of the communication board. |
| 8 | Hardware | Noisy when running | Check the screws and level on the pump head to make sure they are secure. |
| 9 | Software | External control does not work | Check if the pump is in the External Control Mode. |
| 10 | Software | RS485 does not work right | 1. Check if the display shows the communication is ready. 2. Reset the address of the pump. 3. Check whether on the bus there are two pumps using the same address |

If the problem cannot be solved, please contact the manufacturer or distributor.

8 Dimensions


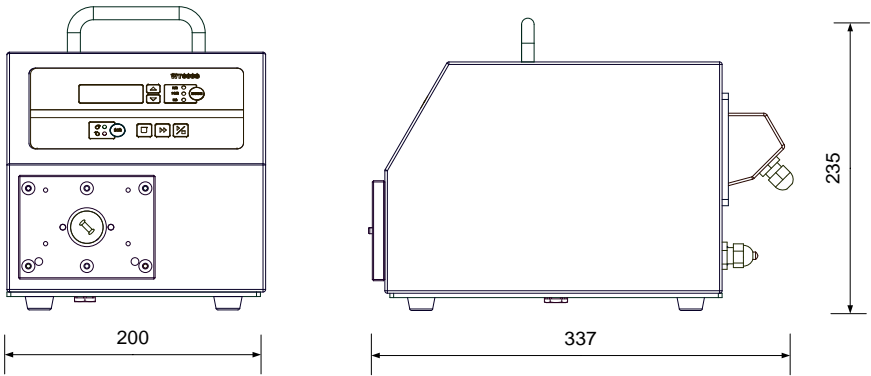
Figure 21. Dimensions (mm)

9 Naming Rule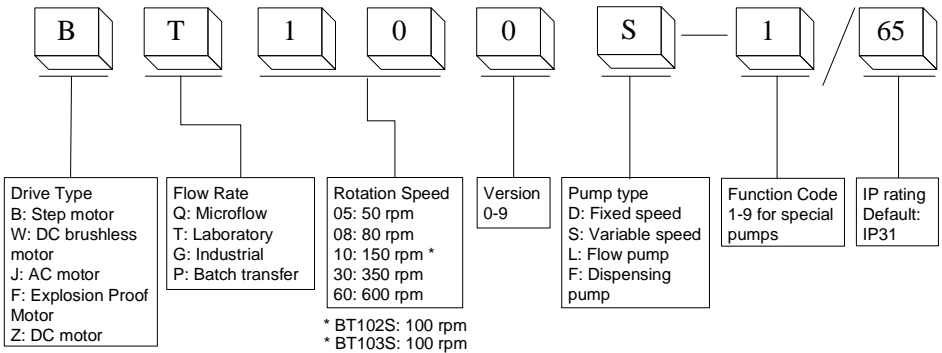

10 Specifications

| | |
|-------------------------------------|--|
| Speed resolution | 1 rpm |
| Speed accuracy | 3% |
| Power supply | AC100-240V 50Hz/60Hz |
| Power consumption | < 200W |
| External logic level control signal | 5V, 12V (standard), 24V (optional) |
| External analog control signal | 0-5V (standard); 0-10V, 4-20mA (optional) |
| Communication interface | RS485 MODBUS |
| Operating condition | Temperature 0-40°C, Relative humidity <80% |
| IP grade | IP31 |
| Display | Four-digit LED |
| Dimensions (LxWxH) | 337x200x235mm (13.27x7.87 x 9.25 inch) |
| Weight | 7.1 kg (15.7 lbs) |

WT600S-65 Applicable Pump Heads and Tubing, Flow Parameters

| Pump head | No. of Channels | Tubing size | Flow rate per channel(mL/min) |
|------------------|------------------------|--------------------------------|--------------------------------------|
| YZ15 | 1 | 13# 14# 16# 19# 25# 17# | 3~2300 |
| YZ25 | 1 | 15# 24# | 50~1700 |
| 2xYZ15 | 2 | 13# 14# 16# 19# 25# 17# | 3~1700 |
| 2xYZ25 | 2 | 15# 24# | 50~1700 |
| KZ25 | 1 | 15# 24# 35# 36# | 80~6000 |
| DT15-24 | 2 | 19# 16# 25# 17# | 14~1600 |
| YT15 | 1 | 13# 14# 16# 19# 25# 17# 18# | 3~2300 |
| YT25 | 1 | 15# 24# 35# 36# | 50~2900 |
| 2xYT15 | 2 | 13# 14# 16# 19# 25# 17# 18# | 3~2300 |
| 2xYT25 | 2 | 15# 24# 35# 36# | 50~2900 |
| KZ 25-L | 1 | 15# 24# 35# 36# | 93~6600 |

Golander LLC

4405 International Blvd
Ste B117, Norcross, GA 30093
USA

Tel: +1 678-587-8806
info@golanderpump.com
www.golanderpump.com

Golander GmbH

Dechant-Heimbach-Str. 29
53177 Bonn
Germany

Tel: +49 228 50446952
info@golander.de
www.golander.de