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(gem. Verordnung (EU) 2016/425, Anhang II, 1.4, CE CAT.III

## Instructions for use (English)

acc. to Regulation (EU) 2016/425, Appendix II, 1.4, CE CAT.III

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## INSTRUCTIONS FOR USE:

EN ENGLISH

Please read carefully before using! You are required to enclose this information leaflet when passing on the personal protective equipment, or to present it personally to the recipient. You may therefore reproduce this leaflet at your own discretion. Marking: Each apron is identified by an inside label indicates the type of protection and some further information for use. (1) ProSafe-2 is the model name for a protective microporous film apron with tie fastening. (2) Manufacturer of the garment. (3) CE-marking - The apron corresponds to the European standards for personal protective equipment of category III.: **Centro Tessile Cottoniero e Abbigliamento S.p.A. (Centrocot), Piazza Sant'Anna 2, 21052 Busto Arsizio VA notified body n. 0624** (Regulation (EU) 2016/425 for Personal Protective Equipment - module C2 and module D). The Notified Body chosen for Conformity to type assessment is: Centrocot NB 0624. (4) The European standards for chemical protective clothing define six types of protection, which are symbolized in one pictogram:

Typ 1 - gas tight clothing  
Typ 2 - non gas tight clothing  
Typ 3 - Liquid tight clothing

Typ 4 - Spray tight clothing  
Typ 5 - Particle tight clothing  
Typ 6 - Limited splash tight clothing

DS products specification correspond to the protection types of the European standards. ProSafe-2 aprons offer protection of type PB and 6B. The size table combines the body measurements with the standard sizes.

(5) The Symbol of the „Open Book“ remembers the wearer to familiarize himself with the „Instructions for Use“. (6) ProSafe-2 aprons are antistatically treated and offer electrostatic protection according to EN1149-1. The person wearing the electrostatic dissipative protective clothing shall be properly earthed. The resistance between the person skin and earth shall be less than  $10^8 \Omega$  e.g. by wearing adequate footwear on dissipative or conductive floors; Electrostatic dissipative protective clothing shall not be open or removed whilst in presence of flammable or explosive atmospheres or while handling flammable or explosive substances; electrostatic dissipative protective clothing is intended to be worn in Zones 1, 2, 20, 21 and 22 (see EN 60079-10-1 [7] and EN 60079-10-2 [8]) in which the minimum ignition energy of any explosive atmosphere is not less than 0,016 mJ; electrostatic dissipative protective clothing shall not be used in oxygen enriched atmospheres, or in Zone 0 (see EN 60079-10-1 [7]) without prior approval of the responsible safety engineer the electrostatic dissipative performance of the electrostatic dissipative protective clothing can be affected by wear and tear, laundering and possible contamination electrostatic dissipative protective clothing shall be worn in such a way that it permanently covers all non-complying materials during normal use (including bending movements). (7) ProSafe-2 aprons are tested against infective agents according to EN 14126:2003. (8) The five care-pictograms indicate:

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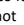
### Performance profile of ProSafe-2 aprons:

Physical Properties	Unit	Result	Test Method	Class	
Abrasion resistance	Cycles	> 1500	EN 530 - method 2	5/6	
Tensile strength	N	> 30	EN ISO 13934-1	1/6	
Flex cracking resistance	Cycles	> 1 000	EN 7854	1/6	
Puncture resistance	N	> 10	EN 863 - EN 13034	2/6	
Trapezoidal tear resistance	N	> 10	EN ISO 9073-4	1/6	
Seams strength	N	> 75	EN ISO 13935-2	3/6	
Resistance to liquid penetration		not applicable	EN ISO 17491-4 met. A EN 13034	not applicable	
Resistance to penetration to liquid	P	R		P	R
Sulphuric acid (H <sub>2</sub> SO <sub>4</sub> )		< 1%	> 95%	3/3	3/3
Sodium hydroxide (NaOH)		< 1%	> 95%	3/3	3/3
o-xylene		< 1%	> 95%	3/3	3/3
Butan-1-ol		< 1%	> 95%	3/3	3/3

P = Penetration index (%penetrated); R = Repellency index (%repelled); Remark: For further information concerning the barrier performances please contact **DS**. EEC-declaration of conformity available under [ProSafe-2 Apron - Lab Coats and Aprons \(dssafetywear.de\)](http://ProSafe-2 Apron - Lab Coats and Aprons (dssafetywear.de))

**PREPARATIONS BEFORE USE:** Do not use incorrect aprons. In case of defect, please contact **DS**. The Correct size combined with correct dressing assures the protective performance of the apron. **LIMITATIONS OF USE:** Please ensure that you have chosen the right garment suitable for your job. For advice, please contact your supplier of the garment **DS**. The user shall be the sole judge for correct combination of full body protective apron and ancillary equipment (gloves, boots, respiratory protective equipment, etc.) and for how long ProSafe-2 aprons can be worn with respect to its protective performance, wear comfort and heat stress. **DS** shall not accept any responsibility whatsoever for improper use of ProSafe-2 aprons. **STORAGE:** The PPE has a storability of 5 years from date of production (9) (see ) unless no damage is visible. ProSafe-2 aprons can be stored according to customers storage practices, store cool and dry, no direct sunlight. **DISPOSAL:** ProSafe-2 aprons without harm to the environment. Restrictions to disposal result only from contaminants introduced during use. (10) LOT NO

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