



LUGOL'S IODINE

VIG-008 – (4 ML CONCENTRATE & 8 OZ. DILUENT)

VIG-016- (8 ML CONCENTRATE & 16 OZ. DILUENT)

VIG-128- (64 ML CONCENTRATE & 1 GAL DILUENT)

Order online at www.volusol.com

WORKING SOLUTION: Add iodine concentrate to diluent just prior to use:

For 1.5% use 1 mL concentrate to 60mL of diluent.

For 3% use 2mL concentrate to 60mL of diluent.

For 6% use 4mL concentrate to 60mL of diluent.

Working solution is stable when tightly covered for 30 days or until visible lightening of solution occurs. Depleted solutions should be replaced not replenished.

Note: Keep working solution out of direct light to prevent lightening of solution and subsequent depletion.