***TriScroll™ 600 Series  
Dry Scroll  
Vacuum Pump***

*INSTALLATION AND  
OPERATION MANUAL*

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*MANUALE  
D'INSTALLAZIONE E D'USO*

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# TriScroll™ 600 Series Dry Scroll Vacuum Pump

*TriScroll is a trademark of Vacuum Technologies.*

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# TriScroll 600 Series Vacuum Pump

---

|   |           |
|---|-----------|
| <b>Technical Information .....</b>                                  | <b>33</b> |
| Electrical Connections.....   | 36        |
| Single Phase Motor Connection.....                                  | 36        |
| Three Phase Motor Connection .....                                  | 37        |
| Mechanical Connections.....   | 38        |
| Pump Location .....   | 38        |
| Rotation of the Pump Inlet<br>with Respect to the Motor Frame ..... | 38        |
| Pump Inlet.....   | 38        |
| Pump Exhaust.....   | 38        |
| Optional Isolation Valve.....                                       | 38        |
| Gas Ballast .....   | 39        |
| Bearing Purge.....  | 39        |
| Purge Kit .....   | 39        |
| Troubleshooting .....   | 40        |
| Maintenance .....   | 41        |
| General Information .....   | 41        |
| Related TriScroll Manuals.....                                      | 41        |
| Maintenance and Tooling Kits .....                                  | 41        |
| Factory Service Options.....  | 42        |
| Accessories.....  | 42        |
| Contacting Vacuum Technologies .....                                | 42        |

## **Request for Return Health and Safety Certification**

# TriScroll 600 Series Vacuum Pump

## List of Figures

| Figure | Caption   | Page |
|--------|---|------|
| 1      | Interface Drawing with Dimensions . . . . .                             | 34   |
| 2      | Pumping Speed Curves . . . . .  | 34   |
| 3      | TriScroll 600 Series Vacuum Pump<br>(Single Phase Pump Shown) . . . . . | 35   |
| 4      | Single Phase Motor Electrical Connections .                             | 36   |
| 5      | Three Phase Motor Electrical Connections .                              | 37   |
| 6      | Isolation Valve Location. . . . .                                       | 38   |

## List of Tables

| Table | Title  | Page |
|-------|--|------|
| 1     | Pump Models and Power Cords.....               | 3    |
| 2     | Circuit Breaker Requirements.....              | 3    |
| 3     | Pumpenmodelle und Versorgungskabel .....       | 7    |
| 4     | Anforderungen an Stromunterbrecher .....       | 7    |
| 5     | Modèles de pompes et câbles .....              | 11   |
| 6     | Exigences des coupe-circuits .....             | 11   |
| 7     | Modelos de bomba y cables de corriente.....    | 15   |
| 8     | Requisitos del disyuntor.....                  | 15   |
| 9     | Modelli di pompe e cavi di alimentazione... .. | 19   |
| 10    | Requisiti dell'interruttore automatico .....   | 19   |
| 11    | ポンプ型式と電源コード .....                              | 23   |
| 12    | 遮断器の要求 .....                                   | 23   |
| 13    | 泵型号和电源电缆 .....                                 | 27   |
| 14    | 断路器要求 .....                                    | 27   |
| 15    | 펌프모델 및 코드 .....                                | 31   |
| 16    | 브레이커 요구 .....                                  | 31   |
| 17    | Specifications .....                           | 33   |
| 18    | Full Load Motor Currents, Amperes.....         | 36   |
| 19    | VPI Valve Installation Data .....              | 39   |
| 20    | Troubleshooting Chart .....                    | 40   |
| 21    | Other Related Manuals .....                    | 41   |
| 22    | Maintenance and Tooling Kits .....             | 41   |
| 23    | Factory Service Options .....                  | 42   |
| 24    | Accessories .....                              | 42   |

*This page intentionally left blank.*

**Declaration of Conformity**  
**Konformitätserklärung**  
**Déclaration de Conformité**  
**Declaración de Conformidad**  
**Verklaring de Overeenstemming**  
**Dichiarazione di Conformità**

We  
Wir  
Nous  
Nosotros  
Wij  
Noi

Varian, Inc.  
Vacuum Technologies  
121 Hartwell Avenue  
Lexington, MA, 02421-3133 USA

declare under our sole responsibility that the product,  
erklären, in alleniniger Verantwortung, daß dieses Produkt,  
déclarons sous notre seule responsabilité que le produit,  
declaramos, bajo nuestra sola responsabilidad, que el producto,  
verklaren onder onze verantwoordelijkheid, dat het product,  
dichiariamo sotto nostra unica responsabilità, che il prodotto,

TriScroll Series Vacuum Pump

to which this declaration relates is in conformity with the following standard(s) or other normative documents.  
auf das sich diese Erklärung bezieht, mit der/den folgenden Norm(en) oder Richtlinie(n) übereinstimmt.  
auquel se réfère cette déclaration est conforme à la (aux) norme(s) ou au(x) document(s) normatif(s).  
al que se refiere esta declaración es conforme a la(s) norma(s) u otro(s) documento(s) normativo(s).  
waarnaar deze verklaring verwijst, aan de volgende norm(en) of richtlijn(en) beantwoordt.  
a cui se riferisce questa dichiarazione è conforme alla/e seguente/i norma/o documento/i normativo/i.

**98/37/EEC, Machinery Directive**

EN 1012-2:1996 Compressors and Vacuum pumps Safety Reqmts; Part 2 Vacuum Pumps  
EN 1050:1996 Safety of machinery - principles for risk assessment  
EN 60204-1 Electrical equipment of industrial machines; general requirements

**73/023/EEC, Low Voltage Directive**

EN 60034 part 1 Rotating electrical machines - Part 1: Rating and performance

**89/336/EEC, Electromagnetic Compatibility Directive**

EN 61000-4-2 Testing and Measurement Techniques - Electrostatic Discharge Immunity Test

A handwritten signature in cursive script that reads "Frederick C. Campbell".

Frederick C. Campbell  
Operations Manager  
Vacuum Technologies  
Varian, Inc.  
Lexington, Massachusetts, USA

March 2003

## Warranty

Products manufactured by Seller are warranted against defects in materials and workmanship for twelve (12) months from date of shipment thereof to Customer, and Seller's liability under valid warranty claims is limited, at the option of Seller, to repair, to replace, or refund of an equitable portion of the purchase price of the Product. Items expendable in normal use are not covered by this warranty. All warranty replacement or repair of parts shall be limited to equipment malfunctions which, in the sole opinion of Seller, are due or traceable to defects in original materials or workmanship. All obligations of Seller under this warranty shall cease in the event of abuse, accident, alteration, misuse, or neglect of the equipment. In-warranty repaired or replaced parts are warranted only for the remaining unexpired portion of the original warranty period applicable to the repaired or replaced parts. After expiration of the applicable warranty period, Customer shall be charged at the then current prices for parts, labor, and transportation.

Reasonable care must be used to avoid hazards. Seller expressly disclaims responsibility for loss or damage caused by use of its Products other than in accordance with proper operating procedures.

Except as stated herein, Seller makes no warranty, express or implied (either in fact or by operation of law), statutory or otherwise; and, except as stated herein, Seller shall have no liability under any warranty, express or implied (either in fact or by operation of law), statutory or otherwise. Statements made by any person, including representatives of Seller, which are inconsistent or in conflict with the terms of this warranty shall not be binding upon Seller unless reduced to writing and approved by an officer of Seller.

## Warranty Replacement and Adjustment

All claims under warranty must be made promptly after occurrence of circumstances giving rise thereto, and must be received within the applicable warranty period by Seller or its authorized representative. Such claims should include the Product serial number, the date of shipment, and a full description of the circumstances giving rise to the claim. Before any Products are returned for repair and/or adjustment, written authorization from Seller or its authorized representative for the return and instructions as to how and where these Products should be returned must be obtained. Any Product returned to Seller for examination shall be prepaid via the means of transportation indicated as acceptable by Seller. Seller reserves the right to reject any warranty claim not promptly reported and any warranty claim on any item that has been altered or has been returned by non-acceptable means of transportation. When any Product is returned for examination and inspection, or for any other reason, Customer shall be responsible for all damage resulting from improper packing or handling, and for loss in transit, notwithstanding any defect or non-conformity in the Product. In all cases, Seller has the sole responsibility for determining the cause and nature of failure, and Seller's determination with regard thereto shall be final.

If it is found that Seller's Product has been returned without cause and is still serviceable, Customer will be notified and the Product returned at its expense; in addition, a charge for testing and examination may be made on Products so returned.

3/1/00

# Instructions for Use

## General Information

This equipment is designed for use by professionals. The user should read this instruction manual and any other additional information supplied by Vacuum Technologies before operating the equipment. Vacuum Technologies will not be held responsible for any events that occur due to non-compliance with these instructions, improper use by untrained persons, non-authorized interference with the equipment, or any action contrary to that provided for by specific national standards.

The TriScroll™ 600 is a dry scroll vacuum pump. This pump is suitable for pumping air or inert gases. The pump is not intended to pump toxic, corrosive, explosive, or particulate-forming gases.

The following paragraphs contain all the information necessary to guarantee the safety of the operator when using the equipment. Detailed information is supplied in “Technical Information” on page 33.

This manual uses the following standard safety protocol:

### WARNING

*The warning messages are for attracting the attention of the operator to a particular procedure or practice which, if not followed correctly, could lead to serious injury.*

### CAUTION

*The caution messages are displayed before procedures, which if not followed, could cause damage to the equipment.*

### NOTE

*The notes contain important information taken from the text.*

## Storage

When transporting and storing the pump, the following environmental requirements should not be exceeded:

Temperature:            -20 °C to 60 °C (-4 °F to 140 °F)  
Relative humidity:    0 to 95% (non-condensing)

## Unpacking and Inspection

The shipping container is a double carton.

1. After opening the outer box, remove the foam packing.
2. Slit open the inner box.
3. Lift the pump with the plywood base out of the inner box.
4. Remove the four bolts securing the pump frame to the plywood base.
5. Locate the NW25 exhaust fitting and set it aside.
6. Inspect the pump for damage.

If there is shipping damage, contact the freight carrier and your local Vacuum Technologies sales office immediately.

7. Save the carton and packing materials.

Total weight of the package, including the pump, is approximately 40 kg (87 lbs).

### WARNING

*When unpacking the pump, be sure not to drop it, and avoid any kind of sudden impact or shock vibration to it.*

### WARNING

*The TriScroll 600 weighs 32 kg (70 lbs). To avoid injury, use proper lifting techniques when moving the pump.*

### NOTE

*Normal exposure to the environment cannot damage the pump. Nevertheless, it is advisable to keep the pump inlet closed until the pump is installed in the system.*

## Installation Requirements

### Safety

Do not remove or modify any safety or insulating equipment from the pump. To do so may create a serious safety hazard and may void the warranty.

#### WARNING

- ❑ *This pump is designed to pump air and inert gases only; it is not designed to pump explosive, flammable, toxic, or corrosive gases. They can cause bodily injury, explosion, or fire.*
- ❑ *Install in an area that is not exposed to rain, steam, or excessive humidity. They can cause electric shock, short circuits, and severe bodily injury.*
- ❑ *Before inspecting or servicing the pump, be sure the electrical supply is disconnected.*
- ❑ *Consult a qualified electrician whenever wiring the pump.*

#### CAUTION

*Although the pump can pump trace particulates normally found in the atmosphere, it is not designed for process solids, chemicals, powders, solvents, condensates, or other particulates. They can damage the equipment, degrade its performance, or shorten its useful life.*

*Single and three phase pumps operate in a clockwise direction when viewed from the motor end. (Note the arrow on the pump frame.) Improper rotation can cause permanent damage to the pump.*

During operation, the following environmental conditions should not be exceeded:

Temperature: +5 °C to +40 °C (41 °F to 104 °F)

Relative humidity: 0 to 95% (non-condensing)

#### CAUTION

*Do not block the fan ducts because the pump can become overheated. A pump surface temperature in excess of 55 °C (131 °F) is potentially damaging. If such conditions are observed, turn pump off and allow to cool. Disassemble, inspect for damage, and repair if necessary.*

#### CAUTION

*As supplied from the factory, the pump is configured as shown in the following table. Verify that the configuration matches the supply voltage.*

*If voltage changeover is required, configure the voltage as described in "Electrical Connections" on page 36 and page 37.*

## TriScroll 600 Series Vacuum Pump

Table 1 lists the available TS-600 pump models and indicates the power cord to be used with each particular pump.

Three phase pumps are shipped without a power cord. A three phase power cable kit, Vacuum Technologies part number S4789001, is available. This kit does not include a plug.

**Table 1 Pump Models and Power Cords**

| Pump Model     | Motor Voltage | Plug Type  | Fusing            | Conductor Size         | Cord Marking | Cord Length |
|----------------|---------------|------------|-------------------|------------------------|--------------|-------------|
| PTS06001UNIV   | 100-115 VAC   | NEMA 5-15P | -                 | 14 Gage                | UL/CSA       | 2.5 m       |
| PTS06001UNIVEU | 200-230 VAC   | CEE 7/7    | -                 | 3x1.50 mm <sup>2</sup> | <HAR>        | 2.5 m       |
| PTS06001UNIVUK | 200-230 VAC   | BS 1363    | BS 1362<br>13 Amp | 3x1.50 mm <sup>2</sup> | <HAR>        | 2.5 m       |
| PTS06003UNIV   | 200-230 VAC   | -          | -                 | -                      | -            | -           |
| PTS06101UNIV   | 100-115 VAC   | NEMA 5-15P | -                 | 14 Gage                | UL/CSA       | 2.5m        |
| PTS06001UVPI   | 100-115 VAC   | NEMA 5-15P | -                 | 14 Gage                | UL/CSA       | 2.5m        |

### Circuit Breakers

**WARNING**

*Protect against short circuits by installing a circuit breaker of the proper capacity.*

Table 2 lists the circuit breaker capacity required for pump motors of various voltages, single and three-phase.

**Table 2 Circuit Breaker Requirements**

| Single Phase Motor | 100-115 VAC<br>50/60 Hz | 200-230 VAC<br>50/60 Hz |
|--------------------|-------------------------|-------------------------|
| Breaker capacity   | 20 Amperes              | 10 Amperes              |
| Three Phase Motor  | 200-230 VAC             | 380-415-460 VAC         |
| Breaker capacity   | 10 Amperes              | 5 Amperes               |

## Operation

In order to reach ultimate vacuum, the pump must be left running for about one hour with the inlet sealed.

Unlike conventional oil-sealed pumps, Vacuum Technologies's dry scroll pumps do not have fluid to cleanse them of accumulated dust and debris. Run the pump periodically at atmosphere for a minute or two to flush out the pump. Flush the pump regularly and adjust this schedule according to your specific conditions.

## Startup Procedure

1. Be sure that the vacuum system isolation valve is closed.
2. Turn on power to the pump.
3. Open the isolation valve.

## Shutdown Procedure

1. Close the vacuum system isolation valve. This prevents debris in pump from being transported into the vacuum system.
2. Turn off power to the pump.

## Maintenance

Personnel responsible for pump operation and maintenance must be well-trained and aware of the accident prevention rules.

### WARNING

- ❑ *Death may result from contact with high voltages. Always take extreme care and observe the accident prevention regulations in force.*
- ❑ *When the machine is powered up, be careful of moving parts and high voltages.*
- ❑ *If you have to perform maintenance on the pump after a considerable time in operation, allow the pump to cool as the temperature of the outer surface may be in excess of 55 °C (131 °F).*
- ❑ *Always disconnect your power supply to the pump before beginning maintenance work.*

### NOTE

*Before returning the pump to the factory for repair, the "Health and Safety" sheet attached to this instruction manual must be completed and sent to the local sales office. A copy of the sheet must be inserted in the pump package before shipping.*

If a pump is to be discarded, it must be disposed of in accordance with specific national and local standards.

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
# Technical Information


**Table 17 Specifications**

|  |  |
|--|--|
| Model                                      | TriScroll™ 600 Series Dry Scroll Vacuum Pump   |
| Interface dimensions                       | See Figure 1   |
| Peak pumping speed                         | 50 Hz: 420 l/m, 25.2 m <sup>3</sup> /hr (14.8 cfm)<br>60 Hz: 500 l/m, 30 m <sup>3</sup> /hr (17.7 cfm)   |
| Media                                      | Clean air. No toxic, corrosive, explosive or particulate forming gases   |
| Ultimate pressure (Torr)                   | 7.0 x 10 <sup>-3</sup> Torr (9.3 x 10 <sup>-3</sup> mbar)  |
| Maximum inlet pressure                     | 1.0 atmosphere (0 psig)  |
| Maximum outlet pressure                    | 1.1 atmosphere (1.5 psig)  |
| Inlet connection                           | NW40   |
| Exhaust connection                         | Female 3/8" National Pipe Thread (NW25 adapter provided)   |
| Gas ballast                                | Female 1/4" National Pipe Thread (40 Micron sintered filter provided)  |
| Ambient operating temperature              | 5 °C to 40 °C (41 °F to 104 °F)  |
| Storage temperature                        | -20 °C to 60 °C (-4 °F to 140 °F)  |
| Motor rating                               | 1.0 HP (0.76 kW)   |
| Operating voltages                         | Single phase models:<br><input type="checkbox"/> 50-60 Hz/100-115:200-230 VAC<br>Three phase models:<br><input type="checkbox"/> 50 Hz/200-230:380-415 VAC<br><input type="checkbox"/> 60 Hz/200-230:460 VAC |
| Motor full load currents                   | See Table 18 on page 36  |
| Motor thermal protection                   | Type U automatic   |
| Operating speed                            | 60 Hz: 1725 RPM, 50 Hz: 1425 RPM   |
| Cooling system                             | Air-cooled   |
| Weight                                     | Pump only: 32 kg (70 lbs)<br>Shipping weight: 40 kg (87 lbs)   |
| Noise level (per ISO 11201)                | 68 dB(A)   |
| Vibration level at inlet (per ISO 10816-1) | 6.3 mm/sec   |

# TriScroll 600 Series Vacuum Pump

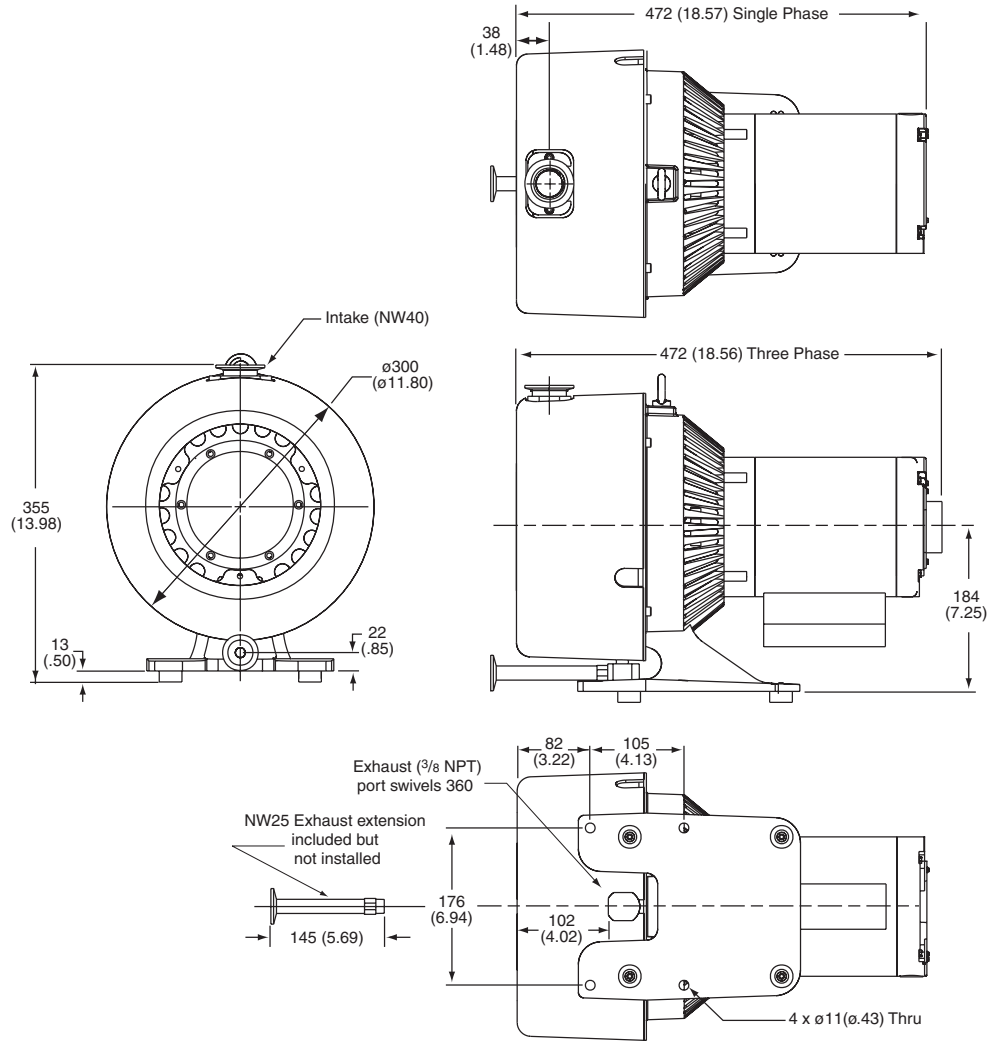
**Figure 1 Interface Drawing with Dimensions**

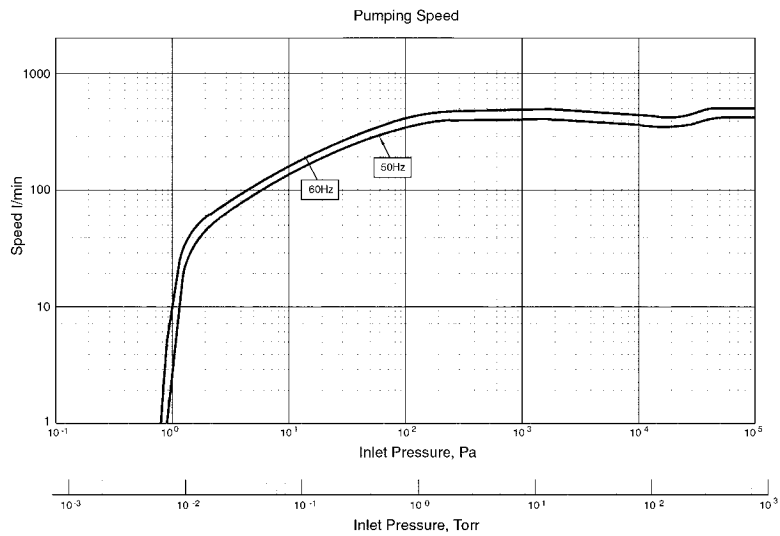**Figure 2 Pumping Speed Curves**

## TriScroll 600 Series Vacuum Pump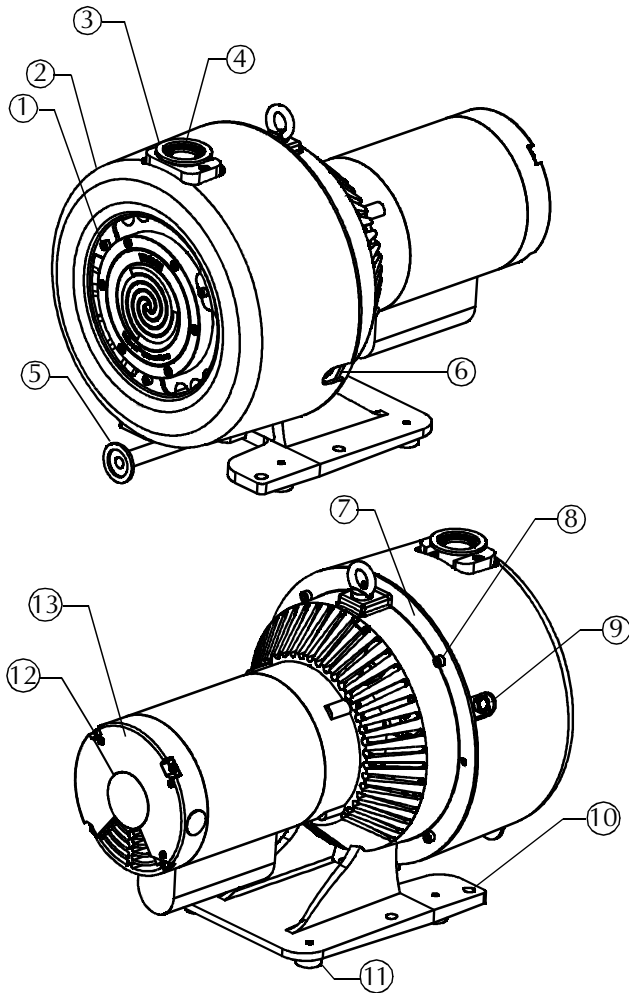

1. Cowling Screws; M5 (3)
2. Cowling
3. Inlet (NW40)
4. Inlet Screen
5. NW25 Exhaust Adapter
6. Bearing Purge Port (1/4" National Pipe Thread)
7. Pump Frame
8. Frame Screws; M6 (4)
9. Gas Ballast Port (1/4" National Pipe Thread)
10. Mounting Holes; 11 mm diameter thru (8)
11. Rubber Feet (4)
12. Motor Cover Screws (3)
13. Motor Electrical Cover

**Figure 3 TriScroll 600 Series Vacuum Pump  
(Single Phase Pump Shown)**

## Electrical Connections


Wire the motor and electrical interlocks (if applicable) in accordance with local electrical codes and the relevant electrical component manufacturer's instructions. Table 18 lists the full load motor currents at various voltages.

**Table 18 Full Load Motor Currents, Amperes**

| 1 phase motor | 100 V | 115 V | 200 V | 230 V |       |  |
|---------------|-------|-------|-------|-------|-------|--|
| 50 Hz         | 14    | 13.6  | 7.0   | 6.8   |       |  |
| 60 Hz         | 15.2  | 12.6  | 7.6   | 6.3   |       |  |
| 3 phase motor | 200 V | 230 V | 380 V | 415 V | 460 V |  |
| 50 Hz         | 4.4   | 4.4   | 2.2   | 2.2   | -     |  |
| 60 Hz         | 4.4   | 4.4   | -     | -     | 2.2   |  |

### Single Phase Motor Connection

The pump can be configured for low voltage, 100 VAC to 115 VAC, or for high voltage, 200 VAC to 230 VAC. Low voltage exists between connection points "4" and "Line 2" whether the pump motor is configured for low or high voltage. Figure 4 shows the electrical connections for a single phase motor.

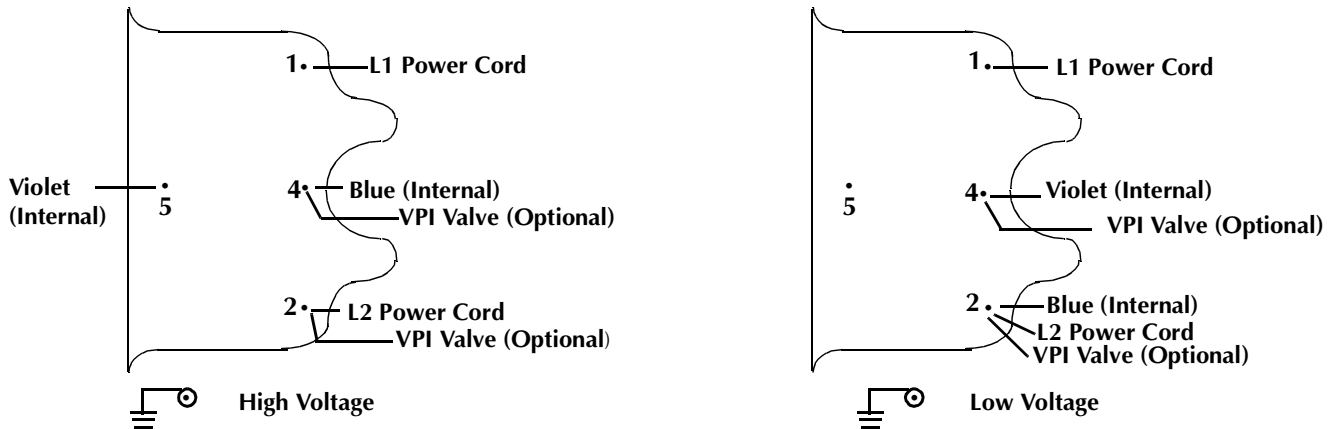**Figure 4 Single Phase Motor Electrical Connections**

1. Verify the electrical supply voltage.
2. Remove the three screws (item 12 in Figure 3 on page 35) that are holding the motor electrical cover (item 13 in Figure 3).
3. Refer to Figure 4 to connect the motor to match your supply voltage.

Two options are available to strain relieve the electrical supply cable.


- The cable can be mechanically held under the electrical cover in the groove provided. The groove is sized for 14 gage cable.
- A 1/2–14 NPSM hole is also provided next to the motor cover.

4. Wire L1 and L2 per Figure 4 using right angle flag connectors or ring connectors. Ensure that no exposed wiring is close to the electrical cover or to other terminals on the board.
5. Secure the ground wire under the ground screw using a ring connector.
6. Replace the motor electrical cover and secure it with the three (3) screws removed in step 2.

# TriScroll 600 Series Vacuum Pump

## Three Phase Motor Connection

The pump can be configured for low voltage, 200 VAC to 230 VAC, or for high voltage, 380 VAC to 415 VAC and 460 VAC. As supplied from the factory, the pump is configured for low voltage. Figure 5 shows the electrical connections for a three phase motor.

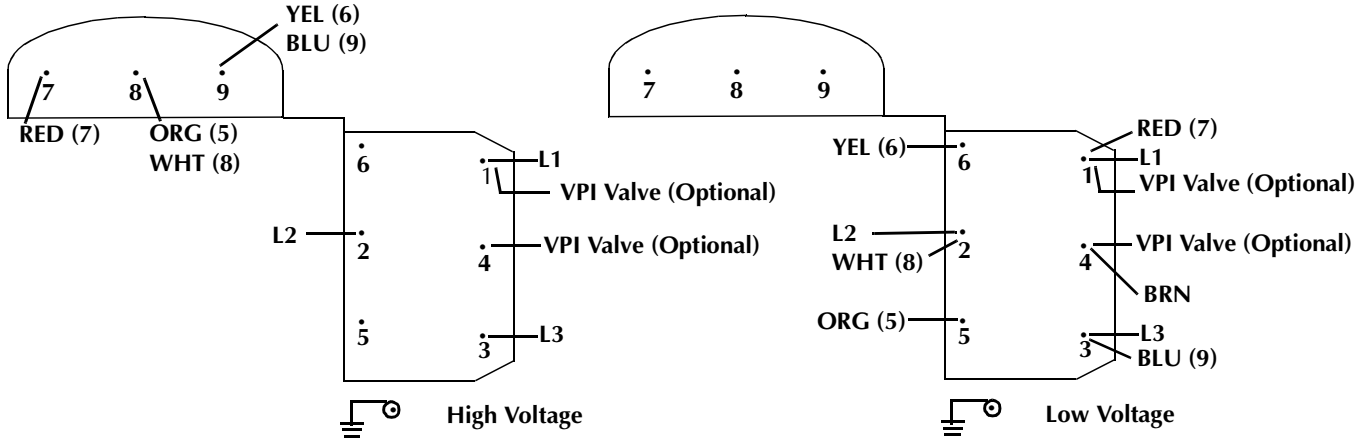**Figure 5 Three Phase Motor Electrical Connections**

1. Verify the electrical supply voltage.
2. Remove the three screws (item 12 in Figure 3 on page 35) that are holding the motor electrical cover (item 13 in Figure 3).
3. Refer to Figure 5 to connect the motor to match your supply voltage.

Two options are available to strain relieve the electrical supply cable.

- ❑ The cable can be mechanically held under the electrical cover in the groove provided. The groove is sized for 14 gage cable.
- ❑ A 1/2–14 NPSM hole is also provided next to the motor cover.

4. L1, L2, and L3 are the power cord leads. Wire L1, L2 and L3 per Figure 5 using right angle flag connectors or ring connectors. Ensure that no exposed wiring is close to the electrical cover or to other terminals on the board.
5. Secure the ground wire under the ground screw using a ring connector.
6. Install the motor electrical cover and secure it with the three (3) screws removed in step 2.
7. Momentarily switch on the pump to verify proper rotation. With the inlet uncovered, a strong air flow should come out of the pump exhaust. If there is no flow out the exhaust, reverse any two of the leads (L1, L2 or L3).
8. Repeat the test to verify correct rotation.

## Mechanical Connections

### Pump Location

Locate the pump on a firm, level surface. Mounting holes provided in the frame can be used to provide permanent attachment. The rubber mounts on the frame can be removed if desired.

### Rotation of the Pump Inlet with Respect to the Motor Frame

The standard pump configuration is with the inlet fitting positioned at the top of the pump as shown in Figure 1 on page 34. The pump inlet has two alternate positions:

- ❑ 90 degrees clockwise
- ❑ 90 degrees counterclockwise

To rotate the inlet:

1. Remove three (3) M5 screws (item 1 on Figure 3 on page 35) that secure the cowling (item 2 on Figure 3) to the scroll module. Remove the cowling.
2. Tilt the pump back so that the rear end of the motor touches the floor.
3. Remove the four (4) M6 screws (item 8 on Figure 3) holding the frame (item 7 on Figure 3) and scroll module together.
4. Axially separate the frame and scroll module.
5. Rotate the module 90 degrees in either direction and realign the two (2) locating pins on the scroll module with the mating frame holes. Ensure that the rubber spider is still on the motor coupling and that the coupling teeth are properly aligned.
6. Rejoin the scroll module with the frame and install and tighten the four (4) M6 screws removed in step 3.
7. Reinstall the cowling using the three (3) M5 screws removed in step 1.

### Pump Inlet

Use NW40, or larger, clean vacuum hardware with as short a length as practical between the pump and the vacuum chamber.

Use a bellows to provide both vibration isolation and strain relief between the pump and the vacuum chamber.

### Pump Exhaust

A female 1/4" National Pipe Thread exhaust fitting is located underneath the scroll module. This fitting swivels 360 degrees. Additionally, an NW25 male adapter with 3/8" National Pipe Thread is provided.


To avoid overheating the pump, do not restrict the exhaust flow with long lengths of small diameter tubing. Use as short as practical lengths of NW25 diameter, or larger, hardware.

## Optional Isolation Valve

Scroll pumps return to atmospheric pressure quickly when shut off, thus the installation of a fast acting, automatic, normally closed isolation valve is strongly recommended to prevent pump debris from being transported back into the vacuum chamber when the pump is turned off.

- ❑ The opening of this valve should occur simultaneously with or after pump startup.
- ❑ Valve closing should occur before, simultaneously with, but no later than 250 ms after pump shut off.

Use an NW40 valve or larger and mount it as close as possible to the pump inlet. Mounting to the pump inlet is ideal. (See Figure 6).

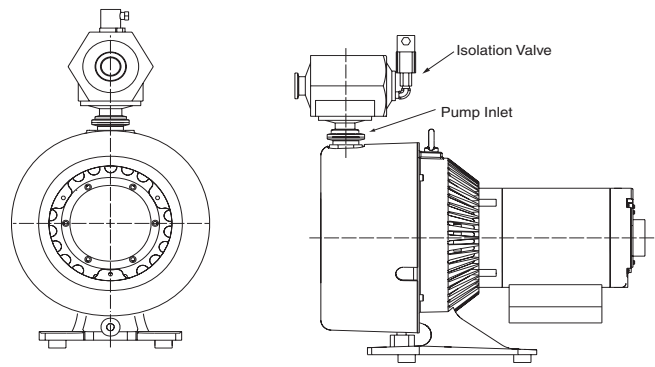**Figure 6 Isolation Valve Location**

Vacuum Technologies offers a variety of manual, electromagnetic and electropneumatic controlled vacuum valves for vacuum applications. The Vacuum Technologies Vacuum Pump Isolation (VPI) Valve is highly recommended for vacuum pump isolation applications. The VPI valve application and installation information for use with TriScroll vacuum pumps has been included below.

Vacuum Technologies also sells the TriScroll 600 pump with the VPI valve preinstalled. The part number is PTS06001UVPI.

### Overview

The status of the VPI Valve is controlled by the TriScroll motor internal thermal switch and its electrical supply. The VPI Valve opens when the pump is running and closes when it is stopped. It will also close when the TriScroll motor thermal overload protection switch shuts down the pump.

### CAUTION

*To prevent damage to the VPI Valve, it must be installed by a qualified electrician and only as specified below.*

## TriScroll 600 Series Vacuum Pump

### Wiring

Using the data in Table 19, verify that the VPI Valve chosen is compatible with the TriScroll supply voltage and frequency. Then, locate the proper wiring diagram to use (Figure 4 on page 36 or Figure 5 on page 37).

1. Remove the three screws (item 12 in Figure 3 on page 35) that are holding the motor electrical cover (item 13 in Figure 3).
2. Connect the VPI Valve solenoid wire leads to the pump motor as shown in Figure 4 or Figure 5.
3. Replace the motor electrical cover and secure it using the three screws removed in step 1. Verify that the valve is properly grounded before applying electrical power.

**Table 19 VPI Valve Installation Data**

| Motor   | TriScroll Vacuum Pump Operating Voltage | VPI Valve Part # NW25 | VPI Valve Part # NW40 | Solenoid Operating Voltage Range | Use the following Wiring Diagram |
|---------|---|-----------------------|-----------------------|----------------------------------|----------------------------------|
| 1 Phase | 100-120 V 50/60 Hz                      | VPI251205060          | VPI401205060          | 90-132 V                         | Figure 4 low voltage             |
| 1 Phase | 200-230 V 50/60 Hz                      | VPI251205060          | VPI401205060          | 90-132 V*                        | Figure 4 high voltage            |
| 3 Phase | 200 V 50/60 Hz                          | VPI251205060          | VPI401205060          | 90-132 V*                        | Figure 5 low voltage             |
| 3 Phase | 230 V 50/60 Hz                          | VPI251335060          | VPI401335060          | 120-146 V*                       | Figure 5 low voltage             |
| 3 Phase | 380 V 50/60 Hz                          | VPI252205060          | VPI402205060          | 180-264 V*                       | Figure 5 high voltage            |
| 3 Phase | 415 V 50/60 Hz                          | VPI252205060          | VPI402205060          | 180-264 V*                       | Figure 5 high voltage            |
| 3 Phase | 460 V 50/60 Hz                          | VPI252665060          | VPI402665060          | 239-293 V*                       | Figure 5 high voltage            |

\* Solenoid operating voltages are lower than the TriScroll vacuum pump operating voltages in order to utilize the TriScroll motor's internal thermal switch to actuate the VPI Valve.

### Gas Ballast

The pump incorporates an automatic gas ballast to prevent water and other condensates from accumulating within the pump. The standard configuration is a sintered filter installed in the 1/4" National Pipe Thread gas ballast port (item 9 on Figure 3 on page 35). This allows enough atmospheric air to enter the pump in order to purge condensates while not affecting pump ultimate pressure or pumping speed.

For applications where the ingress of atmospheric air is undesirable, dry nitrogen at a flow rate of  $\approx 5$  lpm can be bled into the gas ballast port. See "Purge Kit" below. Blocking of the gas ballast port is not recommended.

### Bearing Purge

A 1/4" National Pipe Thread bearing purge port (item 6 on Figure 3 on page 35) protects the main crankshaft bearings in applications where large amounts of water are being pumped. In the standard configuration, this port is sealed.

To enable the bearing purge, dry nitrogen at a flow rate of  $\approx 5$  lpm can be bled into the bearing purge port. This gas supply should be maintained at 2 psig or less and must be kept below 5 psig. See "Purge Kit" below.

#### CAUTION

*The bearing purge port has been disabled on the TriScroll 600 model S4800309. Use of this port could cause internal damage to the pump.*

### Purge Kit

A purge kit (Vacuum Technologies part number PTSPURGEKIT) is available to properly purge either the bearing purge or the gas ballast. This kit contains a flowmeter and all necessary valving and tubing.

## Troubleshooting

Table 20 contains a list of possible problems, their probable causes, and corrective actions.

**Table 20 Troubleshooting Chart**

| <b>Problem</b>             | <b>Probable Cause</b>                             | <b>Corrective Action</b>  |
|----------------------------|---|---|
| Pump won't start           | Circuit breaker open                              | Close breaker.<br>Identify cause of overload.   |
|                            | Motor thermal protector open                      | Allow motor to cool.<br>Identify cause of overload.   |
|                            | Electrical short under the motor electrical cover | Inspect and repair.   |
|                            | Wiring loose or cut                               | Repair or replace.  |
|                            | Excessive voltage drop                            | Check size and length of power supply cable.  |
|                            | Defective motor                                   | Inspect. Contact Vacuum Technologies.   |
| Poor ultimate pressure     | System leak                                       | Locate and repair leak.   |
|                            | Water in pump                                     | Flush pump with air or dry nitrogen.  |
|                            | Gas ballast plugged                               | Replace breather vent. Contact Vacuum Technologies.   |
|                            | Solvent in pump                                   | Flush pump with air or dry nitrogen.<br>Install trap or filter.   |
|                            | Seals worn out                                    | Replace tip seals. (Table 21 and Table 22 on page 41 list maintenance kits and service options.)  |
|                            | Poor conductance to pump                          | Replumb with shorter and/or larger diameter tubing.   |
| Pump makes hammering noise | Pump overheated                                   | Check ambient temperature.<br>Check ventilation to pump.  |
|                            | Debris in pump                                    | Check inlet screen.<br>Flush pump.<br>Disassemble pump and inspect. (Table 21 and Table 22 on page 41 list maintenance kits and service options.) |

## Maintenance

### General Information

Vacuum Technologies TriScroll 600 series pumps are designed to provide years of trouble-free service if maintenance procedures and intervals are observed. Bearing grease replenishment and tip seal replacement is recommended when pump base pressure has risen to an unacceptably high level for your application. Bearings, rotary seals and o-rings should also be replaced if the pump exhibits humming or grinding noises from the bearings. Main bearing life may be shortened if your application requires the pumping of high quantities of water vapor. Use of the bearing purge kit (PTSPURGEKIT) mentioned earlier will keep this water from impacting bearing life.

Maintenance should be performed in accordance with procedures, tooling and materials specified in the manuals listed below.

### Related TriScroll Manuals

Other manuals related to tip seal replacement, pump module replacement, and major maintenance of the TriScroll 600 series pumps are listed in Table 21.

**Table 21 Other Related Manuals**

| Title                          | Applicable TriScroll Model      | Part Number |
|--------------------------------|---------------------------------|-------------|
| Tip Seal Replacement Manual    | All TriScroll 600 Series models | 699904310   |
| Pump Module Replacement Manual | All TriScroll 600 Series models | 699904305   |
| Major Maintenance Manual       | All TriScroll 600 Series models | 699904300   |

### Maintenance and Tooling Kits

Material and tooling required to perform maintenance on TriScroll pumps is provided in kit form. A description of each kit and ordering information is provided in Table 22.

**Table 22 Maintenance and Tooling Kits**

| Description              | Contents   | Applicable TriScroll Model      | Part Number |
|--------------------------|--|---------------------------------|-------------|
| Major Maintenance Kit    | All bearings, bearing seals, bearing lubricant, O-rings, and tip seals required to rebuild TriScroll 600 series pumps.   | All TriScroll 600 Series models | PTSS0600MK  |
| Maintenance Tool Kit     | All fixtures and tools required to perform any maintenance on TriScroll 600 Series pumps.  | All TriScroll 600 Series models | PTSS0600TK  |
| Replacement Tip Seal Set | Replacement tip seals and static O-rings for TriScroll 600 Series pumps.<br><br><i>NOTE:</i> The Maintenance Tool Kit is also required for tip seal replacement. | All TriScroll 600 Series models | PTSS0600TS  |

## TriScroll 600 Series Vacuum Pump

### Factory Service Options

Table 23 lists the factory-rebuild service and advance exchange of complete TriScroll 600 series pumps, as well as factory service options that Vacuum Technologies offers.

**Table 23 Factory Service Options**

| Factory Service Options                                    | Part Number      |
|--|------------------|
| Advance Exchange TriScroll 600 Single Phase (US cord)      | EXPPTS06001      |
| Advance Exchange TriScroll 600 Three Phase                 | EXPPTS06003      |
| Advance Exchange TriScroll 600 Single Phase (EU cord)      | EXPTS06001UNIVEU |
| Advance Exchange TriScroll 600 Single Phase (UK cord)      | EXPTS06001UNIVUK |
| Advance Exchange TriScroll 610 Single Phase                | EXPPTS06101      |
| Advance Exchange TriScroll 610 Three Phase                 | EXPPTS06103      |
| Advance Exchange TriScroll 600 Pump Module Only            | EXPTS0600SC      |
| Advance Exchange TriScroll 610 Pump Module Only            | EXPTS0610SC      |
| Service/Rebuild TriScroll 600 Pump (Single or Three Phase) | PTS0600KMA       |
| Service/Rebuild TriScroll 610 Pump (Single or Three Phase) | PTS0610KMA       |
| Service/Rebuild TriScroll 600 Pump Module Only             | PTS0600SCRP      |
| Service/Rebuild TriScroll 610 Pump Module Only             | PTS0610SCRP      |

### Accessories

The accessories listed in Table 24 are available for use with the TriScroll 600 series pump. Contact your local Vacuum Technologies office to place an order. A list of offices is included on the rear cover of this manual.

**Table 24 Accessories**

|                       |             |
|-----------------------|-------------|
| Purge Kit             | PTSPURGEKIT |
| Three Phase Cable Kit | S4789001    |
| Exhaust Extension     | S4807001    |
| Exhaust Filter Kit    | PTS600EXFIL |

### Contacting Vacuum Technologies

In the United States, you can contact Vacuum Technologies Customer Service at 1-800-8VARIAN. See the back cover of this manual for a listing of our sales and service offices.

Internet users:

- Send email to Customer Service & Technical Support at [vpl.customer.support@varianinc.com](mailto:vpl.customer.support@varianinc.com)
- Visit our web site at [www.varianinc.com/vacuum](http://www.varianinc.com/vacuum)
- Order on line at [www.evarian.com](http://www.evarian.com)

*Request for Return  
Health and Safety Certification*

- Return authorization numbers (RA#) **will not** be issued for any product until this Certificate is completed and returned to a Varian, Inc. Customer Service Representative.
- Pack goods appropriately and drain all oil from rotary vane and diffusion pumps (for exchanges please use the packing material from the replacement unit), making sure shipment documentation and package label clearly shows assigned Return Authorization Number (RA#) VVT cannot accept any return without such reference.
- Return product(s) to the nearest location:

**North and South America**  
 Varian, Inc.  
 Vacuum Technologies  
 121 Hartwell Ave.  
 Lexington, MA 02421  
 Fax: (781) 860-9252

**Europe and Middle East**  
 Varian S.p.A.  
 Via F.lli Varian, 54  
 10040 Leini (TO) – ITALY  
 Fax: (39) 011 997 9350

**Asia and ROW**  
 Varian Vacuum Technologies  
 Local Office

For a complete list of phone/fax numbers see [www.varianinc.com/vacuum](http://www.varianinc.com/vacuum)

- If a product is received at Varian, Inc. in a contaminated condition, **the customer is held responsible** for all costs incurred to ensure the safe handling of the product, and **is liable** for any harm or injury to Varian, Inc. employees occurring as a result of exposure to toxic or hazardous materials present in the product.

|                             |                           |                   |   |
|-----------------------------|---------------------------|-------------------|---|
| <i>CUSTOMER INFORMATION</i> |                           |                   |   |
| Company name: .....         |                           |                   |   |
| Contact person:             | Name: .....               | Tel:.....         | .....   |
|                             | Fax: .....                | E-mail: .....     | .....   |
| Ship method:                | Shipping Collect #: ..... | P.O.#: .....      | .....   |
| Europe only:                | VAT Reg Number: .....     | USA only:         | <input type="checkbox"/> Taxable <input type="checkbox"/> Non-taxable |
| Customer ship to:           | .....                     | Customer bill to: | .....   |
|                             | .....                     |                   | .....   |
|                             | .....                     |                   | .....   |

*PRODUCT IDENTIFICATION*

| Product Description | Varian, Inc. Part Number | Varian, Inc. Serial Number |
|---------------------|--------------------------|----------------------------|
|                     |                          |                            |
|                     |                          |                            |
|                     |                          |                            |

*TYPE OF RETURN (check appropriate box)*

|  |   |  |  |  |
|--|---|--|--|--|
| <input type="checkbox"/> Paid Exchange | <input type="checkbox"/> Paid Repair    | <input type="checkbox"/> Warranty Exchange | <input type="checkbox"/> Warranty Repair | <input type="checkbox"/> Loaner Return |
| <input type="checkbox"/> Credit        | <input type="checkbox"/> Shipping Error | <input type="checkbox"/> Evaluation Return | <input type="checkbox"/> Calibration     | <input type="checkbox"/> Other         |

*HEALTH and SAFETY CERTIFICATION*

|  |                 |            |
|--|-----------------|------------|
| <b>VACUUM TECHNOLOGIES CANNOT ACCEPT ANY BIOLOGICAL HAZARDS, RADIOACTIVE MATERIAL, ORGANIC METALS, OR MERCURY AT ITS FACILITY. CHECK ONE OF THE FOLLOWING:</b>   |                 |            |
| <input type="checkbox"/> I confirm that the above product(s) has (have) <b>NOT</b> pumped or been exposed to any toxic or dangerous materials in a quantity harmful for human contact.                             |                 |            |
| <input type="checkbox"/> I declare that the above product(s) has (have) pumped or been exposed to the following toxic or dangerous materials in a quantity harmful for human contact ( <u>Must be filled in</u> ): |                 |            |
| Print Name .....   | Signature ..... | Date ..... |

**PLEASE FILL IN THE FAILURE REPORT SECTION ON THE NEXT PAGE**

Do not write below this line

Notification (RA) #: ..... Customer ID #: ..... Equipment #: .....

**FAILURE REPORT**

(Please describe in detail the nature of the malfunction to assist us in performing failure analysis):

**TURBO PUMPS AND TURBOCONTROLLERS**

| Claimed Defect   | Position  | Parameters   |
|--|---|--|
| <input type="checkbox"/> Does not start<br><input type="checkbox"/> Does not spin freely<br><input type="checkbox"/> Does not reach full speed<br><input type="checkbox"/> Mechanical Contact<br><input type="checkbox"/> Cooling defective<br><input type="checkbox"/> Noise<br><input type="checkbox"/> Vibrations<br><input type="checkbox"/> Leak<br><input type="checkbox"/> Overtemperature<br><input type="checkbox"/> Clogging | <input type="checkbox"/> Vertical<br><input type="checkbox"/> Horizontal<br><input type="checkbox"/> Upside-down<br><input type="checkbox"/> Other<br>..... | Power:                      Rotational Speed:<br>Current:                    Inlet Pressure:<br>Temp 1:                      Foreline Pressure:<br>Temp 2:                      Purge flow:<br>Operation Time: |
| Describe Failure:  |   |  |
| Turbocontroller Error Message:   |   |  |

**ION PUMPS/CONTROLLERS**

|   |
|---|
| <input type="checkbox"/> Bad feedthrough<br><input type="checkbox"/> Vacuum leak<br><input type="checkbox"/> Error code on display<br><input type="checkbox"/> Poor vacuum<br><input type="checkbox"/> High voltage problem<br><input type="checkbox"/> Other ..... |
| Describe failure:   |
| Customer application:   |

**VALVES/COMPONENTS**

|   |
|---|
| <input type="checkbox"/> Main seal leak<br><input type="checkbox"/> Solenoid failure<br><input type="checkbox"/> Damaged sealing area<br><input type="checkbox"/> Bellows leak<br><input type="checkbox"/> Damaged flange<br><input type="checkbox"/> Other ..... |
| Describe failure:   |
| Customer application:   |

**LEAK DETECTORS**

|   |
|---|
| <input type="checkbox"/> Cannot calibrate<br><input type="checkbox"/> Vacuum system unstable<br><input type="checkbox"/> Failed to start<br><input type="checkbox"/> No zero/high background<br><input type="checkbox"/> Cannot reach test mode<br><input type="checkbox"/> Other ..... |
| Describe failure:   |
| Customer application:   |

**INSTRUMENTS**

|   |
|---|
| <input type="checkbox"/> Gauge tube not working<br><input type="checkbox"/> Communication failure<br><input type="checkbox"/> Error code on display<br><input type="checkbox"/> Display problem<br><input type="checkbox"/> Degas not working<br><input type="checkbox"/> Other ..... |
| Describe failure:   |
| Customer application:   |

**ALL OTHER VARIAN, INC.**

|  |
|--|
| <input type="checkbox"/> Pump doesn't start<br><input type="checkbox"/> Doesn't reach vacuum<br><input type="checkbox"/> Pump seized<br><input type="checkbox"/> Noisy pump (describe)<br><input type="checkbox"/> Overtemperature<br><input type="checkbox"/> Other ..... |
| Describe failure:  |
| Customer application:  |

**DIFFUSION PUMPS**

|   |
|---|
| <input type="checkbox"/> Heater failure<br><input type="checkbox"/> Doesn't reach vacuum<br><input type="checkbox"/> Vacuum leak<br><input type="checkbox"/> Electrical problem<br><input type="checkbox"/> Cooling coil damage<br><input type="checkbox"/> Other ..... |
| Describe failure:   |
| Customer application:   |

## Sales and Service Offices

### Canada

#### Central coordination through:

Varian, Inc.  
121 Hartwell Avenue  
Lexington, MA 02421  
USA  
Tel: (781) 861 7200  
Fax: (781) 860 5437  
Toll Free: (800) 882 7426

### China

#### Varian Technologies - Beijing

Room 1201, Jinyu Mansion  
No. 129A, Xuanwumen Xidajie  
Xicheng District  
Beijing 1000031  
P.R. China  
Tel: (86) 10 6608 1031  
Fax: (86) 10 6608 1541

### France and Benelux

#### Varian s.a.

7 avenue des Tropiques  
Z.A. de Courtaboeuf – B.P. 12  
Les Ulis cedex (Orsay) 91941  
France  
Tel: (33) 1 69 86 38 13  
Fax: (33) 1 69 28 23 08

### Germany and Austria

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Alsfelder Strasse 6  
Postfach 11 14 35  
64289 Darmstadt  
Germany  
Tel: (49) 6151 703 353  
Fax: (49) 6151 703 302

### India

#### Varian India PVT LTD

101-108, 1st Floor  
1010 Competent House  
7, Nangal Raya Business Centre  
New Delhi 110 046  
India  
Tel: (91) 11 5548444  
Fax: (91) 11 5548445

### Italy

#### Varian, Inc.

Via F.lli Varian, 54  
10040 Leini, (Torino)  
Italy  
Tel (39) 011 997 9 111  
Fax (39) 011 997 9 350

### Japan

#### Varian, Inc.

Sumitomo Shibaura Building, 8th Floor  
4-16-36 Shibaura  
Minato-ku, Tokyo 108  
Japan  
Tel: (81) 3 5232 1253  
Fax: (81) 3 5232 1263

### Korea

#### Varian Technologies Korea, Ltd.

Shinsa 2nd Building 2F  
966-5 Daechi-dong  
Kangnam-gu, Seoul  
Korea 135-280  
Tel: (82) 2 3452 2452  
Fax: (82) 2 3452 2451

### Mexico

#### Varian S.A.

Concepcion Beistegui No 109  
Col Del Valle  
C.P. 03100  
Mexico, D.F.  
Tel: (52) 5 523 9465  
Fax: (52) 5 523 9472

### Russia

#### Central coordination through:

Varian, Inc.  
via F.lli Varian 54  
10040 Leini, (Torino)  
Italy  
Tel: (39) 011 997 9 252  
Fax: (39) 011 997 9 316

### Taiwan

#### Varian Technologies Asia Ltd.

18F-13 No.79, Hsin Tai Wu Road  
Sec. 1, Hsi Chih, Taipei Hsien  
Taiwan, R.O.C.  
Tel: (886) 2 2698 9555  
Fax: (886) 2 2698 9678

### UK and Ireland

#### Varian Ltd.

28 Manor Road  
Walton-On-Thames  
Surrey KT 12 2QF  
England  
Tel: (44) 1932 89 8000  
Fax: (44) 1932 22 8769

### United States

#### Varian, Inc.

121 Hartwell Avenue  
Lexington, MA 02421  
USA  
Tel: (781) 861 7200  
Fax: (781) 860 5437

### Other Countries

#### Varian, Inc.

Via F.lli Varian 54  
10040 Leini, (Torino)  
Italy  
Tel: (39) 011 997 9 111  
Fax: (39) 011 997 9 350

### Customer Support and Service:

#### North America

Tel: 1 (800) 882-7426 (toll-free)  
vtl.technical.support@varianinc.com

#### Europe

Tel: 00 (800) 234 234 00 (toll-free)  
vtl.technical.support@varianinc.com

#### Japan

Tel: (81) 3 5232 1253 (dedicated line)  
vtj.technical.support@varianinc.com

#### Korea

Tel (82) 2 3452 2452 (dedicated line)  
vtk.technical.support@varianinc.com

#### Taiwan

Tel: 0 (800) 051 342 (toll-free)  
vtw.technical.support@varianinc.com

### Worldwide Web Site, Catalog and On-line Orders:

[www.varianinc.com](http://www.varianinc.com)

Representatives in most countries

**VARIAN**

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Copyright 2005