


DSW-100 Calibration Manual

1. Power on.
2. Press the CAL switch to start calibration. (*Note: CAL Switch is located underneath of the body.) The display will show "SP on."

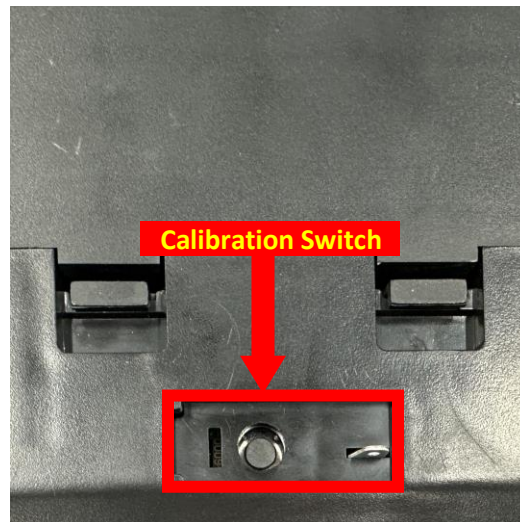

3. While holding the **ZÉRO** button, hit **UNIT** **TARE** **TARE** to enter the calibration mode. The display will show the default calibration weight.
4. Press the **ZÉRO** button to enter zero calibration. (*Note: The platform should be empty.)



5. Press the **ZÉRO** button, and then the display will show "CAL SP."



6. Place the calibration weight and press the **ZÉRO** button.
7. It will return to the main interface.