














8. ANIMAL WEIGHING

MODEL: TBW / TBWS

TBW can use for vibrate loads weigh. This function can use for animal weighing. For settings, see the parameter $P \exists \text{ o}tH \gg \text{An}\bar{n}$

- 1 Press  key during the self checking.
- 2 Display show PN, in succession press ,  and  to enter.
- 3 Display will be show **p 0 chk.**
- 4 Press  key until display show **p 3 oth.**
- 5 Press  key to confirm.
- 6 Press  key to select ON/OFF.
- 7 Press  key to confirm.
- 8 Press  key twice to back to normal weighing mode and scale will start self-checking.
- 9 Press  &  key together to enter animal weighing mode ,the HOLD indicator will be on
- 10 Bring the load on the platform, display will be show the reading
- 11 It can add or remove loads also update the weighing values.
- 12 To complete the current weighing, remove the weight from the platform.
- 13 To exit animal mode ,Press  &  key together again
- 14 To disable the animal weighing function, set the parameter $P \exists \text{ o}tH \gg \text{An}\bar{n} \gg \text{oFF}$

www.VisionTechShop.com

100 TEMPLE AVE, HACKENSACK, NJ 07601