



# International Filter Products

*Revolutionary systems for your industry*

## MKP Series PES Dual Layer Filter Capsules

MKP pharma grade capsules were designed for small-scale filtration, clarification, and filling applications in the pharmaceutical and biotechnology industries worldwide. The PES filter capsules have been specially designed for sterilizing in biopharmaceutical applications. The filters come with pre layer configuration, which increases the filtration volume significantly. The hydrophilic PES membrane does not require pre-wetting agents, thereby eliminating a potential source of contamination. The PES membrane is a low protein binding media. This family of products is particularly suitable for scale-up testing where a full-size capsule or cartridge would be excessive. They are also very suitable for disposable Bio-bag applications. The all-polypropylene construction provides excellent compatibility with a wide range of chemicals. The capsules are thermally sealed, and 100% integrity tested.

### Application

- Buffers and Media
- SVP (Small Volume Parenteral)
- Vaccines
- Scale up processing
- Product Sterilization
- Bio Bags
- Biologics
- Ultrapure Water
- Fermentation Broths
- Pharmaceuticals
- Antibiotics
- Serums

**Part# MKPDS065S020D2H2H-ETO: MKP Filter Capsule with 1/4" Hose Barb Inlet/outlet Connection**





## Technical Data Sheet

### Popular Configuration:

Catalog Number	Micron Rating	Inlet Connection	Outlet Connection	Vent Connection	Sterilization
MKPDS065S020D2H2H-ETO	0.65 µm / 0.2 µm (Dual Layer)	1/4" Hose Barb	1/4" Hose Barb	Luer Loc Cap	ETO Sterilized

❖ Find the complete list of available standard inlet and outlet fittings in page 5.

### Micron Rating:

**Final Membrane:** 0.2 µm PES [Sterilizing Grade Membrane]

**Pre-Filter:** Highly Asymmetric 0.65 µm PES

### Effective Filtration Area:

Mini Capsule; 77.5 in<sup>2</sup> (500 cm<sup>2</sup>)

### Materials of Construction:

Membrane: Pharmaceutical Grade PES (Polyethersulfone) Sterilizing Grade

Membrane Feature: Hydrophilic, Low Protein Binding, High Flow Rate

Media Supports: Polypropylene

Shell, Cage, Core, End Caps: Polypropylene

Valve O-Rings: Silicon (Standard)

Sealing: Thermally Bonded

### Operating Conditions:

Maximum Forward Differential Pressure: **Liquid:** 5.5 bar (80psi) at 77°F/25°C

**Gas:** 4.1 bar (60psi) at 77°F/25°C

Minimum Burst Pressure: 8.3 bar (120psi) at 77°F/25°C

Maximum Reverse Differential Pressure: 3.0 bar (44psi) at 68°F/20°C

Maximum Operating Temperature: 176°F/80°C




# International Filter Products

*Revolutionary systems for your industry*

**Nominal Dimension:**

**Part Number: MKPDS065S020D2H2H-ETO**

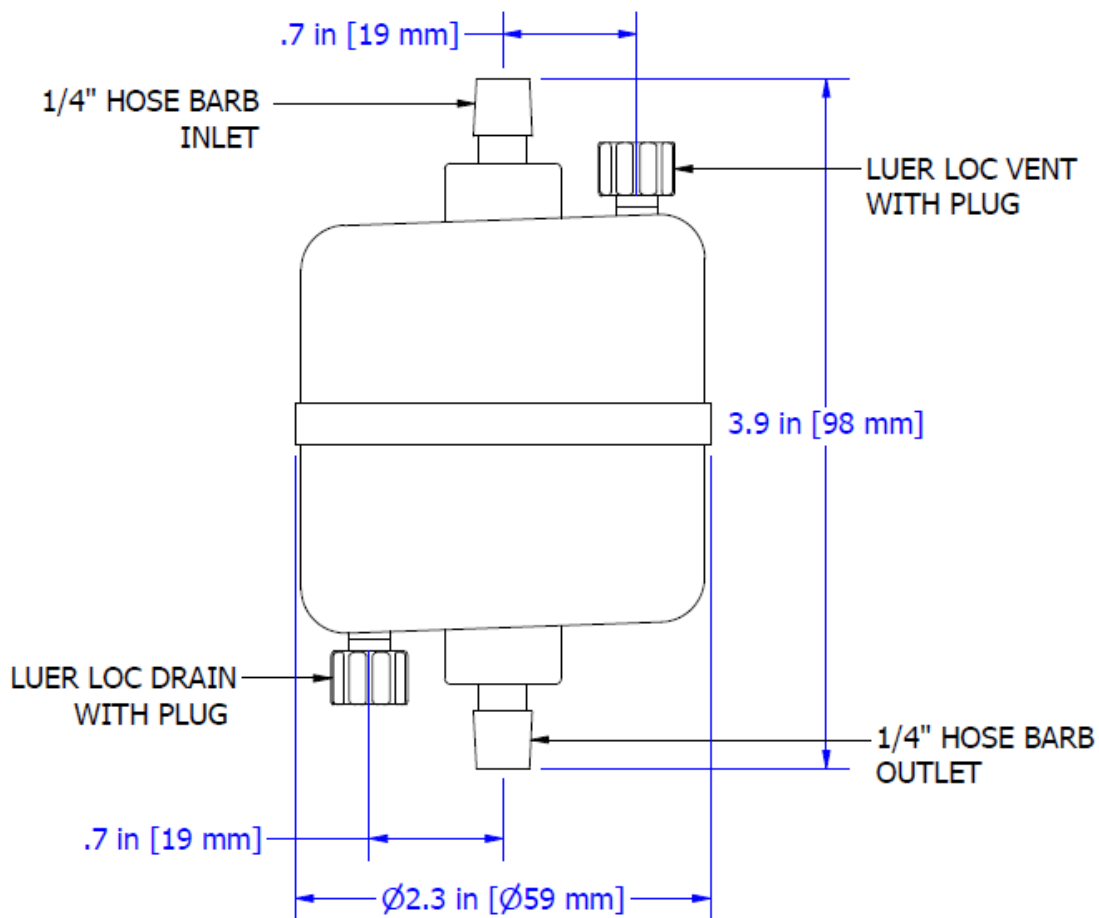

**Hold-Up Volume (Approx.):**

Upstream Volume: 32 ml

Downstream Volume: 11 ml

**Typical Filtered Volume:**

10.0L – 28.0L

**Filter Integrity:**

The finished product was sampled and shown to exhibit a minimum bubble point of  $\geq 18$  psi (1.2 bar) in 60%IPA / 40% water and  $\geq 50$  psi (3.5 bar) in DI Water ( at 22°C ).



# International Filter Products

*Revolutionary systems for your industry*

## **Sterilization:**

Capsules can be autoclaved 25 times at 125 °C (257°F) for 30 minutes or chemically sanitized in situ using common sanitizing agents or hot water at 90 °C (194°F) for a limited time (dependent on time and temperature).

*Warning:* The filters cannot be sterilized by steam-in-place (SIP)

## **Regulatory Compliance 21CFR Part 177 & USP <88>:**

The filters are constructed with polypropylene resins and filtration media in compliance with 21CFR Part 177 of the US Code of Federal Regulations and USP Class VI Biological USP <88> Test for Plastic.

## **USP <85> Bacterial Endotoxins:**

The filters were tested to confirm that an aqueous extraction of this product contains <0.25 EU/ml as determined by the Limulus Amebocyte Lysate (LAL) Test

## **USP <645> Conductivity USP:**

Effluent is tested during the manufacturing process and shown to meet the requirements for USP Sterile Water for Injection for conductivity.

## **Bacterial Retention ASTM F838-05:**

The filters are quantitatively retentive towards *Brevundimonas diminuta* (ATCC #19146) at a minimum challenge level of  $10^7$  CFU per  $cm^2$  of filtration area, consistent with ASTM F838-05

## **USP <87> / ISO 10993-5 Cytotoxicity:**

Extract from this product is non-Cytotoxic

## **Hemolysis ASTM F756-17:**

Extract from this product is non-Hemolytic

## **Human and Veterinarian Use:**

The product is safe for Human and Veterinarian use. CGMP CFR part 210 & 211, additional requirement of 21 CFR part 600 and 21 CFR part 680 are applicable to the aseptic manufacturing process.

## **Animal-Derived Components & TSE/BSE Risk:**

No animal-derived material is intentionally added or used during the manufacture of this product.

## **Shelf Life:**

The MKP capsules have a shelf life of 3 years from the date of sterilization.



# International Filter Products

Revolutionary systems for your industry

## MKP- Series PES Mini Capsules (Pharmaceutical Grade)

### Ordering Guide



[1]	[2]	[3] Inlet and [4] Outlet	[5]
Micron Rating (Pre-Filter/Final Filter)	Filter Size Code [Filtration Area]	Inlet and Outlet Connection	Sterilization
<ul style="list-style-type: none"> <li>S020 : 0.2 <math>\mu\text{m}</math></li> <li>DS065S020 : Highly Asymmetric 0.65/0.2 <math>\mu\text{m}</math></li> <li>DS045S020 : Highly Asymmetric 0.45/0.2 <math>\mu\text{m}</math></li> <li>DS045S010 : Highly Asymmetric 0.45/0.1 <math>\mu\text{m}</math></li> <li>S080020 : 0.8/0.2 <math>\mu\text{m}</math></li> <li>S020010: 0.2/0.1 <math>\mu\text{m}</math></li> </ul> <p>Note: All final membranes are sterilizing grade.</p>	<ul style="list-style-type: none"> <li>B [Filtration Area: 360 <math>\text{cm}^2</math>]</li> <li>C [Filtration Area: 420 <math>\text{cm}^2</math>]</li> <li>D [Filtration Area: 500 <math>\text{cm}^2</math>]</li> </ul>	<ul style="list-style-type: none"> <li>1H : 1/8" Hose Barb</li> <li>1Q : 1/8" Male Quick Coupling with Metal Latch</li> <li>2H : 1/4" Hose Barbs</li> <li>2H-FB : 1/4" Hose barbs with Filling Bell</li> <li>2Q : 1/4" Male Quick Coupling for Metal Latch</li> <li>2N : 1/4" MNPT</li> <li>3H : 3/8" Hose barbs</li> <li>LF : Luer Loc Female</li> <li>LM : Luer Loc Male</li> <li>MT : 1/2" Tri clamps</li> </ul>	<p>-ETO</p> <p>Ethylene Oxide Sterilization</p>

### Example:

**MKPDS065S020D2H2H-ETO**

PureFlo MKP Series Mini Capsule, Highly Asymmetric Dual Layer PES 0.65 $\mu\text{m}$  / 0.2 $\mu\text{m}$  PES membrane (Sterilizing Grade), Filtration Area 500  $\text{cm}^2$ , 1/4" Hose Barb (Inlet/Outlet), ETO Sterilized (Factory Sterilized)

Part Number

Description



# International Filter Products

*Revolutionary systems for your industry*

## Special Configuration:

### Layer Option:

- Single Layer
- Dual Layer

### Micron Rating Option:

- 0.04 Micron
- 0.1 Micron
- 0.2 Micron
- 0.45 Micron
- 0.65 Micron
- 0.8 Micron
- 1.2 Micron

### Inlet Fitting Option:

- 1H : 1/8" Hose Barb
- 1Q : 1/8" Male Quick Coupling with Metal Latch
- 2H : 1/4" Hose Barbs
- 2H-FB : 1/4" Hose barbs with Filling Bell
- 2Q : 1/4" Male Quick Coupling for Metal Latch
- 2N : 1/4" MNPT
- 3H : 3/8" Hose Barbs
- LF : Luer Loc Female
- MT : 1/2" Tri clamps

### Outlet Fitting Option:

- 1H : 1/8" Hose Barb
- 1Q : 1/8" Male Quick Coupling with Metal Latch
- 2H : 1/4" Hose Barbs
- 2H-FB : 1/4" Hose barbs with Filling Bell
- 2Q : 1/4" Male Quick Coupling for Metal Latch
- 2N : 1/4" MNPT
- 3H : 3/8" Hose Barbs
- LM : Luer Loc Male
- MT : 1/2" Tri clamps

