MKP Series PES Single Layer Filter Capsules

MKP pharma grade capsules were designed for small-scale filtration, clarification, and filling applications in the pharmaceutical and biotechnology industries worldwide. The PES filter capsules have been specially designed for sterilizing in biopharmaceutical applications. The filters come with pre layer configuration, which increases the filtration volume significantly. The hydrophilic PES membrane does not require pre-wetting agents, thereby eliminating a potential source of contamination. The PES membrane is a low protein binding media. This family of products is particularly suitable for scale-up testing where a full-size capsule or cartridge would be excessive. They are also very suitable for disposable Bio-bag applications. The all-polypropylene construction provides excellent compatibility with a wide range of chemicals. The capsules are thermally sealed, and 100% integrity tested.

Application

- Buffers and Media
- SVP (Small Volume Parenteral)
- Vaccines
- Scale up processing

- Product Sterilization
- Bio Bags
- Biologics
- Ultrapure Water

- Fermentation Broths
- Pharmaceuticals
- Antibiotics
- Serums

Part# MKPS020D2H2H-ETO: MKP Filter Capsule with 1/4" Hose Barb Inlet/outlet Connection







Technical Data Sheet

Popular Configuration:

Catalog Number	Micron Rating	Inlet Connection	Outlet Connection	Vent Connection	Sterilization
MKPS020D2H2H-ETO	0.2 μm (Single Layer)	1/4" Hose Barb	1/4" Hose Barb	Luer Loc Cap	ETO Sterilized

Find the complete list of available standard inlet and outlet fittings in page 5.

Micron Rating:

Final Membrane: 0.2 µm PES [Sterilizing Grade Membrane]

Effective Filtration Area:

Mini Capsule; 77.5 in² (500 cm²)

Materials of Construction:

Membrane: Pharmaceutical Grade PES (Polyethersulfone) Sterilizing Grade

Membrane Feature: Hydrophilic, Low Protein Binding, High Flow Rate

Media Supports: Polypropylene

Shell, Cage, Core, End Caps: Polypropylene

Valve O-Rings: Silicon (Standard)

Sealing: Thermally Bonded

Operating Conditions:

Maximum Forward Differential Pressure: Liquid: 5.5 bar (80psi) at 77°F/25°C

Gas: 4.1 bar (60psi) at 77°F/25°C

Minimum Burst Pressure: 8.3 bar (120psi) at 77°F/25°C

Maximum Reverse Differential Pressure: 3.0 bar (44psi) at 68°F/20°C


Maximum Operating Temperature: 176°F/80°C





Nominal Dimension:

Part Number: MKPS020D2H2H-ETO

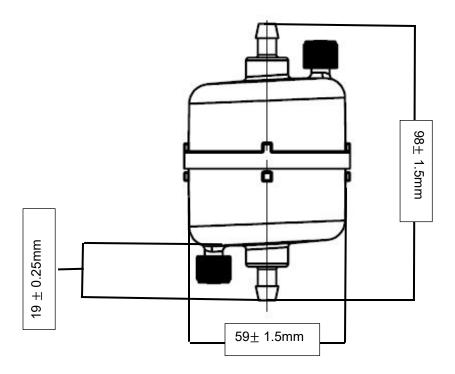

Hold-Up Volume (Approx.):

Upstream Volume: 32 ml Downstream Volume: 11 ml

Typical Filtered Volume:

7.0L - 24.0L

Filter Integrity:

The finished product was sampled and shown to exhibit a minimum bubble point of ≥18 psi (1.2 bar) in 60%IPA / 40% water and ≥ 50 psi (3.5 bar) in DI Water (at 22%).

Sterilization:

Capsules can be autoclaved 25 times at 125 °C (257°F) for 30 minutes or chemically sanitized in situ using common sanitizing agents or hot water at 90 °C (194°F) for a limited time (dependent on time and temperature).

Warning: The filters cannot be sterilized by steam-in-place (SIP)





Regulatory Compliance 21CFR Part 177 & USP <88>:

The filters are constructed with polypropylene resins and filtration media in compliance with 21CFR Part 177 of the US Code of Federal Regulations and USP Class VI Biological USP <88>Test for Plastic.

USP <85> Bacterial Endotoxins:

The filters were tested to confirm that an aqueous extraction of this product contains <0.25 EU/ml as determined by the Limulus Amebocyte Lysate (LAL) Test

USP <645> Conductivity USP:

Effluent is tested during the manufacturing process and shown to meet the requirements for USP Sterile Water for Injection for conductivity.

Bacterial Retention ASTM F838-05:

The filters are quantitatively retentive towards *Brevundimonas diminuta* (ATCC #19146) at a minimum challenge level of 10^7 CFU per cm^2 of filtration area, consistent with ASTM F838-05

USP <87> / ISO 10993-5 Cytotoxicity:

Extract from this product is non-Cytotoxic

Hemolysis ASTM F756-17:

Extract from this product is non-Hemolytic

Human and Veterinarian Use:

The product is safe for Human and Veterinarian use. CGMP CFR part 210 & 211, additional requirement of 21 CFR part 600 and 21 CFR part 680 are applicable to the aseptic manufacturing process.

Animal-Derived Components & TSE/BSE Risk:

No animal-derived material is intentionally added or used during the manufacture of this product.

Shelf Life:

The MKP capsules have a shelf life of 3 years from the date of sterilization.







MKP- Series PES Mini Capsules (Pharmaceutical Grade)

Ordering Guide

M K P [1] [2] [3] [4] [5]						
[1]	[2]	[3] Inlet and [4] Outlet	[5]			
Micron Rating (Pre-Filter/Final Filter]	Filter Size Code [Filtration Area]	Inlet and Outlet Connection	Sterilization			
 S020: 0.2 μm DS065S020: Highly Asymmetric 0.65/0.2 μm DS045S020: Highly Asymmetric 0.45/0.2 μm DS045S010: Highly Asymmetric 0.45/0.1 μm S080020: 0.8/0.2 μm S020010: 0.2/0.1 μm Note: All final membranes are sterilizing grade. 	• B [Filtration Area: 360 cm²] • C [Filtration Area: 420 cm²] • D [Filtration Area: 500 cm²]	 1H: 1/8" Hose Barb 1Q: 1/8" Male Quick Coupling with Metal Latch 2H: 1/4" Hose Barbs 2H-FB: 1/4" Hose barbs with Filling Bell 2Q: 1/4" Male Quick Coupling for Metal Latch 2N: 1/4" MNPT 3H: 3/8" Hose barbs LF: Luer Loc Female LM: Luer Loc Male MT: 1/2" Tri clamps 	-ETO Ethylene Oxide Sterilization			

Example:

MKPS020D3H3H-ETO

PureFlo MKP Series Mini Capsule, Single Layer 0.2µm PES membrane (Sterilizing Grade), Filtration Area 500 cm², 3/8" Hose Barb (Inlet/Outlet), ETO Sterilized (Factory Sterilized)

Part Number Description





Special Configuration:

Layer Option:

- Single Layer
- Dual Layer

Micron Rating Option:

- 0.04 Micron
- 0.1 Micron
- 0.2 Micron
- 0.45 Micron
- 0.65 Micron
- 0.8 Micron
- 1.2 Micron

Inlet Fitting Option:

- 1H: 1/8" Hose Barb
- 1Q: 1/8" Male Quick Coupling with Metal Latch
- 2H : 1/4" Hose Barbs
- 2H-FB: 1/4" Hose barbs with Filling Bell
- 2Q: 1/4" Male Quick Coupling for Metal Latch
- 2N: ¼" MNPT
- 3H : 3/8" Hose Barbs
- LF: Luer Loc Female
- MT : 1/2" Tri clamps

Outlet Fitting Option:

- 1H: 1/8" Hose Barb
- 1Q: 1/8" Male Quick Coupling with Metal Latch
- 2H: 1/4" Hose Barbs
- 2H-FB: 1/4" Hose barbs with Filling Bell
- 2Q: 1/4" Male Quick Coupling for Metal Latch
- 2N: ¼" MNPT
- 3H: 3/8" Hose Barbs
- LM : Luer Loc Male
- MT: 1/2" Tri clamps





