MZP Series PES Dual Layer Filter Capsules

MZP-Series PES capsules are gamma-irradiatable, easy-to-use filtration units, ideally suited for integration into single-use sterile assemblies. The hydrophilic, asymmetric Polyethersulfone (PES) membrane demonstrates exceptional throughput capacity and flow performance in liquid applications. The materials and design were carefully selected to ensure quick and reliable pre- and post-use integrity testing in critical biopharmaceutical applications. The filter capsules are available in a complete spectrum of sizes to simplify scale-up throughout each stage of the drug development process

Application

- **Buffers and Media**
- SVP (Small Volume Parenteral)
- Vaccines
- Scale up processing

- **Product Sterilization**
- **Bio Bags**
- **Biologics**
- **Ultrapure Water**

- Fermentation Broths
- **Pharmaceuticals**
- **Antibiotics**
- Serums

Part# MZPXS020DMTMT: MZP Filter Capsule with 1/2" Tri-Clamp Inlet/outlet Connection







Technical Data Sheet

Popular Configuration:

Catalog Number	Micron	Inlet	Outlet	Vent	Gamma
	Rating	Connection	Connection	Connection	Sterilization
MZPXS020DMTMT	0.65 μm / 0.2 μm (Dual Layer)	½" Tri-Clamp	½" Tri-Clamp	Luer Loc Cap	up to 50kGy

Find the complete list of available standard inlet and outlet fittings in page 5.

Micron Rating:

Final Membrane: 0.2 µm PES [Sterilizing Grade Membrane]

Pre-Filter: Highly Asymmetric 0.65 µm PES

Effective Filtration Area:

Mini Capsule; 77.5 in^2 (500 cm^2)

Materials of Construction:

Membrane: Pharmaceutical Grade PES (Polyethersulfone) Sterilizing Grade

Membrane Feature: Hydrophilic, Low Protein Binding, High Flow Rate

Media Supports: Polyester (PET)

Shell: Gamma Stable Polypropylene

Cage, Core, End Caps: Nylon

Valve O-Rings: Silicon (Standard)

Sealing: Thermally Bonded

Operating Conditions:

Maximum Forward Differential Pressure: Liquid: 5.5 bar (80psi) at 77°F/25°C

Gas: 4.1 bar (60psi) at 77°F/25°C

Minimum Burst Pressure: 8.3 bar (120psi) at 77°F/25°C

Maximum Reverse Differential Pressure: 3.0 bar (44psi) at 68°F/20°C

Maximum Operating Temperature: 176°F/80°C








Nominal Dimension:

Part Number: MZPXS020DMTMT

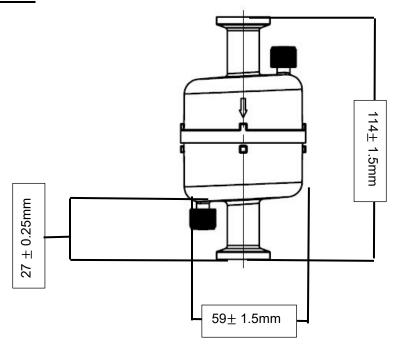

Hold-Up Volume (Approx.):

Upstream Volume: 36 ml Downstream Volume: 13 ml

Typical Filtered Volume:

10.0L - 28.0L

Filter Integrity:

The finished product was sampled and shown to exhibit a minimum bubble point of \geq 18 psi (1.2 bar) in 60%IPA / 40% water and \geq 50 psi (3.5 bar) in DI Water (at 22°C).

Sterilization:

Gamma Irradiation: The filters can be Gamma Sterilized up to 50kGy.

Autoclave: The filters can be autoclaved up to 10 cycles at 125°C (257°F) for 30 minutes.

Warning: The filters cannot be sterilized by steam-in-place (SIP)

Shelf Life:

The MZP capsules have a shelf life of 3 years from the date of manufacture.

Regulatory Compliance:









FP International Filter Products

Revolutionary systems for your industry


Category	Standard or Reference Test		
Physiochemical	 USP Oxidizable Substances USP <645> Conductivity USP <643> Non-Volatile Residue Total Organic Carbon (TOC) Analysis 		
Cleanliness	 USP <85> Bacterial Endotoxin Test USP <788> Particulate Matter 21 CFR 210.72 and 2010.3 (b) (6) Fiber Shedding 		
Biocompatibility	 ASTM F756-17 Hemolysis USP <88> (USP Class VI Plastics) Biological Reactivity Tests, In Vivo USP <87> / ISO 10993-5, In Vitro Cytotoxicity 21 CFR 177 FDA Indirect Food Additive 		
Bacterial Retention	Quantitative retention of 10 ⁷ CFU/cm ² Brevundimonas diminuta per ASTM methodology		
USP Bacterial Endotoxins	A representative sample from the lot is tested to confirm that an aqueous extraction of the product contains <0.25 EU/mL as determined by the Limulus Amebocyte Lysate (LAL) Test.		
Animal-derived Components	No animal-derived material is intentionally added or used during the manufacture of PureFlo Z Series PES Capsules.		
RoHS	Restriction of Hazardous Substances (RoHS 3) Directive 2015/863		
REACH	REACH (Registration, Evaluation, Authorization and Restriction of Chemicals) Article 57, Regulation No. 1907/2006)		
Traceability	 For traceability and easy identification: A Certificate of Conformance is provided with each capsule. Each filter bag and box are labeled with the product part number, lot number, and identifying characteristics. Each capsule is engraved and labeled with the product part number, lot number, and identifying characteristics. 		





MZP- PureFlo® Z Series PES Mini Capsules

Ordering Guide

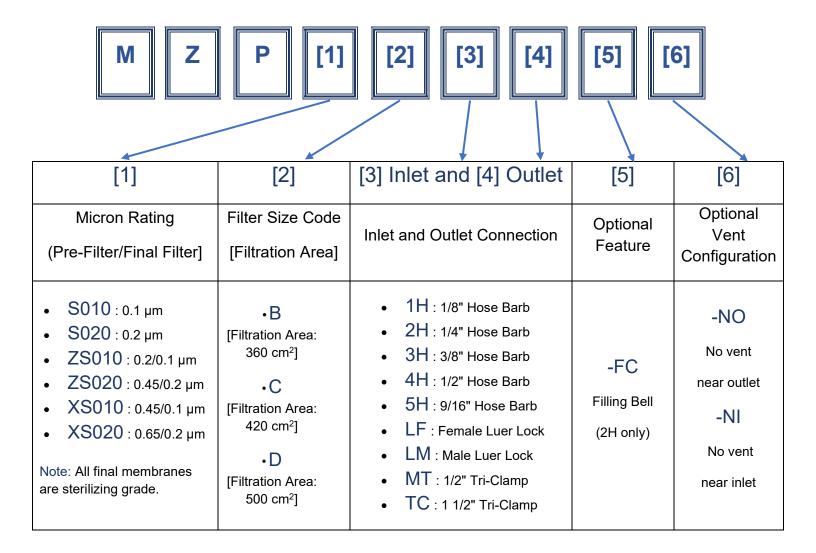

Example:

MZPXS020D3H3H

PureFlo MZP Series Mini Capsule, Highly Asymmetric Dual Layer PES 0.65µm / 0.2µm PES membrane (Sterilizing Grade), Filtration Area 500 cm², 3/8" Hose Barb (Inlet/Outlet)

Part Number Description







Special Configuration:

Layer Option:

- Single Layer
- Dual Layer

Micron Rating Option:

- 0.04 Micron
- 0.1 Micron
- 0.2 Micron
- 0.45 Micron
- 0.65 Micron
- 0.8 Micron
- 1.2 Micron

Inlet Fitting Option:

- 1H : 1/8" Hose Barb
- 1Q: 1/8" Male Quick Coupling with Metal Latch
- 2H: 1/4" Hose Barbs
- 2H-FB : 1/4" Hose barbs with Filling Bell
- 2Q: 1/4" Male Quick Coupling for Metal Latch
- 2N: ¼" MNPT
- LF : Luer Loc Female
- MT : 1/2" Tri clamps

Outlet Fitting Option:

- 1H: 1/8" Hose Barb
- 1Q: 1/8" Male Quick Coupling with Metal Latch
- 2H: 1/4" Hose Barbs
- 2H-FB: 1/4" Hose barbs with Filling Bell
- 2Q: 1/4" Male Quick Coupling for Metal Latch
- 2N: ¼" MNPT
- LM : Luer Loc Male
- MT : 1/2" Tri clamps





