FP International Filter Products

Revolutionary systems for your industry

SZV Series PES – Dual Layer Filter Capsules

SZV-Series PES capsules are gamma-irradiatable, easy-to-use filtration units, ideally suited for integration into single-use sterile assemblies. The hydrophilic, asymmetric Polyethersulfone (PES) membrane demonstrates exceptional throughput capacity and flow performance in liquid applications. The materials and design were carefully selected to ensure quick and reliable pre- and post-use integrity testing in critical biopharmaceutical applications. The filter capsules are available in a complete spectrum of sizes to simplify scale-up throughout each stage of the drug development process.

Application

Buffers and Media	Product Sterilization		Ophthalmic
Final Drug Product	Bio Bags		Pharmaceuticals
Vaccines	Biologics		Antibiotics
 Scale up processing 	Ultra-Pure Water	\blacktriangleright	Serums

Part Number: SZVXS020STCTC













Technical Data Sheet

Popular Configuration:

Catalog Number	Micron Rating	Inlet Connection	Outlet Connection	Vent Connection	Gamma Sterilization	
SZVXS020STCTC	0.65 μm / 0.2 μm (Dual Layer)	1.5" Tri-Clamp	1.5" Tri-Clamp	Bleed Valve with 1/8" Hose Barb	up to 50kGy	
Find the complete list of available standard inlet and outlet fittings in page 5.						

Micron Rating:

Final Media: 0.2 µm PES [Sterilizing Grade Membrane]

Pre-Filter: Highly Asymmetric 0.65 µm PES

Effective Filtration Area:

S-Size; 1.4 *ft*² (1,380 *cm*²)

Materials of Construction:

Membrane: Pharmaceutical Grade PES (Polyethersulfone) Sterilizing Grade

Membrane Feature: Hydrophilic, Low Protein Binding, High Flow Rate

Media Supports: Polyester (PET)

Cage, Core, End Caps: Nylon

Shell: Gamma Stable Polypropylene

Valve O-Rings: Silicon (Standard)

Sealing: Thermally Bonded

Operating Conditions:

Maximum Forward Differential Pressure:

Liquid: 5.5 bar (80psi) at 77°F/25°C

Gas: 4.1 bar (60psi) at 77°F/25°C

Minimum Burst Pressure: 8.3 bar (120psi) at 77°F/25°C

Maximum Reverse Differential Pressure: 3.0 bar (44psi) at 68°F/20°C

Maximum Operating Temperature: 176°F/80°C












Nominal Dimension:

Part Number: SZVXS020STCTC

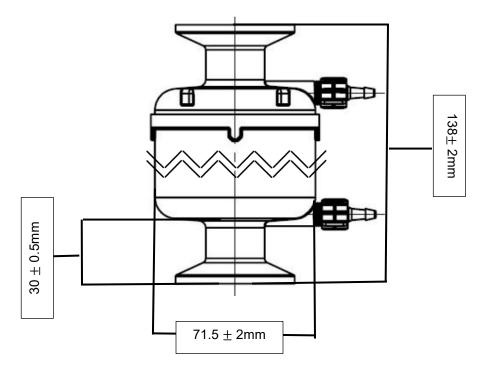

Hold-Up Volume (Approx.):

Upstream Volume: 121 ml

Downstream Volume: 30 ml

Typical Filtered Volume:

28.0L - 52.0L

Filter Integrity:

The finished product was sampled and shown to exhibit a minimum bubble point of \geq 18 psi (1.2 bar) in 60%IPA / 40% water and \geq 50 psi (3.5 bar) in DI Water (at 22°C).

Sterilization:

Gamma Irradiation: The filters can be Gamma Sterilized up to 50kGy.

Autoclave: The filters can be autoclaved up to 10 cycles at 125°C (257°F) for 30 minutes.

Warning: The filters cannot be sterilized by steam-in-place (SIP)

Shelf Life:

The SZV capsules have a shelf life of 3 years from the date of manufacture.









FP International Filter Products

Revolutionary systems for your industry

Regulatory Compliance:

Category	Standard or Reference Test		
Physiochemical	 USP Oxidizable Substances USP <645> Conductivity USP <643> Non-Volatile Residue Total Organic Carbon (TOC) Analysis 		
Cleanliness	 USP <85> Bacterial Endotoxin Test USP <788> Particulate Matter 21 CFR 210.72 and 2010.3 (b) (6) Fiber Shedding 		
Biocompatibility	 ASTM F756-17 Hemolysis USP <88> (USP Class VI Plastics) Biological Reactivity Tests, In Vivo USP <87> / ISO 10993-5, In Vitro Cytotoxicity 21 CFR 177 FDA Indirect Food Additive 		
Bacterial Retention	Quantitative retention of 10 ⁷ CFU/cm ² <i>Brevundimonas diminuta</i> per ASTM methodology		
USP Bacterial Endotoxins	A representative sample from the lot is tested to confirm that an aqueous extraction of the product contains <0.25 EU/mL as determined by the Limulus Amebocyte Lysate (LAL) Test.		
Animal-derived Components	No animal-derived material is intentionally added or used during the manufacture of PureFlo Z Series PES Capsules.		
RoHS	Restriction of Hazardous Substances (RoHS 3) Directive 2015/863		
REACH	REACH (Registration, Evaluation, Authorization and Restriction of Chemicals) Article 57, Regulation No. 1907/2006)		
Traceability	 For traceability and easy identification: A Certificate of Conformance is provided with each capsule. Each filter bag and box are labeled with the product part number, lot number, and identifying characteristics. Each capsule is engraved and labeled with the product part number, number, lot number, and identifying characteristics. 		



www.internationalfilterproducts.com 7914 Ajay Dr. Sun Valley, CA 91352 Tel: (818) 504-8115 - Office Tel: (818) 917-9538 - Cell E-mail: sales@internationalfilterproducts.com



PROUD MEMBER Alliance for Pharmacy Compounding



FP International Filter Products

Revolutionary systems for your industry

SZV- PureFlo® Z Series PES Mid-Size Capsules

Ordering Guide

SZV[1][2][3][4][5][6]						
[1]	[2]	[3] Inlet and [4] Outlet	[5]	[6]		
Micron Rating [Pre-Filter/Final Filter]	Filter Size Code [Filtration Area]	Inlet and Outlet Connection	Optional Feature	Optional Vent Configuration		
 S010 : 0.1 μm S020 : 0.2 μm ZS010 : 0.2/0.1 μm ZS020 : 0.45/0.2 μm XS010 : 0.45/0.1 μm XS020 : 0.65/0.2 μm Note: All final membranes are sterilizing grade.	 H [Filtration Area: 720 cm²] S [Filtration Area: 1380 cm²] L [Filtration Area: 2600 cm²] E [Filtration Area: 3980 cm²] K [Filtration Area: 5200 cm²] 	 1H : 1/8" Hose Barb 2H : 1/4" Hose Barb 3H : 3/8" Hose Barb 4H : 1/2" Hose Barb 5H : 9/16" Hose Barb MT : 1/2" Tri-Clamp TC : 1 1/2" Tri-Clamp 	-FC Filling Bell (2H or 3H only)	-NO No vent near outlet -NI No vent near inlet		

Example:

SZVXS020LTCTC

PureFlo SZV Series Capsule, Highly Asymmetric Dual Layer PES 0.65µm / 0.2µm PES membrane (Sterilizing Grade), 5.0", L-Size, Filtration Area 2,600 cm², 1.5" Tri-Clamps (Inlet/Outlet)

Part Number

Description







International Filter Products

Revolutionary systems for your industry

Special Configuration:

Layer Option:

- Single Layer
- **Dual Layer**

Micron Rating Option:

- 0.1 µm
- 0.2 µm •
- 0.2/0.1 µm
- 0.45/0.2 µm
- 0.45/0.1 µm
- 0.65/0.2 µm

Bleed Valve Seal **Option:**

- Silicone
- EPDM
- Nitrile
- Viton

Inlet Fitting Option:

- 1H = 1/8" Hose Barb •
- 2H = 1/4" Hose Barbs
- 3H = 3/8" Hose Barb
- 4H = 1/2" Hose Barb
- 5H = 5/8" Hose Barb
- MT = 1/2" Tri clamps
- TC = 1-1/2" Tri clamp

Outlet Fitting Option:

- 1H = 1/8" Hose Barb
- 2H = 1/4" Hose Barbs •
- 3H = 3/8" Hose Barb •
- 4H = 1/2" Hose Barb
- 5H = 5/8" Hose Barb
- MT = 1/2" Tri clamps
- TC = 1-1/2" Tri clamp

International Filter Products Revolutionary systems for your industry



