D65R Series Depth Polypropylene Filter Capsules

PureFlo® D65R Filter Capsules (65mm diameter) have been designed for simple, quick, and efficient filtration of fluids and gases used in laboratory, pilot, and small-scale applications. The family of products is particularly suitable for scale up testing. The discs are designed for prefiltration, clarification, and final filtration. The all-polypropylene construction for excellent chemical compatibility.

Application

SVP (Small Volume Parenteral) Fine Chemicals Ink

Clarification Cell Removal Pharmaceuticals

Vaccines Biologics Antibiotics

Scale up processing Water Serums







Technical Data Sheet

Popular Configuration:

Catalog Number	Effective Filtration Area	Inlet Connection	Outlet Connection
D65RDP0452H2H	26 cm ²	¼" Hose Barb	¼" Hose Barb
D65RDP045LFLF	26 cm ²	Luer Loc Female	Luer Loc Female
D65RDP045LFLM	26 cm ²	Luer Loc Female	Luer Loc Male

Find the complete list of available standard inlet and outlet fittings in page 4.

Micron Rating:

Final Media: 4.5 µm Depth Polypropylene

Effective Filtration Area:

 $4 in^2 (26 cm^2)$

Materials of Construction:

Media: **Depth Polypropylene** Media Supports: **Polypropylene**

Shell: Polypropylene

Sealing: Thermally Bonded

Operating Conditions:

Maximum Forward Differential Pressure: <u>Liquid:</u> 5.5 bar (80psi) at 77°F/25°C

Gas: 4.1 bar (60psi) at 77°F/25°C

Minimum Burst Pressure: 8.3 bar (120psi) at 77°F/25°C

Maximum Forward Differential Pressure: 5 bar (72psi) at 68°F/20°C Maximum Reverse Differential Pressure: 3.0 bar (44psi) at 68°F/20°C

Maximum Operating Temperature: 176°F/80°C

Regulatory Compliance:


The filters are constructed with polypropylene resins and filtration media in compliance with 21CFR Part 177 of the US Code of Federal Regulations and USP Class VI Biological Test for Plastic.






Nominal Dimension:


Part Number: D65RDP045LFLM

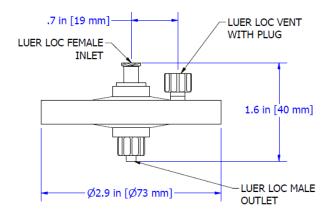

Part Number: D65RDP045LFLF

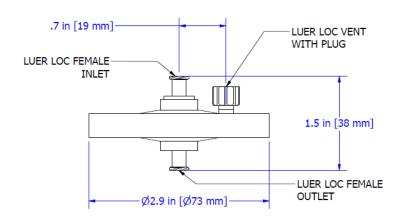

Part Number: D65RDP0452H2H

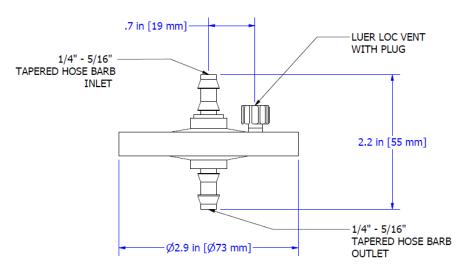







Sterilization:

The filters can be sterilized by autoclaving for 10 cycles at 257°F/125°C for 30 minutes.

Warning: The filters cannot be sterilized by steam-in-place (SIP)

Shelf Life:

The D65R disc filters have a shelf life of 3 years from the date of manufacture.

D65R- PureFlo® Depth Polypropylene Disc filter

Ordering Guide 1 [1] [2] Inlet and [3] Outlet [4] Micron Rating Inlet and Outlet Connection **Optional Feature** [Pre-Filter/Final Filter] 002 : 0.2 μm 1H: 1/8" Hose Barb 005 : 0.5 μm 2H: 1/4" Hose Barb -ETO **010**: 1.0 μm 3H: 3/8" Hose Barb Ethylene Oxide **015** : 1.5 μm LF: Female Luer Lock Sterilization **025**: 2.5 μm LM: Male Luer Lock (Optional) **045**: 4.5 μm MT: 1/2" Tri-Clamp 100 : 10.0 μm



200: 20.0 μm



TC: 1 1/2" Tri-Clamp

Example:

D65RDP045LFLF

PureFlo D65R Series 65mm Disc, Depth PP Media 4.5μm Final Filter, Filtration Area 26 cm², Luer Loc Female Inlet/Outlet.

Part Number

Description

Special Configuration:

Micron Rating Option:

- 0.2 µm
- 0.5 µm
- 1.0 µm
- 1.5 μm
- 2.5 µm
- 4.5 µm
- 10 µm
- 20 µm

Inlet Fitting Option:

- 1H = 1/8" Hose Barb
- 1Q = 1/8" Male Quick Coupling with Metal Latch
- 2H = 1/4" Hose Barbs
- 2HS = 1/4" to 3/8 " Hose Barb
- 2Q = 1/4" Male Quick Coupling for Metal Latch
- 3H = 3/8" Hose Barb
- 4H = 1/2" Hose Barb
- 4Q = 1/2" Male Quick Coupling for Plastic latch
- 5H = 5/8" Hose Barb
- MT = 1/2" Tri clamps
- TC = 1-1/2" Tri clamp

Outlet Fitting Option:

- 1H = 1/8" Hose Barb
- 1Q = 1/8" Male Quick Coupling with Metal Latch
- 2H = 1/4" Hose Barbs
- 2HS = 1/4" to 3/8 " Hose Barb
- 2Q = 1/4" Male Quick Coupling for Metal Latch
- 3H = 3/8" Hose Barb
- 4H = 1/2" Hose Barb
- 4Q = 1/2" Male Quick Coupling for Plastic latch
- 5H = 5/8" Hose Barb
- MT = 1/2" Tri clamps
- TC = 1-1/2" Tri clamp



