



DirecTaq™ 1X Master Mix

Expressed in: *E.coli*

Contents:

DirecTaq 1X Master Mix saves time and cost by enabling direct PCR amplification of unpurified templates. DirecTaq 1X Master Mix is a PCR master mix containing a recombinant, truncated (lacks 5' to 3' exonuclease activity), highly thermostable DNA polymerase from the thermophilic bacterium *Thermus aquaticus*. DirecTaq is thermostable up to 98 °C for PCR assays and is provided as a complete reaction mix consisting of reaction buffer, dNTPs, MgCl₂, and loading dye and only requires the addition of primers and template DNA.

Background:

DirecTaq DNA polymerase is a recombinant, truncated (lacks 5' to 3' exonuclease activity), and highly thermostable DNA polymerase from the thermophilic bacterium *Thermus aquaticus*. The enzyme is thermostable up to 98 °C for PCR assays. It is supplied as a complete reaction mix requiring only primers and template DNA, ideal for colony PCR and genotyping. After the reaction is complete, the PCR products can be loaded directly to an agarose gel for analysis.

Application Notes:

DirecTaq DNA polymerase (exonuclease minus) is resistant to inhibitors and ideal for PCR of GC- rich templates.

**These products are intended for research use only, not for diagnostic use. The safety and efficacy of these products in diagnostic or other clinical uses has not been established.*

Shipping and Storage:

DirecTaq DNA polymerase is supplied in a reaction mix of 10 mM Tris-HCl, 1.5 mM MgCl₂, 50 mM KCl, 0.1% Triton X-100, 15% Sucrose, 0.0268% Orange G, 0.0022% Xylene cyanol FF.

DirecTaq DNA polymerase is shipped on dry or blue ice. On arrival store at -20 °C for optimum stability.

Quality Control:

- DirecTaq DNA polymerase Unit activity: DirecTaq 1X Master Mix contains 3.25 U of DirecTaq. A known polymerase is used to create a standard curve with a real-time qPCR assay against which the activity of this enzyme is measured.
- Purity: >95% as determined by SDS-PAGE analysis
- DirecTaq DNA polymerase is free of detectable RNase and DNase (exo- and endonuclease).
- <0.05 ng contaminating host DNA per 3.25 U of DirecTaq

Reaction set up:

- Thaw all reagents prior to setting up PCR.
- Set up reaction on ice (4 °C).

Component	Stock	~25 μL Rxn	Final Conc.
1X DirecTaq Master Mix	1X	25 μ L	1X
DNA template	Varies	1.0 μ L	
Forward Primer	100 μ M	0.25 μ L	~1.0 μ M
Reverse Primer	100 μ M	0.25 μ L	~1.0 μ M