

## Aqueous-Mount (Low Viscosity)

<b>Description</b>	Aqueous-Mount is formulated for cover slipping sections stained with alcohol soluble chromogens, such as Fast-Red or AEC. No heating is required prior to use. Coverslips may be removed by soaking in water.
<b>Storage</b>	Store at room temperature. Do not use past the expiration date.
<b>Procedure</b>	Apply several drops of Aqueous-Mount and coverslip. To remove coverslips, soak overnight in water.
<b>Precautions</b>	Avoid contact with skin and eyes. Wash after use. Observe all federal, state and local environmental regulations regarding disposal.