

Aqueous-Mount

Description	Aqueous-Mount is formulated for cover slipping sections stained with alcohol soluble chromogens, such as Fast-Red or AEC. No heating is required prior to use. Coverslips may be removed by soaking in water.
Storage	Store at room temperature. Do not use past the expiration date.
Procedure	Apply several drops of Aqueous-Mount and coverslip. To remove coverslips, soak overnight in water.
Precautions	Avoid contact with skin and eyes. Wash after use. Observe all federal, state, and local environmental regulations regarding disposal.