

## PeroxiBlock Basic

<b>Description</b>	PeroxiBlock Basic has been developed to use with immunolabeling techniques for the reduction of nonspecific background staining do to endogenous peroxidase.
<b>Appearance</b>	Clear liquid
<b>Uses</b>	Immunohistochemical procedures.
<b>Limitations</b>	Do not use past expiration date.
<b>Storage</b>	Store PeroxiBlock at 2-8°C. Product is stable for 18 months from date of manufacture.
<b>Procedure</b>	<ol style="list-style-type: none"><li>1. Incubate tissue section for 10 minutes at either room temperature or 37°C prior to application of the primary antibody.</li><li>2. After incubation, rinse once in buffer.</li><li>3. For bulk staining, pour PeroxiBlock Basic in a covered staining tray and dip slides for 5 minutes. Replace with fresh reagent after 5-10 uses.</li></ol>
<b>Precautions</b>	Do not pipette reagent by mouth. Avoid contact with skin and eyes. Wash after use. Observe all federal, state and local environmental regulations regarding disposal.