

FluoroSave Mounting Medium (Hard Set)

Description	FluoroSave Mounting Medium (Hard Set) is a water-soluble medium designed for semi- permanent coverslipping of fluorescent slide preparations. FluoroSave dries to a semi-rigid layer that eliminates tissue damage due to moving coverslips and facilitates long-term storage. The Hard Set formulation provides better drying in humid conditions.
Uses/Limitations	For in vitro diagnostic use only. Histological applications. Do not use past expiration date. Use caution when handling these reagents.
Storage	Store at room temperature.
Procedure	<ol style="list-style-type: none">1. Apply several drops of FluoroSave Mounting Medium (Hard Set) and coverslip.2. Allow FluoroSave to dry for 30 minutes prior to viewing. Product will dry to a semi-solid form overnight.3. FluoroSave is an aqueous based medium and may be removed by soaking in PBS overnight
Precautions	Avoid contact with skin and eyes. Wash after use. Observe all federal, state, and local environmental regulations regarding disposal.