

## FluoroSave Mounting Medium

<b>Description</b>	FluoroSave Mounting Medium is a water-soluble medium designed for semi-permanent coverslipping of fluorescent slide preparations. FluoroSave dries to a semi-rigid layer that eliminates tissue damage due to moving coverslips and facilitates long term storage.
<b>Uses/Limitations</b>	For in vitro diagnostic use only. Histological applications. Do not use past expiration date. Use caution when handling these reagents.
<b>Storage</b>	Store at room temperature.
<b>Procedure</b>	<ol style="list-style-type: none"><li>1. Apply several drops of FluoroSave Mounting Medium and coverslip.</li><li>2. Allow FluoroSave to dry for 30 minutes prior to viewing. Product will dry to a semi-solid form overnight.</li><li>3. FluoroSave is an aqueous based medium and may be removed by soaking in PBS overnight.</li></ol>
<b>Precautions</b>	Avoid contact with skin and eyes. Wash after use. Observe all federal, state and local environmental regulations regarding disposal.