



SilTite®


## Instructions for SilTite® ferrules

### Important

**⚠ Caution** Do not overtighten! Further tightening of the column will cause it to break.

### Instructions

1. For easy installation gradually tighten the nut until the ferrule just begins to hold the fused silica.
2. Then tighten a further 60°; a leak tight seal is guaranteed.

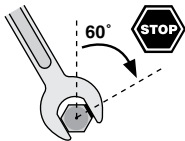

### Information and support

Visit [www.trajanscimed.com](http://www.trajanscimed.com) or  
contact [techsupport@trajanscimed.com](mailto:techsupport@trajanscimed.com)

*Specifications are subject to change without notice.*

MN-0330-A\_RevE © Trajan Scientific Australia Pty Ltd 11/2016