



SilTite[®]

Instructions for Agilent split/splitless injector SilTite[®] kit

Instructions

1. Unscrew the insulation cup fitted around the column nut inside the oven.
2. Remove the column nut (e) and reducing nut (c) from the injector.
3. Remove the base seal from the reducing nut and place the gold SilTite[®] base seal (a) in its place (cross-side up) re-using the original washer (b).
4. Install the column (f) (as above) and tighten the nut until the SilTite ferrule (d) grabs, then tighten a further 60° for a leak tight seal.

Information and support

Visit www.trajanscimed.com or
contact techsupport@trajanscimed.com

Specifications are subject to change without notice.

MN-0331-A_RevC © Trajan Scientific Australia Pty Ltd 11/2016

