



SilTite®


Instructions for SilTite® restrictor seals

Important

⚠ Caution The fused silica restrictor will be broken after approximately 180° of further turning.

Instructions

1. Use fingers (rather than a wrench) to tighten the tee onto the interface (or adapter) until it stops.
2. Use a wrench to tighten to a further 30°. A reliable seal will be achieved. Do not tighten further as the interface may be damaged.

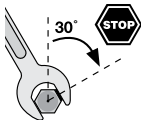

Information and support

Visit www.trajanscimed.com or
contact techsupport@trajanscimed.com

Specifications are subject to change without notice.

MN-0350-A_RevC © Trajan Scientific Australia Pty Ltd 11/2016