

Manufacturing Chemists Since 1953

Fast, Proven, Trusted Since 1963

Professional Strength Biological & Chemical STAIN REMOVER



www.ecamco.com

"There is Simply Nothing Like it "





- Fast-Acting Versatile Readily Biodegradable Surfactants
 - Neutral pH 7-8
 No Bleach, Peroxide or Phosphates

Helps Remove: Most medical, histological & chemical stains including iodine, Wright Stain, Lugol's Stain, fresh blood stains and more.

Erado-Sol helps remove stains from anywhere: hands, glassware, lab equipment, stainless steel, plastic, counters, floors, lab coats, clothing and hospital linen.

Easy to Apply: Simply rub in gently (and like magic): the stain is removed in a matter of seconds. Rinse with water. For more stubborn stains, repeat as needed.

Surfaces and Floors: Simply apply to floors or equipment for difficult stains, let set for up to a couple of hours, then rinse with water. Repeat if necessary.

Colorfast Labcoats & Hospital Linen: Lightly moisten stained area with water. Place small amount of Erado-Sol directly on each stain & massage into fabric. Let sit 15 seconds. Rinse with water. Repeat if needed. For very stubborn stains: let soak, then rinse with water & launder as usual.

Available in the following sizes

207-F-18	18 Wipes	Display Box
207-2	2 fl. oz. / 59 mL	Bottle
207-8	8 fl. oz. / 236 mL	Foamy Dispenser Bottle
207-16	16 fl. oz. / 473 mL	Bottle
207-32	32 fl. oz. / 946 mL	Bottle
207-G	128 fl. oz. / 3.785 L	Bottle
207-D	55 gallon / 208 L	Drum

Liquids also available in fragrance & dye free formula.

*Wipes are offered only in the Fragrance & Dye Free formulation.

Available from your laboratory supply distributor or directly from:

Cambridge Diagnostic Products, Inc. 6880 NW 17th Avenue Fort Lauderdale, FL 33309

For additional information: techinfo@ecamco.com 954.971.4040 or 800.525.6262 www.ecamco.com



Manufacturing Chemists Since 1953