



# Cyclops Metallographic Assembly Guide

## Contents

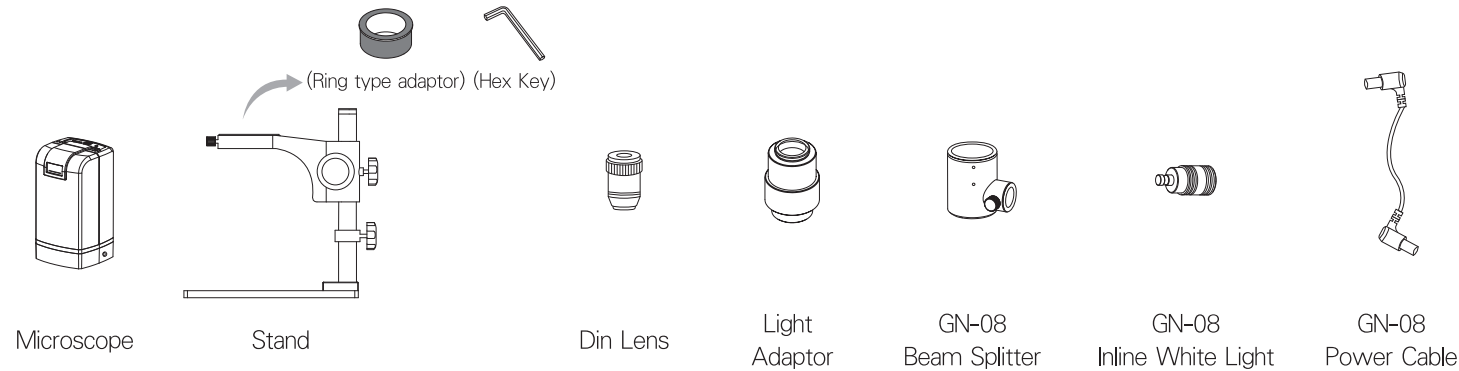

## Ways for Observation

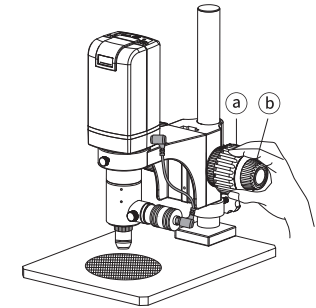


- Place Object on the bottom plate and rotate (a) and (b) to find clear image. (b) for fine tune

## Assembling Steps

- ① Insert the Ring Adaptor from Stand to Light Adaptor, screw tight by hex key.
- ② Screw the Light Adaptor upward to Microscope.
- ③ Insert Microscope into stand, facing front and screw tight.
- ④ Insert Beam Splitter to Light Adaptor and screw tight by hex key. \* Please hold the Beam Splitter until it screwed tight.
- ⑤ Insert Inline White Light.
- ⑥ Screw the Din Lens upright.
- ⑦ Connect Light Control Box and Power Cable.
- ⑧ Assembling Completed. \* Use Remote Control or Software to adjust brightness.

## Adjust Inline White Light Shading

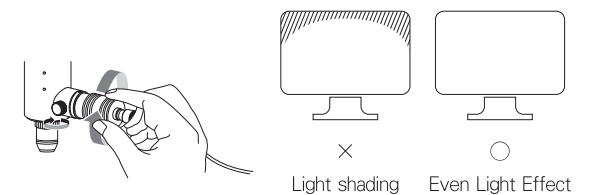

- When seeing Light Shading on the corner of image, rotate the Inline White Light until see clear image. Then screw tight to lock the position.