


## TECHNICAL RELEASE LUZCHEM EXPOSURE STANDARD: LES-LEDGR-14

**General information**: Luzchem Research, Inc. produces and distributes freely these standards as a service to scientists involved in research in photochemistry, photobiology and photostability. These standards are available in Luzchem's website so that references to any of our standards can be used to define conditions of exposure, and should allow other scientists to readily replicate irradiation conditions. Luzchem has measured these exposure standards using a Luzchem SPR-4001 spectroradiometer calibrated against traceable NIST standards within the 6 months preceding the determination. To the best of our knowledge spectral information is accurate within the experimental bandwidth of 2 nm.

Lamp part number	Luzchem LED-GR	Measurement temperature	23 °C
Filter	None, direct exposure	Measurement date	14-August-2014
Filter effect determined by	Not applicable	Spectral range monitored	240 to 880 nm
Photoreactor model	EXPO-LED	Harmonic peak interference	None observed
Number of lamps and	5 overhead lamps,	Lamp type:	T5 Monochromatic
location	approx 10 cm from target	FWHM:	Approx. 31 nm
Cold start wavelength	520 nm <sup>#</sup>	Resolved peaks	520 nm

<sup>&</sup>lt;sup>#</sup> LED lamps start at a shorter wavelength than their stabilized temperature and take about 5 min to equilibrate.

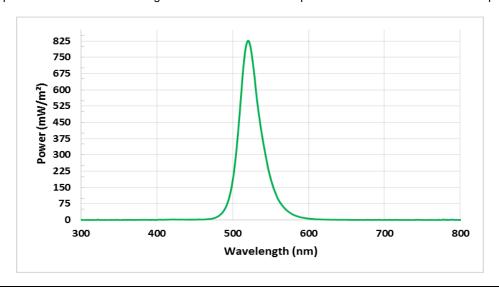

Region	Range nm <sup>a</sup>	Dose mW/m <sup>2</sup>	% Energy <sup>b</sup>
UVA	316-400	56	<1%
Peak			
Wavelength	490-550	26381	>88%
Range			
Visible	400-700	29945	>99%
NIR	701-850	(52) <sup>b</sup>	<1%

Integration range. <sup>b</sup> Within experimental error.

The table to the left shows the energy distribution at the target, expressed as total irradiance and as a percentage of the total energy. The calculations refer to the monitored range indicated.

Toll Free: 1-800-397-0977 Phone: (613)-749-2442 Fax: (613) 749-2393 E-mail: sales@Luzchem.com

Luzchem Research, Inc.