

Quality, Design and Innovation





LIEBHERR

Con	tents	
1	The appliance at a glance	3
1.2 1.3	Overview of appliance and equipmentSpecial features	3
1.4 1.5	Range of appliance use	3
1.6	Conformity	4
2	General safety instructions	4
3 3.1	Start-up Setup conditions	6
3.2	Appliance dimensions	6
3.3 3.4	Transporting the appliance	7
3.5 3.6	Removing the transport lock	7
3.7	Installing the door handle	7
3.8 3.9	Setting up the applianceLeveling out the appliance	8
3.10	Setting up multiple appliances	8
3.11 3.12	After setup Disposal of packaging	9
3.13	Reversing the door	9
3.14 3.15	Aligning the door	16 16
3.16	Switching on the appliance (first use)	17
4 4.1	Use Control and display elements	17 17
4.1.1	Status display	17
4.1.2 4.1.3	Display symbolsAcoustic signals	
4.2	Navigation	18
4.2.1 4.2.2	Navigating with buttons Settings menu	
4.2.3 4.2.4	Customer menu	19
4.2.4	Setting values General rules	19
4.3 4.3.1	Functions Overview of functions	
4.3.2	Switching appliance on and off	. 20
4.3.3 4.3.4	Temperature Temperature recording	21
4.3.5	Settings menu access protection	2]
4.3.6 4.3.7	Access codesMaintenance interval reminder	. 23
4.3.8 4.3.9	Language Temperature unit	
4.3.10	Display brightness	24
4.3.11 4.3.12	Alarm SoundKey Sound	
4.3.13	WiFi connection	24
4.3.14 4.3.15	LAN connectionInfo	
4.3.16 4.3.17	Operating hours Software	
4.3.18	Defrosting	27
4.3.19 4.3.20	Sensor calibration Temperature alarm	
4.3.21	Door alarm	. 28
4.3.22 4.3.23	Alarm simulation	
4.3.24 4.3.25	DemoMode	30
4.4	Factory reset	3]
4.4.1 4.4.2	Warnings Reminders	3] 33
5	Equipment	34
5.1	Safety lock	34
5.2	Sensors	34

6 6.1 6.2 6.3	Maintenance Maintenance schedule Defrosting the appliance Cleaning the appliance	35 35 36 37
7 7.1 7.2 7.3 7.4	Customer support	38 38 38 38 39
8	Shutting down	39
9 9.1 9.2	Disposal Preparing appliance for disposal Disposing of the appliance in an environmentally friendly manner	39 39
10	Additional information	40
11	Quick Start Guide for everyday use	41
	duick Start Guide for everyday use	7.

The manufacturer is constantly working to improve all types and models. Therefore, please be aware that we reserve the right to make changes to the shape, equipment and tech-

nology.			
Symbol	Explanation		
	Read instructions Please read the information in these instructions carefully to understand all of the benefits of your new appliance.		
	You can find detailed instructions on the internet using the QR code on the front of these instruction or by entering the service number at home.liebherr.com/fridgemanuals. The service number can be found on the serial tag:		
	Fig. Example illustration		
	Check appliance Check all parts for transport damage. If you have any complaints, please contact your agent or customer service.		
*	Differences These instructions apply to a range of models, so there may be differences. Sections that apply to certain models only are indicated by		

Videos

Videos about the appliances are available on the YouTube channel of Liebherr-Hausgeräte.

These operating instructions apply to:

an asterisk (*).

Instructions and results

Instructions are marked with a ▶. Results are marked with a \triangleright .

- SRFfg 3501

5.3

- SRFfg 4001
- SRFfg 5501

1 The appliance at a glance

1.1 Included in delivery

Check all parts for transport damage. If you have any issues, please contact your dealer or Customer Service. (see 7.3 Customer service)

The delivery contains the following parts:

- Freestanding appliance
- Equipment *
- Installation materials *
- Operating instructions
- Service brochure
- Quality certificate *
- Power cable
- Handle incl. screws and Torx tool
- Wall fastening kit

1.2 Overview of appliance and equipment

SRFfg 3501 SRFfg 4001

SRFfg 5501

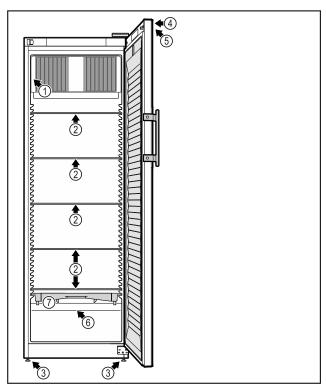


Fig. 1 Example illustration

- (1) Type plate
- (2) Glass shelf
- (3) Adjustable feet
- (4) Lock

Equipment

- (5) Control elements and temperature display
- (6) Safety thermostat sensor
- (7) Defrost water collection tray

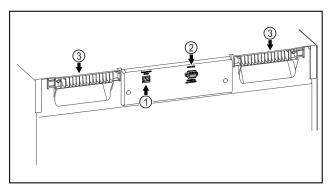


Fig. 2 Example illustration of the rear

Equipment

- (1) LAN/WiFi interface *
- (3) Carrying aid
- (2) Potential-free alarm output

1.3 Special features

Note

Accessories are available via Customer Service (see 7.3 Customer service) and in the Liebherr-Hausgeräte-Shop at https://home.liebherr.com.

SmartModule

The appliance can be equipped with a SmartModule.

This is a WiFi and LAN interface for the connection between the appliance and the Internet.

1.4 Range of appliance use

Intended use

This laboratory refrigerator, with an interior free of ignition sources, is suitable for professional storage of highly flammable products in closed containers, at temperatures between 3 °C and 16 °C.

Explosion-proof interior

The values relating to your appliance are printed on the sticker on the appliance door.



Fig. 3 Example illustration

The interior is certified as an area subject to explosion hazards as per 2014/34/EU (ATEX Directive). Typical products to be stored include research samples, reagents, laboratory inventory etc., that are categorized into the appropriate zone, explosion group, and temperature class.

Refer to the safety data sheet for details of the explosion group of the products to be stored. Contact the supplier of the respective product if there is any uncertainty.

General safety instructions

It is necessary to use an independent alarm system with permanent monitoring if storing valuable or temperature-sensitive substances or products. This alarm system must be designed in such a way that every alarm state is immediately registered by a competent person who then takes appropriate measures.

Foreseeable incorrect use

Do not use the appliance for the following applications:

- Storage and cooling of:
 - · Chemically unstable substances
 - Blood, plasma or other body fluids for the purpose of infusion, application or introduction in the human body.
- Use in potentially explosive atmospheres.
- Use outdoors or in areas exposed to dampness and splashing water.

Any misuse of the appliance may result in damage to or spoilage of stored goods.

Climate classes

The climate class for your appliance is printed on the identification plate.



Fig. 4 Type plate

(X) The climate class states the ambient conditions in which the appliance can be operated safely.

Climate class (X)	max. room temperature in °C	
7	35	75

Note

The minimum permitted room temperature at the setup location is 10 °C.

Slight condensation may form on the glass door on the side walls if the boundary conditions are exceeded.

1.5 Sound emission of the appliance

The A-weighted emission sound pressure level during operation of the appliance is below 70 dB(A) (sound power rel.1 pW).

1.6 Conformity

The refrigerant circuit has been tested for leaks.

This appliance complies with the applicable safety regulations and with EU directives 2014/35/EU, 2014/30/EU, 2014/34/EU and 2011/65/EU.

The full text of the EU Declaration of Conformity is available on the following website: www.Liebherr.com

2 General safety instructions

Please keep these operating instructions in a safe place so you can refer back to them at any time.

If you pass the appliance on, please hand these operating instructions to the new owner.

Read these operating instructions before use in order to use the appliance safely and correctly. Follow the instructions, safety instructions and warning messages included at all times. They are important for ensuring you can operate and install the appliance safely and without any problems.

Dangers for the user:

- This appliance may only be used by specialist and laboratory personnel who have been trained for this purpose and are familiar with all the safety measures for work in a laboratory. Children and persons with impaired physical, sensory or mental abilities or with a lack of experience and knowledge may not commission or operate this appliance.
- The socket must be easily accessible so that the appliance can be disconnected quickly from the electricity in an emergency. It must not be located in the area behind the appliance.
- Always hold the plug of the cable when disconnecting the appliance from the power supply. Do not pull on the cable.
- Remove the plug or disconnect via the fuse if there is a malfunction.
- WARNING: Do not damage the power cable.
 Do not operate the appliance with a faulty power cable.
- WARNING: Multi-sockets/power distributors and other electronic appliances (such as halogen transformers) may not be placed and operated behind appliances.
- WARNING: Do not block the ventilation openings in the appliance housing or in the installation housing.
- WARNING: Dust deposits on fan blades!
 Risk of sparks due to friction!

Do not store dusty objects in the appliance.

The ventilation slots of the recirculation fan must be cleaned each month using a vacuum cleaner.

- Repairs and work on the appliance may only be carried out by Customer Service or other specifically trained qualified personnel.
- Always follow the instructions when assembling, connecting and disposing of the appliance.

Risk of fire:

- The refrigerant contained within the appliance (specifications on the type plate) is environmentally friendly, but flammable. Leaking refrigerant can ignite.
 - WARNING: Do not damage the refrigerant circuit
 - Do not handle ignition sources inside the appliance.
 - WARNING: Do not use electrical appliances inside the food storage compartments of the appliance, unless they are of the type recommended by the manufacturer.
 - If refrigerant leaks: Remove naked flames or ignition sources located near the area of the leak. Ventilate the room well. Contact Customer Service.
- Do not operate the appliance near explosive gases.
- Do not store or use gasoline or other flammable gases and liquids near the appliance.
- Keep lit candles, lamps and other objects with naked flames away from the appliance so they do not cause a fire.

Risk of falling or toppling over:

- WARNING: To avoid a hazard due to instability of the appliance, it must be fixed in accordance with the instructions.
- Do not stand on the base, drawers, doors etc. or use them as improper supports.

Danger of frostbite, feeling of numbness and pain:

 Avoid prolonged skin contact with cold surfaces or chilled/frozen food or take protective measures, e.g. wear gloves.

Risk of injury and damage:

 WARNING: Do not use mechanical devices or other means to accelerate the defrosting process, other than those recommended by the manufacturer.

- WARNING: Risk of injury due to electric shock! There are live electrical parts under the cover
- NOTICE: The appliance must only be operated using original manufacturer accessories or accessories from other providers approved by the manufacturer. The user bears the risk of using accessories which are not approved.

Risk of crushing:

 Do not reach into the hinge when opening and closing the door. Fingers may get trapped.

Specialist personnel qualifications:

The appliance may only be installed, tested, maintained, and commissioned by specialist personnel who are familiar with the installation, commissioning, and operation of the appliance.

Specialist personnel are persons who, on account of their specialist training, knowledge and experience as well as their knowledge of the relevant standards, are able to assess and perform the work assigned to them and identify potential hazards. They must have training, instruction, and authorization to work on the appliance.

Symbols on the device:



The symbol may be located on the compressor. It refers to the oil in the compressor and refers to the following danger: **Can be fatal if swallowed or inhaled.** This notice only applies for recycling. There is no danger during normal operation.



WARNING: Risk of fire / flammable materials. The symbol is located on the compressor and indicates the danger from flammable materials. Do not remove the label.



The symbol is located on the back of the appliance near the alarm relay and indicates the following danger: **Electric shock!** Even if an appliance is disconnected from the mains there may still be extraneous voltage. Do not remove the label.



This label or a similar one may be located on the rear of the appliance. It refers to the foam panels in the door and/or the housing. This notice only applies for recycling. Do not remove the label.

Observe the warning messages and other detailed information in the other sections:

<u> </u>	DANGER	Indicates an immediately hazardous situation, which if not avoided, will result in death or serious injury.
<u>^</u>	WARNING	Indicates a hazardous situation, which if not avoided, could result in death or serious injury.
<u> </u>	CAUTION	Indicates a hazardous situation, which if not avoided, could result in minor or moderate injury.
	NOTICE	Indicates a hazardous situation, which if not avoided, could result in damage to property.
	Note	Indicates useful advice and tips.

3 Start-up

3.1 Setup conditions



WARNING

Risk of fire due to moisture!

If live parts or the power cord get wet, this can cause a short circuit.

► The appliance is designed for use in enclosed spaces. Do not operate the appliance in open space or in damp areas or where there is spray.

3.1.1 Setup location



WARNING

Leaking refrigerant and oil!

Fire. The refrigerant contained within the appliance is environmentally friendly, but flammable. The oil contained within the appliance is flammable. Escaping refrigerant and oil can ignite if they are of high enough concentration and are exposed to an external heat source.

- ▶ Do not damage the pipelines of the coolant circuit and the compressor.
- Do not setup the appliance in direct sunlight, next to an oven, radiator or similar.
- The best place to set up the appliance is a dry and well ventilated room.
- If the appliance is set up in a very humid environment, condensation can form on the outside of the appliance.
 Always ensure sufficient airflow and ventilation in the setup location.
- The more refrigerant there is in the appliance, the larger the room must be in which the appliance is located. If rooms are too small, any leak may create a flammable mixture of gas and air. For every 8 g of refrigerant, the size of the installation space must be at least 1 m³. Specifications regarding the refrigerant contained within the appliance can be found on the type plate inside the appliance.
- The floor of the setup location must be horizontal and even.
- The setup location must be able to withstand the weight of the appliance plus the weight when stocked to maximum capacity. (see 7.1 Technical data)

3.1.2 Electrical connection



WARNING

Danger of fire due to incorrect positioning!

If the power supply cable or plug touches the back of the appliance, the vibration can damage the power supply cable or the plug resulting in a short circuit.

- ► Make sure the power supply cable is not trapped under the appliance when you position the appliance.
- ▶ Install the appliance so that it does not touch any plugs or power cables.
- ▶ Do not connect any appliances to sockets in the area of the back of the appliance.
- ▶ Do not place and operate power strips/power distributors and other electronic devices (such as halogen transformers) at the back of the appliances.

3.2 Appliance dimensions

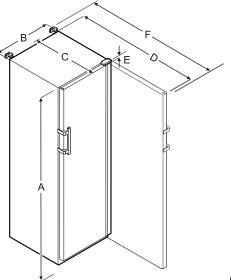


Fig. 5

SRFfg 3501					
A B C		С	D	E	F
1684 mm	597 mm	664 mm	1203 mm	23 mm	1241 mm

SRFfg 4001					
A	С	D	E	F	
1884 mm	597 mm	664 mm	1203 mm	23 mm	1241 mm

SRFfg 5501						
A B C D E F				F		
1684 mm	747 mm	779 mm	1468 mm	23 mm	1506 mm	

- A = appliance height including feet/castors
- **B** = appliance width without handle
- C = appliance depth without handle
- **D** = appliance depth with door open
- **E** = hinge height
- ${f F}$ = appliance depth with protruding handle and anti-tipping device

3.3 Transporting the appliance

- ► Transport the appliance in its packaging.
- ► Transport the appliance upright.

▶ Use two people when transporting the appliance.

3.4 Unpacking the appliance

- ▶ Check the appliance and the packaging for transport damage. Contact the supplier immediately if you suspect any damage. Do not connect the appliance to the power
- ▶ Remove all materials from the back or the side walls of the appliance that may prevent proper installation or prevent air flow and ventilation.

3.5 Removing the transport lock

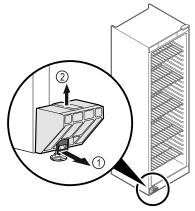


Fig. 6

- Pull tab Fig. 6 (1) forwards.
- Take off transport lock Fig. 6 (2) in upwards direction.

3.6 Installing the door handle


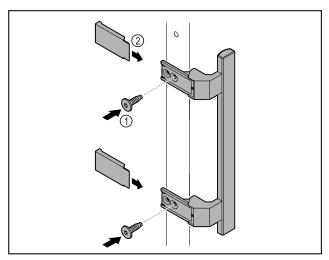


Fig. 7

- ▶ Attach handle from accessory kit to the door using supplied screws Fig. 7 (1).

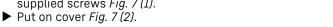


3.7 Mounting the anti-tipping device

Tools





Prevent appliance from tipping over.

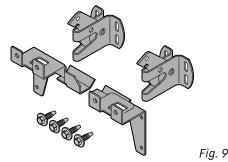

The anti-tipping device, consisting of two retaining parts, two retaining hooks and four screws (4 x 14), is included with the appliance.


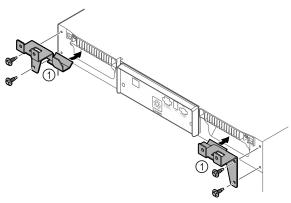


Fig. 10

- Mount retaining parts Fig. 10 (1) on the appliance using the supplied screws.
- Push appliance with retaining parts mounted against the wall.

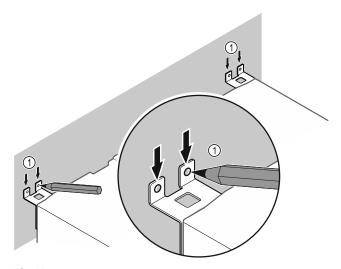

- Make markings on the wall Fig. 11 (1).
- Remove appliance.

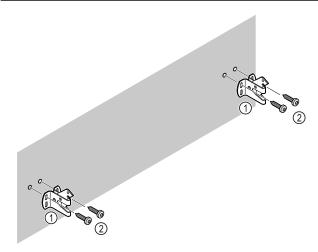


Fig. 12

Use fixing material (e.g. wall anchors) which is appropriate for the nature of the wall (wood, concrete) and sufficient attachment points *Fig. 12 (2)*.

► Attach retaining hooks. Fig. 12 (1)

3.8 Setting up the appliance



CAUTION

Risk of injury and damage.

▶ Use 2 people to set up the appliance.



CAUTION

Risk of injury and damage.

The door can strike against the wall and become damaged as a result. In the case of glass doors, the damaged glass can cause injuries.

- Protect the door from striking against the wall. Attach a door stopper, e.g. felt stopper, to the wall.
- ► Connect all necessary components (e.g. power cable) to the back of the appliance and route to the side.

Note

Cables can be damaged.

▶ Do not crush the cable when pushing the appliance back.

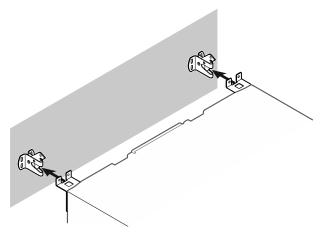


Fig. 13

- ► Slide the appliance up against the wall so that the retaining parts latch into the retaining hooks.
- > The appliance is now secured against tipping.
- > It can be released by folding back the retaining hooks.

3.9 Leveling out the appliance

NOTICE

Appliance body can become deformed and door will not close.

- ► Align appliance horizontally and vertically.
- Compensate for uneven floors using adjustable feet.



WARNING

Incorrect height adjustment of the adjustable foot! Severe or fatal injuries. Incorrect height adjustment can cause the bottom part of the adjustable foot to come loose and the appliance to tip over.

▶ Do not unscrew the adjustable foot too far.



Fig. 14

Raising appliance:

► Turn adjustable foot clockwise.

Lowering appliance:


► Turn adjustable foot counterclockwise.

3.10 Setting up multiple appliances

NOTICE

Risk of damage due to condensation between the side walls.

- ▶ Do not set up the appliance directly next to another refrigeration appliance.
- Set up appliances with a space of 3 cm between appliances.
- ▶ Only set up multiple appliances up to temperatures of 35 °C and 65 % humidity next to one another.
- At higher levels of humidity, increase the space between appliances.

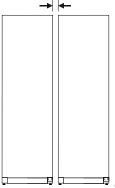

耳 Fig. 15 Side-by-side set up

Note

A side-by-side kit is available as an accessory via Liebherr Customer Service. (see 7.3 Customer service)

3.11 After setup

- ▶ Pull off the protective film from the outside of the housing.
- ► Clean the appliance. (see 6.3 Cleaning the appliance)
- ► If necessary: Disinfect the appliance.
- Keep the invoice so you have the appliance and dealer information available if needed.

3.12 Disposal of packaging



WARNING

Danger of suffocation from packaging materials and films! ► Do not allow children to play with packaging materials.

The packaging is made from recyclable materials:

- Corrugated card/cardboard
- Parts made of foamed polystyrene
- Films and bags from polyethylene
- Packing bands from polypropylene
- Wood frame nailed together with a polyethylene window*
- ▶ Take the packaging material to an official collection point.



Tools











Fig. 16



WARNING

Risk of injury if the door is not reversed correctly!

▶ The door may only be reversed by qualified personnel.



WARNING

Risk of injury and material damage due to heavy door!

- Only perform the conversion if you can carry a weight of 25 kg.
- ▶ Always have someone help you carry out the conversion.

NOTICE

Live parts!

Damage to electrical components.

- ▶ Disconnect the power plug before reversing the door.
- ► Open the door.

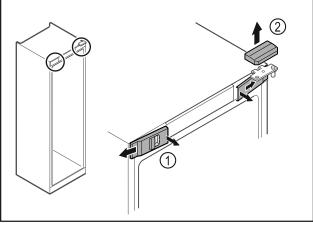


Fig. 17

- ▶ Unlatch the front covers *Fig. 17 (1)* on the inside and remove them sideways.
- Lift off the upper cover Fig. 17 (2).

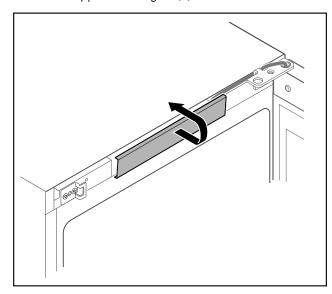


Fig. 18

▶ Unlatch the middle cover and remove it.

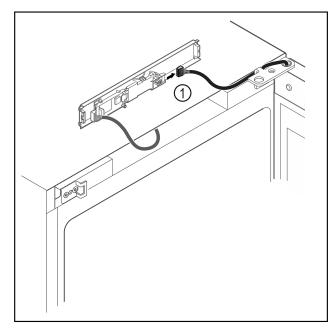


Fig. 19

▶ Disconnect the plug Fig. 19 (1) from the circuit board.

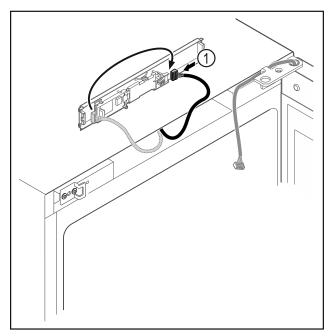


Fig. 20 ► Connect the plug Fig. 20 (1) to the circuit board.

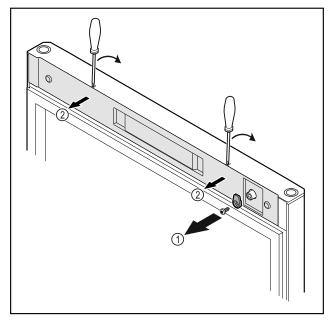


Fig. 21

- Unscrew the door latch Fig. 21 (1).
 Unlatch the cover Fig. 21 (2) with a small screwdriver and remove it.


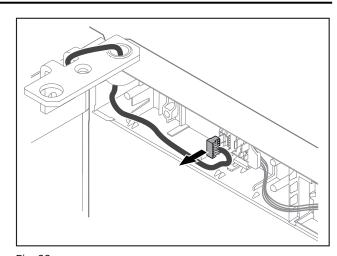


Fig. 22 ► Remove the plug from the plug holder.

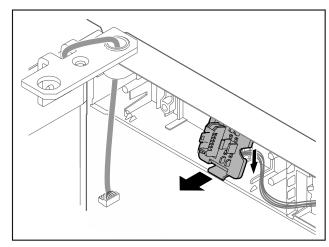

► Unlatch the plug holder downwards.

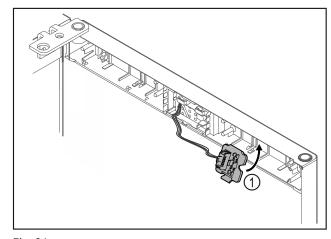


Fig. 24 ▶ Snap in the plug holder Fig. 24 (1) on the opposite side.

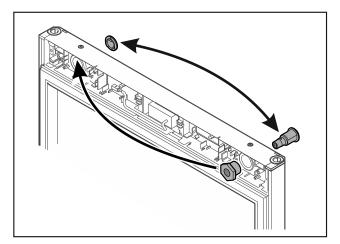


Fig. 25

▶ Put the lock and cover on the opposite side.

Risk of injury if the door tips out!

► Hold the door.

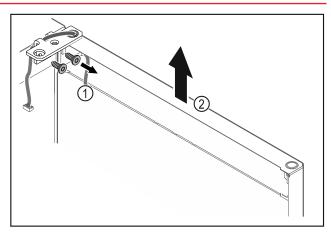


Fig. 26

- Unscrew the hinge Fig. 26 (1).
- Lift the door with the hinge straight up by Fig. 26 (2) roughly 200 mm and take it off.
- ► Carefully place the door on a soft surface.

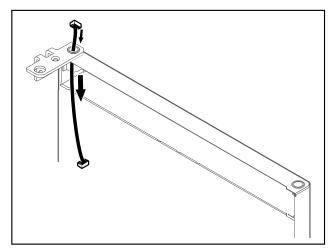


Fig. 27

► Carefully pull out the cable.

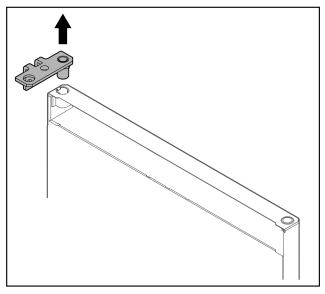


Fig. 28

▶ Pull out the hinge.

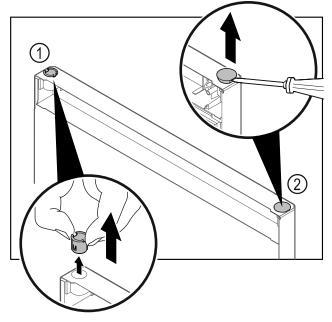


Fig. 29

- Pull out the hinge bushing Fig. 29 (1) with your fingers.
 Carefully lift the cover plug Fig. 29 (2) with a slotted screwdriver and pull it out.


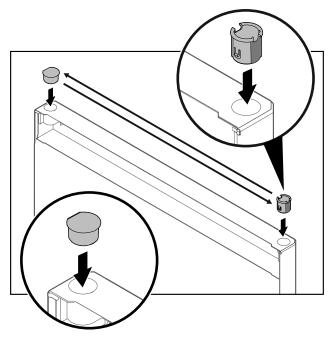


Fig. 30▶ Insert hinge bushing and cover plug on the opposite side (the flattened sides face outwards).

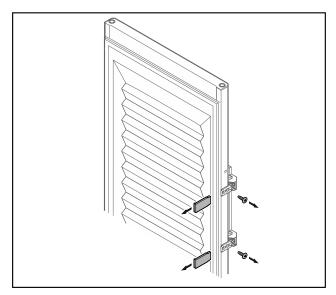

- Fig. 31 Insulated door

 ▶ Remove the panels.
 ▶ Unscrew the handle.

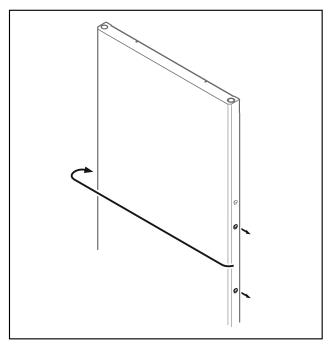


Fig. 32

▶ Put the stopper on the opposite side.

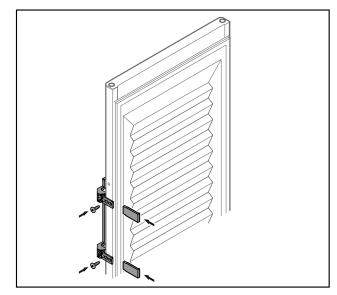


Fig. 33 Insulated door

- Screw the handle onto the opposite side.
 Attach the panels.

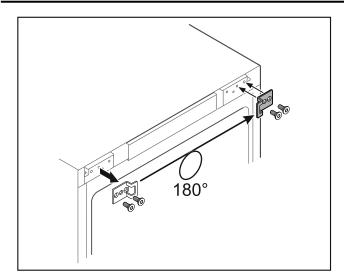


Fig. 34 ▶ Put the closing bracket on the opposite side.

The holes are pre-marked and must be pierced with the selftapping screws.


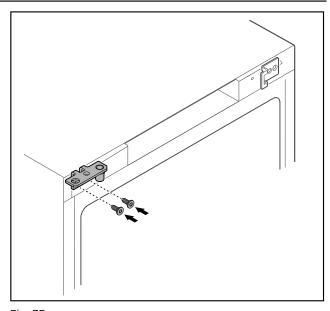


Fig. 35 Move the hinge to the opposite side.



Risk of injury and material damage if the door tips out! ▶ Tighten the bearing pins to the specified torque.

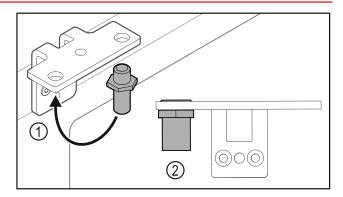

Put the pin back into the hinge Fig. 36 (1).

- ► Tighten the pin Fig. 36 (2) to a torque of 12 Nm.
- ► Unscrew the hinge again.

NOTICE

Risk of injury due to tensioned spring!

▶ Do not disassemble the door closing system Fig. 37 (1).

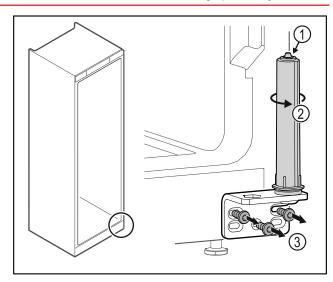


Fig. 37

- Turn the closing system Fig. 37 (2) until it clicks.
- The tension of the closing system is released.
 ▶ Unscrew the hinge Fig. 37 (3).


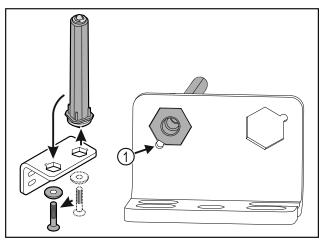


Fig. 38

- Put the door closing system in the hinge. Make sure the pin chamfer Fig. 38 (1) faces the round hole when you put it in.

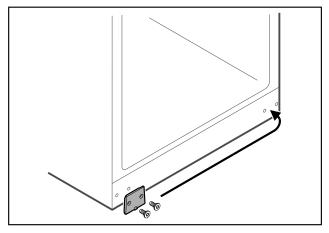

▶ Put the cover plate on the opposite side.

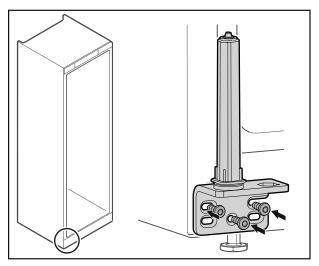


Fig. 40

► Screw the hinge onto the opposite side.

Note

Correct alignment and tension are important for the closing system to work properly.

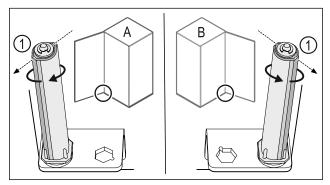


Fig. 41 Left-hinged (A) / Right-hinged (B)

- ▶ Turn the closing system against the resistance until the bar of the closing system *Fig. 41 (1)* points outwards.
- ➤ The closing system automatically stays in this position.
- > The closing system is now aligned and pretensioned.

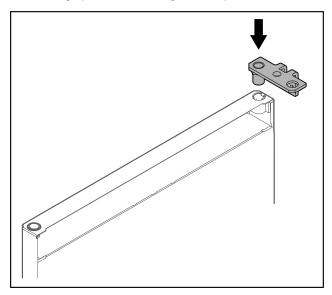


Fig. 42

▶ Put the hinge into the door.

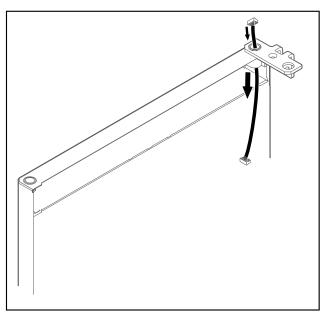


Fig. 43

► Carefully push the cable through.

NOTICE

Risk of injury if the door tips out!

► Hold the door.

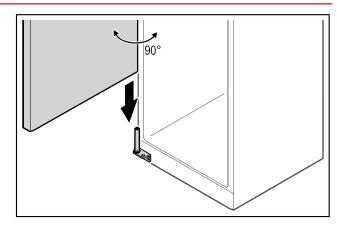



Fig. 44

- ➤ Together with a second person, lift the door from the ground.
- Carefully put the door on the closing system with the door opened at a 90° angle.


NOTICE

Material damage due to incorrect mounting!


▶ Do not pinch the cable when mounting the hinge.

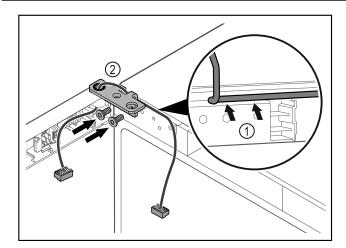

- Fig. 45
 Feed the cable through the recess in the hinge and lay it carefully Fig. 45 (1). Screw on the hinge Fig. 45 (2).

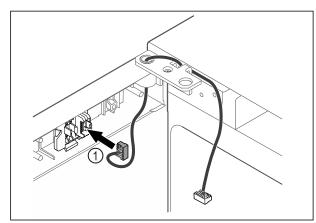

▶ Put the plug Fig. 46 (1) in the plug holder.

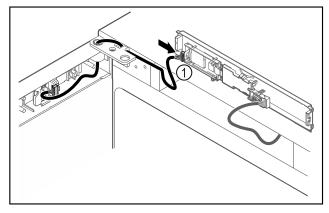

► Connect the plug *Fig. 47 (1)* to the circuit board.

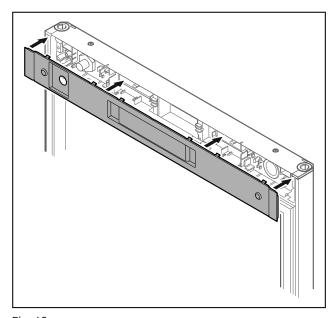


Fig. 48 ▶ Put on the cover.

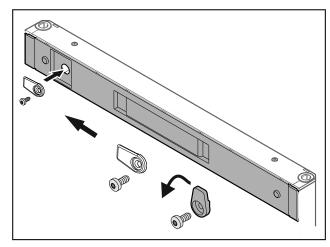


Fig. 49 Screw on the door latch.

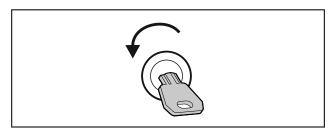


Fig. 50

▶ Open the door.

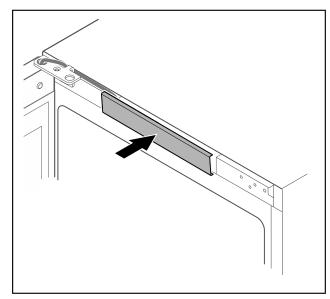


Fig. 51

► Snap in the middle cover.

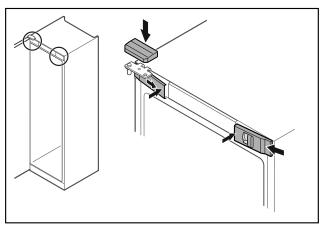


Fig. 52

- ▶ Hook in the front covers Fig. 52 (1) on the side and snap them into place on the inside.
- ▶ Snap on the top cover Fig. 52 (2) from above.
- Close the door.
- The door has now been reversed.

3.14 Aligning the door

If the door is not straight, you can adjust it on the lower hinge.

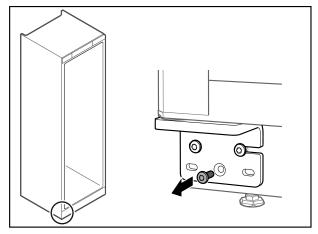


Fig. 53

► Remove the middle screw on the lower hinge.

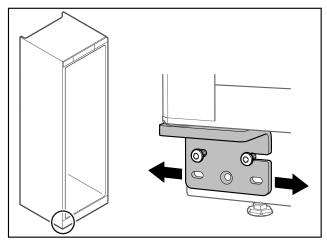


Fig. 54

- ► Slightly undo both screws and move the door with the hinge to the left or right.
- ► Fully tighten the screws (the middle screw is no longer needed).

3.15 Connecting the appliance




WARNING

Incorrect connection!

Risk of fire.

- ▶ Do not use an extension cord.
- ▶ Do not use a multipoint connector strip.

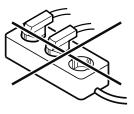

NOTICE

Incorrect connection!

Damage to the electronics.

- ▶ Do not connect the appliance to a stand-alone inverter, e.g. solar power systems and petrol generators.
- Do not use an energy saving plug.

Note

Only use the supplied power cable.

Make sure that the following requirements are fulfilled:

- The appliance is operated using alternating current only.
- The permitted voltage and frequency are printed on the type plate. The type plate location can be seen in the appliance overview chapter. (see 1.2 Overview of appliance and equipment)
- The socket is grounded and fused in accordance with regulations.
- The tripping current for the fuse is between 10 A and 16 $^{\Delta}$
- The socket is easily accessible.
- Check the electrical connection.
- ► Insert the appliance plug on the rear side of the appliance. Ensure that it engages correctly.
- ► Connect the power plug to the power supply.
- ➤ The display switches to the standby symbol.

3.16 Switching on the appliance (first use)

Make sure that the following requirements are fulfilled:

- Appliance is set up and connected.
- All adhesive strips, adhesive and protective films and transport locks are removed from inside and outside the appliance.



Fig. 55 Startup process

The standby symbol flashes until the startup process is complete.

The standby symbol is shown in the display.

If the appliance has been supplied with factory settings, the screen language first needs to be selected when using for the first time.

The following languages are available:

- English
- Spanish
- French
- German
- Dutch
- Danish
- Italian
- Polish
- Czech

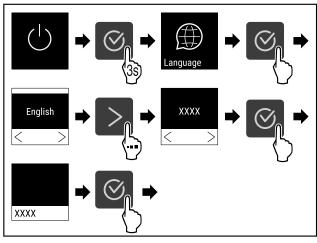


Fig. 56

► Carry out steps (see Fig. 1759).



Fig. 57

▶ The appliance is switched on once the temperature appears on the display.

4 Use

4.1 Control and display elements

The display provides a quick overview of the current appliance status, the temperature setting, the status of functions and settings as well as alarm and error messages.

It is operated using the navigation arrows and confirmation symbol.

Functions can be activated or deactivated and settings values can be changed.



Fig. 58 Display

- (1) Status display
- (2) Backward navigation arrow
- (3) Forward navigation arrow
- (4) Confirm

4.1.1 Status display



Fig. 59 Status display with actual temperature

The status display shows the actual temperature and is the home display. You navigate from here to the other functions and settings. (see 4.2 Navigation) The status display may show a range of display symbols.

4.1.2 Display symbols

The display symbols provide information about the current status of the appliance.

Symbol	Appliance status
	Standby Appliance is switched off.
	Flashing standby symbol Appliance is starting up.
55 °C	Flashing temperature Target temperature not yet reached. Appliance cooling to set temperature.
5,5∘c	Temperature display Displays the current interior temperature.

Symbol	Appliance status
5,5 ⋅c **\(\) Defrost	Display alternating between temperature/defrost symbol with white bar
	Appliance is in manual defrosting mode.
	D in the display Appliance is in DemoMode.
5,5∘c	Appliance is in Bellier load.
Maintenance	Maintenance reminder The time set interval has expired.
n max	Data memory full
min Recording	The data memory (999 hours) is full. From now, the oldest data will be overwritten.
	Flashing symbol A fault is still present.
5,5∘c	A radic is out process.
Error	Error symbol Appliance is still malfunctioning.
	White bar at the bottom Submenu
	White bar at the top Default, active setting or active value.
	Increasing bar Press button for 3 seconds to activate setting.
	Decreasing bar
	Press button for 3 seconds to deactivate setting.

Status display symbols

4.1.3 Acoustic signals

A signal sounds in the following cases:

- If a function or a value is confirmed.
- If a function or a value can neither be activated nor deactivated.
- As soon as a fault occurs.
- If there is an alarm message.

The alarms can be switched on and off in the customer menu.

4.2 Navigation

Navigate in the menu to access the individual functions. The appliance is operated with the buttons next to the display.

4.2.1 Navigating with buttons

9	
Display and button	description
	Forward navigation arrow Advances to next menu (submenu). After the last menu page appears, it goes back to the first menu page.
	Backward navigation arrow Short press: Goes back one menu (submenu). Long press (3 seconds): Returns to the status display.
	Scrolling Pressing multiple times: Scrolls through the menu (submenu).
	Confirm Activate / deactivate function. Open submenu. Confirm selection. The display switches back to the menu.
Back S	Back Display switches back to the function.

4.2.2 Settings menu

The following functions can be accessed in the settings menu:

- Temperature (see 4.3.3 Temperature)
- Temperature alarm (see 4.3.20 Temperature alarm)
- Door alarm (see 4.3.21 Door alarm)
- Alarm simulation (see 4.3.22 Alarm simulation)
- External alarm monitoring (see 4.3.23 Alarm relay)
- Calibration (see 4.3.19 Sensor calibration)
- Maintenance interval reminder (see 4.3.7 Maintenance interval reminder)
- Language (see 4.3.8 Language)
- Temperature unit (see 4.3.9 Temperature unit)
- Display brightness (see 4.3.10 Display brightness)
- LAN connection (see 4.3.14 LAN connection) *
- WiFi connection (see 4.3.13 WiFi connection) *
- Switching off the appliance (see 4.3.2 Switching appliance on and off)

Opening the settings menu

▶ Press the navigation arrow repeatedly until the corresponding function is displayed.



Fig. 60

► Carry out steps (see Fig. 1763).

-or-

- ► Enter access protection PIN code for settings menu. (see 4.3.5 Settings menu access protection)
- > The settings menu is opened. From here you can access all of the above functions.

Settings menu access protection

(see 4.3.5 Settings menu access protection)

4.2.3 Customer menu

The customer menu enables access to the following hidden functions by entering the numerical code **151**:

- Operating hours (see 4.3.16 Operating hours)
- Software version (see 4.3.17 Software)
- Code change for settings menu access protection (see 4.3.6 Access codes)
- Settings menu access protection (see 4.3.5 Settings menu access protection)
- Alarm sounds (see 4.3.11 Alarm Sound)
- Confirmation sounds (see 4.3.12 Key Sound)
- Defrost info (see 4.3.18 Defrosting)
- DemoMode (see 4.3.24 DemoMode)
- Factory reset (see 4.3.25 Factory reset)

Opening the customer menu

- ▶ Open settings menu (see 4.2 Navigation).
- ▶ Press the navigation arrow repeatedly until the corresponding function is displayed.

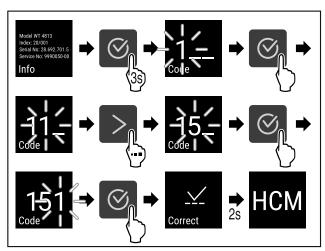



Fig. 61

- ► Carry out steps (see Fig. 1764).
- ➤ The hidden customer menu is opened. From here you can access all the specific menus.

4.2.4 Setting values

Display values can be set in the submenu in combination with the navigation arrow:

Display and button	description
××× >	Increases a value.

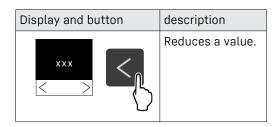

4.2.5 General rules

The following general rules apply in addition to the buttons available:

- Once a value has been selected, the display is shown for 2 seconds before jumping back one level.
- If no selection is made after 10 seconds, the display switches to the status display.
- If a selection is confirmed in a submenu, the display switches back to the menu.

4.3 Functions

4.3.1 Overview of functions



Switching the appliance on and off



Reset logging



Manually starting automatic defrosting



Settings

Access to this function can be blocked using a PIN code. (see 4.3.5 Settings menu access protection)



Info



Temperature



Temperature alarm



Door alarm



Alarm simulation



Alarm forwarding



Sensor calibration



Maintenance reminder interval

Use



Language



Temperature unit



Screen brightness



ΛNI *



WiFi *



Operating hours



Software



Settings lock



Settings lock code



Button sounds



Defrost information



Alarm sounds



DemoMode



Factory reset



4.3.2 Switching appliance on and off

Using this setting the entire appliance can be switched on and off.

Switching on the appliance

Without activated DemoMode:



Fig. 62

► Carry out steps (see Fig. 1765).

With activated DemoMode:

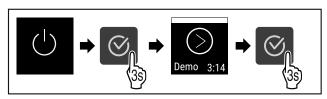


Fig. 63

► Carry out steps (see Fig. 1766).

Note

Deactivate DemoMode before the countdown finishes.



Fig. 64

> The temperature appears on the display.

Switching off the appliance

▶ Press the navigation arrow repeatedly until the corresponding function is displayed.



Fig. 65

- Carry out steps (see Fig. 1768) .
- > The standby symbol appears on the display.

4.3.3 Temperature Temp.

The temperature depends on the following factors:

- How often the door is opened
- How long the door is open for
- The room temperature of the installation site
- The type, temperature and amount of refrigerated items

Note

The temperature may differ from the temperature displayed in some areas of the interior.

At the correct temperature, refrigerated items will keep for longer. This avoids disposing of food unnecessarily.

Setting the temperature

- ► Call up settings menu. (see 4.2 Navigation)
- Press the navigation arrow repeatedly until the corresponding function is displayed.

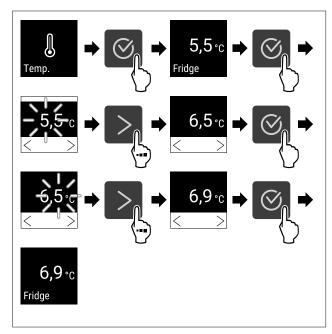


Fig. 66 Changing temperature from 5.5°C to 6.9°C

- ► Carry out steps (see Fig. 1769).



4.3.4 Temperature recording

The appliance displays the minimum and maximum temperatures of the interior. The recording of these temperatures starts automatically after the appliance is switched on; they are recorded at one minute intervals. A note indicating that the data memory is full is displayed after 999 hours (approx. 40 days). The temperature recording should then be reset.

Note

We recommend resetting the temperature recording once after reaching the set temperature when the appliance is commissioned. This ensures that the value displayed for the maximum temperature is a meaningful value.

Displaying temperature recording

The temperature recording displays the length of the recording and the minimum and maximum temperatures measured during this period of time.



Fig. 67

- ▶ Tap navigation arrow repeatedly until corresponding functions are displayed.
- Status screen with the temperature recordings is displayed.

Resetting temperature recording

The displayed minimum and maximum temperatures can be reset at any time. This deletes the displayed values and the recording interval starts again.

 Tap navigation arrow repeatedly until corresponding function is displayed.

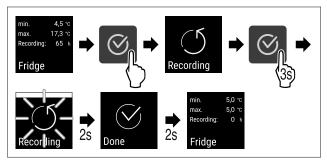


Fig. 68

- ► Complete steps (see Fig. 1771).
- > Temperature recording is reset.

Sett. lock

4.3.5 Settings menu access protection

This setting enables the settings menu access protection via a three-digit PIN code.

Application:

- Prevent settings and functions being changed unintentionally.
- Prevent the appliance from being switched off unintentionally.
- Prevent the temperature from being adjusted unintentionally.

Activating the settings menu access protection

- Open the customer menu. (see 4.2 Navigation)
- ► Tap navigation arrow repeatedly until corresponding function is displayed.

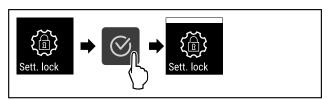


Fig. 69

- Complete steps (see Fig. 1772) .
- > Settings menu access protection is activated.

Deactivating the settings menu access protection

- ▶ Open the customer menu. (see 4.2 Navigation)
- Tap navigation arrow repeatedly until corresponding function is displayed.



Fia. 70

- ► Complete steps (see Fig. 1773).
- > Settings menu access protection is deactivated.

Changing the access protection PIN code for the settings menu

See: (see 4.3.6 Access codes)

Opening protected settings menu

If the access protection is active, it is necessary to enter the PIN code to open the settings menu.

The settings menu remains unlocked for 20 minutes after entering the correct PIN code.

Note

▶ In the following example, the PIN code is: 2 3 4.

Access protection for the settings menu must be active.

► Tap navigation arrow repeatedly until corresponding function is displayed.

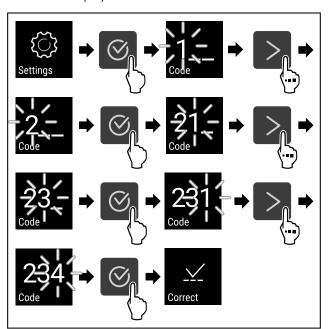


Fig. 71

- ► Complete steps (see Fig. 1774) . Enter PIN code.
- The PIN code is correct.
- ▷ The settings menu opens.

4.3.6 Access codes

Various settings are possible.

Application:

- Changing the settings code.
- Resetting the settings code.



Changing the settings code

This setting allows the settings code for settings menu access protection to be changed.

The setting is made in 3 stages:

- Entering the old settings code
- Entering the new settings code
- Confirming the new settings code

Note

- ▶ In the following example, the default factory-set settings code **1 1 1** is changed.
- ► The new settings code is: 2 3 4
- ➤ Open the customer menu. (see 4.2 Navigation)

 The access protection for the settings menu must be active. (see 4.3.5 Settings menu access protection)
- Press the navigation arrow repeatedly until the corresponding function is displayed.

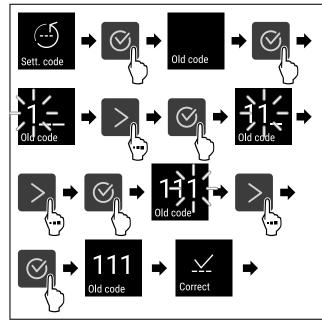


Fig. 72

- Carry out steps (see Fig. 1775).
- Entry of the old settings code successful.

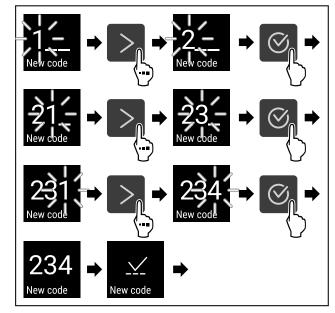


Fig. 73

- ► Carry out steps (see Fig. 1776).
- Entry of the new settings code successful.

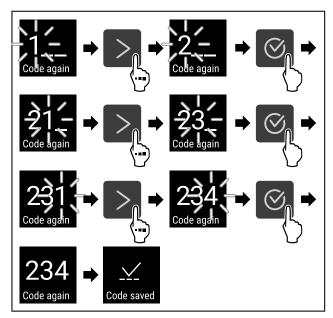


Fig. 74

- ► Carry out steps (see Fig. 1777).
- Confirmation of the new settings code.
- ➤ The settings code has been changed.

Resetting the settings code

The settings code for the settings menu access protection has been forgotten or is not known.

- ► Reset the appliance to factory settings (see 4.3.25 Factory reset).
- > The appliance is reset to the original settings.
- ➤ The factory-set settings code is: 1 1 1



4.3.7 Maintenance interval reminder

Setting for the time interval after which a maintenance reminder is issued.

The following values can be set:

- 7 days
- 14 days
- 30 days
- 60 days
- 90 days
- 180 days
- 360 days
- 720 days
- 1080 days
- Off

Setting maintenance interval reminder

The following steps describe how the maintenance interval is set.

- Call up settings menu. (see 4.2 Navigation)
- ▶ Tap navigation arrow repeatedly until corresponding function is displayed.

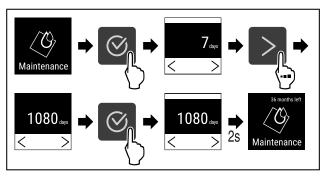


Fig. 75

- Complete steps (see Fig. 1778) .
- The time interval, after which the maintenance reminder is issued, is set.
- > The remaining time is displayed.



This setting allows the display language to be set.

Setting the language

- ▶ Open settings menu (see 4.2 Navigation).
- ▶ Press the navigation arrow repeatedly until the corresponding function is displayed.

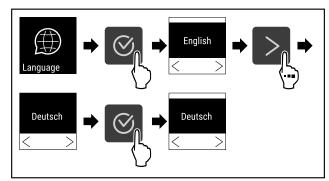


Fig. 76

- ► Complete the steps (see Fig. 1779).
- > The selected language is set.



4.3.9 Temperature unit Temp.

This setting allows the temperature unit to be changed between degrees Celsius and Fahrenheit.

Setting the temperature unit

The example shows how to change the temperature unit from degrees Celsius to Fahrenheit.

- ▶ Open settings menu (see 4.2 Navigation).
- ▶ Press the navigation arrow repeatedly until the corresponding function is displayed.

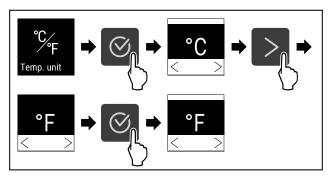


Fig. 77

- ► Complete the steps (see Fig. 1780).
- > The temperature unit is set as Fahrenheit.



4.3.10 Display brightness

This setting allows the display brightness to be adjusted incrementally.

The following brightness values can be set:

- 40%
- 60%
- 80%
- 100% (default)

Setting the display brightness

The following steps describe how to reduce the display brightness, e.g. from . 100 % to 40 %.

- ▶ Open settings menu (see 4.2 Navigation).
- Press the navigation arrow repeatedly until the corresponding function is displayed.

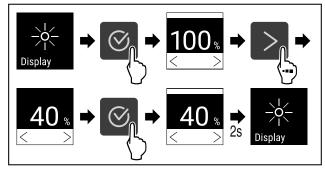


Fig. 78

- ► Complete the steps (see Fig. 1781).
- > The display brightness is changed accordingly.



4.3.11 Alarm Sound

This function enables all alarm sounds, such as the door alarm, to be switched on and off.

Activating Alarm Sound

- ▶ Open the customer menu. (see 4.2 Navigation)
- Press the navigation arrow repeatedly until the corresponding function is displayed.

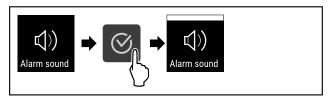


Fig. 79

- ► Complete the steps (see Fig. 1782).

Deactivating Alarm Sound

▶ Open the customer menu. (see 4.2 Navigation)

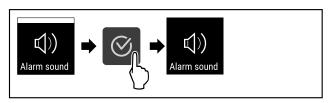


Fig. 80

- ► Complete the steps (see Fig. 1783).
- > Alarm Sound is deactivated.

4.3.12 Key Sound Keys

This function makes it possible to switch all button sounds, confirmation sounds and the Startsound on and off.

Activating Key Sound

- ▶ Open the customer menu. (see 4.2 Navigation)
- Tap navigation arrow repeatedly until corresponding function is displayed.

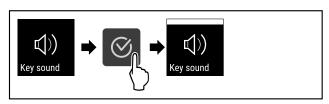


Fig. 81

- ► Complete the steps (see Fig. 1784).

Deactivating Key Sound

▶ Open the customer menu. (see 4.2 Navigation)

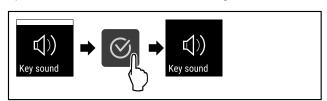


Fig. 82

- ► Complete the steps (see Fig. 1785) .

4.3.13 WiFi connection



Note

Use of the Liebherr SmartMonitoring Dashboard at https://smartmonitoring.liebherr.com requires installation of a SmartModule and a commercial MyLiebherr account. When commissioning online you can register using your login data, or register again and create a company account.

This setting establishes a wireless connection between the appliance and the Internet. The connection is controlled via the SmartModule. The appliance can be integrated via the browser-basedLiebherr SmartMonitoring Dashboard and advanced options and customized methods relating to control, administration and monitoring can be used.

Note

Accessories are available via Customer Service (see 7.3 Customer service) and in the Liebherr-Hausgeräte-Shop at https://home.liebherr.com.

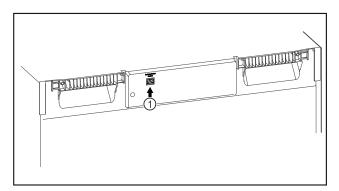


Fig. 83

Make sure that the following requirements are fulfilled:

☐ SmartModule Fig. 83 (1) is in use.

Establishing a connection

You commission and setup your SmartModule online via the Liebherr SmartMonitoring Dashboard on your web-enabled device.

On the Liebherr SmartMonitoring Dashboard information is also available about retrofitting the SmartModule.



Fig. 84

▶ Open the Liebherr SmartMonitoring Dashboard (see Fig. 1787) .

On the fridge or freezer:

- ▶ Open settings menu (see 4.2 Navigation).
- Press the navigation arrow repeatedly until the corresponding function is displayed.

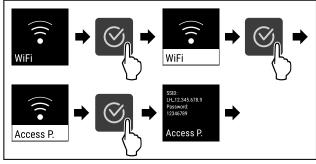


Fig. 85

- ► Carry out steps (see Fig. 1788).
- ► Continue the set-up procedure on your web-enabled device: Liebherr SmartMonitoring Dashboard



Fig. 86

- ► The connection is set up (see Fig. 1789).
- ▶ WiFi connecting appears. The symbol flashes.

- ▶ Follow the Liebherr SmartMonitoring Dashboard instructions.
- > Connection is established.

Disconnecting

- ▶ Open settings menu (see 4.2 Navigation).
- Press the navigation arrow repeatedly until the corresponding function is displayed.

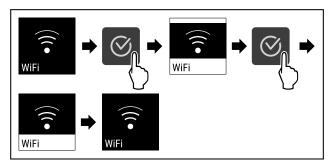


Fig. 87

- ► Carry out steps (see Fig. 1790).
- > There is no longer a connection.

Resetting connection

- ▶ Open settings menu (see 4.2 Navigation).
- Press the navigation arrow repeatedly until the corresponding function is displayed.

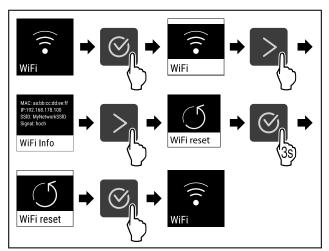


Fig. 88

- ► Carry out steps (see Fig. 1791).
- The WiFi settings are restored to the factory settings.

4.3.14 LAN connection LAN

Note

Use of the Liebherr SmartMonitoring Dashboard at https://smartmonitoring.liebherr.com requires installation of a SmartModule and a commercial MyLiebherr account. When commissioning online you can register using your login data, or register again and create a company account.

This setting establishes a wired connection between the appliance and the Internet. The connection is controlled via the SmartModule. The appliance can be integrated via the browser-based Liebherr SmartMonitoring Dashboard, and advanced options and customized methods relating to control, administration and monitoring can be used.

Note

Accessories are available via Customer Service (see 7.3 Customer service) and in the Liebherr-Hausgeräte-Shop at https://home.liebherr.com.

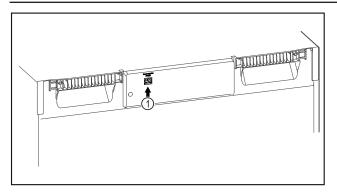


Fig. 89

Make sure that the following requirements are fulfilled:

- ☐ SmartModule Fig. 89 (1) is in use.
- ☐ A network cable is connected.

Establishing a connection

You commission and setup your SmartModule online via the Liebherr SmartMonitoring Dashboard on your web-enabled device.

On the Liebherr SmartMonitoring Dashboard information is also available about retrofitting the SmartModule.



Fig. 90

▶ Open the Liebherr SmartMonitoring Dashboard (see Fig. 1793).

On the fridge or freezer:

- ▶ Open settings menu. (see 4.2 Navigation)
- Press the navigation arrow repeatedly until the corresponding function is displayed.

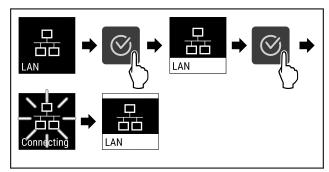


Fig. 91

- ► Carry out steps (see Fig. 1794).
- Connection is established: LAN connecting appears. The symbol flashes.
- ▶ Follow the Liebherr SmartMonitoring Dashboard instructions.
- December Connection is established.

Disconnecting

- ▶ Open settings menu. (see 4.2 Navigation)
- Press the navigation arrow repeatedly until the corresponding function is displayed.

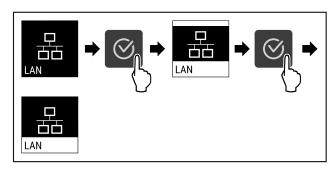


Fig. 92

- ► Carry out steps (see Fig. 1795).
- > There is no longer a connection.



4.3.15 Info

This display shows the model, serial number and service number of the appliance.

Displaying Info

- ▶ Open settings menu (see 4.2 Navigation).
- Press the navigation arrow repeatedly until the corresponding function is displayed.



Fig. 93

> The appliance information appears.



4.3.16 Operating hours

The display shows the appliance operating hours.

Displaying operating hours

► Open the customer menu.



Fig. 94

> The operating hours appear.



4.3.17 Software

The display shows the software version of the appliance.

Displaying the software version

- ▶ Open the customer menu .
- ▶ Press the navigation arrow repeatedly until the corresponding function is displayed.



Fig. 95

> The software version is shown.



The appliance defrosts automatically in normal mode.

Application:

 If there is excessive ice in the interior, the defrosting function can be started manually.

Manually starting automatic defrosting

▶ Tap navigation arrow repeatedly until corresponding function is displayed.

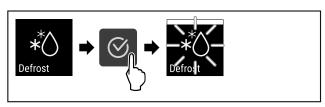


Fig. 96

- ► Complete steps (see Fig. 1799).
- Automatic defrosting has been started. The appliance automatically switches back to normal mode once defrosting has finished.

Canceling manually started defrosting

► Tap navigation arrow repeatedly until corresponding function is displayed.

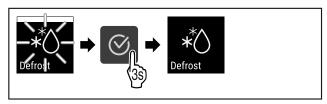


Fig. 97

- ► Complete steps (see Fig. 1800).
- Manually started defrosting is canceled. The appliance automatically switches back to normal mode.



Defrosting info

This setting adjusts the display during an automatic defrosting process.

The following display variants are available:

- Def 1: "Automatic defrosting process" note is displayed in the status screen. The temperature display is not updated and shows the last temperature before the automatic defrosting process was started.
- Def 2: "Automatic defrosting process" note is displayed in the status screen. The temperature display is updated continuously and shows the actual temperature.
- Def 3: The temperature display is not updated and shows the last temperature before the automatic defrosting process was started.
- Def 4: The temperature display is updated continuously and shows the actual temperature.

Defrosting Fig. 98 "Automatic defrosting process" note

- ▶ Open the customer menu. (see 4.2 Navigation)
- ▶ Tap navigation arrow repeatedly until corresponding function is displayed.

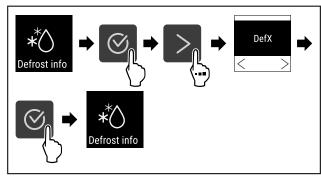


Fig. 99

- ► Complete steps (see Fig. 1802).
- > The desired display variant has been selected.



4.3.19 Sensor calibration

The sensor calibration is used to realign differences between the temperature displayed and the actual temperature in the appliance.

Setting range: +/- 3 Kelvin in 0.1 Kelvin increments.

Calibrating the sensor

- ► Call up settings menu. (see 4.2 Navigation)
- Tap navigation arrow repeatedly until corresponding function is displayed.

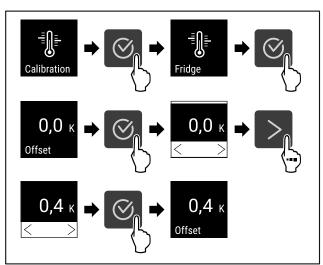


Fig. 100

- ► Complete steps (see Fig. 1803).
- > The sensor is calibrated.

Fridge

4.3.20 Temperature alarm

The temperature alarm is triggered as soon as the interior temperature departs from the settable temperature range.

The following values can be set:

- Lower temperature limit in 0.1 °C increments
- Upper temperature limit in 0.1 °C increments
- Alarm delay time of between 0 and 60 minutes
- Alarm repetition frequency of between 0 and 30 minutes

Setting temperature alarm

- ► Call up settings menu. (see 4.2 Navigation)
- ▶ Tap navigation arrow repeatedly until corresponding function is displayed.



Fig. 101

- ► Complete steps (see Fig. 1804).
- ► Tap navigation arrow repeatedly until corresponding function is displayed.

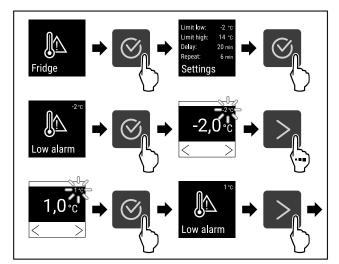


Fig. 102

- ► Complete steps (see Fig. 1805).

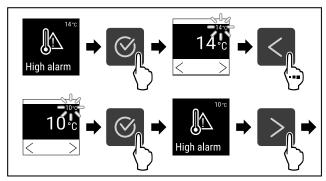


Fig. 103

- ► Complete steps (see Fig. 1806).
- □ Upper temperature limit has been set.

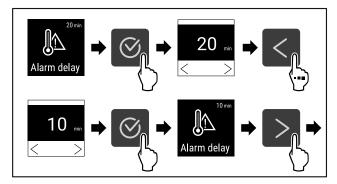


Fig. 104

- ► Complete steps (see Fig. 1807).

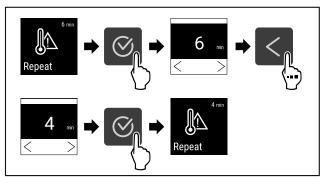


Fig. 105

- ► Complete steps (see Fig. 1808).
- ▷ Temperature alarm has been set.

Displaying set temperature alarm values

- ► Open settings menu (see 4.2 Navigation).
- ► Tap navigation arrow repeatedly until corresponding function is displayed.



Fig. 106

- Complete steps (see Fig. 1809) .
- ► Tap navigation arrow repeatedly until corresponding function is displayed.

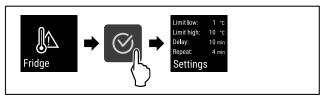


Fig. 107

- ► Complete steps (see Fig. 1810).
- > Set temperature alarm values are displayed.



4.3.21 Door alarm Door alarm

Set the time interval after which the door alarm sounds when the door is opened.

The following values can be set:

- 15 seconds
- 30 seconds
- 60 seconds
- 90 seconds
- 120 seconds
- 150 seconds180 seconds
- Off

Setting the door alarm

The following steps describe how to set the time interval before the door alarm sounds.

- ▶ Open settings menu (see 4.2 Navigation).
- Press the navigation arrow repeatedly until the corresponding function is displayed.



Fig. 108

- Carry out steps (see Fig. 1811) .
- Press the navigation arrow repeatedly until the corresponding function is displayed.

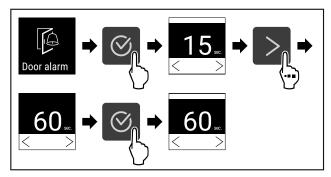


Fig. 109

- ► Carry out steps (see Fig. 1812).
- > The door alarm is set.

Deactivating the door alarm

The following steps describe how to deactivate the door alarm.

- ▶ Open settings menu (see 4.2 Navigation) .
- Press the navigation arrow repeatedly until the corresponding function is displayed.



Fig. 110

- ► Carry out steps (see Fig. 1813).
- Press the navigation arrow repeatedly until the corresponding function is displayed.

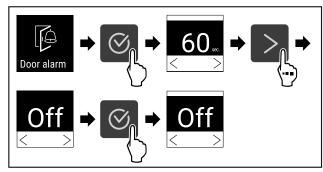


Fig. 111

- ► Carry out steps (see Fig. 1814).
- ➤ The door alarm is deactivated.

4.3.22 Alarm simulation

This setting makes it possible to simulate a change in the appliance temperature to check the function of any safety devices that may be connected.


Starting alarm simulation

- ► Call up settings menu. (see 4.2 Navigation)
- Tap navigation arrow repeatedly until corresponding function is displayed.



Fig. 112

- ► Complete steps (see Fig. 1815).
- ► Tap navigation arrow repeatedly until corresponding function is displayed.

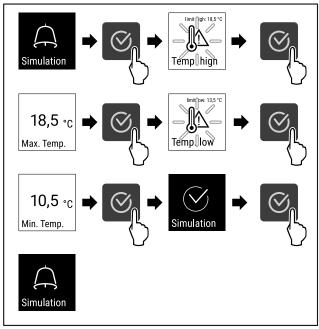

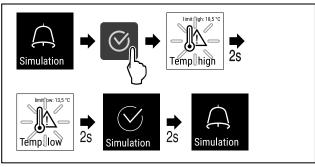


Fig. 113

- ► Complete steps (see Fig. 1816).
- > The simulation has been performed.
- Connected safety devices should have triggered.

4.3.23 Alarm relay

This setting makes it possible to relay one or more alarms to an external alarm source connected to the appliance.

The following settings can be selected:

- Alarm type:
 - Door alarm
 - Temperature alarm
 - Door alarm and temperature alarm
 - All

- Confirm:
 - Yes: Alarm is stopped on the appliance and on the externally connected appliance.
 - No: Alarm is acknowledged on the appliance and remains active on the external appliance until the error is remedied.
- Update (only visible if Yes is selected in response to Confirm):
 - Yes: Alarm is recurring according to the time set in (see 4.3.21 Door alarm)
 - No: Alarm is continuously acknowledged.

Activating alarm relay

- ► Call up settings menu. (see 4.2 Navigation)
- ► Tap navigation arrow repeatedly until corresponding function is displayed.

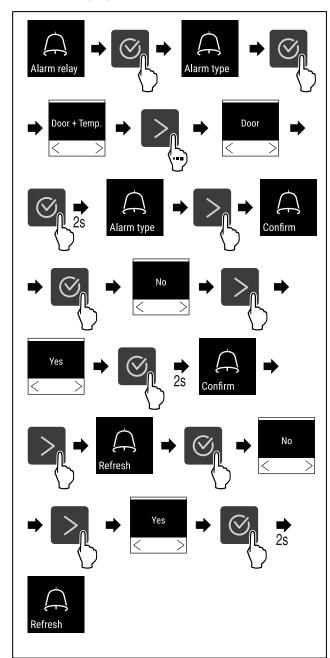


Fig. 114

- ► Complete steps (see Fig. 1817).
- > The alarm relay has been set.

4.3.24 DemoMode



If the appliance is in DemoMode all refrigeration functions are deactivated. This is useful, for example, for product demonstrations.

Activating DemoMode

- ▶ Open the customer menu .
- ▶ Press the navigation arrow repeatedly until the corresponding function is displayed.

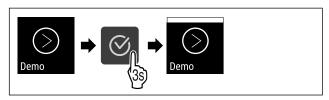


Fig. 115

► Carry out steps (see Fig. 1818).



Fig. 116

- DemoMode is activated.
- > The status display is shown with a "D" in the top left.

Deactivating DemoMode

- ▶ Open the customer menu .
- Press the navigation arrow repeatedly until the corresponding function is displayed.

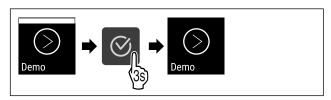


Fig. 117

► Carry out steps (see Fig. 1820).



Fig. 118

- DemoMode is deactivated.
- > The status display appears.

4.3.25 Factory reset Factory res

This function resets the appliance to the factory settings. All settings made up to this point are reset to their original values.

Performing factory reset

- ▶ Open the customer menu.
- Press the navigation arrow repeatedly until the corresponding function is displayed.



Fig. 119

- ► Complete steps (see Fig. 1822) . ▷ Factory reset complete. ▷ The appliance restarts.

4.4 Messages

4.4.1 Warnings

Warnings are issued by means of an audio signal and visually via a symbol on the display. The signal gets louder until the warning is acknowledged.

Message	Cause	Remedy
Door open Door open	The message appears if the door is open for too long.	 Press confirmation button. Alarm stops. Note You can set how long it takes for this message to appear (see 4.3.21 Door
Temperature Temperature alarm	This message appears if the temperature does not match the set temperature. Reasons for the differences in temperature can be: - Warm items for refrigeration have been placed inside. - Too much warm ambient air entering when rearranging and removing food. - The power was cut off for a prolonged period.	Press confirmation button. The highest temperature is shown. Press confirmation button. The current temperature flashes and the display shows the temperature alarm symbol until the set temperature is reached. Check the condition of the refrigerated items.

Message	Cause	Remedy
Error	The appliance is malfunctioning, there is an appliance error or a component of the appliance has a fault.	• Store the refrigerated items elsewhere.
		➤ Press confirmation button. XXXXX ⊗ Error code is displayed. Error code is displayed.
		 Press confirmation button. A further error code is displayed should there be one. or Status screen is displayed with flashing 5,0 °c arrow symbol Pressing the navigation buttons displays the error code(s) again. Make a note of error code(s) and contact Customer Service.
Power Failure	The message appears following an interruption to the electricity supply.	 Press confirmation button. The highest temperature and number of electrical interruptions are displayed. Press confirmation button. Alarm is stopped and current temperature is displayed. Check the condition of the refrigerated items.
*\times_Defrost error Defrost error	The automatic defrosting is not working properly. The previous 5 defrost cycles were not successful.	 ▶ Press confirmation button. ▶ Alarm stops. ▶ Contact Customer Service (see 7.3 Customer service) .

(X) WiFi error WiFi error	WiFi connection is interrupted.	 Check connection. Press confirmation button. Alarm stops.
WWW error WWW error, WiFi	This message appears if there is no Internet connection via WiFi.	 Check connection. Press confirmation button. Alarm stops.
LAN error	LAN connection is interrupted.	 Check connection and network cable. Press confirmation button. Alarm stops.
WWW error WWW error, LAN	This message appears if there is no Internet connection via WiFi.	➤ Check connection. ➤ Press confirmation button. ➤ Alarm stops.

4.4.2 Reminders

Reminders appear when you are prompted to take action. They are issued by means of an audio signal and visually via a symbol on the display. Acknowledge the message by pressing the confirmation button.

Message	Cause	Remedy
Maintenance Maintenance reminder	The message appears when the timer for the set maintenance interval has expired.	Press confirmation button.Alarm stops.
		Note
		The maintenance interval can be set (see 4.3.7 Maintenance interval reminder) .

Equipment



Logging reminder

This message appears when the temperature logging memory (999 hours) is full.



- ▶ Press confirmation button.
- ▶ Alarm stops.
- ▶ The oldest data is overwritten.

Note

The temperature logging can be reset (see 4.3.4 Temperature recording).

5 Equipment

5.1 Safety lock

The lock in the appliance door is equipped with a safety mechanism.

5.1.1 Locking appliance



Fig. 120

- Insert key.
- ► Turn key anticlockwise by 90°.
- \triangleright The appliance is locked.

5.1.2 Unlocking appliance



Fig. 121

- ► Insert key Fig. 121 (1).
- ► Turn key clockwise by 90° Fig. 121 (2).
- \triangleright The appliance is unlocked.

5.2 Sensors

The appliance can be equipped with the following sensors for temperature monitoring:

- Safety thermostat sensor
- C sensor

5.2.1 Safety thermostat sensor

The safety thermostat sensor monitors the temperature drop and thus protects against excessively low temperatures (Freeze Protect). If the temperature drops below the threshold value, the appliance regulates the temperature automatically. The error message FRZ. PROT (Freeze Protect) (see 4.4 Messages) appears.

5.2.2 C sensor (control sensor)

The C sensor is permanently installed and regulates the interior temperature of the appliance. You can calibrate the C sensor if necessary. (see 4.3.19 Sensor calibration)

5.3 Interfaces

There are various connection options on the back of the appliance.

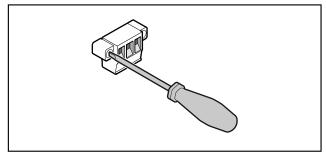


Fig. 122

Note

Screws secure the plugs. To enable disconnection of the plugs, loosen the screws on the left and right.

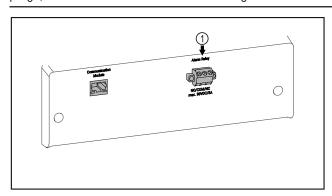


Fig. 123

(1) Potential-free alarm output

5.3.1 Potential-free alarm output



WARNING

Risk of injury due to electric shock!

► The appliance may only be connected to an external alarm device by trained specialist personnel.

These three contacts can be used to connect a visual or an acoustic alarm device. The connection is designed for a maximum of **30 V (DC) - 8 A** from a safety low voltage source SELV (minimum current **150 mA**).

Note

The safety requirements of standard EN 61010 are not met if mains voltage is used at the potential-free alarm contact.

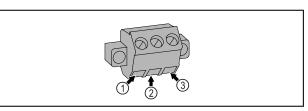


Fig. 124

- (1) NO (normally open) Connection of an alarm indicator lamp or an acoustic alarm signaling device.
- (2) COM (common)
 External voltage
 source
- (3) NC (normally closed)
 Connection of an indicator lamp that displays
 the appliance's normal
 mode.

!\

WARNING

Short-circuit hazard due to live parts! Electric shock or damage to the electronics.

- ► Switch off the appliance.
- ▶ Pull out mains plug or switch off fuse.

Note

A maintenance reminder function can be activated. (see 4.3.7 Maintenance interval reminder)

Note

We recommend creating a maintenance log in which all work (repairs, checks) performed is documented.

Note

Liebherr recommends yearly maintenance. For an individual offer, please contact Customer Service (see 7.3 Customer service) .

6 Maintenance

6.1 Maintenance schedule

Maintenance work is to be performed at regular intervals to ensure proper appliance function.



CAUTION

Danger of injury and damage due to unprofessional maintenance work!

Personal injury and material damage.

► Maintenance work may only be performed by trained specialist personnel.

Component	Activity	Maintenance interval
Sheet metal parts	Check for damage and corrosion. Replacement of damaged parts: (see 7.3 Customer	Yearly
Door, hinges	check alignment (see 3.14 Aligning the door), ease of movement, and tight fit.	Yearly
Lock, door handle	Check for ease of movement and tight fit.	Yearly
Door seal	Check for damage, wear, and tight fit. A replacement door seal is available from customer service.	Yearly
Sensor	Calibrate sensor. (see 4.3.19 Sensor calibration)	Yearly
Sensor feedthrough	Check that the sealing compound is intact.	Yearly
Battery	Replace appliance battery. A replacement battery is available from customer service.	5 to 10 years
Surfaces	Cleaning (see 6.3 Cleaning the appliance)	Recommendation: monthly or as required/specified at installation location

Maintenance

Component	Activity	Maintenance interval
Surfaces	Disinfection Liebherr has tested the following disinfectants: Dismozon pur 1 % Lösung (Bode Chemie), Suma Quicksan (Diversey), Incidin Extra N (Ecolab Healthcare), Acrylan (Antiseptica chempharm. Produkte), Buraton 10 F (Schülke und Mayr), Frankocid N (Franken Chemie), Apesin DSR 50 (Tana), Nüscosept Spray (Dr. Nüsken), Melsept SF (B. Braun Melsungen), Kohrsolin (Bode Chemie), Neoquat S (Dr. Weigert), Indicin Rapid (Ecolab Healthcare), Bacillocid Spezial (Bode Chemie), Neoform K Spray (Dr. Weigert), Apesin Desinf. Spray (Tana), Nüscosept 100 (Dr. Nüsken), Antisept T (Fink Tec), Apesin AP 100 0,50% Lösung (Tana), Perform Pulver 2% Lösung (Schülke und Mayr) Observe the operating instructions of the respective manufacturer. If disinfectants other than those stated are used, test them on a less exposed area first.	Recommendation: monthly or as required/specified at installation location
Installation location, appliance	Decontamination The appliance is suitable for room decontamination with hydrogen peroxide H ₂ O ₂ . In this context, the values of the following application specifications must not be exceeded: Concentration: maximum 250 ppm. Humidity: maximum 85%. Maximum permissible room temperature as per climate class. (see 1.4 Range of appliance use)	Recommendation: as required/specified at installation location, maximum 2x per year
Fan air slots	Cleaning (see 6.3.3 Cleaning fan cover)	Yearly
Alarm relay and sensor	Check plug connections for tight fit.	Yearly
Power cable	Check for damage.	Yearly or if there is a change of location
Tilt angle	Check for proper function and tight fit.	Yearly or if there is a change of location
Adjusting feet/castors	Check for proper function, damage, and tight fit. Replace damaged parts. (see 7.3 Customer service)	Yearly
Defrost water collection tray	Check the fill level of the defrost water collection tray and empty if necessary.	Regular visual inspection

6.2 Defrosting the appliance



WARNING

Appliance incorrectly defrosted! Injuries and damage.

- ▶ Do not use mechanical devices or other means to accelerate the defrosting process, other than those recommended by the manufacturer.
- ▶ Do not use any electrical heating or steam cleaning equipment, naked flames or defrosting sprays.
- ▶ Do not remove ice with sharp objects.

Defrosting is performed automatically. Defrost water drips into the defrost water collection tray and must be emptied at regular intervals.

Note

To avoid having to empty the defrost water collection tray frequently, you can place a container under the drain opening in the tray.

The defrost water collection tray may only be inserted on the lowest support rib.

If the temperature is set very cold and/or the humidity is very high, the rear wall in the interior may ice up.

Manual defrosting:

- ► Switch off appliance. (see 4.3.2 Switching appliance on and off)
- ► Disconnect power plug.
- ► Store refrigerated items elsewhere.
- Leave the appliance door open during defrosting process.
- ▶ Pull out the defrost water collection tray at the front, empty it and clean the appliance.

6.3 Cleaning the appliance

6.3.1 Preparing



WARNING

Short-circuit hazard due to live parts! Electric shock or damage to the electronics.

- ► Switch off the appliance.
- ▶ Pull out mains plug or switch off fuse.



WARNING

Danger of fire

- ▶ Do not damage the refrigerant circuit.
- Empty the appliance.
- ▶ Pull out the power plug.

6.3.2 Cleaning the housing

NOTICE

Improper cleaning!

Damage to the appliance.

- Only use soft cleaning cloths and ph-neutral all-purpose cleaners.
- ▶ Do not use steel wool or sponges that scour or scratch.
- ► Do not use caustic or abrasive cleaning materials or those containing sand, chloride, or acids.



WARNING

Risk of injury or damage due to hot steam.

Hot steam can cause scalding/burns and damage to surfaces.

► Do not use steam cleaners.

NOTICE

Risk of damage due to short circuit.

- When cleaning the appliance, make sure no water gets into the electrical components.
- ▶ Wipe the housing down with soft, clean cloth. If very dirty, use lukewarm water with a neutral cleaner. Glass surfaces can also be cleaned with glass cleaner.
- Clean the condenser coil every year. If the condenser coil is not cleaned, this significantly reduces the efficiency of the appliance.

6.3.3 Cleaning fan cover

NOTICE

Risk of damage due to short circuit.

► When cleaning the appliance, make sure no water gets into the electrical components.

NOTICE

Improper cleaning!

Damage to the appliance.

- Only use soft cleaning cloths and ph-neutral all-purpose cleaners.
- ▶ Do not use steel wool or sponges that scour or scratch.
- ► Do not use caustic or abrasive cleaning materials or those containing sand, chloride, or acids.

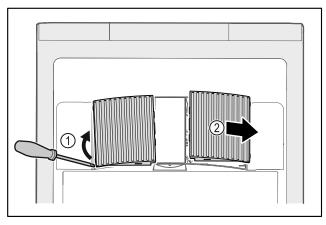


Fig. 125

- ► Carefully lever out fan covers from the side Fig. 125 (1) and remove towards the front Fig. 125 (2).
- Rinse the air slots of the fan covers under running water. If very dirty, use lukewarm water with a neutral cleaner.
- Leave fan covers to air-dry completely.
- Reattach fan covers to the appliance.

6.3.4 Cleaning the interior



WARNING

Short-circuit hazard due to live parts! Electric shock or damage to the electronics.

- Switch off the appliance.
- ▶ Pull out mains plug or switch off fuse.



WARNING

Risk of injury or damage due to hot steam.

Hot steam can cause scalding/burns and damage to surfaces.

▶ Do not use steam cleaners.

NOTICE

Improper cleaning!

Damage to the appliance.

- Only use soft cleaning cloths and ph-neutral all-purpose cleaners.
- ▶ Do not use steel wool or sponges that scour or scratch.
- ▶ Do not use caustic or abrasive cleaning materials or those containing sand, chloride, or acids.

Note

Defrost water drain (see 1.2 Overview of appliance and equipment): Remove deposits using a thin object (e.g. a cotton bud).

Note

Do not damage or remove the type plate on the inside of the appliance. The type plate is important for after sales service.

- ▶ Open the door.
- ► Empty the appliance.
- ► Clean the interior and equipment parts with lukewarm water and a little dish detergent. Do not use any gritty or acidic cleaning agent or any chemical solvent.

6.3.5 After cleaning

- ▶ Wipe the appliance and equipment parts dry.
- Connect and switch on appliance.

When the temperature is sufficiently cold:

Place in items for refrigeration.

Customer support

- Clean regularly.
- ► Clean and dust appliances with a heat exchanger (metal grill on rear of appliance) once per year.

7 Customer support

7.1 Technical data

Temperature range		
Refrigeration	+3 to +16 °C	

Maximum load per shelf ^{A)}			
Appliance width (see 3.2 Appliance dimensions)	Grid shelves	Glass shelves	Wooden grid shelf
600 mm	45 kg	40 kg	45 kg
750 mm	60 kg	40 kg	60 kg

 $^{^{\}mbox{\scriptsize A}\mbox{\scriptsize)}}$ Heavier loads may result in a slight bending of the shelves.

Maximum net capacity

See net capacity on the type plate (see 1.2 Overview of appliance and equipment).

Setup conditions			
Degrees (surroundi		pollution appliance)	

Setup conditions	
Ambient temperature	+10 °C to +35 °C
Maximum relative ambient moisture	75%, not condensing
Maximum operating elevation (meters above sea level)	2000 m / 1500 m (glass door)

Electrical values	
Nominal voltage	220 - 240 V ~
Maximum power supply fluctuation	+/-10%
Overvoltage categories	II (OVC II)

7.2 Technical malfunction

Your appliance is designed and built to ensure it works reliably and has a long service life. If a malfunction nonetheless occurs during operation, please check whether the malfunction is due to an operating error. If this is the case, you will be charged for the cost incurred even if this falls within the warranty period.

You can rectify the following malfunctions yourself.

7.2.1 Appliance function

Defect	Cause	Remedy
The appliance is	→ The appliance is not switched on.	► Switch on the appliance.
not working.	→ The power plug is not properly inserted in the socket.	► Check the power plug.
	→ There is a problem with the wall socket breaker.	► Check the breaker.
	→ Power failure	 ▶ Keep the appliance closed. ▶ If necessary move the refrigerated/frozen items to another fridge or freezer if the power is off for a prolonged period.*
	→ The appliance power plug is not correctly inserted in the appliance.	► Check the appliance power plug.
The temperature is not cold enough.	→ The appliance door is not closed properly.	► Close the appliance door.
	→ Ventilation is not sufficient.	▶ Unclog the ventilation grill and clean it.
	→ The ambient temperature is too high.	Observe the suitable ambient conditions:
	→ The appliance was opened too many times or for too long.	► Wait to see if the required temperature is achieved once again on its own. If not, contact Customer Service (see 7.3 Customer service) .
	→ The temperature is set incorrectly.	▶ Set a colder temperature and check after 24 hours.
	→ The appliance is too close to a heat source (stove, heater etc).	► Move either the appliance or the heat source.

7.3 Customer service

First check whether you can remedy the fault yourself . If this is not the case, please contact Customer Service.

Please refer to the enclosed "Liebherr Service" brochure for the address.



WARNING

Unprofessional repair! Injuries.

- ▶ Repairs and work on the appliance and the power supply cable not described in the Manual (see 6 Maintenance) should only be carried out by a qualified service provider.
- A damaged power cable may only be replaced by the manufacturer, the manufacturer's Customer Service or a similarly qualified person.
- For appliances with an IEC connector, the change can be made by the customer.

7.3.1 Contacting customer service

Make sure you have the following appliance information ready:

- Appliance name (model and index)
- ☐ Service no. (service)
- Serial no. (S no.)
- Access the appliance information via the display (see 4.3.15 Info) .

-or-

- Refer to the type plate for appliance information. (see 7.4 Type plate)
- Note appliance information.
- Notify Customer Service: Report faults and the appliance
- This will help us to provide fast and focused service.
- Follow any further instructions provided by Customer Service.

7.4 Type plate

The type plate is inside the appliance. See appliance overview

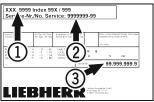


Fig. 126

- (1) Appliance designation
- (3) Serial No.
- (2) Service No.

8 Shutting down

- Empty the appliance.
- Switch off appliance.
- Pull out the power plug.
- Remove appliance connector if required: remove and at the same time move from left to right.
- Clean the appliance. (see 6.3 Cleaning the appliance)
- ▶ Leave the door open so that no bad odors form.

Damage to the appliance and malfunctions!

After shutting down, only store the appliance at the permitted room temperature of between -10 °C and +50 °C.

9 Disposal

9.1 Preparing appliance for disposal



Liebherr uses batteries in some of its appliances. In the EU, for environmental reasons, legislation obliges the end user to remove these batteries before disposing of the old appliances. If your appliance contains batteries, a corresponding notice is attached to the appliance.

Lamps

If you can remove the lamps yourself without destroying them, also remove these before disposal.

- Take the appliance out of operation.
- Appliance with batteries: Remove the batteries. For description, see Maintenance chapter.
- If possible: Remove lamps without destroying them.

9.2 Disposing of the appliance in an environmentally friendly manner



The appliance still contains valuable materials and must be collected separately from unsorted municipal waste so it can be recycled.





Dispose of batteries separately from the old appliance. You can by this returning batteries free of charge to retailers or taking them to recycling centers and collection centers for recyclable materials.

Lamps

For Germany:

Dispose of removed lamps via respective collection systems.

You can dispose of the appliance free of charge via the class 1 collection containers at the local recycling and recyclable material collection centers. When purchasing a new fridge/ freezer, if the sales area > 400 m², retailers will also take back the old appliance free of charge.

WARNING

Leaking refrigerant and oil!

Fire. The refrigerant contained within the appliance is environmentally friendly, but flammable. The oil contained within the appliance is flammable. Escaping refrigerant and oil can ignite if they are of high enough concentration and are exposed to an external heat source.

- ▶ Do not damage the pipelines of the coolant circuit and the compressor.
- Transport the appliance away without damaging it.
- Dispose of batteries, lamps, and the appliance according to the above specifications.

Additional information

10 Additional information

Further information on warranty conditions can be found in the enclosed service brochure or on the Internet at https://home.liebherr.com.

If there is a fault with the appliance, inform:	
Model ID /	
Inventory number:	

Quick Start Guide for everyday use

Information about using this Quick Start Guide:

- This Quick Start Guide is not a substitute for the full operating instructions nor is it a substitute for appliance training.
- This Quick Start Guide provides support for everyday use and handling the appliance.
- Read the full operating instructions and the safety instructions contained therein.
- This Quick Start Guide is for: SRFfg 35../40../55..
- Scanning the QR code will open the full operating instructions:





Intended use

This laboratory refrigerator, with an interior free of ignition sources, is suitable for professional storage of highly flammable products in closed containers, at temperatures between 3 °C and 16 °C.

Foreseeable incorrect use

Do not use the appliance for the following applications:

- Storage and cooling of:
 - Chemically unstable substances
 - Blood, plasma or other body fluids for the purpose of infusion, application or introduction in the human body
- Use in areas with risk of explosion
- Outdoor use or in areas where it is exposed to splash water or damp conditions

When stocking, observe the following points:

- ☐ Position grid shelves according to the height required.
- ☐ Observe maximum load.
- When the storage temperature has been reached, load the appliance (compliance with cold chain).
- Keep the ventilation slots of the recirculation fan free in the interior.
- ☐ Items for refrigeration must not come into contact with the evaporator on the rear wall.
- Keep liquids in closed containers.
- Leave space when storing items for refrigeration to ensure adequate air circulation.

Status display



Fig. 127 Status display with actual temperature

Navigating with buttons

	,		
Display and button	Description		
	Forward navigation arrow Advances to next menu (submenu). The first menu appears again after the last menu.		
	Backward navigation arrow		
	Short press: Goes back one menu (submenu). Long press (3 seconds): Returns to the status display.		
	Confirm Activates/deactivates function. Opens submenu. Confirms selection. The display switches back to the menu.		
Back	Back Display switches back to the function.		

Displaying temperature recording

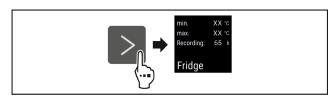


Fig. 128

- ► Complete steps (see Fig. 1831).
- ▷ Temperature recording displays minimum temperature, maximum temperature and length of recording.

Resetting temperature recording

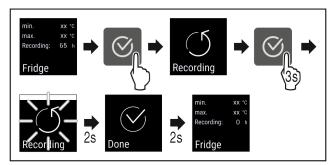


Fig. 129

- ► Complete steps (see Fig. 1832).
- > Temperature recording is reset.
- Recording interval starts again.

Quick Start Guide for everyday use

Unlocking appliance



Fig. 130

- ► Insert the key. *Fig. 130 (1)*
- ► Turn the key clockwise by 90°. Fig. 130 (2)
- \triangleright The appliance is unlocked.

Locking appliance



Fig. 131

- ► Insert the key. *Fig. 131 (1)*
- ► Turn the key anticlockwise by 90°. Fig. 131 (2)
- > The appliance is locked.

Cleaning appliance

- ► Empty the appliance.
- ▶ Pull out the power plug.
- ▶ Observe the information in the operating instructions.
- ► Dry the appliance and equipment thoroughly after cleaning.
- ► Connect the appliance.
- ► Switch on the appliance.
- ► Load the appliance once the storage temperature is reached.



What should I do if there is a fault?

A message appears in the display if there is a fault. Messages have different meanings:

Symbol color in the display	Meaning	Remedy
YELLOW	Reminder Reminder of general procedures. You can carry out these procedures to eliminate the message.	Consult the operating instructions and perform the troubleshooting steps.
RED	Warning Appears when there are malfunctions. You can eliminate simple malfunctions yourself.	Consult the operating instructions and perform the troubleshooting steps.

If you cannot remedy the fault yourself, contact a competent member of staff or your nearest customer service outlet.

Cut out the page





home.liebherr.com/fridge-manuals

