



UPLC[®] and CE Methods for Orthogonal Analysis of N-Glycans

A summary of Ultra-Performance Liquid Chromatography (UPLC) and Capillary Electrophoresis (CE) methods used in the ProZyme poster, Orthogonal Methods for Glycoanalysis: CE and UPLC (CE Pharm 2013)¹

UPLC Methods

Analytical System

LC System: Waters ACQUITY UPLC[®] (H-class) with FLR detector, Empower[™] software

Columns

- Waters ACQUITY UPLC BEH Glycan 1.7 μm (Glycan Separation Technology, GST column)

Method	Column Size	Waters Part Number
5-Minute Gradient	2.1 x 50 mm	186004740
10-Minute Gradient	2.1 x 100 mm	186004741
60-Minute Gradient A ^{2,3}	2.1 x 150 mm	186004742
60-Minute Gradient B	2.1 x 150 mm	186004742

- No guard column was used
- Column Temperature: 60°C

Sample Preparation Recommendations

- Prepare samples following protocols set forth in corresponding manual for GlykoPrep[®] Rapid N-Glycan Preparation with 2-AB (ProZyme product code GP96NG-AB).
- Inject 1 μl of aqueous sample per UPLC run.
- Remaining samples may be stored frozen in the dark. Use sealing film if storing in 96-well plates.

3832 Bay Center Place
Hayward, CA 94545-3619

TOLL FREE (800) 457-9444
PHONE (510) 638-6900
FAX (510) 638-6919

E-MAIL info@prozyme.com
WEB www.prozyme.com

These suggestions and data are based on information we believe to be reliable. They are offered in good faith, but without guarantee, as conditions and methods of use of our products are beyond our control. We recommend that the prospective user determine the suitability of our materials and suggestions before adopting them on a commercial scale. Suggestions for use of our products or the inclusion of descriptive material from patents and the citation of specific patents in this publication should not be understood as recommending the use of our products in violation of any patent or as permission to license to use any patents of ProZyme, Inc.

ProZyme and GlykoPrep are trademarks or registered trademarks of ProZyme, Inc. in the United States and other countries; Waters[®], UPLC[®], ACQUITY[®] and Empower[™] are registered trademarks of Waters Corporation, Milford, MA, USA; Beckman Coulter[®] and 32 Karat[™] are trademarks of Beckman Coulter, Inc., Brea, CA, USA.

Solvents

- Solvent A: Acetonitrile
- Solvent B: 100 mM Ammonium Formate pH 4.4

Excitation and Emission

Excitation: 360 nm

Emission: 428 nm

5-Minute Gradient

Method Details

Time (min)	Flow Rate (ml/min)	% A	% B
0	1.5	73	27
4	1.5	55	40
4.05	1.5	30	60
4.25	1.5	30	60
4.3	1	73	27
4.7	1.5	73	27
5	1.5	73	27

10-Minute Gradient

Method Details

Time (min)	Flow Rate (ml/min)	% A	% B
0	1	75	25
8	1	60	40
8.1	0.5	40	60
8.5	0.5	40	60
8.6	1	40	60
8.8	1	75	25
10.0	1	75	25

60-Minute Gradient A^{2,3}

Method Details

Time (min)	Flow Rate (ml/min)	% A	% B
0	0.5	78	22
38.5	0.5	55.9	44.1
39.5	0.25	20	80
44.5	0.25	20	80
46.5	0.5	78	22
60	0.5	78	22

60-Minute Gradient B

Method Details

Time (min)	Flow Rate (ml/min)	% A	% B
0	0.75	75	25
54	0.75	62.5	37.5
55	0.25	40	60
56	0.25	40	60
57	0.5	75	25
58	0.75	75	25
60	0.75	75	25

CE Methods

Analytical System

CE System: Beckman Coulter® PA800 *plus* with LIF detector, 32-Karat™ software

Capillaries

- Beckman Coulter N-CHO and eCAP (developmental)

Method	Capillary	Beckman Part Number
10-Minute Separation ⁴	eCAP	NA
15-Minute Separation ⁶	N-CHO	477601
20-Minute Separation ²	N-CHO	477601
35-Minute Separation ⁵	N-CHO	477601

Sample Preparation Recommendations

- Prepare samples following protocols set forth in corresponding manual for GP96NG-APTS
- Load each sample at 2 psi for 10 seconds
- Remaining samples may be stored frozen in the dark.

Separation Buffers

Method-specific; see individual methods

Excitation and Emission

Excitation: 360 nm

Emission: 428 nm

10-Minute Separation⁴

Method Details

Length	60 cm (effective: 50 cm)
I.D.	30 μ m
Separation Buffer	25 mM boric acid, 8 mM ammonium acetate; pH 9.0
Temperature	35°C
Voltage	30 kV for 10 minutes
Polarity	Reversed

15-Minute Separation⁶

Method Details

Length	60 cm (effective: 50 cm)
I.D.	50 μ m
Separation Buffer	40 mM ϵ -aminocaproic acid, 40 mM acetic acid, 0.2% hydroxymethylcellulose
Temperature	20°C
Voltage	30 kV for 15 minutes
Polarity	Reversed

20-Minute Separation²

Method Details

Length	60 cm (effective: 50 cm)
I.D.	50 μ m
Separation Buffer	N-linked Carbohydrate Separation Gel Buffer (Part No. 477623)
Temperature	20°C
Voltage	30 kV for 20 minutes
Polarity	Reversed

35-Minute Separation⁵

Method Details

Length	60 cm (effective: 50 cm)
I.D.	50 μ m
Separation Buffer	1:1 Mixture of N-linked Carbohydrate Separation Gel Buffer (Part No. 477623) and eCAP ds DNA 1000 Gel Buffer (Part No. 477628)
Temperature	20°C
Voltage	20 kV for 35 minutes
Polarity	Reversed

References

1. Szabo *et al.* Orthogonal Methods for Glycoanalysis: CE and UPLC. Poster presented at: CE in Biotechnology & Pharmaceutical Industries: 15th Symposium on the Practical Applications for the Analysis of Proteins, Nucleotides and Small Molecules (CE Pharm 2013), October 6-10, 2013, Arlington, VA, USA.
2. N-Glycan Mapping Study Using Orthogonal Methods: CE and UPLC. Workshop at CE in Biotechnology & Pharmaceutical Industries: 15th Symposium on the Practical Applications for the Analysis of Proteins, Nucleotides and Small Molecules (CE Pharm 2013), October 6-10, 2013, Arlington, VA, USA.
3. Ahn, J. *et al.* UPLC-FLR Method Development of 2-AB-Labeled Glycan Separation in Hydrophilic Interaction Chromatography (HILIC). Application Note. Waters Corporation, Milford, MA, USA.
4. Szabo, Z. *et al.* Rapid High-Resolution Characterization of Functionally Important Monoclonal Antibody N-Glycans by Capillary Electrophoresis. *Analytical Chemistry*, 2011, 83, (13), 5329-5336.
5. Rampal, S. *et al.* Separation of Fucosylated, non-Fucosylated, and Complex Carbohydrates Associated with Monoclonal Antibodies using Capillary Electrophoresis. *Chromatography Today* 2011, 3, 38-40.
6. Hamm, M. *et al.* Characterization of N-Linked Glycosylation in a Monoclonal Antibody Produced in NS0 Cells Using Capillary Electrophoresis with Laser-Induced Fluorescence Detection. *Pharmaceuticals*, 2013, 6, 393-406.