

Use:

NON-Sterile – Disinfect before each and every use.

Ensure all fittings are secure and leak-free.

Tubing connection requires 1/4" OD, 1/8" ID

Insert animal into mask until tight seal is achieved.

Face mask can be trimmed. Do Not Remove Ports.

Left Port = Gas Supply (≤ 1.0 Liter/minute)

Right Port = Exhaust to appropriate filter

Ports can be rotated 360°

To remove mask from tubing, cut a slit in tubing over port barb.

CAUTION:

Misuse of this product can lead to leakage of Isoflurane

For Animal Research Purposes Only

Authorized Personnel Must Supervise and Ensure Safe Use

Manufacturer:

Xenotec, Inc. d/b/a Ocuscience

202 West Ninth Street

Suite 301

Rolla, Missouri 65401

816.529.5712

www.xenotec.us

Rodent Isoflurane Facemask Kit



CONTENTS:

- (1) Small Facemask with ports (Mouse)**
- (1) Medium Facemask with ports (Mouse/Rat)**
- (1) Large Facemask with ports (Rat)**
- (2) 6 feet medical grade PVC tubing**
- (2) 1/4 Inch Adapters**
- (2) 3/8 Inch Adapters**



Xenotec, Inc.

Medical Device Development